


making a difference

Download our
latest catalogue:



BIOSCIENCE CATALOGUE

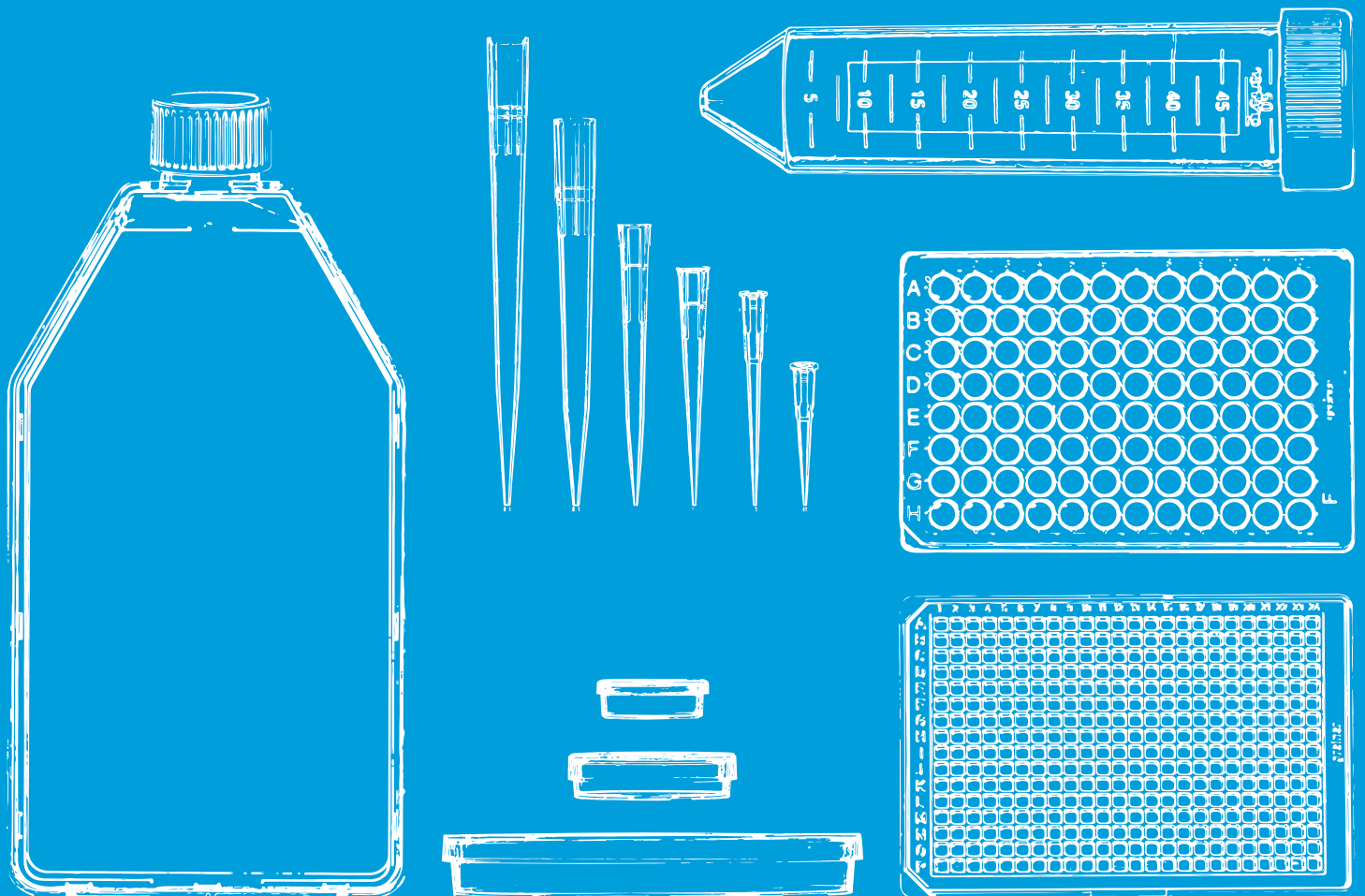

EDITION 2026

www.gbo.com


greiner
BIO-ONE

making a difference



To us, health is the highest personal priority. This inspires our daily mission of serving life sciences and healthcare professionals with advanced and sustainable products and solutions.

It is the passion and commitment of our employees employees that makes the difference. With a targeted approach, our global team continuously strives for excellence and works closely with our customers.

We empower people to make a difference in health.

CONTENT


6



PRODUCTS

Here you can find our product overview


10

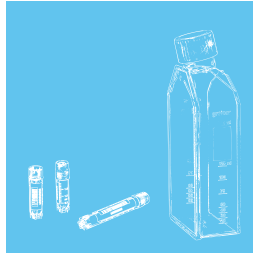

BIOSCIENCE

Making a difference in research and science


12

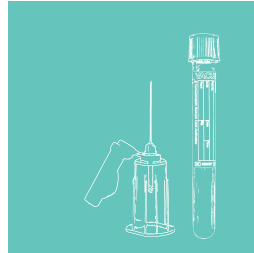

PREANALYTICS

Preanalytical specimen collection systems and safety solutions

14

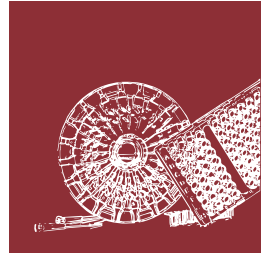

OEM

Tailor made solutions

17



SERVICE

E-Commerce Solutions from Greiner Bio-One

18



VALUE PROPOSITIONS


About our value propositions

251



TECHNICAL APPENDIX

284

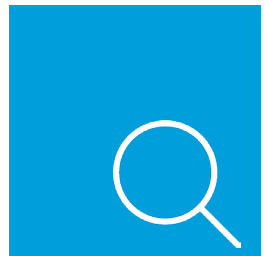

NUMERICAL INDEX

PRODUCTS

37

CELL CULTURE

Cell Culture Flasks	38
CELLSTAR® Cell Culture Dishes	44
CELLSTAR® Cell Culture Multiwell Plates	48
ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates	52
ThinCert® 96 Well HTS Insert	56
CELLSTAR® Cell Culture Microplates	60
CELLSTAR® Cell Culture Tubes	68
EASYstrainer Cell Strainers	70
Media Bottles	72

75

MASS CELL CULTURE

CELLMASTER Cell Culture Roller Bottles	76
CELLdisc Cell Culture Device	79

87

3D CELL CULTURE

CELLSTAR® Cell Culture Vessels Cell-Repellent Surface	88
Magnetic 3D Cell Culture	92
ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates	97

103

PRODUCTS FOR MICROSCOPY

CELLview Dish Cell Culture Dish with Glass Bottom	104
CELLview Slide Cell Culture Slide with Glass Bottom	106
CELLview Plate Cell Culture Plate with Glass Bottom	108
SCREENSTAR Microplates	110
SensoPlate Glass Bottom Microplates	112



shop.gbo.com

Online Shop

115 HTS-MICROPLATES

96 Well Microplates	116
384 Well Microplates	120
1536 Well Microplates	124
Polypropylene Storage Plates	126
Microplates for Compound Storage	129
Non-binding Microplates	131
Streptavidin-coated Microplates	134
UV-Star® Microplates	136
Lids / Sealers / CapMats	138
Protein Crystallisation Plates	141

145 IMMUNOLOGY / HLA

ELISA Microplates	146
ELISA Strip Plates	148
Immuno Tubes	151
Terasaki Plates	153

157 MICROBIOLOGY

Dishes / Plates / Others	158
--------------------------	-----

165 TUBES / MULTIPURPOSE BEAKERS

Tubes without closures	166
Tubes with closures	170
Leucosep Separation Tubes	174
Multipurpose Containers / Beakers	176

PRODUCTS

181 CRYOTECHNICS

Cryo.s Cryo Tubes	182
Cryo.s with Barcode	188
Cryo.s Biobanking-Tubes	191

201 REACTION TUBES / ANALYSER CUPS

Reaction Tubes / Analyser Cups	202
Semi-micro / Macro Cuvette	206

209 MOLECULAR BIOLOGY

Sapphire PCR Tubes	210
Sapphire PCR Microplates	214

217 LIQUID HANDLING

Sapphire Pipettes	218
Sapphire Pipette Tips	221
CELLSTAR® Serological Pipettes	230

235 LAB EQUIPMENT

Lab Equipment	236
---------------	-----



241 TECHNICAL APPENDIX



shop.gbo.com

Online Shop

BIOSCIENCE

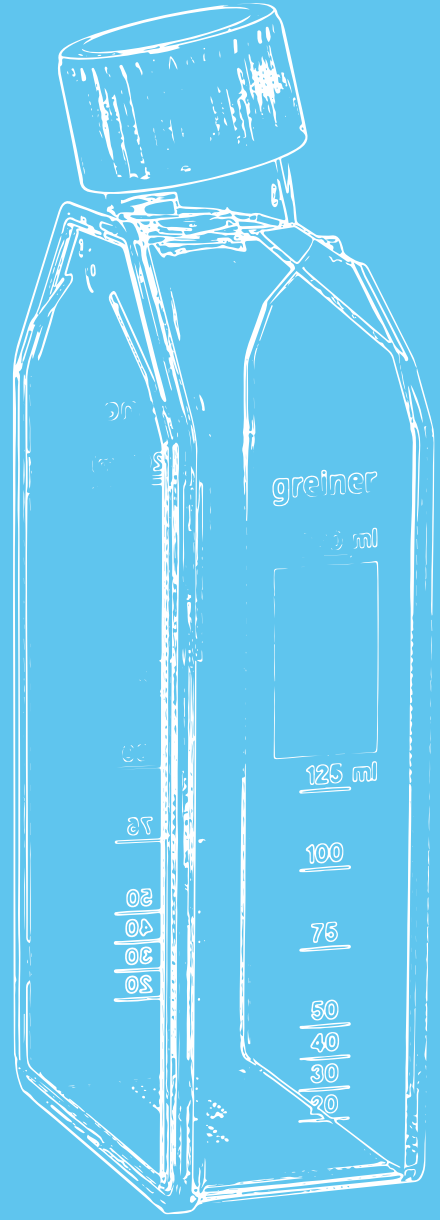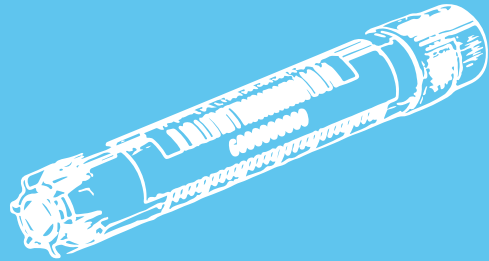

MAKING A DIFFERENCE IN RESEARCH AND SCIENCE

Greiner Bio-One is one of the leading suppliers of special products for the cultivation and analysis of cell and tissue cultures. Under the brand name CELLSTAR®, we offer cell culture flasks, dishes and plates in a wide variety of formats and surface modifications, so that you will always find exactly the right product for every application and all cell types.


In addition, we have been developing and producing microplates for high-throughput screening for more than 60 years, enabling

industry and research to carry out fast and efficient drug testing. Among other things, Greiner Bio-One was the first manufacturer to introduce a microplate in 1536 well format to the market, setting new standards in terms of automation, performance and cost reduction.

Based on decades of experience in the cryogenic storage of samples, we also offer comprehensive solutions for automated storage systems in biobanks.

- / **High-quality products for your lab**
- / **Special products for cell and tissue culture**
- / **Microplates for industry and research**
- / **Cryogenic storage systems**

PREANALYTICS

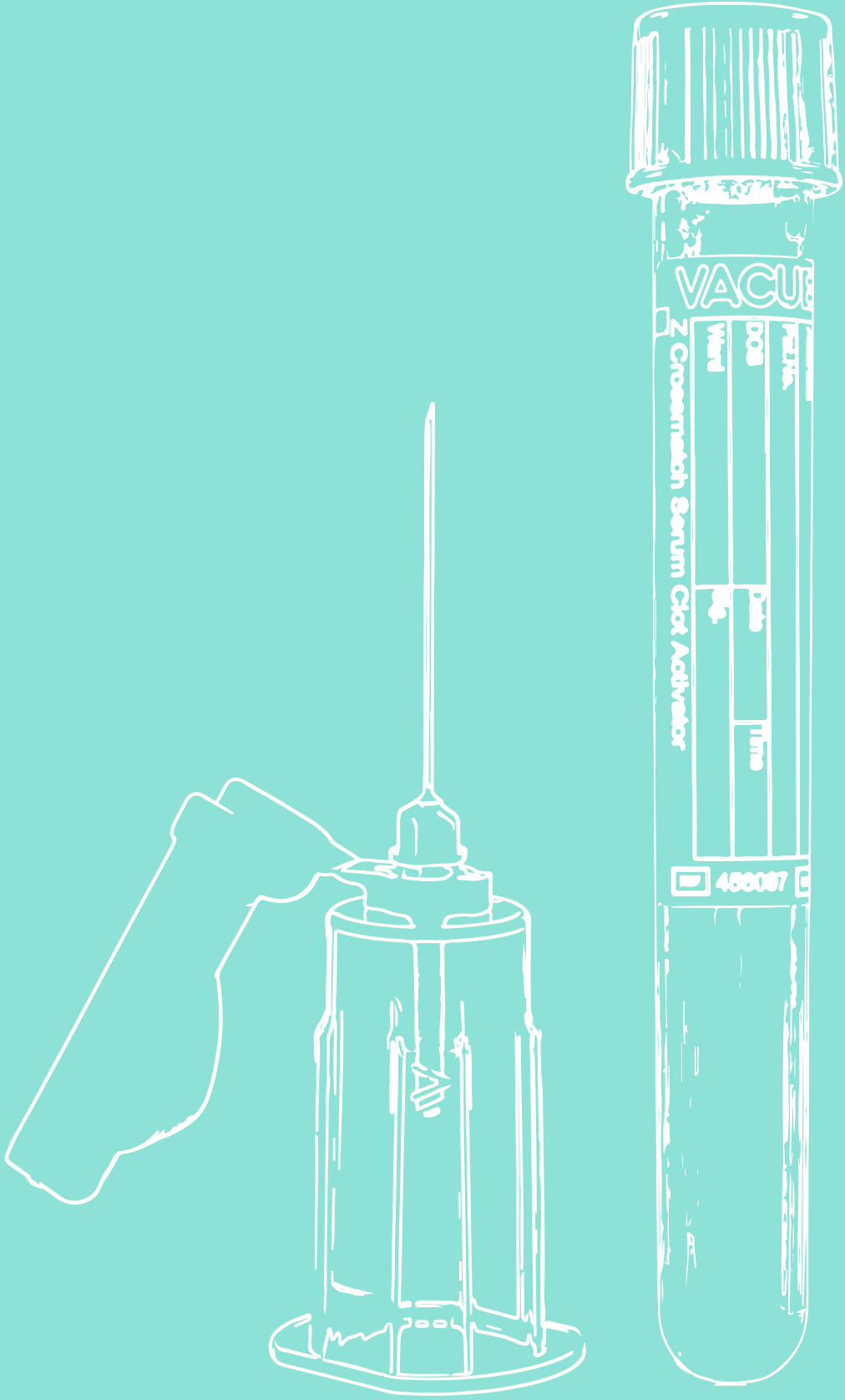

PREANALYTICAL SPECIMEN COLLECTION SYSTEMS AND SAFETY SOLUTIONS

Greiner Bio-One was the first company in the world to successfully launch a blood collection system made out of virtually unbreakable PET plastic including a safety cap. Many components of the VACUETTE® System have often been copied, but never reached our high quality standards.


As a trendsetter in the field

of sample collection, we offer solutions that meet the highest demands of our customers. Because your work is an important contribution to health protection.

Continuously new and improved product solutions, perfect functionality, maximum safety and high product quality are the result of decades of experience.

- / **Provide safety**
- / **Simplify handling**
- / **Provide support**
- / **Increase efficiency**
- / **Save costs**

OEM

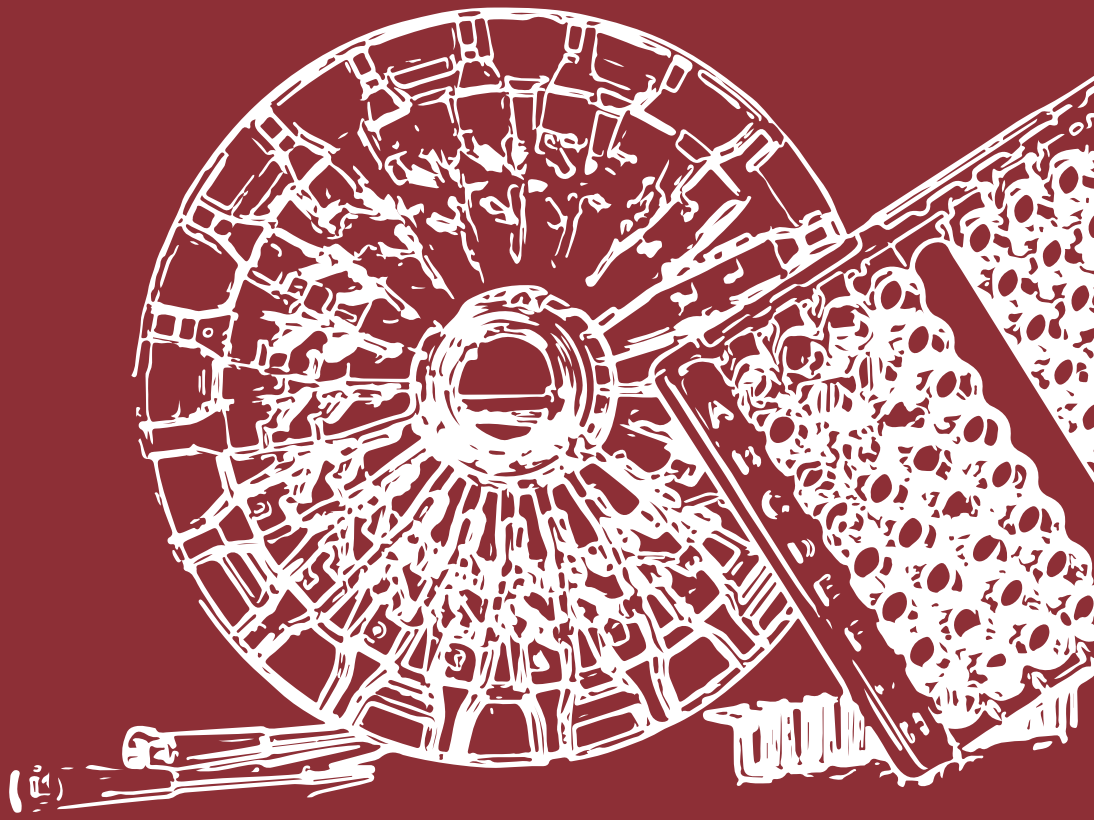

TAILOR MADE SOLUTIONS

Greiner Bio-One as an Original Equipment Manufacturer (OEM) is a long-term partner of the pharmaceutical industry, biotechnology, diagnostic and medical technology industries. The company manufactures numerous products through injection moulding and offers the entire product development and production process. From design,

via prototyping, through fully automated manufacturing, Greiner Bio-One delivers the complete solution with support of experienced scientists, engineers and specialists.

With a global sales network specifically established for the field of OEM, Greiner Bio-One offers custom-made solutions with personal service.

- / Research & development with diverse application and research laboratories**
- / Construction of moulds, planning of plants and assembly lines**
- / Modern, fully automated production**
- / Post-production according to customer requirements**
- / Comprehensive quality assurance and management system**
- / Worldwide networked warehousing and customer service**



FAST. SECURE. EFFICIENT. E-COMMERCE SOLUTIONS FROM GREINER BIO-ONE

As your strong partner in the fields of biotechnology, diagnostics, medical devices and research, Greiner Bio-One already provides you with solutions on a daily basis. In addition, we can offer you a wide selection of e-business solutions to ensure the optimal integration of our products and services into your purchasing process. That in turn will significantly reduce your process-related costs and warehouse costs.

Take advantage of e-business connectivity and make your processes faster, more secure and more efficient.

Greiner Bio-One offers you numerous solutions and the necessary expertise to support your processes, regardless of which ERP system you use. Would you like to purchase via our online shop or are you interested in one of our eProcurement solutions such as EDI or OCI?

In order to stay informed, you can subscribe to our Greiner Bio-One Newsletter and never miss any news about our products and events.




- / Online-Shop
- / EDI (Electronic Data Interchange)
- / OCI (Open Catalogue Interface)
- / Electronic catalogues



Subscribe to our newsletter
and never miss
the latest news!



A male scientist in a white lab coat and hairnet is smiling and gesturing with his hand in a laboratory. The lab coat has a 'greiner BIO-ONE' logo on the pocket. In the background, there is a piece of laboratory equipment with a yellow emergency stop button.

As scientists, we
share your passion for
your research success.
We are happy to support you
wherever and whenever
we are needed.

TRUE TRUST

Take advantage of our experience and passion for your projects and tasks



WE KEEP OUR PROMISES

For us, real trust means keeping promises. This is one of the core values that we are committed to as a company, both towards you as our customers and towards our employees.

As a family-owned company, we believe that trusting, active partnership is the key to sustainable and reliable processes. We


understand your requirements, speak your language and are happy to support you in advancing your scientific progress. We are at your side at every stage – our competent team of experts is always on hand, from the first contact and flawless delivery of our products, to all forms of after-sales service.

PUT YOUR TRUST IN US:

- / Partner for sustainable and reliable processes
- / More than 60 years of experience in science and healthcare
- / Innovation with tradition
- / Support during every work step



DISCOVER MORE:

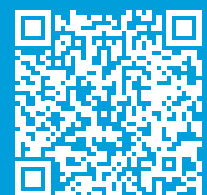

With state-of-the-art technology,
we develop high-performance
solutions for consistent,
reproducible results that
you can always rely on.



QUALITY

Reproducible results
you can rely on

CONSISTENTLY HIGH QUALITY FOR RELIABLE RESULTS

Quality has been our top priority for more than 60 years. This applies all over the world, as all of our production facilities are committed to the same high standard.

Our products are made from strictly controlled, certified raw materials. With state-of-the-art production technology and

continuous in-process checks by specially trained employees, we ensure consistent quality of the products. Quality promise not only relates to the manufacturing process and therefore to product quality, but to every process and every interaction with you. We are ISO 9001 and ISO 13485 certified.




CONSISTENT HIGH QUALITY:

- / Location-independent
- / State-of-the-art manufacturing technologies
- / Controlled and certified raw materials
- / ISO 9001 and ISO 13485 certified



DISCOVER MORE:





Thanks to our presence in over 100 countries, you can rely on us to provide you with consistent high quality, availability and support. Wherever you are.

GLOBAL PRESENCE

There to support you
- wherever you are

ON-SITE FOR YOU WORLDWIDE

By global presence we mean being there for you whenever and wherever you need us.

Our seven production sites in Europe, Asia and North and South America meet the same high quality standards and

ensure low-risk delivery routes and a secure product supply. The support teams of our 28 offices and distributors in more than 100 countries are helpful and competent, and are always on hand when you need help or support.



WE PLACE IMPORTANCE ON:

- / **Competent local support**
- / **Optimum product availability**
- / **Consistent high quality worldwide**



DISCOVER MORE:





Our innovative spirit
supports the results of
your pioneering work.
In the past, present
and future.

STATE- OF-THE-ART TECHNOLOGY

Pioneers for
progress

IDEAS BECOME SOLUTIONS

For more than 60 years, our passion for plastic technology and manufacturing processes has driven us to constantly innovate to support you optimally in your daily work and ensure your success.

We have been instrumental in the development of consumables in the diagnostics and life sciences industries, as well as in the medical market, since 1963.

We are proud of our rich pioneering history – from the production of the first plastic petri dishes, to Europe’s first plastic vacuum sampling system or the invention of the world’s first 1536-well microplates. We will continue to strive for new and innovative product solutions that are precisely tailored to your requirements and optimally support you in your daily work.



PROGRESS NEEDS PIONEERS:

- / Product solutions for current and emerging trends
- / Innovative production technologies



DISCOVER MORE:



We work hand in hand with you to
develop tailored solutions to suit
your requirements.



CUSTOMER FOCUS

We focus on you and your goals

WE SUPPORT YOU SO WE CAN MAKE BIG THINGS HAPPEN TOGETHER

Everything we do and the way we do it is designed to make your work easier. By understanding your wishes and needs, as well as your requirements and workflows, we are able to offer you real added value and thus trust in long-term and sustainable customer relationships.

We are here to provide you with the best products and solutions

to suit your needs. Our dedicated team ensures that processes run smoothly and is always on hand when needed. Our multidisciplinary teams have many years of practical experience in science and laboratory work and understand the challenges of your daily work first hand. We are happy to use this expert knowledge to create training materials tailored to your specific needs.




WE OFFER:

- / Long-term customer relationship
- / Solutions to suit your requirements
- / Training and training materials



DISCOVER MORE:

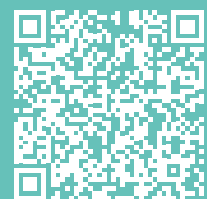

We use over 60 years of experience
in the laboratory and healthcare
sectors, and our passion
for innovation, to meet your
requirements and grow with you.



EXPERTISE

In all relevant market trends

WE SPEAK YOUR LANGUAGE

We understand your tasks and their workflows with all their requirements. Our experts are at your side with more than 60 years of experience – from the development of products and solutions to support with regulatory issues or individual training.

More than 80% of our sales and application specialists have a background in natural sciences

and practical experience in the laboratory. We would be happy to advise you on site – at the place where your requirements arise. You can rely on our expertise to quickly integrate our products into your processes, because we know and understand your workflow. We are happy to support you in every step of your individual workflow, for example with product training, support documentation and service.



WHAT WE OFFER:

- / **Comprehensive portfolio for your workflow needs**
- / **Experienced team of industry experts**
- / **Best practices for optimal efficiency**



DISCOVER
MORE:



We help you to optimize your operations and processes so that you can always achieve the best possible results.



EFFICIENCY

We are your partner
for increasing
efficiency – inside
and outside the
laboratory

WE INCREASE YOUR EFFICIENCY

In the laboratory and day-to-day work, increasing efficiency is no longer just an option, but an essential requirement. As a strategic cooperation partner for day-to-day routine activities in and outside the laboratory, we support you in overcoming your current and future challenges.

We offer significant added value by supporting the efficien-

cy of your workflows. For example, our digital solutions create a connection between pre-bar-coded test tubes and patient data from the laboratory information system. This reliably rules out mix-ups and minimizes sampling errors. Your workflows can thus be measurably optimized, better results can be achieved with minimum time, resources, and work.



YOUR STRATEGIC PARTNER:

- / Increased efficiency
- / Digital solutions
- / Best practice recommendations



DISCOVER MORE:



We are committed to climate protection and limiting global warming to 1.5°C.



SUSTAINABILITY

The principle behind our processes for many years

BECAUSE IT'S REALLY IMPORTANT TO US

Sustainable action is the basic prerequisite for a future worth living in. As a plastics processing company, we strive to develop sustainable solutions that provide answers to the major challenges of our time.

Greiner AG's overarching "Blue-Plan" strategy focuses on the three pillars of climate, circular economy and people. Greiner


Bio-One is also working intensively on these sub-areas.

As a producer of plastic medical products, it is our declared long-term goal to combine ecological, economic and social aspects. We attach particular importance not only to your excellent supply of high-quality products, but also to energy-efficient and sustainable operation as a plastics producer.




OUR GOALS UNTIL 2030:

- / **Comprehensive circular approach**
- / **Preparation of employees for the challenges of the future**

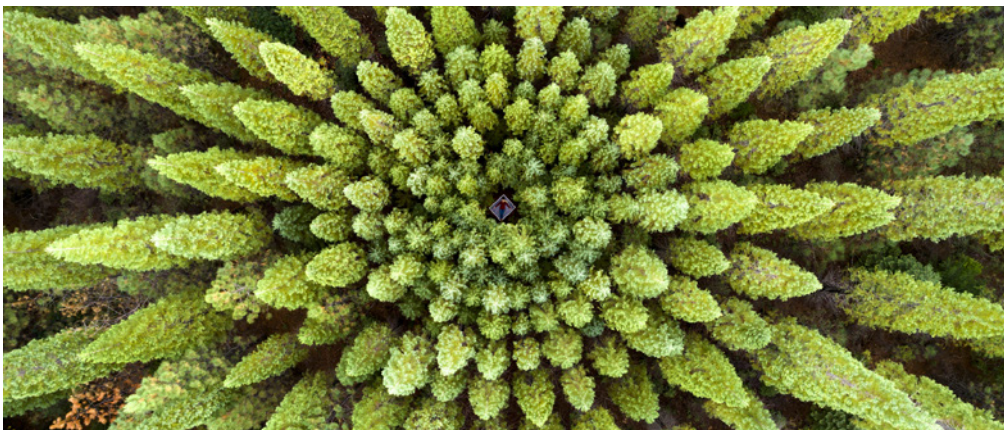

DISCOVER MORE:

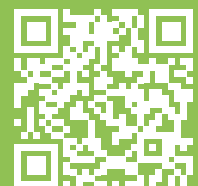

Learn more about the
ACT[®] Ecolabel.



my green lab.
ACT Ecolabel



my green lab.
ACT Ecolabel

Labware certified with
ACT® Ecolabel

MAKE AN ECO-CONSCIOUS CHOICE

At Greiner Bio-One we believe in taking meaningful actions to create sustainability and transparency in the life sciences industry. This commitment is reflected in the variety of products that have already received ACT® Ecolabel certification from My Green Lab®.

By investing in the independently audited ACT® Ecolabel, we aim to provide our customers with the information they need to make eco-conscious choices. The ACT® Ecolabel offers a clear and reliable assessment of the environmental impact of our products, helping you to make decisions that support a greener future.


The potential to minimise the environmental footprint of laboratories through smarter pur-

chasing decisions is immense. The ACT® Environmental Impact Factor Label was created to meet the needs of both scientists and procurement professionals by providing clear, third-party verified data on the environmental impact of lab products. By focusing on Accountability, Consistency, and Transparency (ACT®), it simplifies the process of selecting more sustainable products.


The ACT® Ecolabel uses a science-based, 100-point weighted score, with details on product impacts such as energy use, materials, chemical hazards, and end-of-life options - plus an evaluation of the manufacturing facility and company commitments to reduce greenhouse gas emissions.

WHY ACT® ECOLABEL?

- / Transparent 3rd-party verified sustainability data
- / Includes carbon footprint
- / Improved communication of sustainability attributes

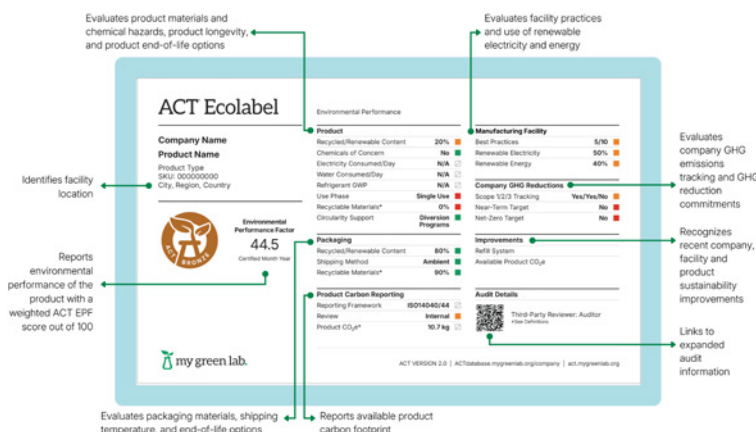

OUR CERTIFIED PRODUCTS:

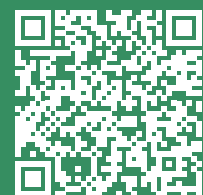



The secret of the continuing success of the CELLSTAR® product line from Greiner Bio-One lies in its high quality and reliability. For more than 60 years, Greiner Bio-One has been setting standards for future-oriented research and new technologies in the field of cell culture.

CELL CULTURE

/ Cell Culture Flasks	38	ThinCert® 96 Well HTS Insert (Membrane and Receiver Plates).....	57
CELLSTAR® Standard Cell Culture Flasks	39	ThinCert® 96 Well HTS Insert (Membrane, Receiver and Feeder Plates).....	57
Cell Culture Flasks - Advanced TC	40	Receiver plates for ThinCert® 96 Well HTS Insert	58
Cell Culture Flasks - CELLCOAT®	40	Feeder Plates for ThinCert® 96 Well HTS Insert	59
Suspension Culture Flasks	41		
Cell Culture Flasks - Cell-Repellent Surface	42		
AutoFlask - Cell Culture Flask For Automated Systems	42		
Cell Scraper	43		
/ CELLSTAR® Cell Culture Dishes	44	/ CELLSTAR® Cell Culture Microplates	60
Cell Culture Dishes	45	Cell Culture Microplates 96 Well.....	61
Cell Culture Dishes - Advanced TC	45	Cell Culture Microplates 96 Well - µClear®	62
Cell Culture Dishes - CELLCOAT®	46	Cell Culture Microplates 96 Well - Half Area.....	62
Cell Culture Dishes - Cell-Repellent Surface.....	46	Cell Culture Microplates 96 Well - CELLCOAT®	63
CELLview Dish	47	Cell Culture Microplates 96 Well - Advanced TC	64
OneWell Plate / FourWell Plate.....	47	Suspension Culture Microplates 96 Well	64
		Cell Culture Microplates 384 Well.....	65
		Cell Culture Microplates 384 Well - Small Volume	66
		Cell Culture Microplates 1536 Well.....	66
		Lids	67
/ CELLSTAR® Cell Culture Multiwell Plates	48	/ CELLSTAR® Cell Culture Tubes	68
Multiwell Plates 6 / 12 / 24 / 48 Well Format	49	CELLreactor	69
Multiwell Plates Advanced TC	49		
Multiwell Plates CELLCOAT®	50		
Multiwell Plates Cell-Repellent Surface	51		
/ ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates	52	/ EASYstrainer Cell Strainers	70
ThinCert® Cell Culture Inserts 6 Well.....	53	EASYstrainer	71
ThinCert® Cell Culture Inserts 12 Well	54		
ThinCert® Cell Culture Inserts 24 Well.....	55		
ThinCert® Plate 6 / 12 Well.....	55	/ Media Bottles	72
		Media Bottles	73
/ ThinCert® 96 Well HTS Insert	56		



/ Learn more about our cell culture flasks:



CELL CULTURE FLASKS

Greiner Bio-One offers standard and filter cap cell culture flasks with different surfaces. All Greiner Bio-One cell culture flasks are made of high-grade polystyrene and are free of detectable DNases, RNases, human DNA, endotoxins and are non-cytotoxic.

The specific design facilitates optimal access to the flask when using a cell scraper or pipette. The top stacking rim stabilizes

the flask and facilitates easy flask stacking in the incubator. Both sides of the flask have graduation marks to aid liquid fill. Filter or standard screw caps options are available.

Greiner Bio-One flasks offer a specific physical surface treatment to ensure adhesion and proliferation of adherent cells. In addition, the portfolio includes flasks with a hydrophobic surface, ideal for suspension cultures and hybridoma cells.

- / Standard or filter screw cap
- / Canted neck
- / User-friendly packaging
- / Growth area for adherent cells: 25, 75 and 175 cm²
- / Total volume: 50, 250, 550 and 650 ml



CELLSTAR® Standard Cell Culture Flasks

For the maintenance of adherent cells Greiner Bio-One offers a specific physical surface treatment which ensures adhesion and proliferation of these cells.

/ Item No. 690170 with measuring grid



Raw material: PS, Surface treatment: TC, Sterile: +

Item no.	Flask design	Growth area	Cap color	Working volume	Total volume	Cap design	Triple-packed	Sterile	Qty. inner / outer
690170		25 cm ²	● red	5 ml - 10 ml	50 ml	screw cap		+	10 / 200
690175		25 cm ²	● red	5 ml - 10 ml	50 ml	filter screw cap		+	10 / 200
690175-TRI		25 cm ²	● red	5 ml - 10 ml	50 ml	filter screw cap	yes	+	10 / 130
658170		75 cm ²	● red	15 ml - 40 ml	250 ml	screw cap		+	5 / 120
658175		75 cm ²	● red	15 ml - 40 ml	250 ml	filter screw cap		+	5 / 120
658175-TRI		75 cm ²	● red	15 ml - 40 ml	250 ml	filter screw cap	yes	+	5 / 90
660160	flat	175 cm ²	● red	20 ml - 75 ml	550 ml	screw cap		+	5 / 50
660175	flat	175 cm ²	● red	20 ml - 75 ml	550 ml	filter screw cap		+	5 / 50
660175-TRI	flat	175 cm ²	● red	20 ml - 75 ml	550 ml	filter screw cap	yes	+	5 / 40
661160	high	175 cm ²	● red	20 ml - 85 ml	650 ml	screw cap		+	4 / 40
661175	high	175 cm ²	● red	20 ml - 85 ml	650 ml	filter screw cap		+	4 / 40

/ The item no. of triple-packed products is composed of the standard item no. plus -TRI (e.g. 604160-TRI). In general, all Greiner Bio-One cell culture products can be produced triple-packed. Please contact your Greiner Bio-One sales representative regarding customised triple-packed products.



Cell Culture Flasks - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.


- / Consistent and even cell attachment
- / Homogeneous and optimised cell growth

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Raw material: PS, Surface treatment: Advanced TC, Sterile: +

Item no.	Flask design	Growth area	Cap color	Working volume	Total volume	Cap design	Sterile	Qty. inner / outer
658970		75 cm ²	● blue	15 ml - 40 ml	250 ml	screw cap	+	5 / 120
658975		75 cm ²	● blue	15 ml - 40 ml	250 ml	filter screw cap	+	5 / 120
660960	flat	175 cm ²	● blue	20 ml - 75 ml	550 ml	screw cap	+	5 / 50
660975	flat	175 cm ²	● blue	20 ml - 75 ml	550 ml	filter screw cap	+	5 / 50

/ Application Report „Advanced TC for improving the cultivation / differentiation of embryonic stem cells“(F076036)

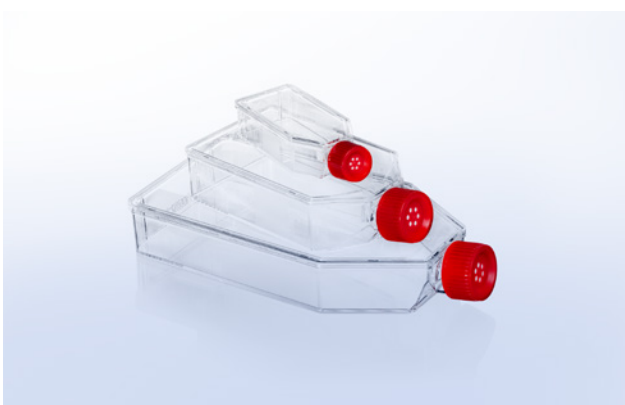

Cell Culture Flasks - CELLCOAT®

- / Canted neck for optimal access
- / Graduation on both sides
- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

ASEPTIC

Raw material: PS, Surface treatment: CELLCOAT®, Cap design: filter screw cap

Item no.	Flask design	Growth area	Protein coating	Cap color	Working volume	Total volume	Aseptic	Qty. inner / outer
690950		25 cm ²	Collagen Type I	● red	5 ml - 10 ml	50 ml	+	10 / 50
658950		75 cm ²	Collagen Type I	● red	15 ml - 40 ml	250 ml	+	5 / 50
661950	high	175 cm ²	Collagen Type I	● red	20 ml - 85 ml	650 ml	+	5 / 40
690940		25 cm ²	Poly-D-Lysine	● red	5 ml - 10 ml	50 ml	+	10 / 50

Item no.	Flask design	Growth area	Protein coating	Cap color	Working volume	Total volume	Aseptic	Qty. inner / outer
658940		75 cm ²	Poly-D-Lysine	● red	15 ml - 40 ml	250 ml	+	5 / 50
661940	high	175 cm ²	Poly-D-Lysine	● red	20 ml - 85 ml	650 ml	+	5 / 40
690920		25 cm ²	Fibronectin	● red	5 ml - 10 ml	50 ml	+	10 / 10
658920		75 cm ²	Fibronectin	● red	15 ml - 40 ml	250 ml	+	10 / 10
661920	high	175 cm ²	Fibronectin	● red	20 ml - 85 ml	650 ml	+	5 / 5
690910		25 cm ²	Laminin	● red	5 ml - 10 ml	50 ml	+	5 / 10
658910		75 cm ²	Laminin	● red	15 ml - 40 ml	250 ml	+	5 / 10
661910	high	175 cm ²	Laminin	● red	20 ml - 85 ml	650 ml	+	5 / 5




Suspension Culture Flasks

- / Hydrophobic surface, ideal for suspension cultures, hybridoma and embryonic stem cells
- / Graduation on both sides
- / Sterile and user-friendly packaging
- / Canted neck for optimal access

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Raw material: PS, Surface treatment: suspension, Sterile: +

Item no.	Flask design	Cap color	Total volume	Cap design	Triple-packed	Sterile	Qty. inner / outer
690190		○ white	50 ml	screw cap		+	10 / 200
690195		○ white	50 ml	filter screw cap		+	10 / 200
690195-TRI		○ white	50 ml	filter screw cap	yes	+	10 / 130
658190		○ white	250 ml	screw cap		+	5 / 120
658195		○ white	250 ml	filter screw cap		+	5 / 120
658195-TRI		○ white	250 ml	filter screw cap	yes	+	5 / 90
660190	flat	○ white	550 ml	screw cap		+	5 / 50
661190	high	○ white	650 ml	screw cap		+	4 / 40
661195	high	○ white	650 ml	filter screw cap		+	4 / 40
661195-TRI	high	○ white	650 ml	filter screw cap	yes	+	4 / 28

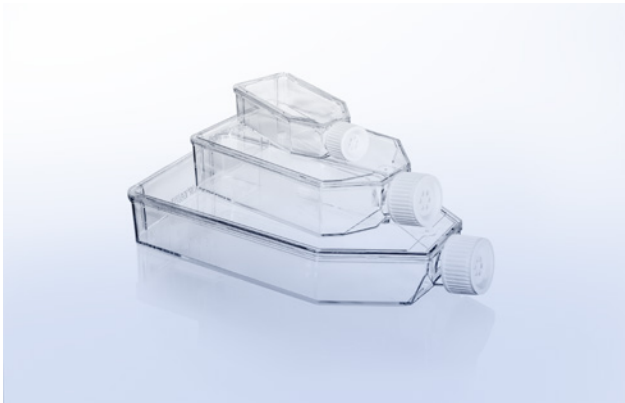

**Cell Culture Flasks -
Cell-Repellent Surface**

A cell-repellent surface reliably prevents cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Flask design: flat, Raw material: PS, Surface treatment: cell-repellent, Cap design: filter screw cap, Sterile: +

Item no.	Cap color	Total volume	Sterile	Qty. inner / outer
690985	○ white	50 ml	+	10 / 20
658985	○ white	250 ml	+	5 / 15
660985	○ white	550 ml	+	5 / 5



**AutoFlask - Cell Culture Flask
For Automated Systems**

- / Standard microplate footprint
- / Compatible with a wide range of cell culture and liquid handling systems
- / Hydrophobic filter membrane
- / Different surface treatments
- / Customised barcoding on request

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Flask design: AutoFlask, Growth area: 83.6 cm², Raw material: PS, Total volume: 110 ml, Sterile: +

Item no.	Barcode	Surface treatment	Product color	Color code	Working volume	Sterile	Qty. inner / outer
779160	yes	TC	○ clear	● red	20 ml - 40 ml	+	10 / 100
779190	yes	suspension	○ clear	○ white	60 ml - 80 ml	+	10 / 100




STERILE

Cell Scraper

- / For gentle mechanical removal of adherent cells
- / Optimised blade design for maximum cell harvest
- / Blade length: 1.8 cm
- / Pivot angle 60°
- / Sterile individual packaging

Description: Cell Scraper, Feature: blade length 1.8 cm, Sterile: +

Item no.	Length	Sterile	Qty. inner / outer
541070	28 cm	+	50 / 100
541080	40 cm	+	100 / 100

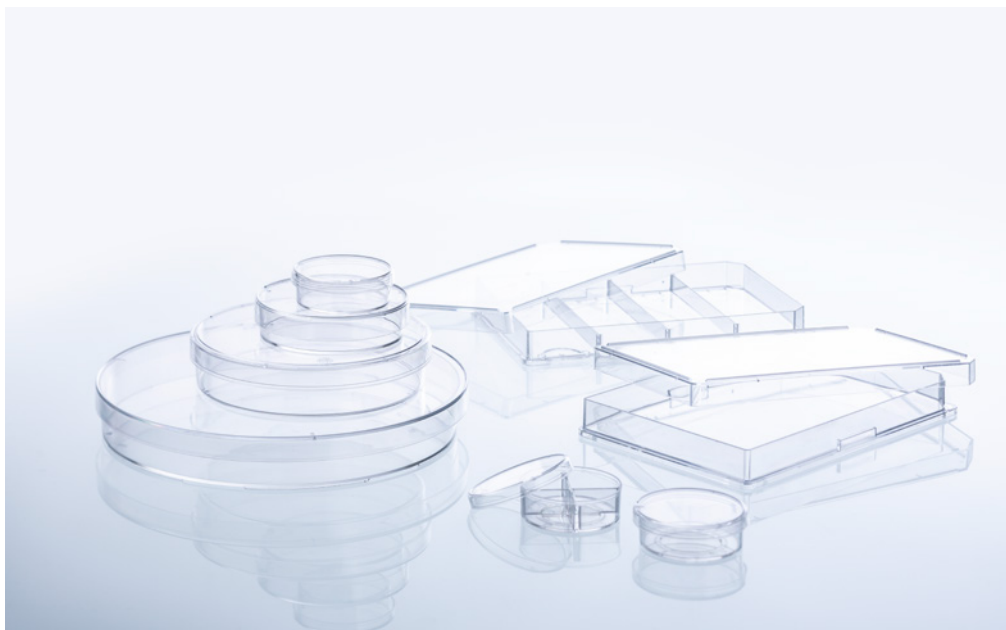

- / For exact dimensions of our cell culture dishes, please refer to the product data sheets on our website.



CELLSTAR® CELL CULTURE DISHES

Our CELLSTAR® cell culture dishes are manufactured to meet the highest quality standards at Greiner Bio-One.

This assortment of dishes offers a range of growth sizes from 8.7 to 143 cm². These dimensions are provided in nominal sizes of 35, 60, 100, and 145 mm. For culture dishes associated with growth areas of either 58 or 143 cm², we supply variants with side heights of 20 mm (for special applications). Our standard cell culture dishes have been treated in order to optimise growth and adhesion.

However, we are also capable of providing sterile cell culture trays without any type of coating for suspension culture. Similar to all of our cell culture laboratory products, our range of dishes is packed user-friendly. This allows technicians to safely handle their samples with a wide variety of cell culture methods without the risk of cell culture contamination.

- / Easy stacking
- / Maximal transparency for excellent microscopic analysis
- / Improved cell adhesion through physical surface treatment



Cell Culture Dishes

- / Improved cell adhesion through physical surface treatment
- / Maximal transparency for excellent microscopic analysis
- / Sterile and user-friendly packaging
- / Easy stacking

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Surface treatment: TC, Vent nock: yes, Sterile: +

Item no.	Height	Ø nominal size	Growth area	Working volume	Total volume	Triple-packed	Sterile	Qty. inner / outer
627160	10 mm	35 mm	8.7 cm ²	≥3 ml	10 ml		+	10 / 740
628160	15 mm	60 mm	21 cm ²	6 ml - 7 ml	28 ml		+	10 / 600
628160-TRI	15 mm	60 mm	21 cm ²	6 ml - 7 ml	28 ml	yes	+	10 / 300
664160	20 mm	100 mm	58 cm ²	16 ml - 17 ml	100 ml		+	15 / 360
664160-TRI	20 mm	100 mm	58 cm ²	16 ml - 17 ml	100 ml	yes	+	15 / 180
639160	20 mm	145 mm	143 cm ²	25 ml - 27 ml	240 ml		+	5 / 120
639160-TRI	20 mm	145 mm	143 cm ²	25 ml - 27 ml	240 ml	yes	+	5 / 120

/ For exact dimensions of our cell culture dishes, please refer to the product data sheets on our website



Cell Culture Dishes - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.


- / Consistent and even cell attachment
- / Homogeneous and optimised cell growth

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Height: 20 mm, Surface treatment: Advanced TC, Vent nock: yes, Sterile: +

Item no.	Ø nominal size	Growth area	Working volume	Total volume	Sterile	Qty. inner / outer
664960	100 mm	58 cm ²	16 ml - 17 ml	100 ml	+	15 / 360
639960	145 mm	143 cm ²	25 ml - 27 ml	240 ml	+	5 / 120

/ Application Report „Advanced TC for improving the cultivation / differentiation of embryonic stem cells“(F076036)

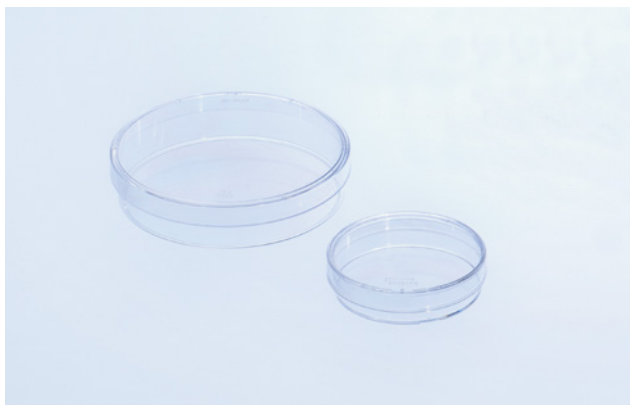


Cell Culture Dishes - CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Surface treatment: CELLCOAT®, Vent neck: yes

Item no.	Height	Ø nominal size	Growth area	Protein coating	Working volume	Total volume	Aseptic	Qty. inner / outer
628950	15 mm	60 mm	21 cm ²	Collagen Type I	6 ml - 7 ml	28 ml	+	20 / 100
664950	20 mm	100 mm	58 cm ²	Collagen Type I	16 ml - 17 ml	100 ml	+	10 / 40
628940	15 mm	60 mm	21 cm ²	Poly-D-Lysine	6 ml - 7 ml	28 ml	+	20 / 100
664940	20 mm	100 mm	58 cm ²	Poly-D-Lysine	16 ml - 17 ml	100 ml	+	10 / 40
628930	15 mm	60 mm	21 cm ²	Poly-L-Lysine	6 ml - 7 ml	28 ml	+	20 / 100
628920	15 mm	60 mm	21 cm ²	Fibronectin	6 ml - 7 ml	28 ml	+	5 / 20
664920	20 mm	100 mm	58 cm ²	Fibronectin	16 ml - 17 ml	100 ml	+	5 / 10
628910	15 mm	60 mm	21 cm ²	Laminin	6 ml - 7 ml	28 ml	+	5 / 20
664910	20 mm	100 mm	58 cm ²	Laminin	16 ml - 17 ml	100 ml	+	10 / 10

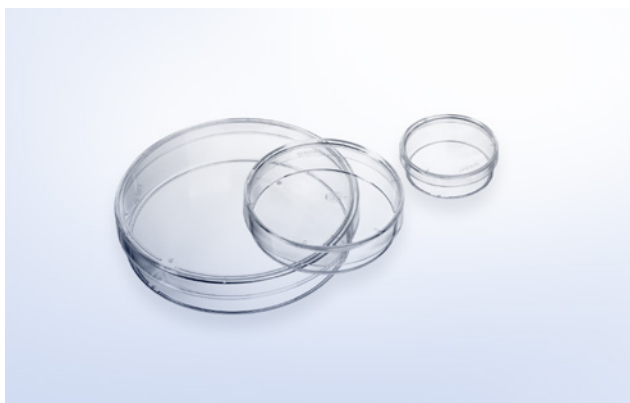


**Cell Culture Dishes -
Cell-Repellent Surface**

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Raw material: PS, Surface treatment: cell-repellent, Vent neck: yes, Sterile: +

Item no.	Height	Ø nominal size	Working volume	Total volume	Sterile	Qty. inner / outer
627979	10 mm	35 mm	≤3 ml	10 ml	+	10 / 40
628979	15 mm	60 mm	6 ml - 7 ml	28 ml	+	10 / 20
664970	20 mm	100 mm	16 ml - 17 ml	100 ml	+	1 / 5

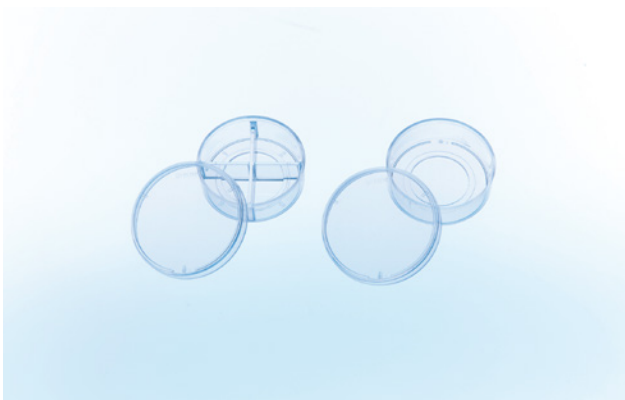

CELLview Dish

The CELLview cell culture dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of *in-vitro* cultivated cultures.

/ Embedded glass bottom for maximal planarity


STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic

 non-
pyrogenic

Ø: 35 mm, Bottom: glass, Vent nock: yes, Sterile: +

Item no.	Compartment	Growth area	Growth area / unit	Surface treatment	Working volume	Total volume	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
627860	1	8.7 cm ²		TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627965	1	8.7 cm ²		Advanced TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627870	4		1.9 cm ²	TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40
627975	4		1.9 cm ²	Advanced TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40

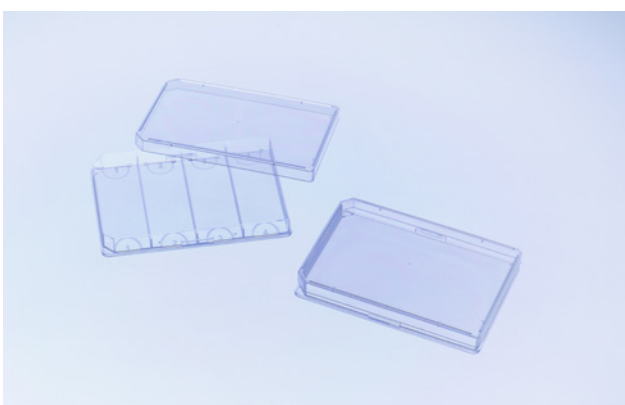


OneWell Plate / FourWell Plate

- / For the cultivation of a large number of cells the OneWell Plate with ANSI standard dimensions offers the best possible conditions.
- / The FourWell Plates allows users to culture cells on slides while remaining in accordance with an HTS-compatible plate. This four-chambered plate is particularly suitable for parallel experiments.

Height: 14.4 mm, Length: 127.8 mm, Width: 85.5 mm, Lid: yes, Vent nock: yes

Item no.	Well format	Growth area	Surface treatment	Product color	Total volume (well)	Sterile	Qty. inner / outer
96077307	4		suspension	○ clear	18.6 ml	+	8 / 32
670190	1	95 cm ²	suspension	○ clear	113.7 ml		8 / 32
670180	1	95 cm ²	TC	○ clear	113.7 ml	+	8 / 32

/ Samples are available on request.

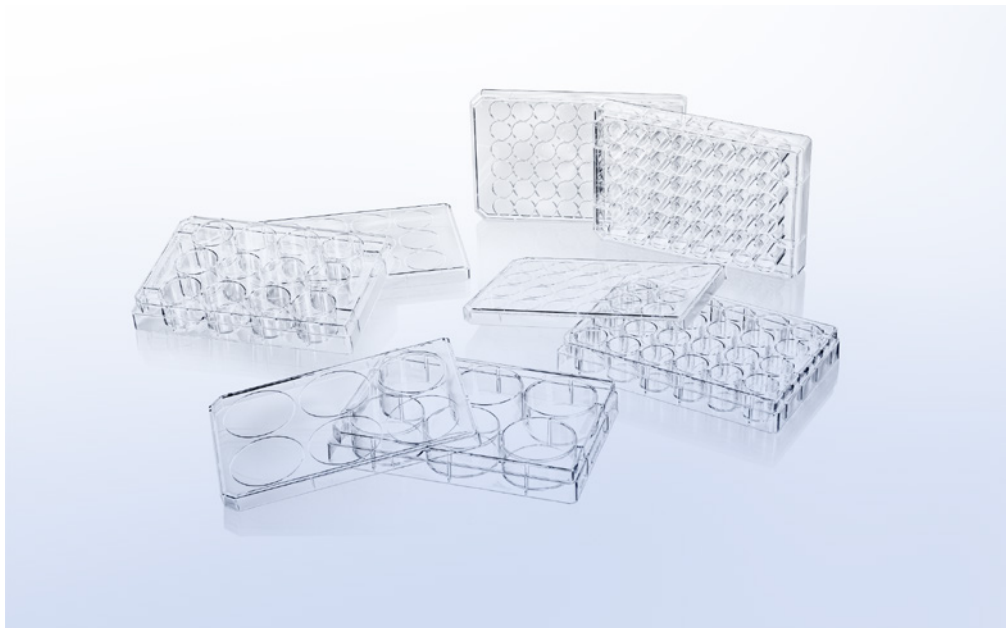

/ Learn more about our Cell Culture Multiwell Plates:




CELLSTAR® CELL CULTURE MULTIWELL PLATES

For both adherent and suspension cell culture, choose Greiner Bio-One's multiwell plates in either 6, 12, 24 or 48 well format.

The special chimney wells in conjunction with low evaporation lids significantly reduce the risk of cross-contamination between the wells. The plates are easily stacked, both with and without lids. Due to their advanced design, the pla-

tes still exhibit important features like constant heat and gas exchange even when stacked. Uniform plate dimensions (well depth, diameter and distance between the wells) comply with ANSI standards, enabling easy manual and automated handling. Each plate is sterile, free of detectable DNase, RNase, human DNA and is non-pyrogenic.

- / With hydrophilic surface (TC surface treatment) for improved cell adhesion
- / With hydrophobic surface for suspension cultures and hybridoma cells
- / High clarity and low autofluorescence
- / Alphanumeric well coding

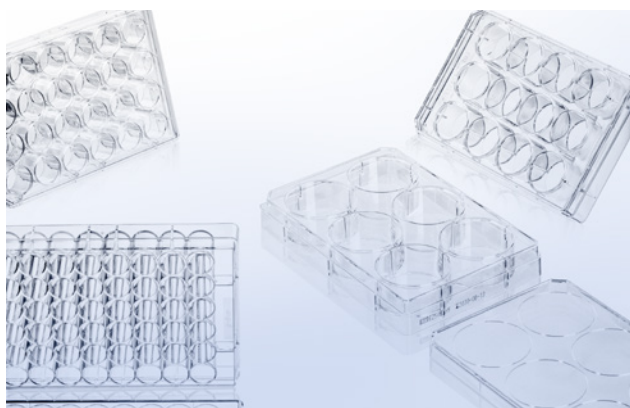

Multiwell Plates

6 / 12 / 24 / 48 Well Format

- / High transparency
- / Alphanumeric well coding
- / With hydrophilic surface (TC surface treatment) for improved cell adhesion
- / With hydrophobic surface for suspension cultures and hybridoma cells



Bottom: solid, Raw material: PS, Lid: yes, condensation rings, Sterile: +

Item no.	Well format	Growth area / unit	Surface treatment	Product color	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657160	6	9.6 cm ²	TC	○ clear	16 ml	2 ml - 5 ml	+	1 / 100
665180	12	3.9 cm ²	TC	○ clear	6.5 ml	2 ml - 4 ml	+	1 / 100
662160	24	1.9 cm ²	TC	○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1 / 100
677180	48	1 cm ²	TC	○ clear	1.7 ml	0.5 ml - 1 ml	+	1 / 100
657185	6		suspension	○ clear	16 ml	2 ml - 5 ml	+	1 / 100
665102	12		suspension	○ clear	6.5 ml	2 ml - 4 ml	+	1 / 100
662102	24		suspension	○ clear	3.3 ml		+	1 / 100
677102	48		suspension	○ clear	1.7 ml	0.5 ml - 1 ml	+	1 / 100



Multiwell Plates - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

- / Consistent and even cell attachment




Bottom: solid, Raw material: PS, Surface treatment: Advanced TC, Lid: yes, condensation rings, Sterile: +

Item no.	Well format	Growth area / unit	Product color	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657960	6	9.6 cm ²	○ clear	16 ml	2 ml - 5 ml	+	1 / 100
665980	12	3.9 cm ²	○ clear	6.5 ml	2 ml - 4 ml	+	1 / 100

Item no.	Well format	Growth area / unit	Product color	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
662960	24	1.9 cm ²	○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1 / 100

/ Application Report „Advanced TC for improving the cultivation / differentiation of embryonic stem cells“(F076036)

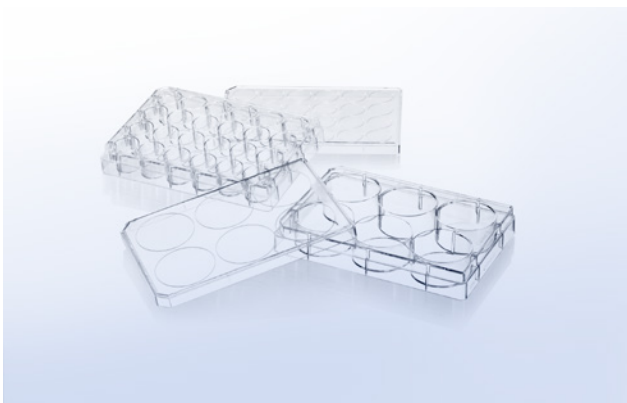


Multiwell Plates - CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells



Bottom: solid, Raw material: PS, Surface treatment: CELLCOAT®, Lid: yes, condensation rings

Item no.	Well format	Growth area / unit	Protein coating	Product color	Total volume (well)	Working volume (well)	Aseptic	Qty. inner / outer
657950	6	9.6 cm ²	Collagen Type I	○ clear	16.1 ml	2 ml - 5 ml	+	5 / 50
662950	24	1.9 cm ²	Collagen Type I	○ clear	3.3 ml	0.5 ml - 1 ml	+	5 / 50
657940	6	9.6 cm ²	Poly-D-Lysine	○ clear	16.1 ml	2 ml - 5 ml	+	5 / 50
662940	24	1.9 cm ²	Poly-D-Lysine	○ clear	3.3 ml	0.5 ml - 1 ml	+	5 / 50
657930	6	9.6 cm ²	Poly-L-Lysine	○ clear	16.1 ml	2 ml - 5 ml	+	5 / 50
662930	24	1.9 cm ²	Poly-L-Lysine	○ clear	3.3 ml	0.5 ml - 1 ml	+	5 / 50

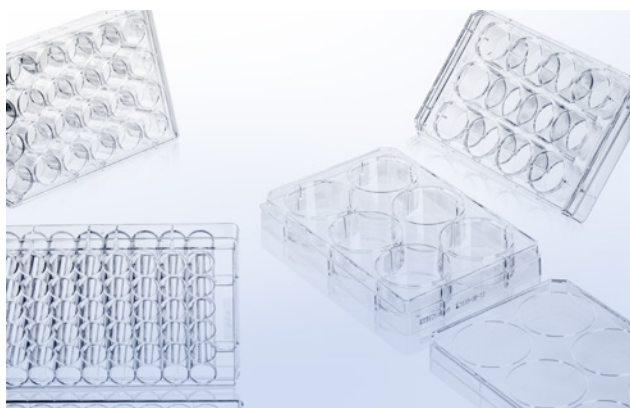

Multiwell Plates - Cell-Repellent Surface

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic

 non-
pyrogenic

Well profile: F-bottom, Bottom: solid, Raw material: PS, Surface treatment: cell-repellent, Lid: yes, condensation rings,
Sterile: +

Item no.	Well format	Product color	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
657970	6	○ clear	16.1 ml	2 ml - 5 ml	+	1 / 5
665970	12	○ clear	6.5 ml	2 ml - 4 ml	+	1 / 5
662970	24	○ clear	3.3 ml	0.5 ml - 1.5 ml	+	1 / 5
677970	48	○ clear	1.7 ml	0.5 ml - 1 ml	+	1 / 5

/ Forum No. 17: CELLSTAR® Cell Culture Vessels with Cell-Repellent Surface (F073777)

/ Application Report „Advantage of CELLSTAR® Cell Culture Vessels with Cell-Repellent Surface for 3-D Cell Culture in Hydrogels“ (F073792)



- / Forum No. 8: ThinCert® cell culture products - Overview (F073017)
- / Application Note „Immuno cytochemistry“(F073100)
- / For further information and application notes please refer to our website: www.gbo.com

THINCERT® CELL CULTURE INSERTS FOR 6, 12 AND 24 WELL MULTIWELL PLATES

For advanced cell and tissue culture applications, Greiner Bio-One offers a broad range of membrane supports - ThinCert®.

Combining 6 different membrane specifications (pore size and density) in geometries to fit 6, 12 and 24 well plates, the ThinCert® cell culture inserts are suitable for a wide range of applications including transport, secretion and diffusion studies, migrational experiments, cytotoxicity testing, co-cultures, trans epithelial electric resistance (TEER) measurements, as well as pri-

mary cell culture.

ThinCert® cell culture inserts are compatible with standard CELLSTAR® multiwell plates from Greiner Bio-One, and are pre-packed together with the requisite number of plates. The automated production process includes double optical control of each insert produced, ensuring that any biological contamination is avoided. The sterility of the single blisterpacked inserts and multiwell plates is ensured by irradiation.

- / Stable clear polystyrene housing
- / Sealed PET capillary pore membrane
- / Pre-configured multiwell plates with ThinCert® cell culture inserts available on request



ThinCert® Cell Culture Inserts

6 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic

 non-
pyrogenic

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 24.85 mm, Ø external: 27.85 mm, Cultural surface: 452.4 mm², Surface treatment: TC, Working volume (ThinCert®): 1 ml - 3.6 ml, Working volume (well): 2 ml - 4.15 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
657640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 24
657641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 24
657610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 24
657630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 24
657631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 24
657638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 24



ThinCert® Cell Culture Inserts

12 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

STERILE

FREE OF
detectable
DNase

FREE OF
detectable
human DNA


FREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 13.85 mm, Ø external: 15.85 mm, Cultural surface: 113.1 mm², Surface treatment: TC, Working volume (ThinCert®): 0.3 ml - 1 ml, Working volume (well): 1 ml - 2 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
665640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 48
665641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 48
665610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 48
665630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 48
665631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 48
665638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 48

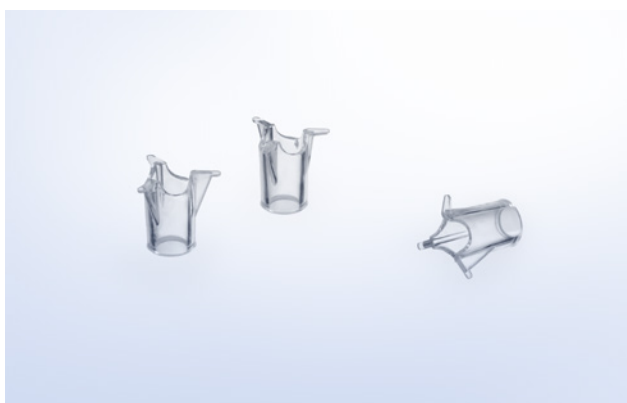**ThinCert® Cell Culture Inserts****24 Well**

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange


STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic

 non-
pyrogenic

Feature: 2 multiwell plates / box, Height: 16.25 mm, Ø internal: 8.4 mm, Ø external: 10.4 mm, Cultural surface: 33.6 mm², Surface treatment: TC, Working volume (ThinCert®): 0.1 ml - 0.35 ml, Working volume (well): 0.4 ml - 1.2 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
662640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 48
662641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 48
662610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 48
662630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 48
662631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 48
662638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 48

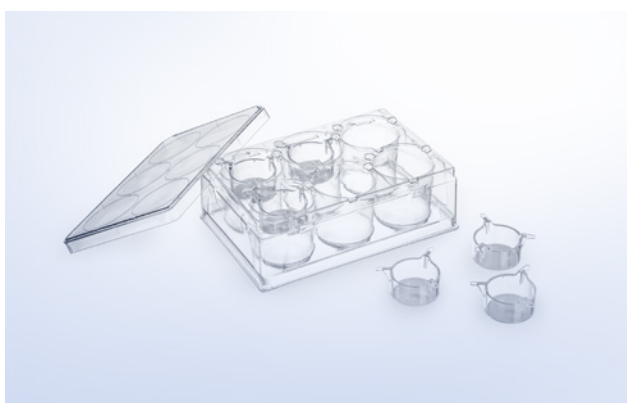**ThinCert® Plate****6 / 12 Well**

- / Optimised for use with ThinCert® cell culture inserts
- / Deep wells for an increased volume of medium in air-lift culture
- / Notches for fixed insert position
- / Available in 6 and 12 well format

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic

 non-
pyrogenic

Height: 39.5 mm, Length: 129.5 mm, Width: 86.6 mm, Lid: yes, condensation rings, Sterile: +

Item no.	Well format	Working volume (well)	Sterile	Qty. inner / outer
657110	6	≤20 ml	+	1 / 50
665110	12	≤4 ml	+	1 / 60



- / Application Note "High-throughput in-vitro airway modelling with ThinCert® 96 Well HTS Insert" (F073140)
- / Brochure «ThinCert® 96 Well HTS Insert» (F073149)



THINCERT® 96 WELL HTS INSERT

ThinCert® 96 Well HTS Insert is the latest high-throughput addition to our ThinCert® family which opens doors to automation and offers an efficient solution to scale up your membrane-based cell culture application. Take advantage of the assay miniaturisation to maximise both time and cost savings when compared to single-insert solutions.

- / 96 well system for high-throughput applications
- / Automation-friendly geometry
- / Pore sizes of 0.4 µm, 3 µm and 8 µm covering a range of different applications
- / High permeability and transparency for optimal diffusion rates and optical features

ThinCert® 96 Well HTS Inserts offer a diverse selection of membrane specifications (pore size and density) within a 96 well format. The automation-friendly inserts are highly versatile and can be utilised for various applications such as transport, secretion, and diffusion studies, migration/invasion experiments, co-cultures, cytotoxicity tests, TEER measurements and

the generation of tissue models. The two-component system comprises a 96 well porous membrane plate manufactured entirely from polycarbonate (PC) and a corresponding polystyrene (PS) receiver plate. The tissue culture treated PC membrane allows for an ideal exchange of nutrients and substances, creating in vivo-like cultivation conditions to facilitate optimal cell growth, monolayer formation, and tissue differentiation. For preparatory cell culture, particularly for transport and permeability studies, a single-well feeder plate is available. Using the two feeder ports in the membrane plate allows for a one-step media change, effectively minimising both handling effort and the potential for contamination.



ThinCert® 96 Well HTS Insert (Membrane and Receiver Plates)

- / Pore sizes of 0.4 µm, 3 µm and 8 µm covering a range of different applications
- / High permeability and transparency for optimal diffusion rates and optical features
- / Working volume per well (membrane plate): 15 - 160 µl
- / Working volume per well (receiver plate): 120 - 300 µl
- / ACT® Ecolabel certification

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic


 non-
pyrogenic

 my green lab.
ACT Ecolabel
NEW
+

Well format: 96, Surface treatment: TC, Lid: yes, condensation rings, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
655630	2 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 5
655640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 5
655641	2 x 10 ⁷ / cm ²	0.4 µm	optimized transparency	+	1 / 5
655680	1 x 10 ⁵ / cm ²	8 µm	clear	+	1 / 5

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

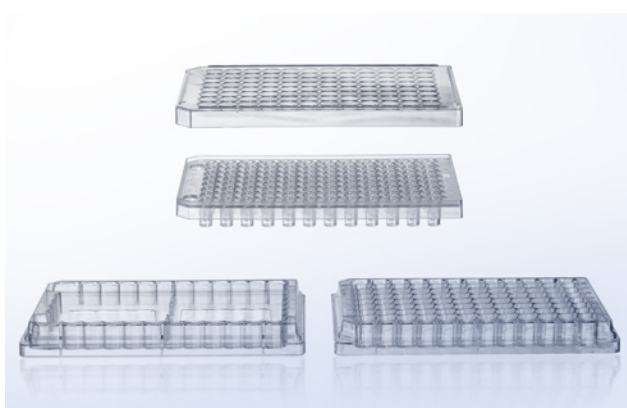

ThinCert® 96 Well HTS Insert (Membrane, Receiver and Feeder Plates)

- / Pore sizes of 0.4 µm, 3 µm and 8 µm covering a range of different applications
- / Working volume per well (membrane plate): 15 - 160 µl
- / Working volume per well (receiver plate): 120 - 300 µl
- / Additional Feeder Plate for simultaneous feeding and homogeneous conditions for all wells (Working volume: 20 - 30 ml)
- / ACT® Ecolabel certification

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
cytotoxic


 non-
pyrogenic

 my green lab.
ACT Ecolabel
NEW
+

Well format: 96, Surface treatment: TC, Lid: yes, condensation rings, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
655642	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 3
655643	2 x 10 ⁷ / cm ²	0.4 µm	optimised transparency	+	1 / 3
655644	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 3

- / Item no. 655642 and 655643 include a Feeder Plate with cross-shaped media stabiliser.
- / Item no. 655644 includes a Feeder Plate without media stabiliser.
- / All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

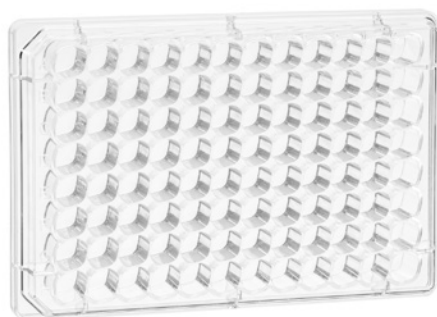

**Receiver Plates for
ThinCert® 96 Well HTS Insert**

- / Reduced wicking effect
- / High permeability and transparency for optimal diffusion rates and optical features
- / Working volume per well (receiver plate): 120 - 300 µl

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
NEW +					

Lid: yes, condensation rings, Sterile: +

Item no.	Growth area	Surface treatment	Sterile	Qty. inner / outer
655169		untreated	+	8 / 32
655167	53 mm ²	TC	+	8 / 32



**Feeder Plates for
ThinCert® 96 Well HTS Insert**

- / Simultaneous feeding of all wells
- / Homogeneous conditions for all wells
- / Reduced risk of wicking during cultivation

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic

Working volume: 20 - 30 ml, Lid: yes, condensation rings, Sterile: +

Item no.	Sterile	Qty. inner / outer
670640	+	1 / 5
670641	+	1 / 5

- / Item no. 670640 with cross-shaped media stabiliser.
- / Item no. 670641 without media stabiliser.



- / Selected cell culture microplates are certified with ACT® Ecolabel. Learn more:



CELLSTAR®

CELL CULTURE MICROPLATES


Cell culture treated microplates are available in the following versions: 96, 384, 1536 well format.

Depending on the application, the well profile is a key feature in a 96 well cell culture microplate. The chimney well cell culture microplate has the same well profile as the standard F-bottom plate. The difference to the standard plate is the chimney-like arrangement of the wells i.e. each well stands on its own. Therefore the risk of contamination from sample material being carried over is minimised. Clear-bottom microplates have pig-

mented walls and a transparent thin film bottom, the so-called μ Clear® bottom. In contrast to our standard microplates with a solid polystyrene bottom, they are ideal for cell culture and microscopic applications using fluorescence or luminescence detection methods.

For many applications, a reduction of the sample volume is an important feature. 96 well half area microplates offer an interesting alternative here. They can be pipetted automatically as well as manually without any problem and allow a reduction of the sample volume up to 50 %.

- / Available with different cell culture surfaces for optimal cell culture conditions
- / Footprint compatible with automated systems
- / Alphanumeric well coding

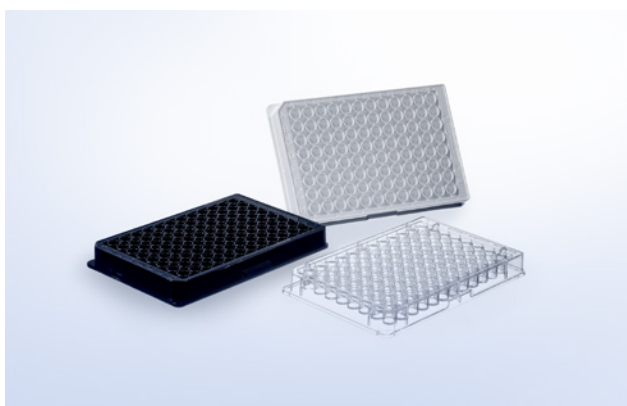

Cell Culture Microplates

96 Well

- / With U-bottom, V-bottom or F-bottom
- / Clear / black / white
- / Chimney well design, raised wells and condensation rings in lids prevent cross-contamination
- / Improved cell adhesion through physical surface treatment
- / Selected product versions certified with ACT® Ecolabel (*)

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

 my green lab.
ACT Ecolabel


Well format: 96, Bottom: solid, Raw material: PS, Surface treatment: TC, Sterile: +

Item no.	Growth area / unit	Well profile	Product color	Working volume (well)	Lid	Sterile	Qty. inner / outer
650160	35 mm ²	U-bottom	○ clear	40 µl - 280 µl	no	+	1 / 100
650180*	35 mm ²	U-bottom	○ clear	40 µl - 280 µl	yes	+	1 / 100
651160	28 mm ²	V-bottom	○ clear	40 µl - 200 µl	no	+	1 / 100
651180*	28 mm ²	V-bottom	○ clear	40 µl - 200 µl	yes	+	1 / 100
655160	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	no	+	1 / 100
655162	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	no	+	5 / 100
655180*	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	yes, condensation rings	+	1 / 100
655182*	34 mm ²	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	yes, condensation rings	+	10 / 160
655073	34 mm ²	F-bottom / Chimney Well	○ white	25 µl - 340 µl	no	+	10 / 40
655083*	34 mm ²	F-bottom / Chimney Well	○ white	25 µl - 340 µl	yes, condensation rings	+	8 / 32
655079	34 mm ²	F-bottom / Chimney Well	● black	25 µl - 340 µl	no	+	10 / 40
655086*	34 mm ²	F-bottom / Chimney Well	● black	25 µl - 340 µl	yes, condensation rings	+	8 / 32

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

/ For selected products Greiner Bio-One also offers user-friendly bulk packaging.

/ Barcode labelling on request

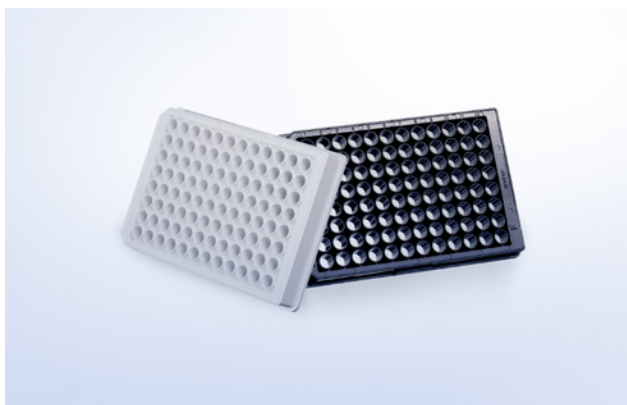

Cell Culture Microplates

96 Well - µClear®

Clear-bottom microplates have pigmented walls and a transparent thin film bottom, the so-called µClear® bottom. They are ideal for cell culture and microscopic applications using fluorescence or luminescence detection methods. The plates are certified with ACT® Ecolabel.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------

Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Bottom: µClear®, Raw material: PS, Surface treatment: TC, Working volume (well): 25 µl - 340 µl, Sterile: +

Item no.	Product color	Lid	Sterile	Qty. inner / outer
655088	○ white	no	+	10 / 40
655098	○ white	yes, condensation rings	+	8 / 32
655087	● black	no	+	10 / 40
655090	● black	yes, condensation rings	+	8 / 32

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



Cell Culture Microplates

96 Well - Half Area

For many applications, a reduction of the sample volume is an important feature. 96 well half area microplates offer an interesting alternative here. They can be pipetted automatically as well as manually without any problem and allow a reduction of the sample volume up to 50 %.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------


Well format: 96, Growth area / unit: 15 mm², Well profile: F-bottom, Raw material: PS, Surface treatment: TC, Plate design: half area, Working volume (well): 15 µl - 175 µl, Lid: yes, Sterile: +

Item no.	Bottom	Product color	Sterile	Qty. inner / outer
675180	solid	○ clear	+	8 / 32
675083	solid	○ white	+	8 / 32
675086	solid	● black	+	8 / 32

Item no.	Bottom	Product color	Sterile	Qty. inner / outer
675090*	µClear®	● black	+	8 / 32

/ For selected products Greiner Bio-One also offers user-friendly bulk packaging

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>.

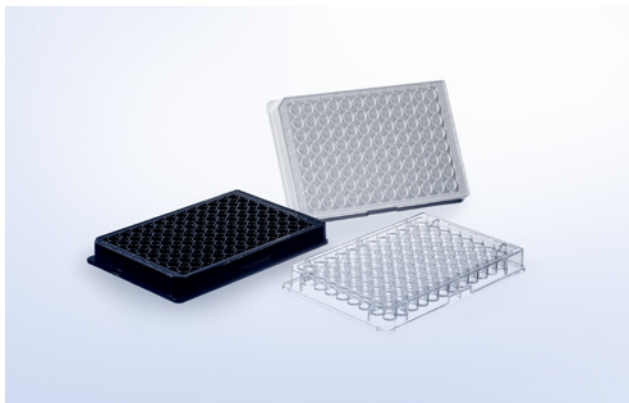

Cell Culture Microplates

96 Well - CELLCOAT®

- / Improved adhesion and cell proliferation
- / Reduced-serum or serum-free cultivation
- / Improved growth of primary cells

Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Raw material: PS, Surface treatment: CELLCOAT®, Working volume (well): 25 µl - 340 µl, Lid: yes, condensation rings

Item no.	Bottom	Protein coating	Product color	Aseptic	Qty. inner / outer
655950	solid	Collagen Type I	○ clear	+	5 / 20
655956	µClear®	Collagen Type I	● black	+	5 / 20
655940	solid	Poly-D-Lysine	○ clear	+	5 / 20
655944	µClear®	Poly-D-Lysine	○ white	+	5 / 20
655946	µClear®	Poly-D-Lysine	● black	+	5 / 20
655948	µClear®	Poly-D-Lysine	● black	+	20 / 120
655930	solid	Poly-L-Lysine	○ clear	+	5 / 20
655936	µClear®	Poly-L-Lysine	● black	+	5 / 20



Cell Culture Microplates

96 Well - Advanced TC

The Advanced TC surface provides optimal conditions for the cultivation of sensitive and fastidious cells or the usage of restricted growth conditions.

- / Consistent and even cell attachment


STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Well format: 96, Growth area / unit: 34 mm², Well profile: F-bottom / Chimney Well, Raw material: PS, Surface treatment: Advanced TC, Working volume (well): 25 µl - 340 µl, Lid: yes, condensation rings, Sterile: +

Item no.	Bottom	Product color	Sterile	Qty. inner / outer
655980	solid	○ clear	+	1 / 100
655983	µClear®	○ white	+	8 / 32
655986	µClear®	● black	+	8 / 32

Suspension Culture Microplates

96 Well

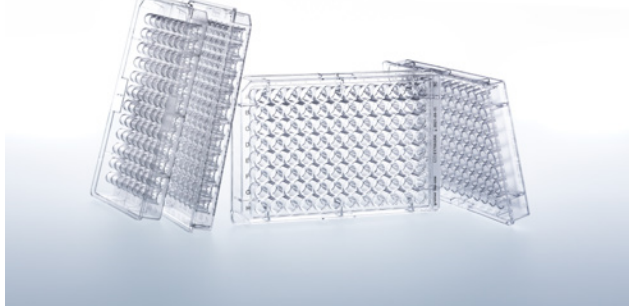


- / Hydrophobic surface, ideal for suspension cultures, hybridoma and embryonic stem cells
- / Solid bottom
- / ACT® Ecolabel certification

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------

Well format: 96, Bottom: solid, Raw material: PS, Surface treatment: suspension, Sterile: +

Item no.	Well profile	Product color	Working volume (well)	Lid	Sterile	Qty. inner / outer
650185	U-bottom	○ clear	40 µl - 280 µl	yes	+	60 / 240
655185	F-bottom / Chimney Well	○ clear	25 µl - 340 µl	yes, condensation rings	+	60 / 240

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

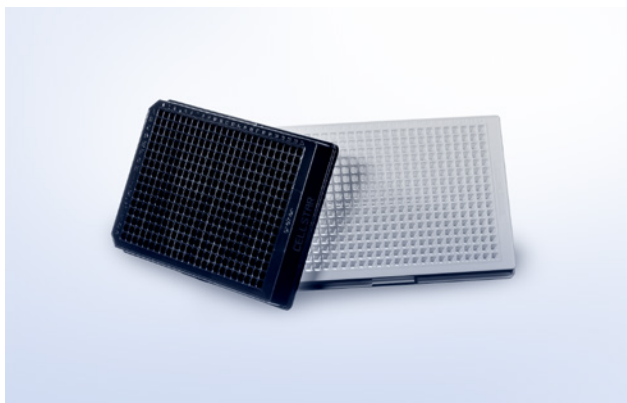

Cell Culture Microplates

384 Well

- / Clear / black / white
- / Solid bottom or μ Clear® film bottom
- / Barcode labelling on request
- / Alphanumeric well coding
- / Selected product versions certified with ACT® Ecolabel (*)



Well format: 384, Growth area / unit: 10 mm², Well profile: F-bottom, Raw material: PS, Working volume (well): 15 μ l - 110 μ l

Item no.	Bottom	Surface treatment	Protein coating	Product color	Total volume (well)	Lid	Sterile	Aseptic	Qty. inner / outer
781165	solid	TC		○ clear		no	+		10 / 40
781182	solid	TC		○ clear		yes	+		8 / 32
781073	solid	TC		○ white		no	+		10 / 40
781080	solid	TC		○ white		yes	+		8 / 32
781079	solid	TC		● black		no	+		10 / 40
781086	solid	TC		● black		yes	+		8 / 32
781093*	μ Clear®	TC		○ white		no	+		10 / 40
781098*	μ Clear®	TC		○ white		yes	+		8 / 32
781092*	μ Clear®	TC		● black		no	+		10 / 40
781091*	μ Clear®	TC		● black		yes	+		8 / 32
781090*	μ Clear®	TC		● black		yes	+		20 / 120
781950	solid	CELLCOAT®	Collagen Type I	○ clear	131 μ l	yes		+	5 / 20
781956	μ Clear®	CELLCOAT®	Collagen Type I	● black	131 μ l	yes		+	5 / 20
781940	solid	CELLCOAT®	Poly-D-Lysine	○ clear	131 μ l	yes		+	5 / 20
781944	μ Clear®	CELLCOAT®	Poly-D-Lysine	○ white	131 μ l	yes		+	5 / 20
781946	μ Clear®	CELLCOAT®	Poly-D-Lysine	● black	131 μ l	yes		+	5 / 20
781948	μ Clear®	CELLCOAT®	Poly-D-Lysine	● black	131 μ l	yes		+	20 / 120
781936	μ Clear®	CELLCOAT®	Poly-L-Lysine	● black	131 μ l	yes		+	5 / 20
781983	μ Clear®	Advanced TC		○ white	131 μ l	yes	+		8 / 32
781986	μ Clear®	Advanced TC		● black	131 μ l	yes	+		8 / 32

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>




Cell Culture Microplates
384 Well - Small Volume

HiBase

- / Perfect for top reading even at low working volumes
- / Savings in reagent similar to 1536 well microplates
- / Made of black / white polystyrene for fluorescence or luminescence measurements

Well format: 384, Growth area / unit: 2.7 mm², Well profile: F-bottom, Bottom: solid, Raw material: PS, Plate geometry: HiBase, Plate design: Small Volume, Working volume (well): 4 µl - 25 µl, Lid: yes

Item no.	Surface treatment	Protein coating	Product color	Sterile	Aseptic	Qty. inner / outer
784080	TC		○ white	+		8 / 32
784946	CELLCOAT®	Poly-D-Lysine	● black		+	5 / 30

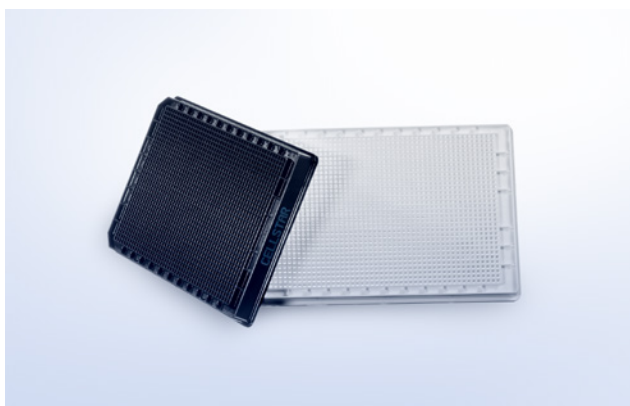

Cell Culture Microplates
1536 Well


- / Clear / black / white
- / Solid bottom or µClear® film bottom
- / Barcode labelling on request
- / Selected product versions certified with ACT® Ecolabel (*)



Well format: 1536, Growth area / unit: 2.3 mm², Well profile: F-bottom, Raw material: PS, Plate geometry: HiBase, Working volume (well): 3 µl - 10 µl

Item no.	Bottom	Surface treatment	Protein coating	Product color	Lid	Sterile	Aseptic	Qty. inner / outer
782180	solid	TC		○ clear	yes	+		1 / 32
782073	solid	TC		○ white	no	+		15 / 60
782080	solid	TC		○ white	yes	+		10 / 40
782078	solid	TC		● black	no	+		15 / 60
782086	solid	TC		● black	yes	+		10 / 40
782093*	µClear®	TC		○ white	no	+		15 / 60
782092*	µClear®	TC		● black	no	+		15 / 60
782946	µClear®	CELLCOAT®	Poly-D-Lysine	● black	yes		+	5 / 20

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

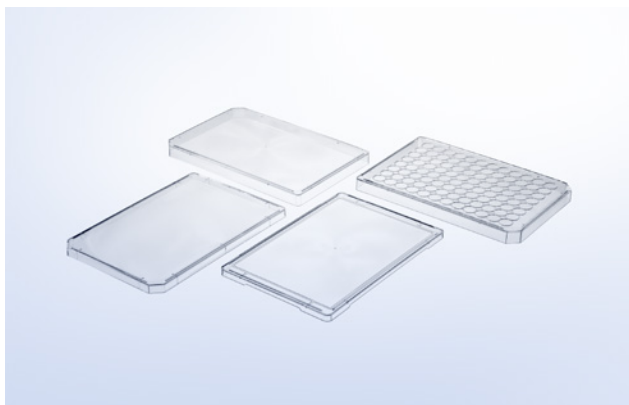

Lids

All sterile lids are non-cytotoxic.

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	 non-pyrogenic
--------------------------	------------------------------	--------------------------	---

Description: Lid, Raw material: PS

Item no.	Height	Product color	Condensation rings	Lid type	Sterile	Qty. inner / outer
656101	9 mm	○ clear	no	high		1 / 100
656161	9 mm	○ clear	no	high	+	1 / 100
656170	9 mm	○ clear	yes	high		1 / 100
656171	9 mm	○ clear	yes	high	+	1 / 100
656190	6 mm	○ clear	no	flat		20 / 200
656191	6 mm	○ clear	no	flat	+	20 / 200
691101	4.8 mm	○ clear	no	ultra low		25 / 100
691161	4.8 mm	○ clear	no	ultra low	+	25 / 100



/ Learn more about our Cell Culture Tubes:



CELLSTAR® CELL CULTURE TUBES

- / Bioreactor for suspension and spheroid cells
- / Expansion of aerobic bacteria, yeast and microorganisms
- / Storage of components and liquids requiring gas exchange

CELLSTAR® CELLreactor tube can be used as small bioreactor for suspension and spheroid cell culture, facilitating miniaturisation of large-scale setups and maximising the number of parallel experiments. Each CELLreactor tube cap features several holes and a membrane with a pore size of 0.2 µm to guarantee maximal sterility while providing excellent gas exchange. In case the aeration has to be reduced, individual openings can be sealed.

Agitation of internal liquids is achieved with standard shaking lab equipment minimising foam formation and shearing forces

induced by integrated mixing devices. Compared to cell culture and spinner flask as well as other cultivation disposables, no transfer for cell harvest is required. Based on the conical design, the tubes fit in standard 15 ml / 50 ml centrifuge rotors and cells can be spun down in the same tube. In addition to cell culture applications, the CELLSTAR® CELLreactor tube can also be applied for the expansion of aerobic bacteria, yeast or other microorganisms in shaken cultures as well as storage of components and liquids requiring gas exchange.

**CELLreactor**

15 ml and 50 ml polypropylene tube with filter screw cap

- / For cultivation of suspension cells and expansion of aerobic microorganisms
- / Facilitates a high number of parallel experiments
- / Maximal sterility and excellent gas exchange
- / Conical tube design and in-tube harvest

STERILE

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Cap design: filter screw cap, Sterile: +

Item no.	Height	Ø	Cap color	Working volume	Nominal volume	Sterile	Qty. inner / outer
188241	120 mm	17 mm	● blue	1 ml - 5 ml	15 ml	+	20 / 500
227245	115 mm	30 mm	● blue	1 ml - 35 ml	50 ml	+	20 / 500



/ Watch our video:



EASYSTRAINER CELL STRAINERS

With EASYstrainer cell strainers, Greiner Bio-One offers an innovative and user-friendly solution for the filtration of cell suspensions.

EASYstrainer can be used, for example, after enzymatic tissue digestion for primary cell isolation or for cell preparation prior to flow cytometry. The large cell strainers fit on all conical 50 ml tubes and are available with mesh sizes of 40, 70 and 100 μm . EASYstrainer Small fits 15 ml tubes as well as smaller tubes and reaction vessels. It is available with mesh sizes of 20, 40, 70 and 100 μm .

The stackability of EASYstrainer with different mesh sizes allows for the separation of cells with different sizes in one step. In addition, the upper part of EASYstrainer Small can be inverted. This allows retained cells to be rinsed out for further use. EASYstrainer makes working aseptically much easier: It can be held via a surrounding rim or a handle in order to avoid accidental contact with the sterile filter material. Additional safety is provided by the transparent blister packaging from which the strainers can be conveniently and aseptically removed.

- / Flexible fit, suitable for tubes from 1.5 ml to 50 ml
- / Available Mesh Sizes: 20, 40, 70 und 100 μm
- / Innovative ventilation of cell strainers prevents liquid overspill



EASYstrainer

- / Flexible fit, suitable for tubes from 1.5 ml to 50 ml
- / Handle and ridged skirt for improved aseptic handling
- / Venting slot enables fast filtration and prevents liquid overspill
- / Single packaging reduces risk of contamination

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
 recommended					

Sterile: +

Item no.	Description	Mesh size	Product color	Sterile	Qty. inner / outer
542040	EASYstrainer for 50 ml tubes	40 µm	● green	+	1 / 50
542070	EASYstrainer for 50 ml tubes	70 µm	● blue	+	1 / 50
542000	EASYstrainer for 50 ml tubes	100 µm	● yellow	+	1 / 50
542120	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	20 µm	● red	+	1 / 50
542140	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	40 µm	● green	+	1 / 50
542170	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	70 µm	● blue	+	1 / 50
542100	EASYstrainer Small for tubes: 1.5 / 5 / 15 ml and 12x75 mm	100 µm	● yellow	+	1 / 50



/ Learn more about our media bottles:



MEDIA BOTTLES

- / In three sizes available (100 ml, 500 ml, 1,000 ml)
- / Screw cap with or without original ring
- / Graduation on both sides
- / Triple-packed without box tray for GMP compliant workflow

Media bottles from Greiner Bio-One have been developed for storing a wide variety of solutions and, thanks to their square design, enable space-saving use of the available storage area.

The use of high-quality PET as a

raw material offers exceptional transparency and high resilience, even at low storage temperatures. The optional tamper-evident screw cap eliminates the need for shrink wrap after filling and offers additional convenience and safety.



Media Bottles

- / USP Class VI certified raw material
- / Sterility Assurance Level SAL 10⁻⁶
- / Temperature range -80 °C to +60 °C
- / Leak proofness according to IATA 95 kPa
- / Manufactured in the EU

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
NEW					

Flask design: tetragonal, Raw material: PET, Sterile: +

Item no.	Cap color	Total volume	Cap design	Sterile	Qty. inner / outer
950710	○ white	500 ml	screw cap	+	25 / 50
951710	○ white	100 ml	screw cap	+	49 / 98
952710	○ white	1,000 ml	screw cap	+	12 / 24
950720	○ white	500 ml	tamper-evident	+	25 / 50
951720	○ white	100 ml	tamper-evident	+	49 / 98
952720	○ white	1,000 ml	tamper-evident	+	12 / 24



The cultivation of cells as mass cultures has become increasingly important over the past few decades. Mass cell cultures are mainly used for the production of vaccines or recombinant proteins for therapeutic approaches.

MASS CELL CULTURE

/	CELLMASTER Cell Culture Roller Bottles	76
	Roller Bottles Polystyrene	77
	Roller Bottles Polyethylene Terephthalate(PET).....	78
/	CELLdisc Cell Culture Device	79
	CELLdisc 1 / 4 / 8 / 12 / 16 / 24 / 40 Layers	80
	CELLdisc External Filter.....	81
	CELLdisc Closed Filling Caps.....	82
	CELLstage Filling Accessory.....	83
	CELLlevator	83
	CELLring	84



/ For further information, please refer to our website: www.gbo.com



CELLMASTER CELL CULTURE ROLLER BOTTLES

Roller bottles made of polystyrene or PET range in size from 850 cm² to 4250 cm² growth area and are available either with standard or filter screw caps. The standard screw cap enables a tight closing and contamination-free cultivation. The filter screw cap contains a membrane with a pore size of 0.2 µm guarantees optimal protection against contamination. It provides a high gas exchange rate which enables a mass cultivation of cells using e.g. bicarbonate

buffer and CO₂.

The product range includes a short and a long form, labelled as X and XL, accordingly. Both sizes are available with a smooth or radially ribbed surface. The wave-like design increases the growth area of the roller bottle without changing space requirements. All roller bottles are sterilized by irradiation according to validated procedures (ISO 11137). Endotoxin testing is conducted in accordance with USP 85 with a tolerance level of 0.03 EU/ml.

- / PS or PET roller bottles
- / Different sizes with or without radially ribbed surface
- / Graduations from 200 to 2000 ml
- / Seamless production technique rules out leaking seams



Roller Bottles
Polystyrene

- / For adherent cell culture
- / Manufactured from crystal clear polystyrene
- / Complies with USP Class VI test requirements
- / Lot number and expiry date on each bottle
- / Filter screw cap with 0.2 µm pore size

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Ø: 122 mm, Graduation: yes, Raw material: PS, Surface treatment: TC, Sterile: +

Item no.	Flask design	Height	Growth area	Cap colour	Surface	Total volume	Cap design	Sterile	Qty. inner / outer
680660	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	screw cap	+	2 / 24
680665	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	screw cap	+	24 / 24
680648	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	screw cap	+	24 / 48
680658	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	filter screw cap	+	2 / 24
680668	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	filter screw cap	+	24 / 24
680645	1 X	271 mm	850 cm ²	● blue	smooth	2,520 ml	filter screw cap	+	24 / 48
681670	2.5 X	271 mm	2,125 cm ²	● blue	ribbed	2,300 ml	screw cap	+	2 / 24
681675	2.5 X	271 mm	2,125 cm ²	● blue	ribbed	2,300 ml	screw cap	+	24 / 24
681672	2.5 X	271 mm	2,125 cm ²	● blue	ribbed	2,300 ml	filter screw cap	+	2 / 24
682660	1 XL	500 mm	1,700 cm ²	● blue	smooth	4,970 ml	screw cap	+	1 / 12
682612	1 XL	500 mm	1,700 cm ²	● blue	smooth	4,970 ml	screw cap	+	12 / 12
682624	1 XL	500 mm	1,700 cm ²	● blue	smooth	4,970 ml	screw cap	+	12 / 24
682615	1 XL	500 mm	1,700 cm ²	● blue	smooth	4,970 ml	filter screw cap	+	12 / 12
682625	1 XL	500 mm	1,700 cm ²	● blue	smooth	4,970 ml	filter screw cap	+	12 / 24
682670	5 XL	500 mm	4,250 cm ²	● blue	ribbed	4,640 ml	screw cap	+	1 / 12
682672	5 XL	500 mm	4,250 cm ²	● blue	ribbed	4,640 ml	screw cap	+	12 / 24
682678	5 XL	500 mm	4,250 cm ²	● blue	ribbed	4,640 ml	filter screw cap	+	1 / 12

/ Thread enables quick opening with a 2/3 turn. Screw caps with larger knurls for improved grip and ease of opening / closing.

/ Sterile, individually packed screw caps: standard (item no. 383361) and filter (item no. 383382).



Roller Bottles

Polyethylene Terephthalate (PET)

- / For adherent cell culture
- / Complies with USP Class VI test requirements
- / With standard screw cap

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Flask design: 1 X, Height: 271 mm, Ø: 116 mm, Growth area: 850 cm², Graduation: yes, Raw material: PET, Surface: smooth, Total volume: 2,300 ml, Cap design: screw cap, Sterile: +

Item no.	Cap colour	Sterile	Qty. inner / outer
680190	● blue	+	1 / 18
680195	● blue	+	30 / 30



/ Watch our video:



CELLDISC CELL CULTURE DEVICE

CELLdisc is a multilayer cell culture device covering a range of 250 cm² to 10,000 cm² growth area. The innovative ergonomic design provides a versatile system from research scale to industrial batches with a 40 % higher surface/volume ratio than conventional multilayer systems.

A centrally located ventilation channel assures uniform distribution of gas throughout the device whereas the wide opening port simplifies manual filling. The compact and robust cylindrical device is ideally suited for automation and upscaling of cell cultures.

To guarantee an ideal cell culture environment CELLdisc is provided with two surface treatments: Beside the **TC surface** for standard cells and applications CELLdisc is also available with the **Advanced TC surface**, improving cultivation of sensitive cells, even under restricted growth conditions and increasing cellular adhesion, proliferation and transfection rates.

- / Mass Cell Culture
- / Antibody, virus and vaccine production
- / Production of recombinant or therapeutic proteins



CELLdisc

1 / 4 / 8 / 12 / 16 / 24 / 40 Layers

- / 40 % higher surface/volume ratio than comparable systems
- / Easy operation and minimal space occupation
- / Media exchange without contact to cell layers

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Ø: 200 mm, Raw material: PS, Cap design: screw cap, Sterile: +

Item no.	Layers	Height	Growth area	Surface treatment	Cap colour	Working volume	Sterile	Qty. inner / outer
678101	1	61 mm	250 cm ²	TC	● red	15 ml - 50 ml	+	1 / 8
678104	4	93 mm	1,000 cm ²	TC	● red	60 ml - 200 ml	+	1 / 4
678108	8	135 mm	2,000 cm ²	TC	● red	120 ml - 400 ml	+	1 / 3
678112	12	177 mm	3,000 cm ²	TC	● red	180 ml - 600 ml	+	1 / 2
678116	16	220 mm	4,000 cm ²	TC	● red	240 ml - 800 ml	+	1 / 2
678124	24	304 mm	6,000 cm ²	TC	● red	360 ml - 1,200 ml	+	1 / 2
678140	40	474 mm	10,000 cm ²	TC	● red	600 ml - 2,000 ml	+	1 / 1
678904	4	93 mm	1,000 cm ²	Advanced TC	● blue	60 ml - 200 ml	+	1 / 4
678908	8	135 mm	2,000 cm ²	Advanced TC	● blue	120 ml - 400 ml	+	1 / 3
678912	12	177 mm	3,000 cm ²	Advanced TC	● blue	180 ml - 600 ml	+	1 / 2
678916	16	220 mm	4,000 cm ²	Advanced TC	● blue	240 ml - 800 ml	+	1 / 2
678924	24	304 mm	6,000 cm ²	Advanced TC	● blue	360 ml - 1,200 ml	+	1 / 2
678940	40	474 mm	10,000 cm ²	Advanced TC	● blue	600 ml - 2,000 ml	+	1 / 1



CELLdisc
External Filter

- / Triple-packed CELLdiscs, pre-assembled with external filters
- / USP VI certified venting filters, manufactured with reinforced hydrophobic membranes
- / A defined pore size of 0.2 µm prevents contamination and ensures sufficient airflow to the cells

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Feature: External Filter, Ø: 200 mm, Raw material: PS, Surface treatment: TC, Cap design: screw cap, Sterile: +

Item no.	Layers	Height	Growth area	Cap colour	Working volume	Sterile	Qty. inner / outer
678101-EXF	1	61 mm	250 cm ²	● red	15 ml - 50 ml	+	1 / 4
678104-EXF	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1 / 3
678108-EXF	8	135 mm	2,000 cm ²	● red	120 ml - 400 ml	+	1 / 2
678112-EXF	12	177 mm	3,000 cm ²	● red	180 ml - 600 ml	+	1 / 2
678116-EXF	16	220 mm	4,000 cm ²	● red	240 ml - 800 ml	+	1 / 2
678124-EXF	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1 / 2
678140-EXF	40	474 mm	10,000 cm ²	● red	600 ml - 2,000 ml	+	1 / 1

/ Assembled product is sterilized (SAL 10⁻⁶) to reduce risk of contamination



CELLdisc Closed Filling Caps

- / Triple-packed CELLdiscs, pre-assembled with external filters and closed filling cap
- / Choice of single tubing (CF1) or double tubing (CF2) depending on desired emptying process
- / Tubing with MPC type connector for quick and safe connection /disconnection

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Ø: 200 mm, Raw material: PS, Surface treatment: TC, Cap design: screw cap, Sterile: +

Item no.	Feature	Layers	Height	Growth area	Cap colour	Working volume	Sterile	Qty. inner / outer
678101-CF1	hose connector without Dip-In	1	61 mm	250 cm ²	● red	15 ml - 50 ml	+	1 / 4
678104-CF1	hose connector without Dip-In	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1 / 3
678104-CF2	hose connector with Dip-In	4	93 mm	1,000 cm ²	● red	60 ml - 200 ml	+	1 / 3
678108-CF1	hose connector without Dip-In	8	135 mm	2,000 cm ²	● red	120 ml - 400 ml	+	1 / 2
678108-CF2	hose connector with Dip-In	8	135 mm	2,000 cm ²	● red	120 ml - 400 ml	+	1 / 2
678112-CF1	hose connector without Dip-In	12	177 mm	3,000 cm ²	● red	180 ml - 600 ml	+	1 / 2
678112-CF2	hose connector with Dip-In	12	177 mm	3,000 cm ²	● red	180 ml - 600 ml	+	1 / 2
678116-CF1	hose connector without Dip-In	16	220 mm	4,000 cm ²	● red	240 ml - 800 ml	+	1 / 2
678116-CF2	hose connector with Dip-In	16	220 mm	4,000 cm ²	● red	240 ml - 800 ml	+	1 / 2
678124-CF1	hose connector without Dip-In	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1 / 2
678124-CF2	hose connector with Dip-In	24	304 mm	6,000 cm ²	● red	360 ml - 1,200 ml	+	1 / 2
678140-CF1	hose connector without Dip-In	40	474 mm	10,000 cm ²	● red	600 ml - 2,000 ml	+	1 / 1
678140-CF2	hose connector with Dip-In	40	474 mm	10,000 cm ²	● red	600 ml - 2,000 ml	+	1 / 1

/ Assembled product is sterilized (SAL 10⁻⁶) to reduce risk of contamination



CELLstage

Filling Accessory

- / Available for CELLdisc 4-24 layers and CELLdisc 40 layers
- / Creates the optimum angle and position for CELLdisc filling
- / Stainless steel allows multiple sterilization methods
- / Suitable for left- and right handed users

Item no.	Description	Qty. inner / outer
878072	for CD4 - CD24	1 / 1
878073	for CD40	1 / 1



CELLlevator

Easy and secured CELLdisc stacking

- / Maximum loading capacity 8 kg
- / Space saving storage
- / Autoclavable (120 °C, 2 bar)

Item no.	Qty. inner / outer
878071	1 / 9

/ Autoclaving the CELLlevator more than 3 times is not recommended. After each autoclave cycle the integrity of the segments has to be analysed by the user.



CELLring

Levelling ring

/ Ensures exact planar positioning of CELLdisc


Item no.	Qty. inner / outer
878075	1 / 3



In 3D cell culture, cells form in a spatial orientation - similar to the body. This enables, for example, the replacement of animal experiments and - usually as high-throughput screening - research on drugs, stem cells and tumor development.

3D CELL CULTURE

/ CELLSTAR® Cell Culture Vessels Cell-Repellent Surface	88
Cell Culture Dishes - Cell-Repellent Surface	89
Cell Culture Flasks - Cell-Repellent Surface	89
Multiwell Plates / Microplates - Cell-Repellent Surface.....	90
SCREENSTAR Microplates - Cell-Repellent Surface.....	91
/ Magnetic 3D Cell Culture	92
Spheroid Bioprinting 96 Well.....	93
Spheroid Bioprinting 384 Well.....	93
Magnetic Levitation 6 / 24 Well	94
Screening 96 / 384 Well	95
MagPen Single / 24 Well / 96 Well	95
Consumables / Accessories Magnetic 3D Cell Culture.....	96
/ ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates	97
ThinCert® Cell Culture Inserts 6 Well.....	98
ThinCert® Cell Culture Inserts 12 Well	99
ThinCert® Cell Culture Inserts 24 Well.....	100
ThinCert® Plate 6 / 12 Well.....	100

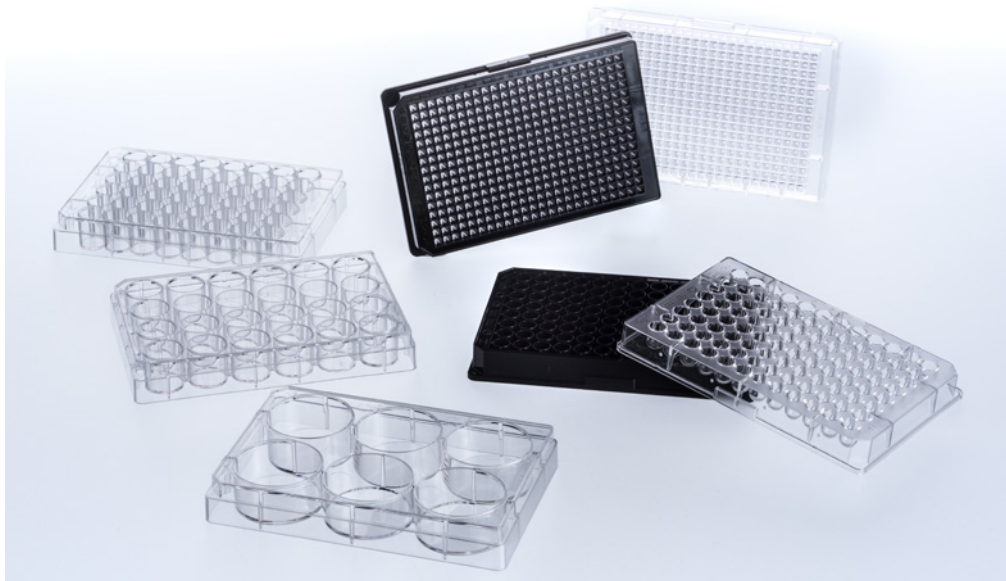

/ Forum No. 17:
CELLSTAR® Cell
Culture Vessels with
Cell-Repellent Surface
(F073777)

CELLSTAR® CELL CULTURE VESSELS CELL-REPELLENT SURFACE


Cell culture is an essential tool in drug discovery, tissue engineering, toxicology testing, stem cell research, as well as in basic research.

Beside conventional two-dimensional (2D) monolayer cell culture, 3D cell culture models are becoming a routine tool which enables the expression of extracellular matrix (ECM) components as well as the formation of cell-cell and cell-matrix interactions. These characteristics are important for replicating *in-*

vivo cell differentiation, proliferation, and function *in vitro*.

Greiner Bio-One developed CELLSTAR® cell culture vessels with cell-repellent surface specifically for culturing in 3D. The cell-repellent surface effectively prevents cell adherence and therefore can promote the spontaneous formation of three-dimensional spheroids. Cell culture vessels with cell-repellent surface are also ideal platforms for long-term cultivation in hydrogels.

- / Spheroid and organoid culture
- / Aggregation of stem cells
- / Suspension culture of semi-adherent/adherent cell lines
- / 3D culture in hydrogels

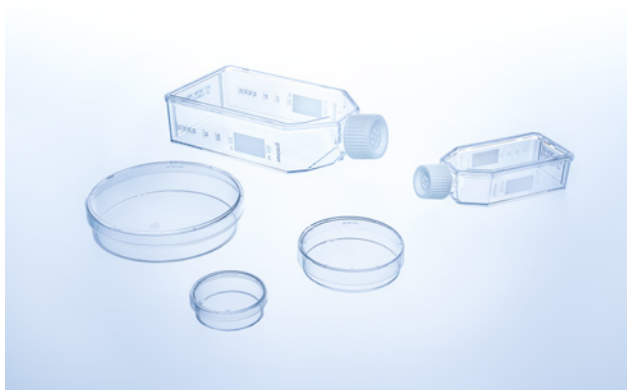


Cell Culture Dishes - Cell-Repellent Surface

Cell culture vessels with cell-repellent surface reliably prevent cell attachment in suspension cultures of semi adherent / adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.



Raw material: PS, Surface treatment: cell-repellent, Vent neck: yes, Sterile: +

Item no.	Height	Ø nominal size	Working volume	Total volume	Sterile	Qty. inner / outer
627979	10 mm	35 mm	≤3 ml	10 ml	+	10 / 40
628979	15 mm	60 mm	6 ml - 7 ml	28 ml	+	10 / 20
664970	20 mm	100 mm	16 ml - 17 ml	100 ml	+	1 / 5

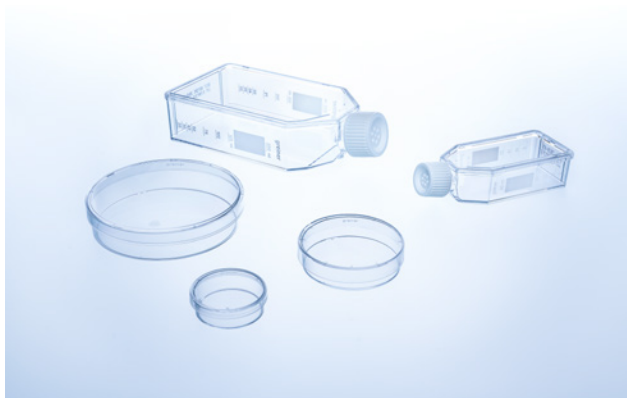


Cell Culture Flasks - Cell-Repellent Surface

A cell-repellent surface reliably prevent cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.



Flask design: flat, Raw material: PS, Surface treatment: cell-repellent, Cap design: filter screw cap, Sterile: +

Item no.	Cap colour	Total volume	Sterile	Qty. inner / outer
690985	○ white	50 ml	+	10 / 20
658985	○ white	250 ml	+	5 / 15
660985	○ white	550 ml	+	5 / 5

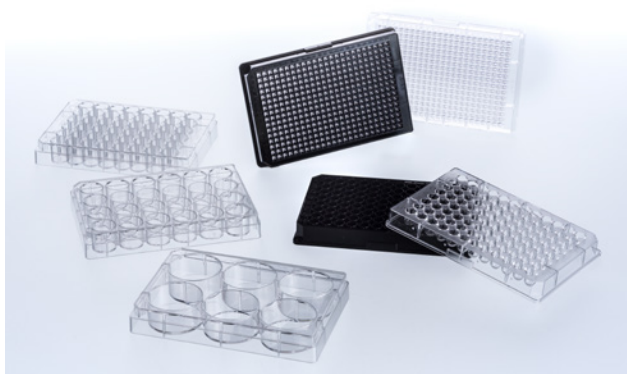

Multiwell Plates / Microplates - Cell-Repellent Surface

- / 6 / 12 / 24 / 48 well multiwell plates available
- / 96 / 384 / 1536 well plates with various well geometries and optional μ Clear® film bottom


STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Raw material: PS, Surface treatment: cell-repellent, Sterile: +

Item no.	Well format	Well profile	Bottom	Product colour	Total volume (well)	Working volume (well)	Lid	Sterile	Qty. inner / outer
657970	6	F-bottom	solid	○ clear	16.1 ml	2 ml - 5 ml	yes, condensation rings	+	1 / 5
665970	12	F-bottom	solid	○ clear	6.5 ml	2 ml - 4 ml	yes, condensation rings	+	1 / 5
662970	24	F-bottom	solid	○ clear	3.3 ml	0.5 ml - 1.5 ml	yes, condensation rings	+	1 / 5
677970	48	F-bottom	solid	○ clear	1.7 ml	0.5 ml - 1 ml	yes, condensation rings	+	1 / 5
650970	96	U-bottom	solid	○ clear	323 μ l	40 μ l - 280 μ l	yes, condensation rings	+	1 / 6
650979	96	U-bottom	solid	○ clear	323 μ l	40 μ l - 280 μ l	yes, condensation rings	+	8 / 32
651970	96	V-bottom	solid	○ clear	234 μ l	40 μ l - 200 μ l	yes, condensation rings	+	1 / 6
655970	96	F-bottom / Chimney Well	solid	○ clear	392 μ l	25 μ l - 340 μ l	yes, condensation rings	+	1 / 6
655976	96	F-bottom / Chimney Well	μ Clear®	● black	392 μ l	25 μ l - 340 μ l	yes, condensation rings	+	8 / 32
655976-SIN	96	F-bottom / Chimney Well	μ Clear®	● black	392 μ l	25 μ l - 340 μ l	yes, condensation rings	+	1 / 32
781970	384	F-bottom	solid	○ clear	131 μ l	15 μ l - 110 μ l	yes	+	1 / 60
781974	384	F-bottom	μ Clear®	○ white	131 μ l	15 μ l - 110 μ l	yes	+	8 / 32
781976	384	F-bottom	μ Clear®	● black	131 μ l	15 μ l - 110 μ l	yes	+	8 / 32
781976-SIN	384	F-bottom	μ Clear®	● black	131 μ l	15 μ l - 110 μ l	yes	+	1 / 32

Item no.	Well format	Well profile	Bottom	Product colour	Total volume (well)	Working volume (well)	Lid	Sterile	Qty. inner / outer
787979	384	U-bottom	solid	○ clear	122 µl	10 µl - 90 µl	yes	+	8 / 32
782974*	1536	F-bottom	µClear®	○ white	12.6 µl	3 µl - 10 µl	yes	+	10 / 40

/ *Plate geometry: HiBase

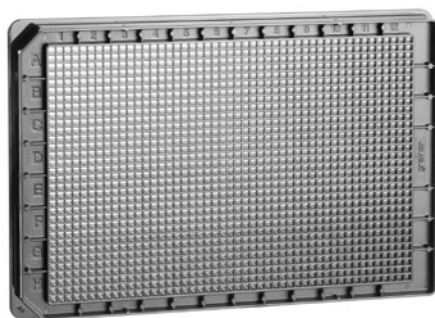

SCREENSTAR Microplates - Cell-Repellent Surface

- / For sophisticated microscopic applications and high content screening
- / Entirely made of cycloolefin
- / Universal compatibility with microscope objectives

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Well format: 1536, Well profile: F-bottom, Bottom: Cycloolefin film, Raw material: COP, Surface treatment: cell-repellent, Plate geometry: LoBase, Lid: yes, Sterile: +

Item no.	Product colour	Sterile	Qty. inner / outer
789979	● black	+	17 / 68



/ Brochure "3D Cell Culture" (F071076)


MAGNETIC 3D CELL CULTURE

The core technology of Greiner Bio-One's magnetic 3D cell culture is the magnetisation of cells with NanoShuttle-PL. The cells can be aggregated with magnetic forces, either by levitation or Bioprinting, to form structurally and biologically representative 3D models in vitro. NanoShuttle-PL consists of gold, iron oxide, and Poly-L-Lysine. These nanoparticles ($\emptyset < 50$ nm) magnetise cells by electrostatically attaching to cell membranes during an overnight static incubation. Magnetised cells will ap-

pear peppered with dark nanoparticles after incubation.

NanoShuttle-PL is biocompatible, having no effect on metabolism, proliferation and inflammatory stress. Additionally, it does not interfere with experimental techniques, such as fluorescence or Western blotting. With magnetised spheroids, solution addition and removal are made easy by using magnetic force to hold them in a stationary position during aspiration, thereby limiting spheroid loss.

- / 3D in a 2D workflow
- / Easy handling
- / Fast 3D tissue assembly
- / No sample loss
- / Scalable
- / Ready for automation

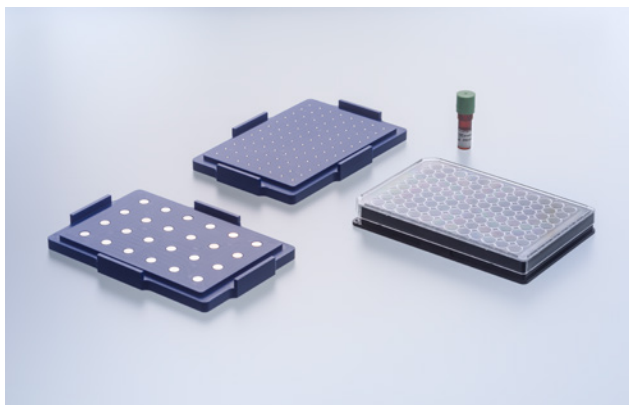


Spheroid Bioprinting

96 Well

Magnetized cells are brought together to form spheroids at the well bottom using weak magnetic forces.

Item no.	Description	Content kit	Qty. inner / outer
655840	96 Well Bioprinting Kit, clear	NanoShuttle-PL (3 vials), Spheroid Drive, Holding Drive, 96 well cell culture microplates (clear) with cell-repellent surface (2 x 655970)	1 / 1
655841	96 Well Bioprinting Kit, black, μ Clear [®]	NanoShuttle-PL (3 vials), Spheroid Drive, Holding Drive, 96 well cell culture microplates (black, μ clear [®]) with cell-repellent surface (2 x 655976-SIN)	1 / 1
655850	96 Well Ring Drive	96 Well Ring Drive for the formation of 3D ring structures	1 / 1
655830	96 Well Spheroid and Holding Drive	Spheroid Drive, Holding Drive	- / 1

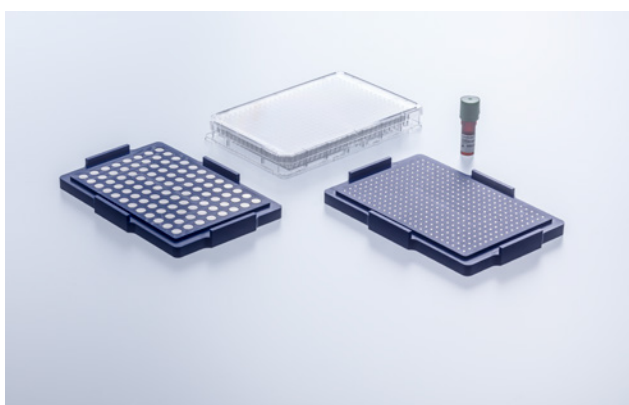


Spheroid Bioprinting

384 Well

Magnetized cells are brought together to form spheroids at the well bottom using weak magnetic forces.

Item no.	Description	Content kit	Qty. inner / outer
781840	384 Well Bioprinting Kit, clear	NanoShuttle-PL (2 vials), Spheroid Drive, Holding Drive, 384 well cell culture microplates (clear) with cell-repellent surface (2 x 781970)	1 / 1
781841	384 Well Bioprinting Kit, black, μ Clear [®]	NanoShuttle-PL (2 vials), Spheroid Drive, Holding Drive, 384 well cell culture microplates (black, μ Clear [®]) with cell-repellent surface (2x 781976-SIN)	1 / 1
781850	384 Well Ring Drive	384 Well Ring Drive for the formation of 3D ring structures	1 / 1

Item no.	Description	Content kit	Qty. inner / outer
781830	384 well spheroid and holding drive	Spheroid Drive, Holding Drive	- / 1

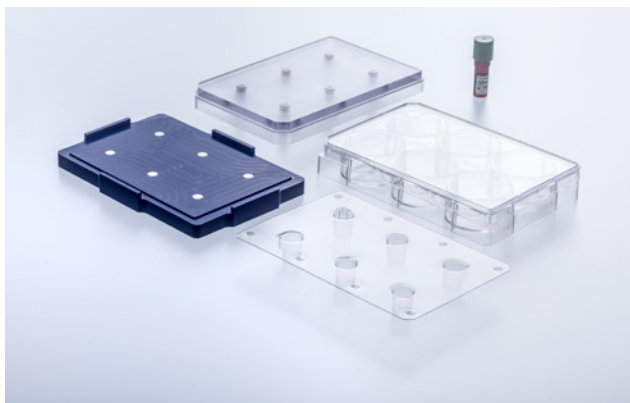


Magnetic Levitation

6 / 24 Well

Magnetised cells are levitated off the plate bottom by a magnet and rapidly form aggregates.


Item no.	Description	Content kit	Sterile	Qty. inner / outer
657840	6 Well Bio-Assembler Kit	Levitation Drive, Holding Drive, NanoShuttle-PL (2 Vials), 6 well cell culture multiwell plates (2 x 657970) and 6 Well Intermediate lid (2 x 657825) with cell-repellent surface		1 / 1
657825	6 Well Intermediate lid	Intermediate lid with cell-repellent surface	+	2 / 10
657830	6 Well Levitation and Holding Drive	Levitation Drive (1), Holding Drive (1)		- / 1
662840	24 Well Bio-Assembler Kit	Levitation Drive, Holding Drive, NanoShuttle-PL (2 Vials), 24 well cell culture multiwell plates (2 x 662970) and 24 Well Intermediate lid (2 x 662825) with cell-repellent surface		1 / 1
662825	24 Well Intermediate lid	Intermediate lid with cell-repellent surface	+	2 / 10
662830	24 well Levitation and Holding Drive	Levitation Drive (1), Holding Drive (1)		- / 1

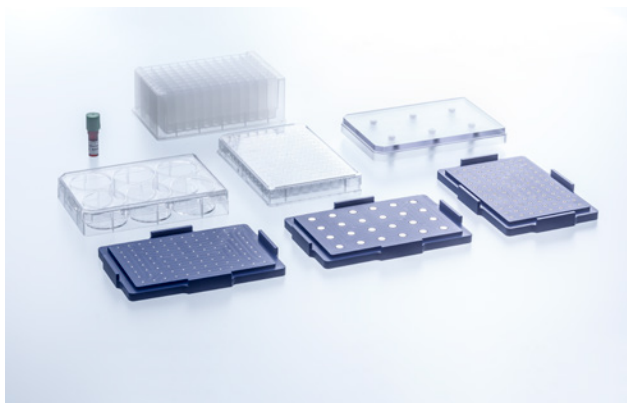

Screening
96 / 384 Well

Ideal kits for combining applications and examination methods.

Item no.	Description	Content kit	Qty. inner / outer
655846	96 Well BiO Assay Kit	NanoShuttle-PL (3 vials), 6 Well Levitation Drive, 96 Well Spheroid, Holding and Ring Drive, 96 Well Deep Well Plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 96 Well cell culture microplates (clear) with cell-repellent surface (2 x 655970), 6 Well Intermediate Lid with cell-repellent surface (2 x 657825)	1 / 1
781846	384 Well BiO Assay Kit	NanoShuttle-PL (2 vials), 6 Well Levitation Drive, 384 Well Spheroid and Holding Drive, 96 Well Deep Well plate, 6 Well cell culture multiwell plates with cell-repellent surface (2 x 657970), 384 Well cell culture microplates (clear) with cell-repellent surface (2 x 781970), 6 Well Intermediate Lid with cell-repellent surface (2 x 657825)	1 / 1

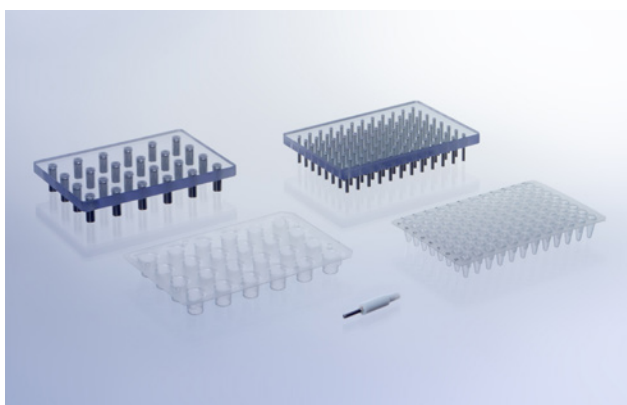

MagPen
Single / 24 Well / 96 Well

MagPen is a smart assistant for easy and fast transfer and collection of magnetized cell cultures by a simple "pick up-and-drop"-step.

Item no.	Description	Content kit	Sterile	Qty. inner / outer
657850	MagPen 3-pack	Teflon caps (3), magnets (3)		- / 3
657824	24 Well Multi-MagPen Kit	24 Well Multi-MagPen Drive and 24 Well Multi-MagPen Sleeve (2 x 651524) with cell-repellent surface		- / 1
651524	24 Well Multi-MagPen Sleeve	Multi-MagPen Sleeve with cell-repellent surface	+	1 / 10
657896	96 Well Multi-MagPen Kit	96 Well Multi-MagPen Drive und 96 Well Multi-MagPen Sleeve (2 x 61596) with cell-repellent surface		- / 1

Item no.	Description	Content kit	Sterile	Qty. inner / outer
651596	96 Well Multi-MagPen Sleeve	Multi-MagPen Sleeve with cell-repellent surface	+	1 / 10



Consumables / Accessories

Magnetic 3D Cell Culture

- / NanoShuttle-PL consists of gold, iron oxide and Poly-L-Lysine



Item no.	Description	Content kit	Aseptic	Qty. inner / outer
657841	NanoShuttle-PL	600 µl vials of NanoShuttle-PL (1)	+	- / 1
657843	NanoShuttle-PL 3-pack	600 µl vials of NanoShuttle-PL (3)	+	- / 3
657846	NanoShuttle-PL 6-pack	600 µl vials of NanoShuttle-PL (6)	+	- / 6
657852	NanoShuttle-PL 12-pack	600 µl vials of NanoShuttle-PL (12)	+	- / 12



- / Forum No. 8: ThinCert® cell culture products – Overview (F073017)
- / Application Note „Immuno cytochemistry“(F073100)
- / For further information and application notes please refer to our website: www.gbo.com



THINCERT® CELL CULTURE INSERTS FOR 6, 12 AND 24 WELL MULTIWELL PLATES

For advanced cell and tissue culture applications, Greiner Bio-One offers a broad range of membrane supports - ThinCert®.

Combining 6 different membrane specifications (pore size and density) in geometries to fit 6, 12 and 24 well plates, the ThinCert® cell culture inserts are suitable for a wide range of applications including transport, secretion and diffusion studies, migrational experiments, cytotoxicity testing, co-cultures, trans epithelial electric resistance (TEER) measurements, as well as pri-

mary cell culture.

ThinCert® cell culture inserts are compatible with standard CELLSTAR® multiwell plates from Greiner Bio-One, and are pre-packed together with the requisite number of plates. The automated production process includes double optical control of each insert produced, ensuring that any biological contamination is avoided. The sterility of the single blisterpacked inserts and multiwell plates is ensured by irradiation.

- / Stable clear polystyrene housing
- / Sealed PET capillary pore membrane
- / Pre-configured multiwell plates with ThinCert® cell culture inserts available on request

3D Cell Culture

ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates



ThinCert® Cell Culture Inserts

6 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

STERILE

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 24.85 mm, Ø external: 27.85 mm, Cultural surface: 452.4 mm², Surface treatment: TC, Working volume (ThinCert®): 1 ml - 3.6 ml, Working volume (well): 2 ml - 4.15 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
657640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 24
657641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 24
657610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 24
657630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 24
657631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 24
657638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 24



ThinCert® Cell Culture Inserts

12 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

STERILE

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

Feature: 4 multiwell plates / box, Height: 16.25 mm, Ø internal: 13.85 mm, Ø external: 15.85 mm, Cultural surface: 113.1 mm², Surface treatment: TC, Working volume (ThinCert®): 0.3 ml - 1 ml, Working volume (well): 1 ml - 2 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
665640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 48
665641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 48
665610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 48
665630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 48
665631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 48
665638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 48

3D Cell Culture

ThinCert® Cell Culture Inserts for 6, 12 and 24 Well Multiwell Plates



ThinCert® Cell Culture Inserts

24 Well

- / Hanging geometry
- / Improved cell adhesion through physical surface treatment
- / Simplified pipetting due to self-lift geometry
- / Enhanced pipetting access and gas exchange

STERILE

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

non-
cytotoxic

non-
pyrogenic

Feature: 2 multiwell plates / box, Height: 16.25 mm, Ø internal: 8.4 mm, Ø external: 10.4 mm, Cultural surface: 33.6 mm², Surface treatment: TC, Working volume (ThinCert®): 0.1 ml - 0.35 ml, Working volume (well): 0.4 ml - 1.2 ml, Sterile: +

Item no.	Pore density	Ø pore	Optical membrane properties	Sterile	Qty. inner / outer
662640	1 x 10 ⁸ / cm ²	0.4 µm	translucent	+	1 / 48
662641	2 x 10 ⁶ / cm ²	0.4 µm	clear	+	1 / 48
662610	2 x 10 ⁶ / cm ²	1 µm	clear	+	1 / 48
662630	0,6 x 10 ⁶ / cm ²	3 µm	clear	+	1 / 48
662631	2 x 10 ⁶ / cm ²	3 µm	translucent	+	1 / 48
662638	0,15 x 10 ⁶ / cm ²	8 µm	translucent	+	1 / 48



ThinCert® Plate

6 / 12 Well

- / Optimised for use with ThinCert® cell culture inserts
- / Deep wells for an increased volume of medium in air-lift culture
- / Notches for fixed insert position
- / Available in 6 and 12 well format

STERILE

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

non-
cytotoxic

non-
pyrogenic

Height: 39.5 mm, Length: 129.5 mm, Width: 86.6 mm, Lid: yes, condensation rings, Sterile: +

Item no.	Well format	Working volume (well)	Sterile	Qty. inner / outer
657110	6	≤20 ml	+	1 / 50

Item no.	Well format	Working volume (well)	Sterile	Qty. inner / outer
665110	12	≤4 ml	+	1 / 60



Microscopy is a basic and important method, which is often used in research as well as in medical diagnostics. For this purpose, Greiner Bio-One offers a variety of tailor-made solutions, which guarantee optimal basic conditions for microscopic experiments.

PRODUCTS FOR MICROSCOPY

- / CELLview Dish Cell Culture Dish with
Glass Bottom 104
 CELLview Dish 105

- / CELLview Slide Cell Culture Slide with
Glass Bottom 106
 CELLview Slide 107

- / CELLview Plate Cell Culture Plate with
Glass Bottom 108
 CELLview Plate 109

- / SCREENSTAR Microplates 110
 SCREENSTAR Microplates 96 / 384 / 1536 Well. 111

- / SensoPlate Glass Bottom Microplates .. 112
 SensoPlate Glass Bottom Plates 24 / 96 / 384 /
 1536 Well..... 113



/ Instructions for use
for CELLview Cell
Imaging Consumables
(F074004)

CELLVIEW DISH

CELL CULTURE DISH WITH GLASS BOTTOM


CELLview Dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of in-vitro cultivated cultures.

It is made from high-grade polystyrene combined with an integrated glass bottom. The innovative design of the dish provides a single-plane, flat bottom with a consistent working distance and maximal planarity. Moreover, the dish bottom configuration facilitates optimal thermal conductivity and avoids thermal variations in heated platforms

used for live cell imaging.

The subdivided version of CELLview Dish enables simultaneous multiplex analyses of different cell lines, various stimulations or diverse transfections. Quartering the cell culture dish provides four individual compartments with a growth area of approximately 1.9 cm², allowing minimisation of cells and reagents required per individual assay. In addition to the untreated glass surface, Greiner Bio-One provides a tissue culture surface as well as the Advanced TC surface modification to enhance the attachment of adherent cells, thus eliminating the need for protein coating.

- / TC and Advanced TC surface modification available
- / Maximal spectral transmission
- / No autofluorescence, no depolarisation of light

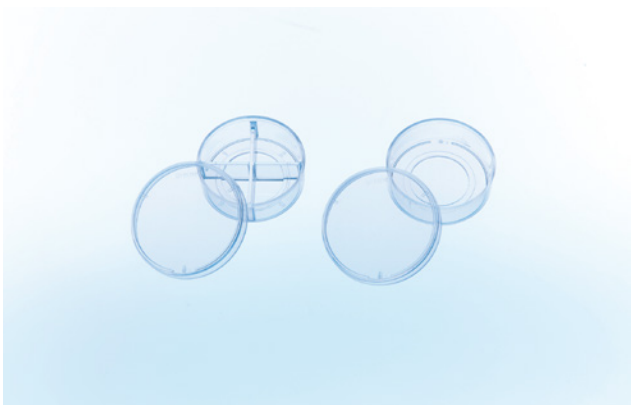

CELLview Dish

The CELLview cell culture dish combines the convenience of a standard size 35 mm disposable plastic cell culture dish with the optical quality of glass, providing superior high-resolution microscopic images of *in-vitro* cultivated cultures.

/ Embedded glass bottom for maximal planarity

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Ø: 35 mm, Bottom: glass, Vent nock: yes, Sterile: +

Item no.	Compartment	Growth area	Growth area / unit	Surface treatment	Working volume	Total volume	Total volume (well)	Working volume (well)	Sterile	Qty. inner / outer
627860	1	8.7 cm ²		TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627965	1	8.7 cm ²		Advanced TC	2.5 ml - 5 ml	10 ml			+	10 / 40
627870	4		1.9 cm ²	TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40
627975	4		1.9 cm ²	Advanced TC			1.5 ml	0.1 ml - 0.5 ml	+	10 / 40

/ Samples are available on request.



/ Instructions for use
for CELLview Cell
Imaging Consumables
(F074004)

CELLVIEW SLIDE CELL CULTURE SLIDE WITH GLASS BOTTOM

CELLview Slide consists of a transparent slide with a black upper housing that effectively subdivides the slide into 10 compartments, which have been designed to mimic the size and layout of a standard 96 well microplate. Because of this standard layout, the slide is compatible with multichannel pipettes making it simple and efficient to use. Furthermore, the round well design helps to reduce meniscus effects for optimum results in cell culture and microscopic analysis. The slide has a 0.17 mm thin cover glass embed-

ded in its bottom for improved optical clarity and imaging. Embedding the cover glass guarantees an even focal plane which is a prerequisite for all high-speed and high-resolution microscopy applications. Furthermore, the black upper housing reduces cross talk between adjacent wells during fluorescence microscopy and the optical glass, which exhibits virtually no autofluorescence, allows for maximum spectral transmission without depolarisation of transmitted light.

- / 10 wells with alphanumeric coding
- / Black detachable compartmentalization
- / Well distance is equal to a 96 well microplate
- / Positioning notch for automated microscopy
- / Highly transparent achromatic borosilicate glass bottom



CELLview Slide

- / 10 wells with alphanumeric coding
- / Positioning notch for automated microscopy
- / Black detachable compartmentalization
- / Reduced meniscus effect due to round well design
- / Glass thickness: 0.17 mm

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Well format: 10, Height: 12.6 mm, Length: 75 mm, Width: 25 mm, Growth area / unit: 34 mm², Bottom: glass, Total volume (well): 440 µl, Sterile: +

Item no.	Surface treatment	Sterile	Qty. inner / outer
543078	TC	+	1 / 20
543079	TC	+	5 / 45
543978	Advanced TC	+	1 / 20
543979	Advanced TC	+	5 / 45

/ Samples are available on request.



- / For plates with an optical cycloolefin film bottom, please refer to our SCREENSTAR plates in this chapter.

CELLVIEW PLATE

CELL CULTURE PLATE WITH GLASS BOTTOM

CELLview glass bottom microplates are designed for demanding and high-resolution microscopic applications.

They consist of a cycloolefin-based black frame with a 0.17 mm thin borosilicate glass bottom providing superior images of in-vitro cultures. The optimised microplate geometry and

the recessed bottom facilitate imaging of all peripheral wells even with immersion objectives. The round conical well design reduces the meniscus effect in order to assure equal cellular distribution and constant imaging results. An appropriate surface treatment improves cellular attachment and growth.

- / Outstanding image quality and resolution
- / Cycloolefin-based frame with 0.17 mm glass bottom
- / Excellent optical transparency
- / Recessed well bottom facilitating the use of lenses with low working distance and high aperture
- / Compatible with advanced confocal microscopic systems



CELLview Plate

- / For outstanding image quality and resolution
- / Recessed well bottom facilitating the use of objectives with low working distance
- / Ditch at the perimeter can be filled with liquid to minimise edge effects and evaporation
- / Compatible with advanced automated microscopic systems

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Well format: 96, Growth area / unit: 33 mm², Well profile: F-bottom / Chimney Well, Bottom: glass, Raw material: COP, Surface treatment: TC, Working volume (well): 25 µl - 440 µl, Lid: yes, Sterile: +

Item no.	Product colour	Sterile	Qty. inner / outer
655891	● black	+	1 / 16



/ Forum No. 15:
SCREENSTAR: A 1536
Well Microplate for
High-Content and High-
Throughput Screening
(F073120)


SCREENSTAR MICROPLATES

SCREENSTAR are specialised microplates for sophisticated microscopic applications, in high content screening (HCS) or high-resolution microscopy with water and oil immersion objectives.

They combine outstanding glass-like optical properties with an excellent surface for adherent cell culture. Moreover, the plates display ex-

cellent optical properties with reduced autofluorescence in the lower UV range, low birefringence and a refractive index of 1.53 comparable to glass. SCREENSTAR microplates enable complete periphery access for high magnification objectives. They are entirely manufactured out of cycloolefin with a black pigmented cycloolefin frame and a 190 µm ultraclear cycloolefin film bottom.

- / 96 / 384 / 1536 well format
- / For sophisticated microscopic applications and high content screening
- / Highly transparent cycloolefin film bottom
- / Adherent TC surface treatment

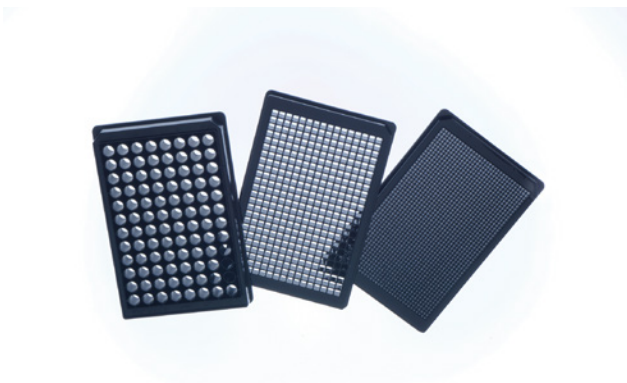

SCREENSTAR Microplates

96 / 384 / 1536 Well

- / For sophisticated microscopic applications and high content screening
- / Highly transparent cycloolefin film bottom
- / Adherent TC surface treatment

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Bottom: Cycloolefin film, Raw material: COP, Surface treatment: TC, Sterile: +

Item no.	Well format	Growth area / unit	Well profile	Product colour	Working volume (well)	Lid	Sterile	Qty. inner / outer
655866	96	33 mm ²	F-bottom / Chimney Well	● black	25 µl - 440 µl	yes	+	1 / 16
781866	384	7.0 mm ²	F-bottom	● black	10 µl - 110 µl	yes	+	8 / 32
789866	1536	2.1 mm ²	F-bottom	● black	3 µl - 15 µl	no	+	17 / 68

/ Forum No. 15: SCREENSTAR: A 1536 Well Microplate for High-Content and High-Throughput Screening (F073120)



/ For further information, please refer to our website: www.gbo.com



SENSOPLATE GLASS BOTTOM MICROPLATES

The research of biomolecular processes on the level of single molecules and in volume ranges equivalent to the size of a single bacterium is of immense importance, both in basic research and in industrial high-throughput screening. The combination of modern confocal optics, new fluorescent dyes, sensitive photomultipliers and improved data processing has revolutionised the technique of fluorescence correlation spectroscopy (FCS).

Over the past few years this has led to its widespread application,

and alongside the technological advances in hardware development, Greiner Bio-One worked hand-in-hand with customers and instrument suppliers to develop the glass bottom microplates. These better satisfy the requirements of fluorescence correlation spectroscopy with regards to optical clarity and deformation when compared to standard polystyrene plates. The SensoPlate family was developed in a complete product line consisting of 24, 96, 384 and 1536 well glass bottom formats.

- / For fluorescence correlation spectroscopy and microscopic applications
- / 24 / 96 / 384 / 1536 well format
- / Black frame with highly transparent glass bottom
- / Glass bottom thickness of 175 μm is equivalent to the light path of standard coverslips
- / With lid and single-packed

SensoPlate Glass Bottom Plates


24 / 96 / 384 / 1536 Well



- / For fluorescence correlation spectroscopy and microscopic applications
- / 24 / 96 / 384 / 1536 well format
- / Black frame with highly transparent glass bottom
- / Glass bottom thickness of 175 µm is equivalent to the light path of standard coverslips
- / With lid and single-packed

Well profile: F-bottom, Bottom: glass, Raw material: PS, Surface treatment: untreated, Lid: yes


Item no.	Well format	Product colour	Plate geometry	Working volume (well)	Sterile	Qty. inner / outer
662892	24	● black		0.5 ml - 1.5 ml	+	1 / 12
655892	96	● black		25 µl - 340 µl	+	1 / 16
781892	384	● black		10 µl - 130 µl	+	1 / 16
781856	384	● black	LoBase	10 µl - 130 µl		1 / 16
783892	1536	● black	LoBase	3 µl - 10 µl	+	1 / 16

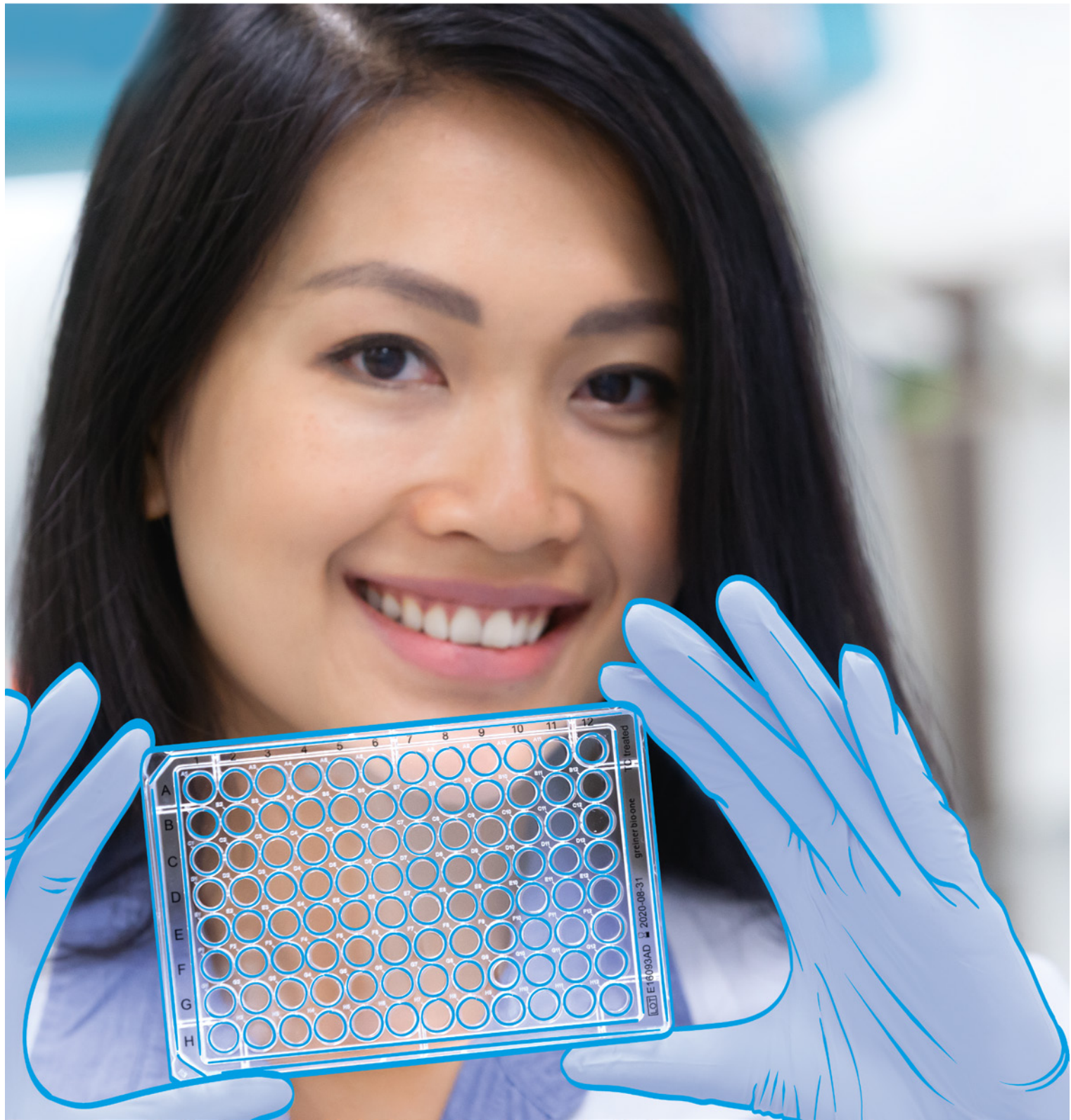

Continued progress in research and related technologies, such as microscopy, imaging, detection and liquid handling, has given rise to a wide variety of platforms used in basic science, biotechnology and pharmaceutical drug development. Today, researchers can select application-specific microplates among a broad range of products that differ in format, design, base material, colour and surface properties.

HTS-MICROPLATES

/ 96 Well Microplates	116	/ UV-Star® Microplates	136
96 Well Microplates Polystyrene	117	UV-Star® Microplates	137
96 Well Microplates Polystyrene - Half Area....	118		
96 Well Microplates Polypropylene.....	119	/ Lids / Sealers / CapMats	138
		Lids	139
/ 384 Well Microplates	120	Sealers	139
384 Well Microplates Polystyrene	121	CapMats	140
384 Well Microplates Polystyrene - Small Volume		
.....	122	/ Protein Crystallisation Plates	141
384 Well Microplates Polypropylene.....	122	96 Well CrystalQuick Plates	142
/ 1536 Well Microplates	124		
1536 Well Microplates	125		
/ Polypropylene Storage Plates	126		
96 Well MASTERBLOCK® Polypropylene.....	127		
384 Deep Well MASTERBLOCK®		
Polypropylene.....	127		
1536 Deep Well Microplates Polypropylene	128		
/ Microplates for Compound Storage	129		
Microplates for Compound Storage 384 and 1536		
Well	130		
/ Non-binding Microplates	131		
Non-binding Microplates 96 Well	132		
Non-binding Microplates 384 and 1536 Well ...	132		
/ Streptavidin-coated Microplates.....	134		
Streptavidin-coated Microplates 96 and 384 Well		
.....	135		

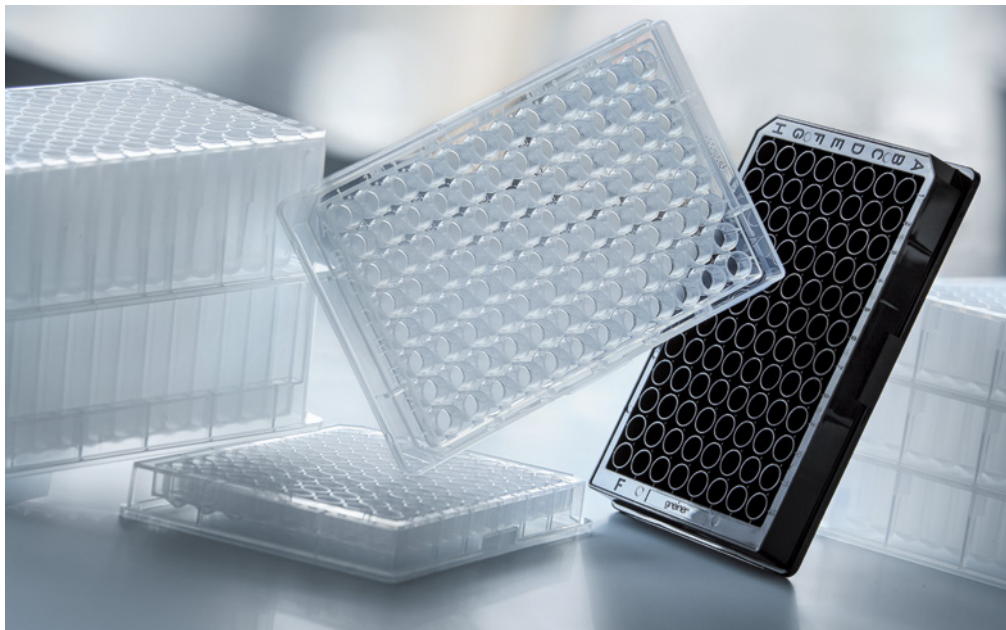

/ Selected microplates are certified with ACT® Ecolabel. Learn more:




96 WELL MICROPLATES

Since its introduction in the 1960's applications for the 96 well microplate have continually increased to the extent that it is impossible to envisage modern research and industry without it today. Greiner Bio-One has been manufacturing microplates and strip microplates for diagnostics

and immunological research for over 40 years. A large number of different 96 well microplates is available in a wide variety of surface treatments. The spectrum ranges from clear bottom microplates and completely black or white microplates to UV-Star® products.

- / Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- / With U-bottom, V-bottom or F-bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality
- / In non-binding quality

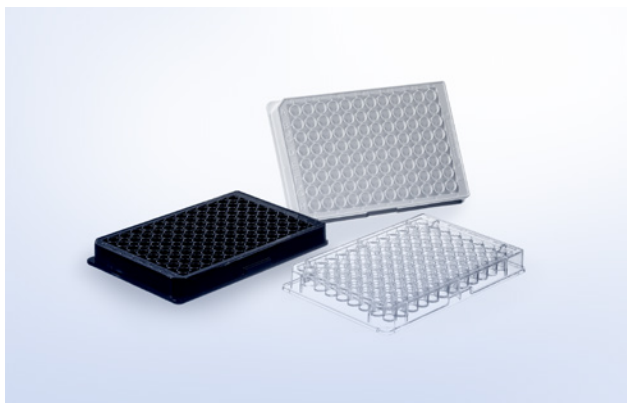

96 Well Microplates
Polystyrene

- / With U-bottom, V-bottom or F-bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality
- / With solid bottom or μ Clear[®] film bottom
- / Selected product versions certified with ACT[®] Ecolabel (*)

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	my green lab. ACT Ecolabel
--------------------------	------------------------------	--------------------------	---------------	----------------------------

Well format: 96, Raw material: PS, Lid: no

Item no.	Well profile	Bottom	Binding character- istic	Binding brand name	Surface treatment	Product colour	Working volume (well)	Sterile	Qty. inner / outer
650101*	U-bottom	solid			untreated	○ clear	40 μ l - 280 μ l		10 / 100
650161	U-bottom	solid			untreated	○ clear	40 μ l - 280 μ l	+	2 / 100
651101*	V-bottom	solid			untreated	○ clear	40 μ l - 200 μ l		10 / 100
651161	V-bottom	solid			untreated	○ clear	40 μ l - 200 μ l	+	2 / 100
655101*	F-bottom	solid			untreated	○ clear	25 μ l - 340 μ l		10 / 100
655161	F-bottom	solid			untreated	○ clear	25 μ l - 340 μ l	+	2 / 100
655075*	F-bottom / Chimney Well	solid			untreated	○ white	25 μ l - 340 μ l		10 / 40
655074*	F-bottom / Chimney Well	solid	high- binding	LUMITRAC 600		○ white	25 μ l - 340 μ l	+	10 / 40
655076*	F-bottom / Chimney Well	solid			untreated	● black	25 μ l - 340 μ l		10 / 40
655077*	F-bottom / Chimney Well	solid	high- binding	FLUOTRAC 600		● black	25 μ l - 340 μ l	+	10 / 40
655095	F-bottom / Chimney Well	μ Clear [®]			untreated	○ white	25 μ l - 340 μ l		10 / 40
655094	F-bottom / Chimney Well	μ Clear [®]	high- binding			○ white	25 μ l - 340 μ l	+	10 / 40
655097	F-bottom / Chimney Well	μ Clear [®]	high- binding			● black	25 μ l - 340 μ l	+	10 / 40
655096	F-bottom / Chimney Well	μ Clear [®]			untreated	● black	25 μ l - 340 μ l		10 / 40

/ Item numbers marked with * are certified with ACT[®] Ecolabel. The ACT[®] registered trademark is owned and controlled by My Green Lab Inc.
More details at <https://act.mygreenlab.org>



96 Well Microplates

Polystyrene - Half Area

- / Sterile / non-sterile
- / Reduction of sample volume by up to 50 %
- / Standardised pathlength (1 cm=170 µl, 0.5 cm=80 µl)
- / Non-treated or in high-binding quality
- / With solid bottom or µClear® film bottom

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic

Well format: 96, Well profile: F-bottom, Raw material: PS, Plate design: half area, Working volume (well): 15 µl - 175 µl, Lid: no

Item no.	Bottom	Binding characterisitc	Surface treatment	Product colour	Sterile	Qty. inner / outer
675161	solid		untreated	○ clear	+	10 / 40
675101	solid		untreated	○ clear		10 / 40
675074	solid	high-binding		○ white	+	10 / 40
675075	solid		untreated	○ white		10 / 40
675077	solid	high-binding		● black	+	10 / 40
675076	solid		untreated	● black		10 / 40
675096	µClear®		untreated	● black		10 / 40

/ Further information on Half Area Microplates: Forum No. 16: 96 Well Half Area Microplates and their Application in Fluorescence, Luminescence and Transmission Measurements (F073121)



96 Well Microplates Polypropylene

- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance
- / Black plates for fluorescence polarisation
- / Selected product versions certified with ACT[®] Ecolabel(*)

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase


 non-
pyrogenic

 my green lab.
ACT Ecolabel

Well format: 96, Bottom: solid, Raw material: PP, Surface treatment: untreated, Lid: no

Item no.	Well profile	Product colour	Working volume (well)	Sterile	Qty. inner / outer
650201*	U-bottom / Chimney Well	○ natural	50 µl - 300 µl		10 / 100
650261*	U-bottom / Chimney Well	○ natural	50 µl - 300 µl	+	10 / 100
650209	U-bottom / Chimney Well	● black	50 µl - 300 µl		10 / 100
651201*	V-bottom/Chimney Well	○ natural	50 µl - 335 µl		10 / 100
651209	V-bottom/Chimney Well	● black	50 µl - 335 µl		10 / 100
655201*	F-bottom / Chimney Well	○ natural	25 µl - 370 µl		10 / 100
655209*	F-bottom / Chimney Well	● black	25 µl - 370 µl		10 / 100

/ Item numbers marked with * are certified with ACT[®] Ecolabel. The ACT[®] registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

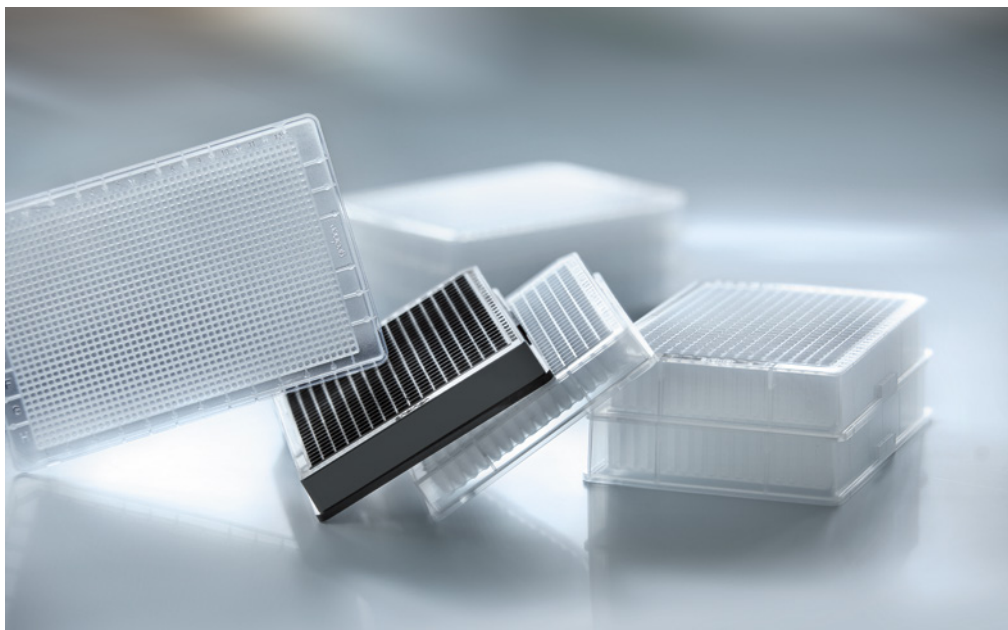

/ Selected microplates are certified with ACT® Ecolabel. Learn more:




384 WELL MICROPLATES

Drug screening has undergone rapid development over the past years.

The number of tests with new targets and the number of active agents to be tested is constantly increasing. Volume reduction, simple testing and cost savings are some of the highest priorities and high-format microplates with a low well volume are one of the most important tools in achieving this. One of the first

higher-format microplates was the 384 well plate, launched by Greiner Bio-One in 1994/1995. Compared with the 96 well microplate, the number of wells is quadrupled in this microplate, combined with a volume reduction from 382 μ l to 131 μ l. The well-to-well spacing is 4.5 mm (96 well plate: 9 mm). The external dimensions of the 384 well microplates are compatible with automated systems.

- / Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- / With F-bottom, V-bottom or as Small Volume option
- / Solid bottom or μ Clear® film bottom
- / Non-treated or in high-binding quality
- / In non-binding quality

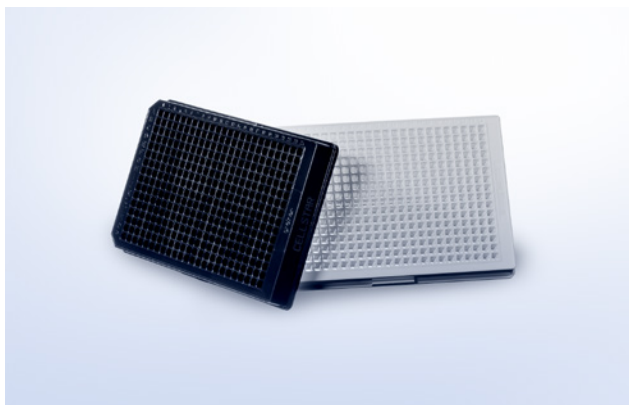


384 Well Microplates
Polystyrene

- / Solid bottom or μ Clear® film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	 non- pyrogenic
--------------------------------	------------------------------------	--------------------------------	---

Well format: 384, Well profile: F-bottom, Raw material: PS, Working volume (well): 15 μ l - 110 μ l

Item no.	Bottom	Binding characterisic	Binding brand name	Surface treatment	Product colour	Lid	Sterile	Qty. inner / outer
781101	solid			untreated	○ clear	no		10 / 100
781061	solid	high-binding			○ clear	no	+	10 / 40
781162	solid			untreated	○ clear	no	+	10 / 100
781185	solid			untreated	○ clear	yes	+	1 / 32
781186	solid			untreated	○ clear	yes	+	8 / 32
781074	solid	high-binding	LUMITRAC 600		○ white	no	+	10 / 40
781075	solid			untreated	○ white	no		10 / 40
781077	solid	high-binding	FLUOTRAC 600		● black	no	+	10 / 40
781076	solid			untreated	● black	no		10 / 40
781095	μ Clear®			untreated	○ white	no		10 / 40
781097	μ Clear®	high-binding			● black	no	+	10 / 40
781096	μ Clear®			untreated	● black	no		10 / 40

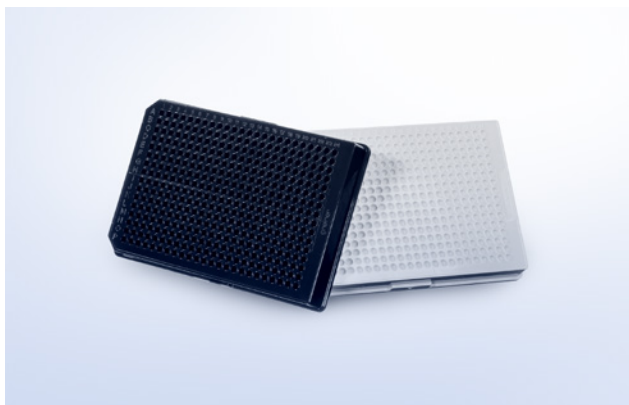

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic


384 Well Microplates

Polystyrene - Small Volume

- / Perfect for top reading even at low working volumes
- / Savings in reagent similar to 1536 well microplates
- / Made of clear / black / white polystyrene for transmission, fluorescence or luminescence measurements

Well format: 384, Well profile: F-bottom, Bottom: solid, Raw material: PS, Surface treatment: untreated, Plate geometry: HiBase, Plate design: Small Volume, Working volume (well): 4 µl - 25 µl, Lid: no

Item no.	Product colour	Qty. inner / outer
784101	○ clear	10 / 40
784075	○ white	10 / 40
784075-25	○ white	25 / 150
784076	● black	10 / 40
784076-25	● black	25 / 150

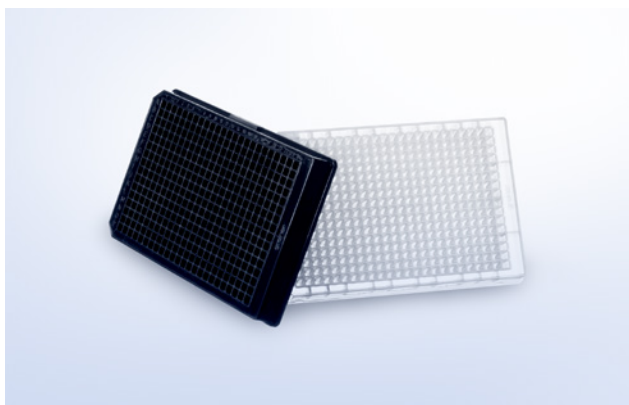

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic

 my green lab.
ACT Ecolabel

384 Well Microplates

Polypropylene

- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance
- / Available natural or black
- / Item No. 781201-906 for Acoustic Liquid Handling
- / Selected product versions certified with ACT® Ecolabel (*)

Well format: 384, Bottom: solid, Raw material: PP, Surface treatment: untreated, Lid: no

Item no.	Well profile	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781201-906	F-bottom	○ natural			15 µl - 145 µl	10 / 100
781201*	F-bottom	○ natural			15 µl - 145 µl	10 / 100
784201*	V-bottom	○ natural	Deep Well	Small Volume	1 µl - 90 µl	10 / 100

Item no.	Well profile	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781209	F-bottom	● black			15 μ l - 145 μ l	10 / 100
781280*	V-bottom	○ natural			13 μ l - 120 μ l	10 / 100

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc.
More details at <https://act.mygreenlab.org>



- / For cell culture treated microplates, please refer to chapter cell culture

1536 WELL MICROPLATES


The highest possible degree of automation, optimal performance and cost savings continue to be the requirements placed on microplates for high-throughput screening.

In 1997, shortly after the launch of the 384 well microplates, Greiner Bio-One was the first manufacturer to introduce another innovative microplate format – the 1536 well microplate. The external dimensions were

the same as those used in the 96 well and 384 well microplates. However, to utilise the available space most efficiently, the number of wells was increased fourfold from 384 to 1536.

Close cooperation with numerous users has now led to the development of a broad product range, and the constant drive towards improvements in quality has, for example, led to a reduction in curvature of the plates to < 100 µm.

- / Available in polystyrene, polypropylene and cycloolefin
- / Clear / black / white
- / HiBase for top-reading applications
- / Solid bottom or µClear® film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

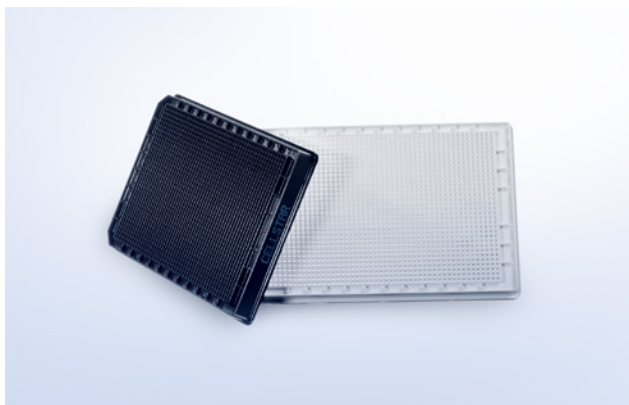

1536 Well Microplates

- / Solid bottom or μ Clear® film bottom
- / Sterile / non-sterile
- / Non-treated or in high-binding quality

FREE OF
detectable
DNase


FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic

Well format: 1536, Well profile: F-bottom, Raw material: PS, Plate geometry: HiBase, Working volume (well): 3 μ l - 10 μ l, Lid: no

Item no.	Bottom	Binding characterisitc	Binding brand name	Surface treatment	Product colour	Sterile	Qty. inner / outer
782101	solid			untreated	○ clear		15 / 60
782061	solid	high-binding	MICROLON 600		○ clear	+	15 / 60
782075	solid			untreated	○ white		15 / 60
782074	solid	high-binding	LUMITRAC 600		○ white	+	15 / 60
782076	solid			untreated	● black		15 / 60
782095	μ Clear®			untreated	○ white		15 / 60
782096	μ Clear®			untreated	● black		15 / 60

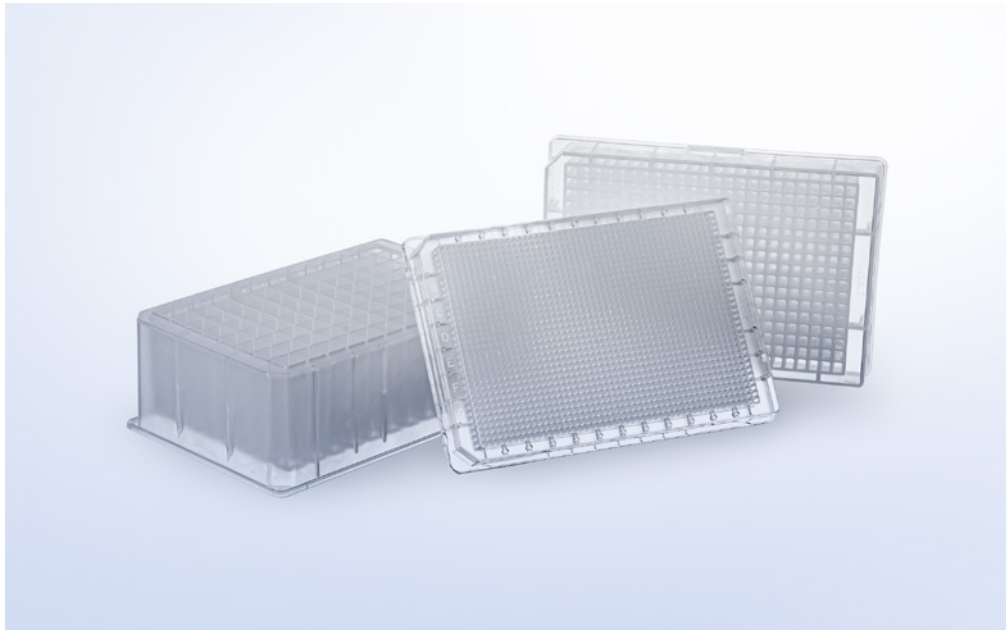

/ Selected microplates are certified with ACT® Ecolabel. Learn more:



POLYPROPYLENE STORAGE PLATES


Greiner Bio-One polypropylene microplates are perfect storage plates for active agents, patient samples or biomolecules.

Their most important properties are biological inertness, resistance to numerous solvents commonly used in the laboratory, such as DMSO and high temperature resistance. The footprint is compatible with

automated systems.

The microplates are also characterised by elevated well walls which make it possible to easily seal them. The portfolio ranges from the 96 well MASTERBLOCK® with volumes of 0.5 ml, 1 ml or 2 ml to a MASTERBLOCK® with 384 wells or 1536 wells.

- / Ideally suited for the storage of active agents, DNA/RNA or stock cultures
- / High chemical resistance and temperature tolerance

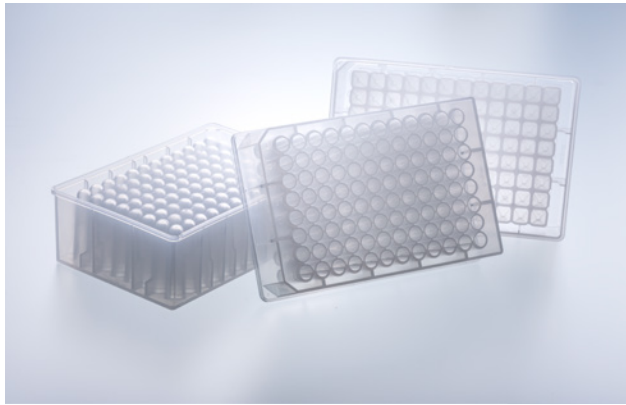

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	my green lab. ACT Ecolabel
--------------------------	------------------------------	--------------------------	---------------	----------------------------


96 Well MASTERBLOCK® Polypropylene

- / Ideal for storing non-human sample material and cultivating bacteria / yeast
- / Uniform external dimensions and tolerances
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers
- / Sealable with CapMats
- / ACT® Ecolabel certification

Well format: 96, Bottom: solid, Raw material: PP

Item no.	Well profile	Product colour	Total volume (well)	Lid	Sterile	Qty. inner / outer
780201	U-bottom	○ natural	1 ml	CapMat 381070, 381061		1 / 50
780215	U-bottom	○ natural	1 ml	CapMat 381070, 381061		5 / 50
780261	U-bottom	○ natural	1 ml	CapMat 381070, 381061	+	1 / 50
786201	V-bottom	○ natural	0.5 ml	CapMat 381070, 381061		8 / 80
786261	V-bottom	○ natural	0.5 ml	CapMat 381070, 381061	+	1 / 80
780270	V-bottom	○ natural	2 ml	CapMat 381080, 381081		1 / 50
780271	V-bottom	○ natural	2 ml	CapMat 381080, 381081	+	1 / 50
780285	V-bottom	○ natural	2 ml	CapMat 381080, 381081		5 / 50

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

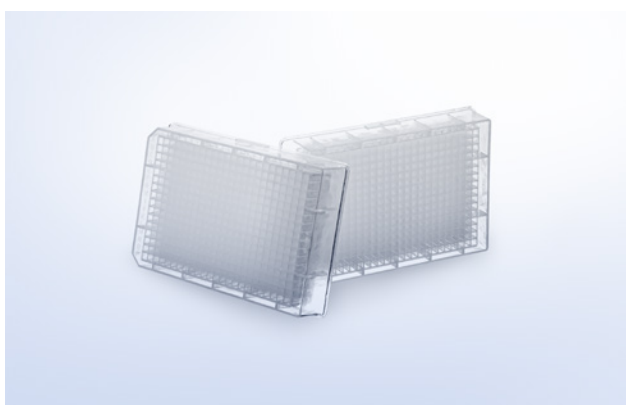

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	my green lab. ACT Ecolabel
--------------------------	------------------------------	--------------------------	---------------	----------------------------


384 Deep Well MASTERBLOCK® Polypropylene

- / Ideal for compound libraries and applications with larger volumes
- / Conical well shape for precise pipetting
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers
- / ACT® Ecolabel certification

Well format: 384, Well profile: V-bottom, Bottom: solid, Raw material: PP, Plate geometry: Deep Well, Lid: no

Item no.	Product colour	Sterile	Qty. inner / outer
781270	○ natural		6 / 60
781271	○ natural	+	6 / 60

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

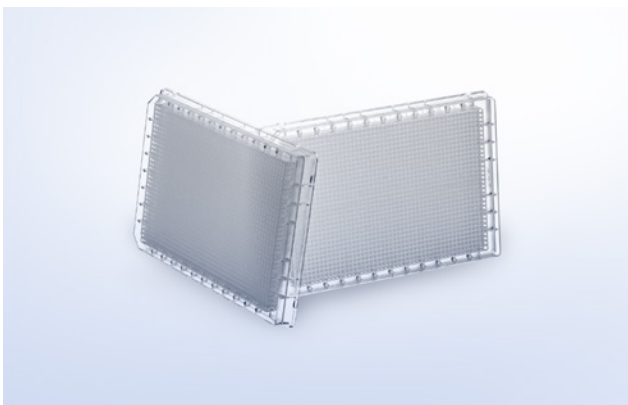

1536 Deep Well Microplates

Polypropylene

- / Uniform external dimensions and tolerances
- / Total volume of 18 μ l
- / Alphanumeric well coding
- / Sealable with adhesive films and heat sealers

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase


 non-
pyrogenic

Well format: 1536, Well profile: V-bottom, Bottom: solid, Raw material: PP, Plate geometry: Deep Well, Working volume: 3 μ l - 15 μ l, Lid: no

Item no.	Product colour	Sterile	Qty. inner / outer
782261	<input type="radio"/> natural	+	15 / 60
782270	<input type="radio"/> natural		15 / 60



- / You can find more information about this product in our web-shop.

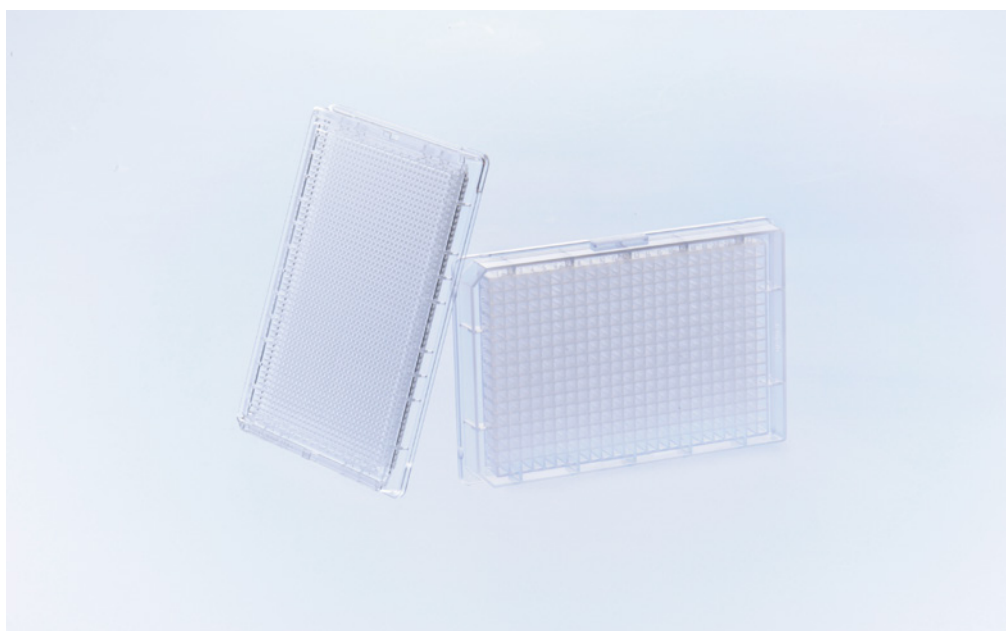


MICROPLATES FOR COMPOUND STORAGE

- / Suited for compound storage
- / Resistant against polar solvents
- / Excellent water and vapour barrier function
- / Very low level of leachables
- / Low biomolecule binding
- / Glass-like optical properties

Polypropylene is still the material of choice for storage plates, but the material class of cycloolefins is becoming more routinely used because of its unsurpassed performance for a wide range of applications.

In compound storage, plates made from cycloolefins offer the

best combination of chemical resistance to polar solvents, like DMSO, and optical clarity. In addition, the dimensional stability and glass-like optical properties make this material ideally suited for plates in fully automated systems.

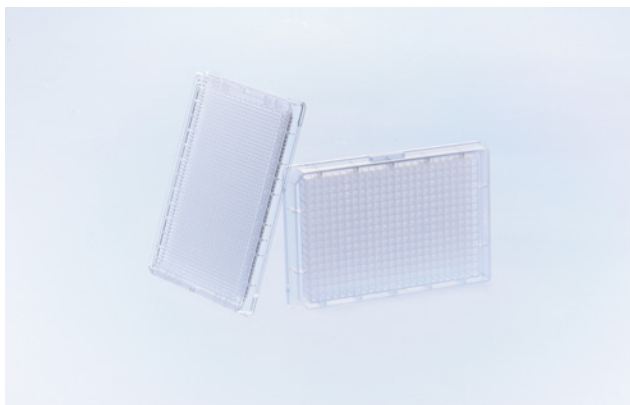

Microplates for Compound Storage

384 and 1536 Well

Microplates for Acoustic Liquid Handling

- / Stringent production specifications for a constant bottom quality
- / Microplates are deionised and packed in antistatic bags

Item no.	Well format	Well profile	Bottom	Raw material	Surface treatment	Product colour	Plate geometry	Plate design	Working volume (well)	Lid	Sterile	Qty. inner / outer
781201-906	384	F-bottom	solid	PP	untreated	○ natural			15 µl - 145 µl	no		10 / 100
788860-906	384	F-bottom	solid	cycloolefin		○ clear	LoBase	Small Volume	4 µl - 25 µl	no		10 / 80
793855	384	F-bottom	solid	cycloolefin		○ clear	HiBase	Small Volume	1 µl - 25 µl	no		15 / 60
782855	1536	F-bottom	solid	cycloolefin		○ clear	HiBase		1 µl - 10 µl	no		15 / 60
782865	1536	F-bottom	solid	cycloolefin	TC	○ clear	HiBase		1 µl - 10 µl	no	+	15 / 60
792870-906	1536	F-bottom	solid	cycloolefin		○ clear			1 µl - 14 µl	no		15 / 60



/ For more detailed information on well profiles and technical details please refer to our data sheets on our website: www.gbo.com




NON-BINDING MICROPLATES

- / 96 / 384 / 1536 well format
- / Clear / black / white
- / Solid bottom or μ Clear® film bottom
- / Ultra low non-specific biomolecular binding properties
- / Long-term surface stability without degradation or leaching

Polystyrene microplates with non-treated surfaces are commonly used for homogeneous biochemical HTS assays and demonstrate low and reproducible biomolecule binding.

However, even low amounts of biomolecular binding (e. g. DNA, RNA, proteins, peptides) can cause an undesirable increase in background, resulting in decreased signal-to-noise ratio.

Greiner Bio-One's non-binding microplate surfaces prevent unwanted non-specific binding, especially advantageous for sensitive biochemical assays. Characterised by low protein, DNA, RNA and peptide binding properties the non-binding surfaces significantly increase assay sensitivity by reducing background and improving signal-to-noise ratio.

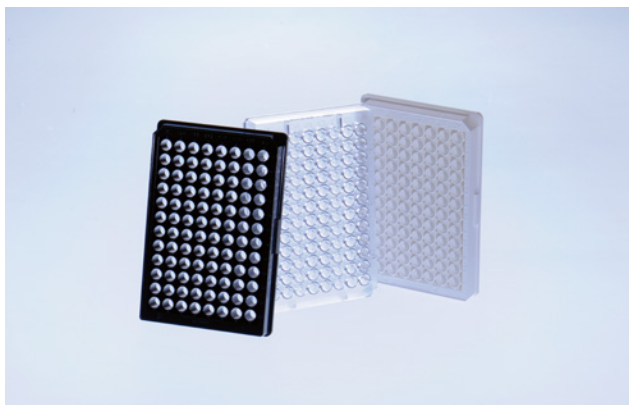

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
--------------------------	------------------------------	--------------------------	---------------


Non-binding Microplates

96 Well

- / Clear / black / white
- / U-bottom, V-bottom or F-bottom/chimney well
- / Solid bottom or μ Clear® film bottom

Well format: 96, Raw material: PS, Binding characterisitc: Non-binding, Lid: no

Item no.	Well profile	Bottom	Product colour	Working volume (well)	Qty. inner / outer
650901	U-bottom	solid	○ clear	40 μ l - 280 μ l	10 / 40
651901	V-bottom	solid	○ clear	40 μ l - 200 μ l	10 / 40
655901	F-bottom / Chimney Well	solid	○ clear	25 μ l - 340 μ l	10 / 40
655904	F-bottom / Chimney Well	solid	○ white	25 μ l - 340 μ l	10 / 40
655900	F-bottom / Chimney Well	solid	● black	25 μ l - 340 μ l	10 / 40
655903	F-bottom / Chimney Well	μ Clear®	○ white	25 μ l - 340 μ l	10 / 40
655906	F-bottom / Chimney Well	μ Clear®	● black	25 μ l - 340 μ l	10 / 40

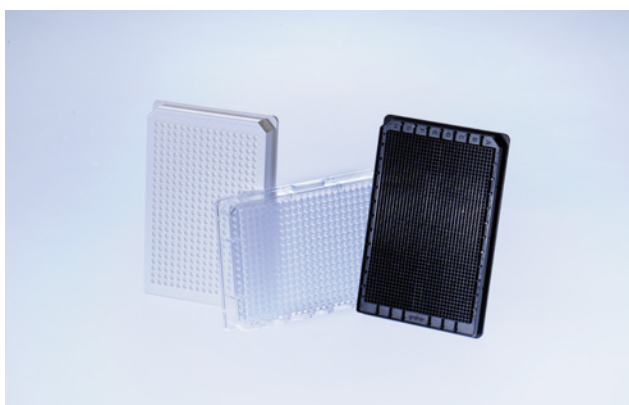

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
--------------------------	------------------------------	--------------------------	---------------

Non-binding Microplates


384 and 1536 Well

- / Clear / black / white
- / With F-bottom or as Small Volume option
- / Solid bottom or μ Clear® film bottom

Well profile: F-bottom, Raw material: PS, Binding characterisitc: Non-binding, Lid: no

Item no.	Well format	Bottom	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781901	384	solid	○ clear			15 μ l - 110 μ l	10 / 40
781904	384	solid	○ white			15 μ l - 110 μ l	10 / 40

Item no.	Well format	Bottom	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
781900	384	solid	● black			15 µl - 110 µl	10 / 40
781903	384	µClear®	○ white			15 µl - 110 µl	10 / 40
781906	384	µClear®	● black			15 µl - 110 µl	10 / 40
784904	384	solid	○ white	HiBase	Small Volume	4 µl - 25 µl	10 / 40
784900	384	solid	● black	HiBase	Small Volume	4 µl - 25 µl	10 / 40
782904	1536	solid	○ white	HiBase		3 µl - 10 µl	15 / 60
782900	1536	solid	● black	HiBase		3 µl - 10 µl	15 / 60

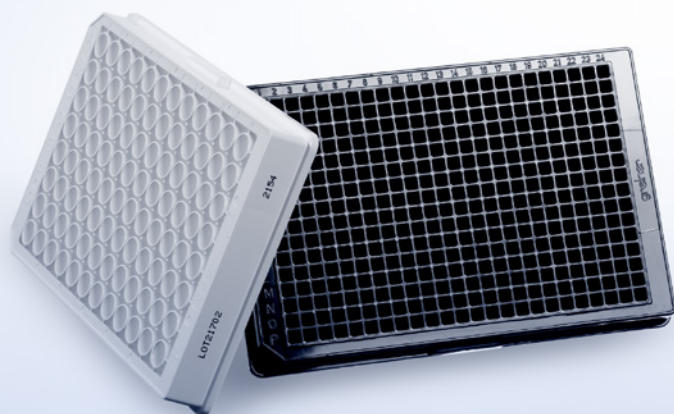

/ For more detailed information on well profiles and technical details please refer to our data sheets on our website: www.gbo.com

STREPTAVIDIN-COATED MICROPLATES


Streptavidin-coated solid phases serve as reliable binding surfaces for all types of biotinylated molecules.

Numerous ligands can be biotinylated simply and due to the low molecular weight of biotin (244 Da) the functionality of the molecules is normally not impaired. Thus streptavidin-coated solid phases make it possible to rapidly isolate, determine and quantify components from a reaction mixture. By immobilising

the biotinylated substance, it is also possible to reproduce complete reaction chains on a streptavidin solid phase, e. g. enzyme immunoassays, enzyme activity assays, DNA hybridisation techniques and quantification of PCR products.

The high-purity streptavidin is bound to the plate surface in a uniform and stable layer. The coefficient of variation from well to well is under 5 % for 96 well microplates and under 8 % for 384 well microplates.

- / Available in 96 well and 384 well format
- / Clear / black / white
- / 3-year shelf life at room temperature
- / All plates pre-blocked and ready-to-use

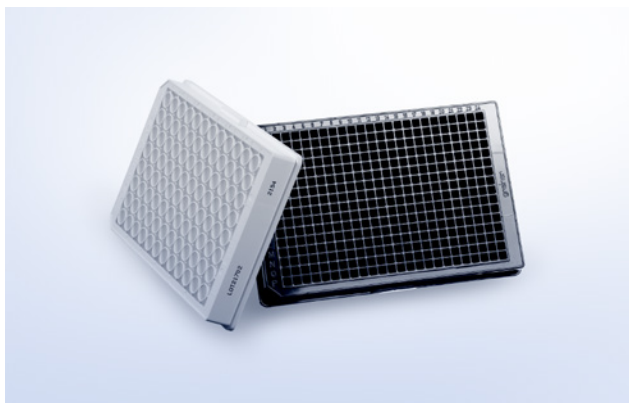


**Streptavidin-coated
Microplates 96 and 384 Well**

- / Long shelf life at room temperature
- / Lot number on each pack
- / Pre-blocked with BSA

Raw material: PS

Item no.	Well format	Well profile	Bottom	Surface treatment	Product colour	Lid	Qty. inner / outer
655990	96	C-bottom	solid	Streptavidin	○ clear	no	5 / 40
655997	96	C-bottom	solid	Streptavidin	● black	no	5 / 40
655994	96	F-bottom	solid	Streptavidin	○ white	no	5 / 40
655996	96	F-bottom	μClear®	Streptavidin	● black	no	5 / 40
781990	384	F-bottom	solid	Streptavidin	○ clear	no	5 / 40
781995	384	F-bottom	solid	Streptavidin	○ white	no	5 / 40
781997	384	F-bottom	solid	Streptavidin	● black	no	5 / 40

/ Further streptavidin-coated microplates are available on request.

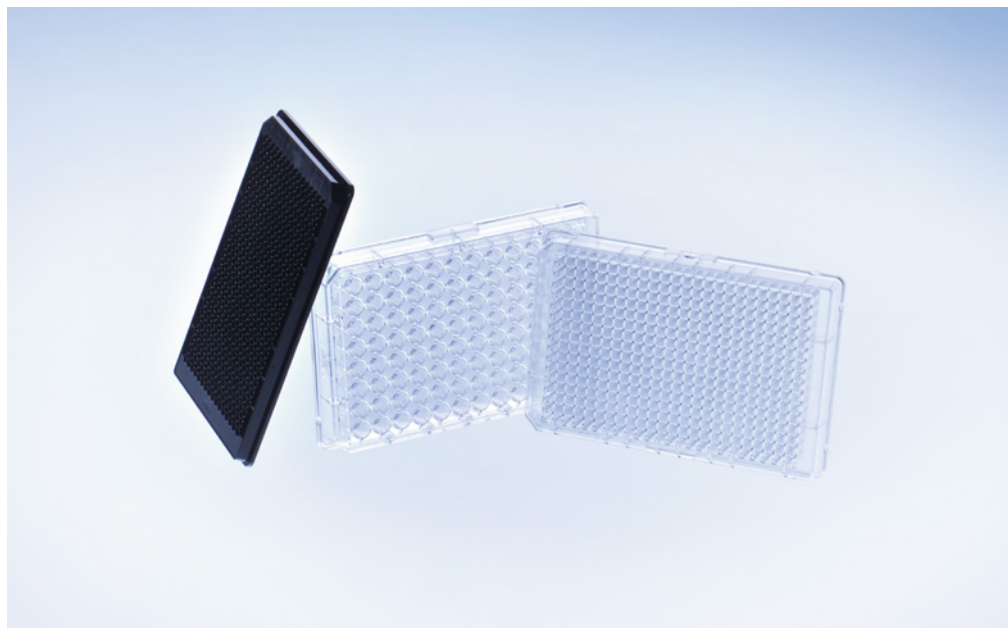

/ Application Note: UV/
VIS Spectroscopy in
Microplates UV-Star®,
µClear®, MICROLON and
CELLSTAR® (F073041)

UV-STAR® MICROPLATES

UV / VIS spectroscopy is a classical analytical method for determining the chemical constitution of a substance and its concentration in aqueous solution.


UV / VIS spectroscopy is usually conducted in quartz glass cuvettes. However, cuvettes do not provide sufficient throughput when dealing with large amounts of samples, and microplates can be used to speed up work.

Standard polystyrene microplates are only partially suitable for transmission measurements in the UV. Polystyrene absorbs UV especially in the short-wavelength range (< 320 nm). µClear® microplates with a thin polystyrene film base already

have much lower background values and can be used up to 340 nm without any problem. The adaptation of the patented µClear® process technology to a new, innovative UV-transparent material has made it possible to produce microplates that extend the transmission range up to 230 nm.

For the determination of nucleic acid and protein concentrations at 260 nm or 280 nm without background interference UV-Star® microplates are the ideal alternative to expensive and fragile quartz glass plates or cuvettes. UV-Star® plates are also DMSO-resistant and can be stored at -20 °C without any problem.

- / Available in 96 well and 384 well format
- / With cycloolefin film bottom
- / Optical window down to 230 nm ideal for nucleic acid determinations at 260 nm/280 nm
- / For measurements of protein concentration at 280 nm

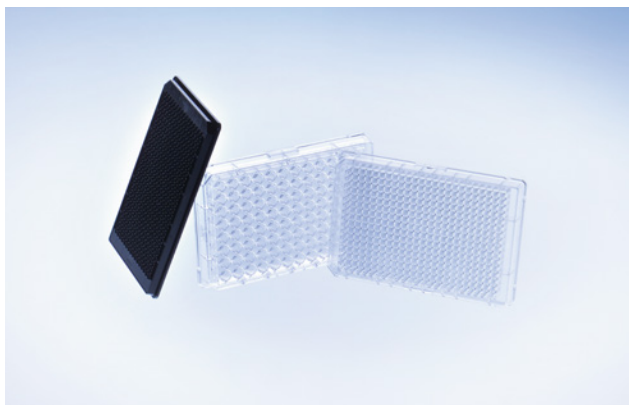

UV-Star® Microplates

- / Available in 96 well and 384 well format
- / With cycloolefin film bottom
- / Optical window down to 230 nm ideal for nucleic acid determinations at 260 nm/280 nm
- / For measurements of protein concentration at 280 nm

FREE OF
detectable
DNase


FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic

Bottom: Cycloolefin film, Raw material: COC, Surface treatment: untreated, Lid: no

Item no.	Well format	Well profile	Product colour	Plate geometry	Plate design	Working volume (well)	Qty. inner / outer
655801	96	F-bottom / Chimney Well	○ clear			25 µl - 340 µl	10 / 40
655809	96	F-bottom / Chimney Well	● black			25 µl - 340 µl	10 / 40
675801	96	F-bottom	○ clear		half area	15 µl - 175 µl	10 / 40
781801	384	F-bottom	○ clear			15 µl - 110 µl	10 / 40
788876	384	F-bottom	● black	LoBase	Small Volume	4 µl - 25 µl	10 / 80

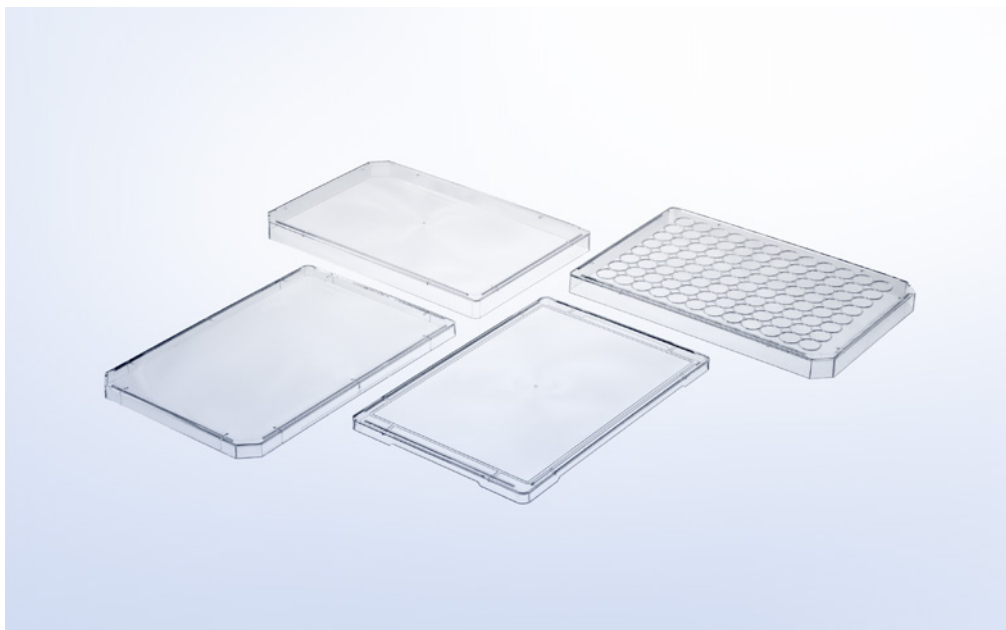

- / Forum No 6: Sealers for microplates and their areas of application in molecular biology and cell culture (F073013)

LIDS / SEALERS / CAPMATS


Lids offer protection against contamination and evaporation during sample storage and cell cultivation.

In addition, Greiner Bio-One offers a full set of adhesive sealers for a wide range of routine and specialised applications, e.g. high-throughput screening, im-

munology, microbiology, molecular biology and PCR.

An alternative method for sealing 96 well plates are CapMats. They are made of ethyl vinyl acetate (EVA), are resistant to DMSO and can be used in a temperature range between -20°C and $+60^{\circ}\text{C}$. CapMats are not pierceable.

- / Polystyrene lids in three different profile heights with and without condensation rings
- / Classical adhesive films coated with acrylate adhesive as well as advanced adhesive films with pressure-sensitive silicone adhesive
- / EVA CapMats for sealing of 96 well plates

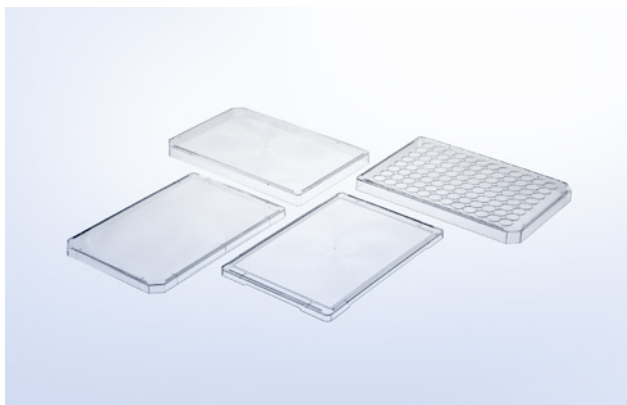


Lids

All sterile lids are non-cytotoxic.



Description: Lid, Raw material: PS

Item no.	Height	Product colour	Condensation rings	Lid type	Sterile	Qty. inner / outer
656101	9 mm	○ clear	no	high		1 / 100
656161	9 mm	○ clear	no	high	+	1 / 100
656170	9 mm	○ clear	yes	high		1 / 100
656171	9 mm	○ clear	yes	high	+	1 / 100
656190	6 mm	○ clear	no	flat		20 / 200
656191	6 mm	○ clear	no	flat	+	20 / 200
691101	4.8 mm	○ clear	no	ultra low		25 / 100
691161	4.8 mm	○ clear	no	ultra low	+	25 / 100

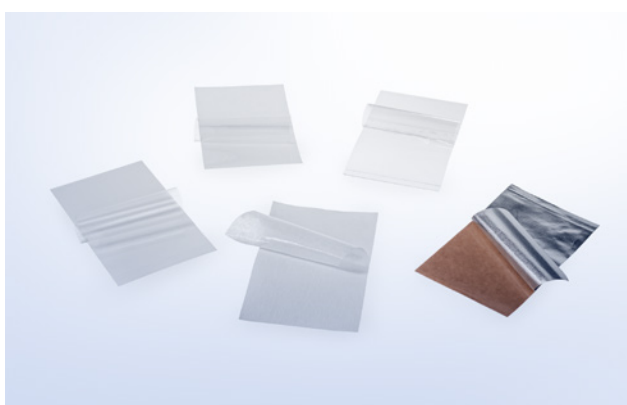


Sealers

The classical sealers such as EASYseal, AMPLIseal, SILVERseal and BREATHseal are coated with an acrylate adhesive. The advanced sealer VIEWseal is coated with a pressure-sensitive silicone adhesive.

Item no.	Description	Feature	Pierceable	Sterile	Qty. inner / outer
676001	EASYseal	clear			100 / 2,000
676090	SILVERseal	aluminium foil	yes		100 / 1,200
676070	VIEWseal	clear			100 / 1,200
676040	AMPLIseal	clear			100 / 2,000
676050	BREATHseal	Gas-permeable			50 / 2,500

Item no.	Description	Feature	Pierceable	Sterile	Qty. inner / outer
676051	BREATHseal	Gas-permeable		+	50 / 500

/ Forum No. 6: Sealers for microplates and their areas of application in molecular biology and cell culture (F073013)

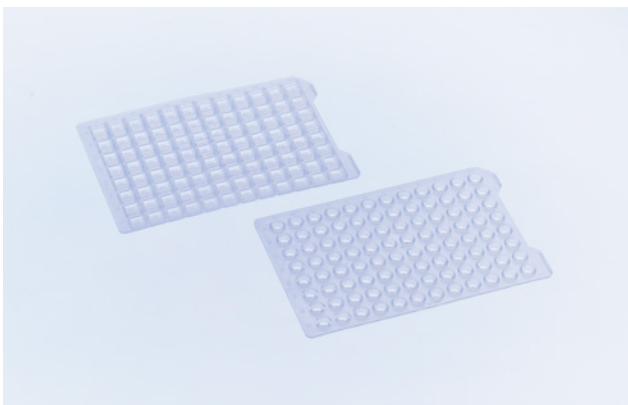

CapMats


- / Made of ethyl vinyl acetate (EVA)
- / Available for 96 well microplates and MASTER-BLOCK®
- / Sterile / non-sterile

Description: 96 Well CapMat, Pierceable: no, Raw material: EVA

Item no.	Nap shape	Sterile	Qty. inner / outer
381070	round		10 / 50
381061	round	+	1 / 50
381080	square		10 / 50
381081	square	+	1 / 50



/ For further information, please refer to our website: www.gbo.com

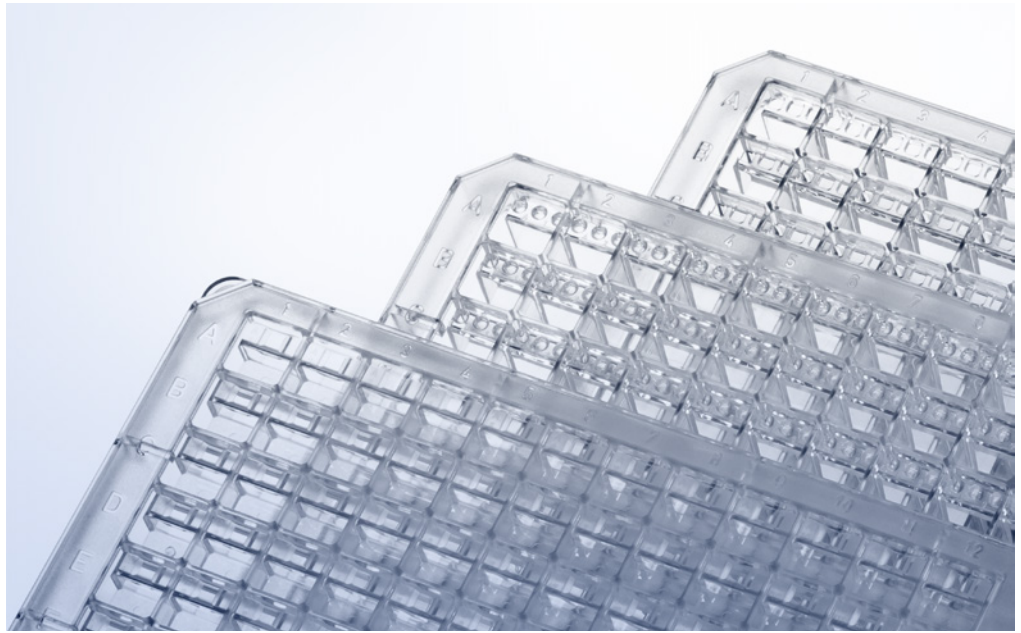

PROTEIN CRYSTALLISATION PLATES


Detailed knowledge of interactions between the structure and function of biomolecules is fundamental for biological, medical and pharmaceutical research. A reliable method is the X-ray structure analysis of protein crystals.

In order to show three-dimensional molecular structures, the proteins must first be crystallised. Since every protein is different and a variety of factors influence protein crystallisa-

tion, the optimal conditions are usually tested with high-throughput technologies (pipetting robots). Greiner Bio-One has developed specific microplates to meet the requirements of high-throughput crystallisation in a short time and with relatively small amounts of protein. Techniques such as vapour diffusion (sitting drop) can be carried out effectively.

LBR (low birefringence) plates are specifically designed for the use of polarised light.

- / For vapour diffusion
- / Made of polystyrene or polyolefin (LBR plates)
- / Footprint conforming to ANSI 1-2004
- / Barcode labelling on request
- / CrystalQuick plates with alphanumeric well coding

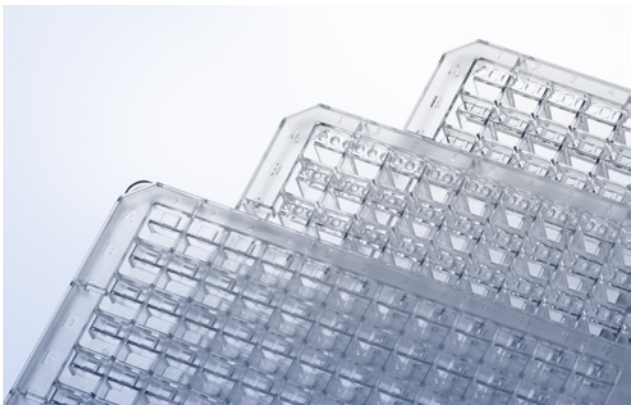

96 Well CrystalQuick Plates

- / Crystallisation plates for sitting drop applications with different well profiles and material properties
- / Optimised for sealing with VIEWseal or AMPLIseal
- / Alphanumeric well coding

Well format: 96


Item no.	Height	Material properties	Well profile	Plate design	Well / reservoir	Well volume, max	Volume per reservoir	Qty. inner / outer
609171	8 mm	standard	flat	square (LP)	1	3.9 μ l	140 μ l	20 / 80
609871	8 mm	LBR	flat	square (LP)	1	3.9 μ l	140 μ l	20 / 80
609120	14.4 mm	standard	concave	round (RW)	3	1.9 μ l	320 μ l	10 / 40
609820	14.4 mm	LBR	concave	round (RW)	3	1.9 μ l	320 μ l	10 / 40



The ELISA technique (Enzyme-Linked Immunosorbent Assay) is one of the most widely used biochemical methods in analytical laboratories and in diagnostics. With this method, analytes such as peptides, proteins, antibodies and hormones are detected and quantified selectively and in low concentrations. In addition, ELISAs can be automated and carried out with a high sample throughput, are relatively inexpensive and the test results are available promptly.

IMMUNOLOGY / HLA

/ ELISA Microplates	146
96 Well ELISA Microplates U-Bottom / V-Bottom / F-Bottom / Half Area	147
/ ELISA Strip Plates	148
96 Well ELISA Strip Plates	149
Single-break Strip Plates	149
/ Immuno Tubes	151
Immuno Tubes	152
/ Terasaki Plates	153
Terasaki Plates	154

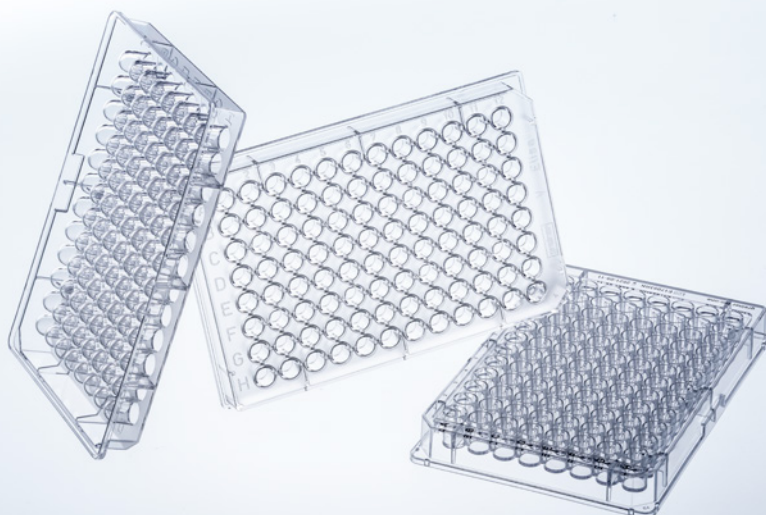

/ You can find more information about this product in our web-shop.



ELISA MICROPLATES


Greiner Bio-One has been manufacturing microplates for diagnostics and immunological research for over 30 years.

The microplate footprint is compatible with automated systems. A key step in ELISA is the binding of one assay component – antigen or antibody – to the solid surface by passive adsorption. Therefore, the features of this surface are crucial for the

performance of the assay. All ELISA microplates from Greiner Bio-One are made out of high-quality virgin polystyrene. The resin is highly transparent and therefore ideally suited for optical measurements.

The quality of our immunological products is constantly controlled in our quality assurance laboratory by means of ELISA.

- / With U-Bottom / V-Bottom / F-Bottom standard (ST) / F-bottom chimney well / half area well geometry
- / With medium and high-binding surfaces

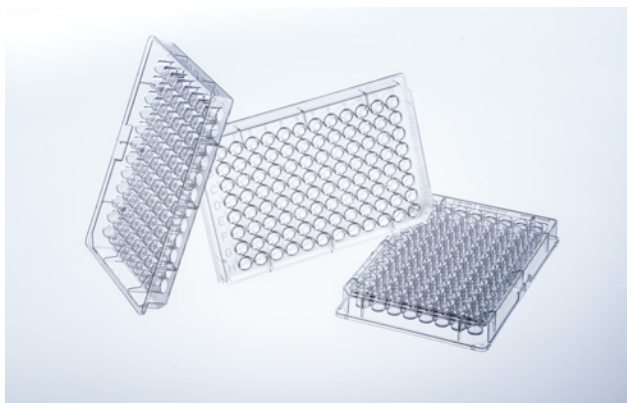

96 Well ELISA Microplates

U-Bottom / V-Bottom / F-Bottom / Half Area

/ Manufactured from crystal clear polystyrene

FREE OF
detectable
DNase


FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
pyrogenic

Well format: 96, Bottom: solid, Raw material: PS, Lid: no

Item no.	Well profile	Binding characterisitic	Binding brand name	Product colour	Plate design	Working volume (well)	Qty. inner / outer
650001	U-bottom	med. binding	MICROLON 200	○ clear		40 µl - 280 µl	10 / 40
650061	U-bottom	high-binding	MICROLON 600	○ clear		40 µl - 280 µl	10 / 40
651001	V-bottom	med. binding	MICROLON 200	○ clear		40 µl - 200 µl	10 / 40
651061	V-bottom	high-binding	MICROLON 600	○ clear		40 µl - 200 µl	10 / 40
655001	F-bottom/ST	med. binding	MICROLON 200	○ clear		25 µl - 340 µl	10 / 40
655061	F-bottom/ST	high-binding	MICROLON 600	○ clear		25 µl - 340 µl	10 / 40
655080	F-bottom / Chimney Well	med. binding	MICROLON 200	○ clear		25 µl - 340 µl	10 / 40
655081	F-bottom / Chimney Well	high-binding	MICROLON 600	○ clear		25 µl - 340 µl	10 / 40
675001	F-bottom	med. binding	MICROLON 200	○ clear	half area	15 µl - 175 µl	10 / 40
675061	F-bottom	high-binding	MICROLON 600	○ clear	half area	15 µl - 175 µl	10 / 40

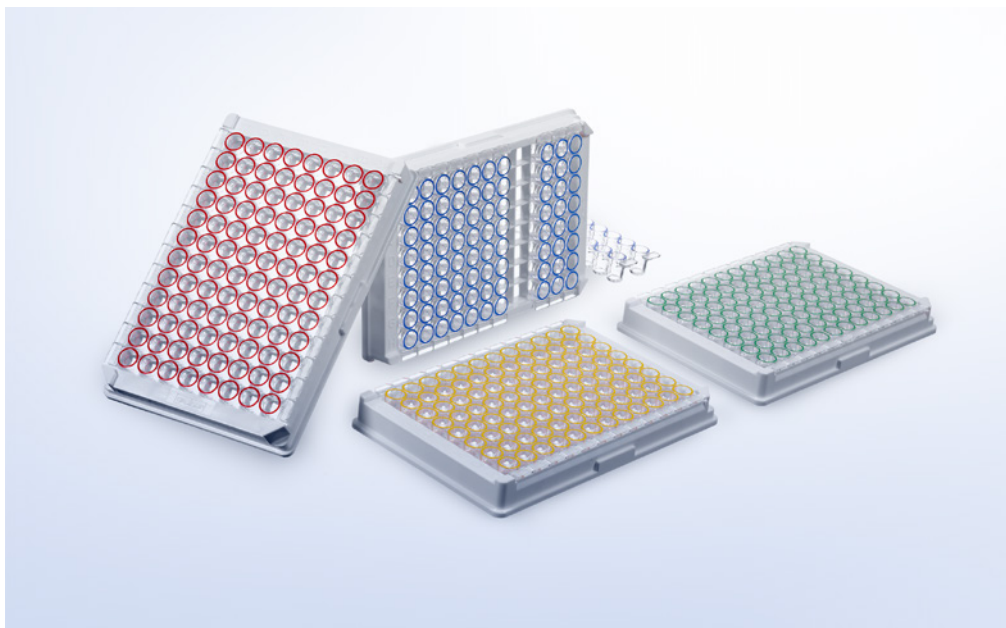

/ You can find more information about this product in our web-shop.




ELISA STRIP PLATES

Greiner Bio-One is offering a large variety of 96 well strip microplates for diagnostic and immunological research applications.

Microplates in strip format offer the advantage of greater flexibility in diagnostics. Individual strips can be removed from the support frame so that the number of tests to be performed can be adjusted to the number of samples and is not predetermined by the microplate format used. In addition, the individual strips can be subjected to a wide variety of different test conditions. The quality of our immunological products is constantly

controlled in our quality assurance laboratory by means of ELISA. Clear MICROLON products are tested in a colourimetric immunoassay, black FLUOTRAC plates in a fluorescence immunoassay, white LUMITRAC plates in a luminescence immunoassay. The portfolio is completed by C8 single-break strip plates for immunological applications. These are supplied as twelve 8-well strips in a support frame with 96 spaces (12 x 8 matrix). The individual wells can be broken off separately ("single-break" option) and the number of tests performed can thus be precisely adjusted to the number of samples.

96 Well ELISA Strip Plates

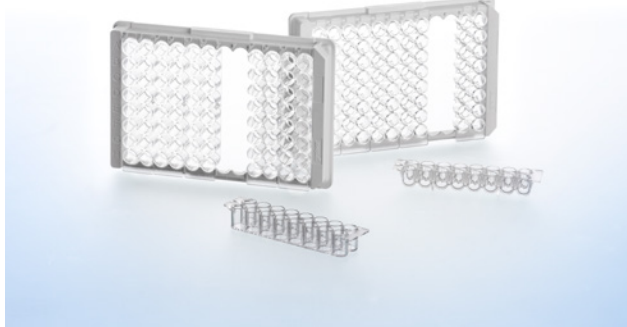


- / Strips mounted in frame
- / Black and white F16 strip plates on request

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
--------------------------	------------------------------	--------------------------	---------------

Well format: 96, Well profile: F-bottom, Bottom: solid, Raw material: PS, Working volume (well): 20 µl – 350 µl, Lid: no

Item no.	Binding characterisitc	Binding brand name	Product colour	Plate design	Qty. inner / outer
762070	med. binding	MICROLON 200	○ clear	strip plate 12x8	10 / 100
762071	high-binding	MICROLON 600	○ clear	strip plate 12x8	10 / 100
762075	med. binding	LUMITRAC 200	○ white	strip plate 12x8	10 / 100
762074	high-binding	LUMITRAC 600	○ white	strip plate 12x8	10 / 100
762076	med. binding	FLUOTRAC 200	● black	strip plate 12x8	10 / 100
762077	high-binding	FLUOTRAC 600	● black	strip plate 12x8	10 / 100
756070	med. binding	MICROLON 200	○ clear	strip plate 6x16	10 / 100
756071	high-binding	MICROLON 600	○ clear	strip plate 6x16	10 / 100

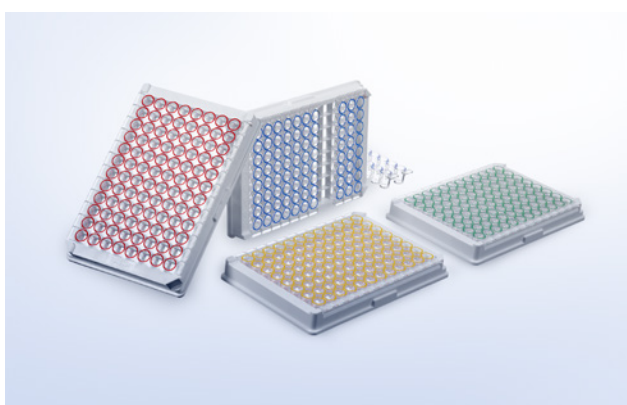

Single-break Strip Plates

- / Clear with / without colour coding
- / Strips mounted in frame
- / C8 Single-break strip plates without colour coding cell culture treated on request

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
--------------------------	------------------------------	--------------------------	---------------

Well format: 96, Well profile: C-bottom, Bottom: solid, Raw material: PS, Plate design: strip plate 12x8, Working volume (well): 20 µl – 300 µl, Lid: no

Item no.	Binding characterisitc	Binding brand name	Colour stripe	Colour coding well rim	Qty. inner / outer
705070	med. binding	MICROLON 200	○ clear		10 / 100

Item no.	Binding characterisitc	Binding brand name	Colour stripe	Colour coding well rim	Qty. inner / outer
705071	high-binding	MICROLON 600	○ clear		10 / 100
705063	med. binding	MICROLON 200	○ clear	● red	10 / 100
705073	high-binding	MICROLON 600	○ clear	● red	10 / 100
705074	high-binding	MICROLON 600	○ clear	● blue	10 / 100
705065	med. binding	MICROLON 200	○ clear	● green	10 / 100
705075	high-binding	MICROLON 600	○ clear	● green	10 / 100
705076	high-binding	MICROLON 600	○ clear	● yellow	10 / 100



/ More tubes can be found in the chapter Tubes / Multipurpose containers



IMMUNO TUBES

Immuno tubes are often used for determining hormone levels, for example TSH (thyroid stimulating hormone). With a length of 75 mm and a diameter of 12 mm, they are available in both MICROLON 200 and MICROLON 600 quality.

The quality of our immunological products is constantly controlled in our quality assurance laboratory by means of ELISA.



Immuno Tubes


/ More tubes can be found in the chapter Tubes / Multipurpose containers

Height: 75 mm, Ø: 12 mm, Raw material: PS, Nominal volume: 5 ml

Item no.	Binding characterisitc	Binding brand name	Bottom shape	Qty. inner / outer
115001	med. binding	MICROLON 200	round	250 / 2,000
115061	high-binding	MICROLON 600	round	250 / 2,000



/ We provide Terasaki plates with one years' stability on the surface treatment.

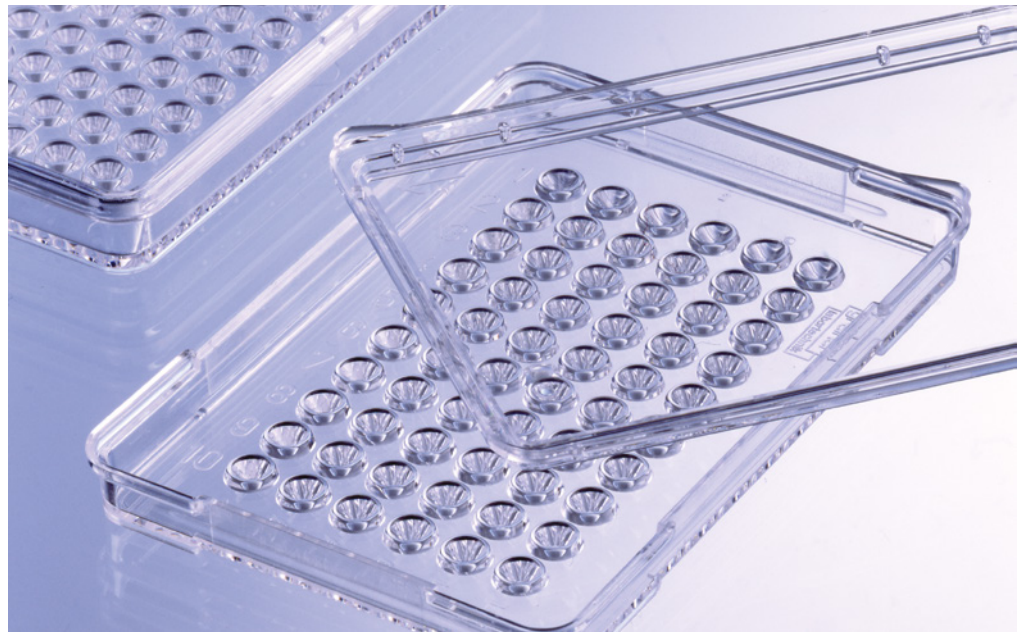

TERASAKI PLATES

Greiner Bio-One Terasaki plates are suitable for all applications for serological determination of HLA antigens. The plates are supplied with 60 wells.

The human leucocyte antigen (HLA) system is the major histocompatibility complex (MHC) of humans and is composed of the two polymorphic classes HLA-I (A, B, and C) and HLA-II (DR, DQ, and DP). Basically, four different areas of indication can be distinguished for HLA typing: transplantation, transfusion, disease association and forensic.

The serological determination of HLA proteins of the HLA-A, -B, -C and -DR genetic positions is primarily performed with the complement-dependent microlymphocytotoxicity test (LCT) or Terasaki test, which has been standardised since 1964. The basis for this test method is the cytolysis of the lymphocytes to be tested, which is caused by the antibody-antigen mediated activation of the complement system. Permeabilised lymphocytes are generally stained with chromophores or fluorophores and evaluated microscopically.


Terasaki Plates

- / 60 well Terasaki plates
- / Manufactured from crystal clear polystyrene



Well format: 60, Stackable: no, Surface treatment: TC, Well volume, max: 11.5 μ l, Working volume (well): \leq 10 μ l, Lid: yes

Item no.	Qty. inner / outer
653180	10 / 270
653190	120 / 480

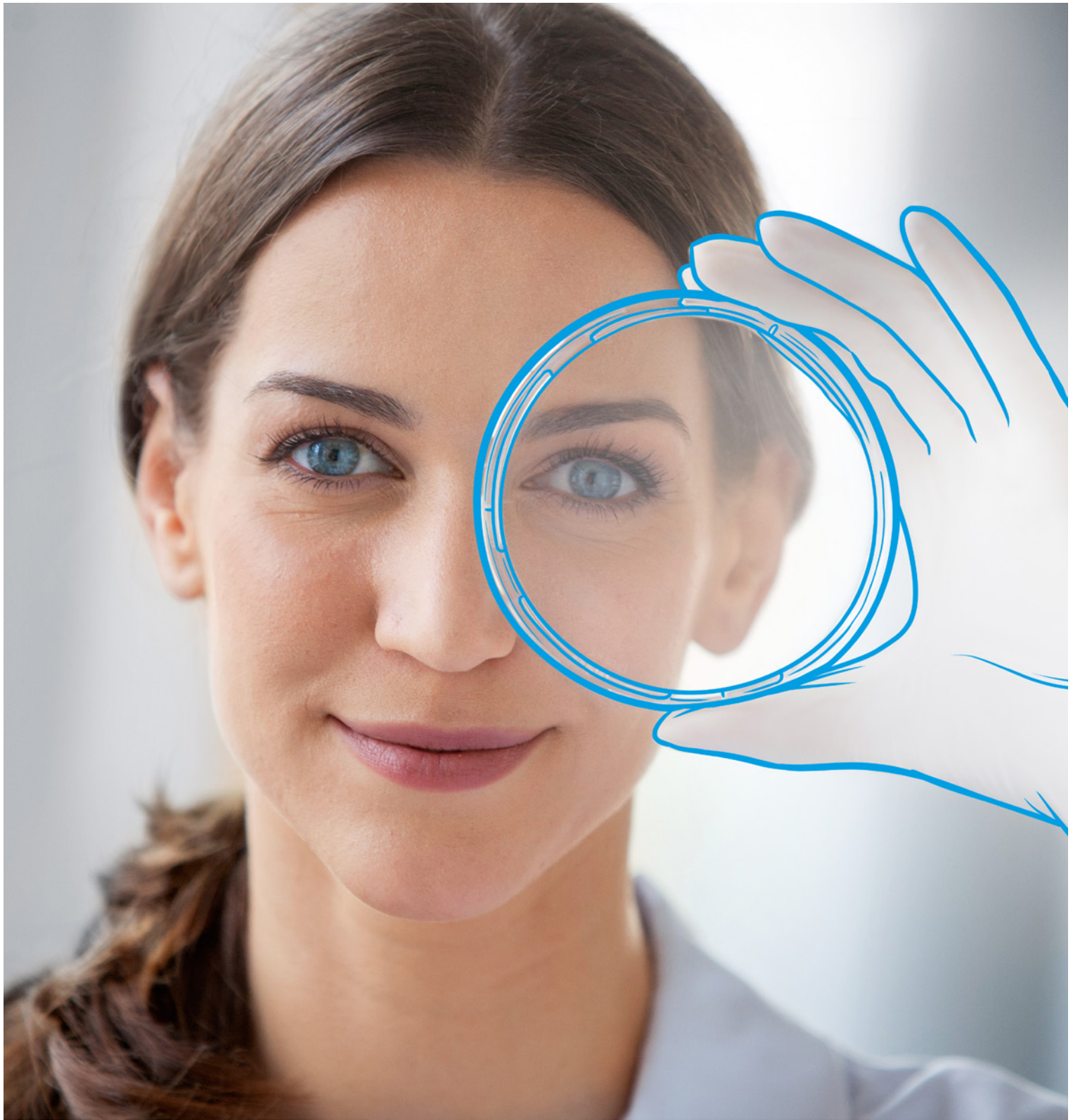

The study of microorganisms, their metabolism and their effects on other (macro-)organisms generally refers to the field of microbiology. The various detection methods are used in many research areas and in industry. For example, tests for the detection of microorganisms and for quality assurance are also applied in the production of pharmaceuticals, cosmetics, food and beverages.

MICROBIOLOGY

/ Dishes / Plates / Others	158
Petri Dishes	159
Divided Petri Dish	160
Square Petri Dish / OneWell Plate	160
Swab Tube with Cotton Swab	161
Disposable Inoculation Loops / Needles	161
Disposal Bags	162



/ For exact dimensions of our petri dishes, please refer to the product data sheets on our website.

DISHES / PLATES / OTHERS

Greiner Bio-One offers petri dishes with diameters of 35, 60, 94, 100 and 145 mm. Manufactured from highly transparent polystyrene (PS), they are heat-resistant up to 60 °C for use with hot agar. Depending on the application vented and non-vented design can be selected. Besides the standard round design,

a square-profile dish (120 mm x 120 mm x 17 mm) and the One-Well Plate in SBS format are available.


The product range is completed by a special model with two compartments which enables parallel experiments with different samples or parameters in one dish.



Petri Dishes

- / Available in different sizes
- / Easy stacking
- / With vents
- / Manufactured from crystal clear polystyrene

Item no.	Height	Ø nominal size	Ventnock	Sterile	Qty. inner / outer
627102	10 mm	35 mm	yes		10 / 740
628102	15 mm	60 mm	yes		20 / 600
628161	15 mm	60 mm	yes	+	20 / 600
632180	16 mm	94 mm	no		20 / 480
633180	16 mm	94 mm	yes		20 / 480
633181	16 mm	94 mm	yes	+	20 / 480
664102	20 mm	100 mm	yes		15 / 360
664161	20 mm	100 mm	yes	+	15 / 360
639102	20 mm	145 mm	yes		15 / 120
639161	20 mm	145 mm	yes	+	15 / 120

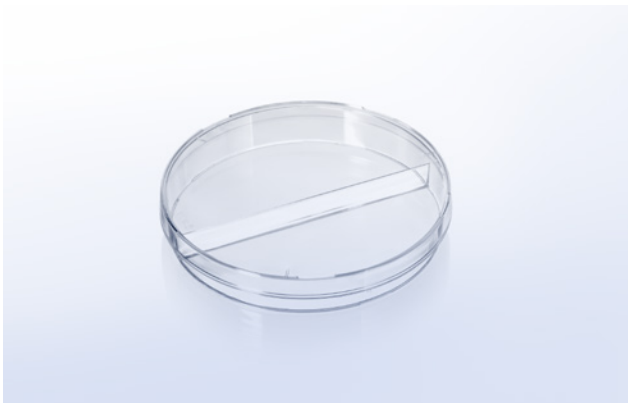


Divided Petri Dish

- / With vents
- / Manufactured from crystal clear polystyrene

Compartments: 2, Height: 15 mm, Ø Nominal size: 94 mm, Total volume (well): 20 ml, Vent nock: yes

Item no.	Qty. inner / outer
635102	20 / 480

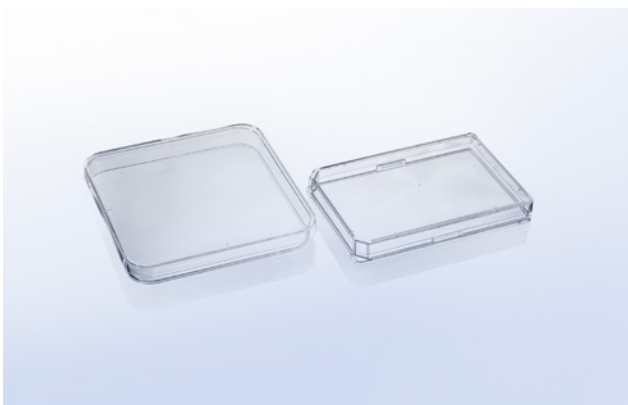


Square Petri Dish / OneWell Plate

- / With vents
- / Manufactured from crystal clear polystyrene
- / Optimised space requirement compared to round dishes

Lid: yes, Vent nock: yes

Item no.	Feature	Well format	Length	Width	Product colour	Total volume (well)	Qty. inner / outer
670102	microplate format	1	127.8 mm	85.5 mm	○ clear	113.7 ml	8 / 32
688102	square		120 mm	120 mm	○ clear		10 / 240

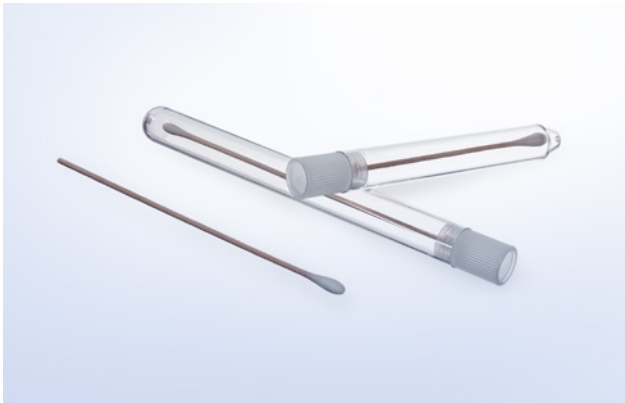


STERILE

Swab Tube with Cotton Swab

- / Suitable for taking non-human bacteriological, serological or cytological samples in veterinary research
- / For hygienic controls in food industry as well as in environmental sampling

Feature: swab tube, polystyrene, Ø: 16 mm, Sterile: +

Item no.	Height	Sterile	Qty. inner / outer
420161	110 mm	+	1,300 / 1,300
421161	152 mm	+	- / 1,000

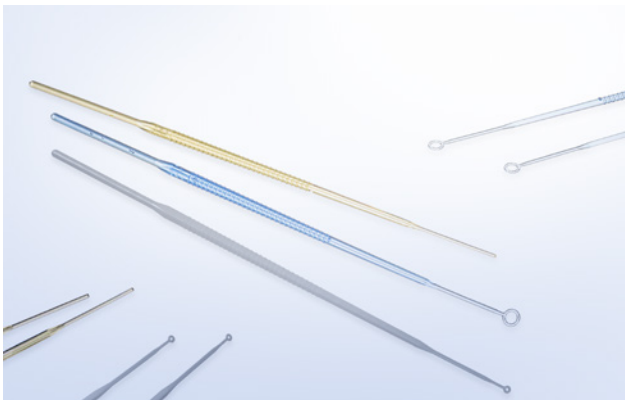

STERILE

Disposable Inoculation Loops / Needles

- / Needles ideal for picking single colonies
- / Flexible loops for easier collection and inoculation
- / Colour coding for volume differentiation

Length: 200 mm, Sterile: +

Item no.	Description	Product colour	Volume	Sterile	Qty. inner / outer
731101	inoculation loop	○ white	1 µl	+	50 / 2,000
731161	inoculation loop	○ white	1 µl	+	1 / 600
731165	inoculation loop	○ white	1 µl	+	10 / 3,000
731170	inoculation loop	● blue	10 µl	+	50 / 2,000
731171	inoculation loop	● blue	10 µl	+	1 / 600
731175	inoculation loop	● blue	10 µl	+	10 / 3,000
731180	inoculation needle	● yellow		+	50 / 2,000
731181	inoculation needle	● yellow		+	1 / 600

Item no.	Description	Product colour	Volume	Sterile	Qty. inner / outer
731185	inoculation needle	● yellow		+	1 / 3,000



Disposal Bags

Disposal bags made of polypropylene for sterilisation in the steam autoclave. For users of hot-air sterilisers, disposal bags made of polyamide are suitable for use up to +160 °C.

Foil thickness: 0.05 mm


Item no.	Feature	Length	Width	Nominal capacity	Suitable for steam autoclaves	Suitable for hot air sterilizers	Material	Qty. inner / outer
643201		500 mm	300 mm	10 l	yes		PP	500 / 500
643203	Imprint "Biohazard"	500 mm	300 mm	10 l	yes		PP	500 / 500
643401		500 mm	300 mm	10 l		yes	PA	500 / 500
644201		780 mm	400 mm	30 l	yes		PP	500 / 500
644203	Imprint "Biohazard"	780 mm	400 mm	30 l	yes		PP	500 / 500
644401		780 mm	400 mm	30 l		yes	PA	500 / 500
646201		780 mm	600 mm	65 l	yes		PP	500 / 500
646203	Imprint "Biohazard"	780 mm	600 mm	65 l	yes		PP	500 / 500
646401		780 mm	600 mm	65 l		yes	PA	300 / 300
649201		1,100 mm	700 mm	130 l	yes		PP	350 / 350
649203	Imprint "Biohazard"	1,100 mm	700 mm	130 l	yes		PP	350 / 350
649401		1,100 mm	700 mm	130 l		yes	PA	200 / 200



Greiner Bio-One offers a comprehensive range of tubes and multipurpose beakers for various applications. The vessels are available in different materials, with and without caps and with labelling options.

TUBES / MULTIPURPOSE BEAKERS

/ Tubes without closures	166
Tube Polystyrene	167
Tube Polypropylene	167
Grip Stoppers	168
Tube Polypropylene for Storage Box	168
96 Well Storage Box	169
/ Tubes with closures	170
Centrifuge Tube Polypropylene - 15 ml	171
Centrifuge Tube Polypropylene - 50 ml	171
CELLreactor	172
Tubes with closures Polystyrene	173
Tubes with Two-position Cap	173
/ Leucosep Separation Tubes	174
Leucosep 12 ml and 50 ml	175
/ Multipurpose Containers / Beakers	176
Multipurpose Containers Polystyrene	177
Multipurpose Beaker Polypropylene	177
Drosophila Containers	178
Ceaprene Stoppers for Drosophila Containers	178

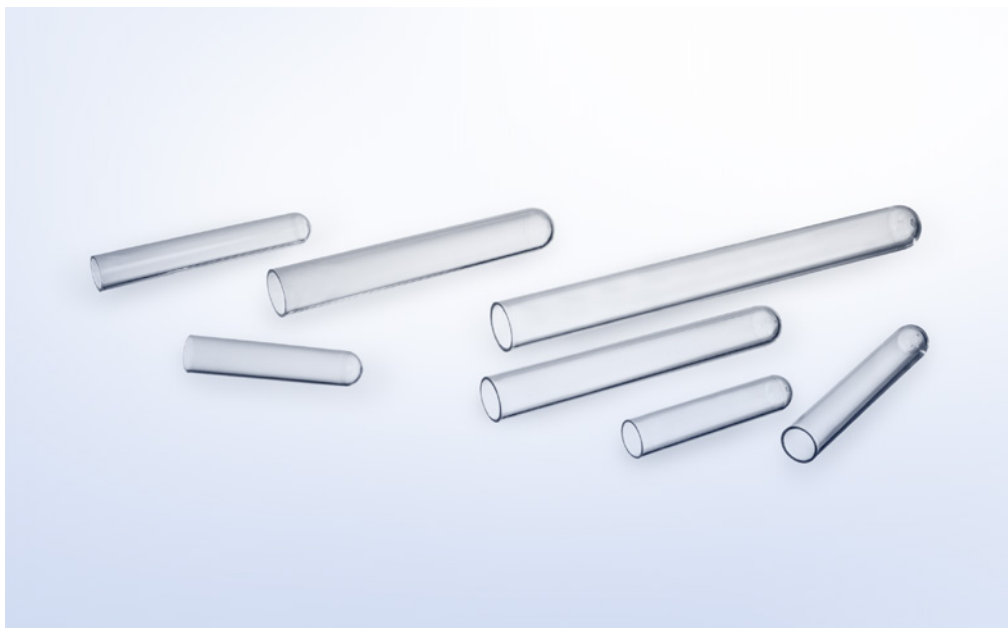

- / For the max. relative centrifugal force (RCF) and the chemical/thermal resistance of our tubes please refer to the technical appendix.


TUBES WITHOUT CLOSURES

The range of Greiner Bio-One tubes is very versatile and meets a wide variety of different demands.

Our tubes are made of different materials: **Polystyrene (PS)** is ideally suited for optical measurements as a result of its high clarity. **Polypropylene (PP)** is re-

commended for the storage of chemical and biological samples due to its thermal, mechanical and chemical resistance. Greiner Bio-One tubes without closure are available with round bottom. In addition, it is possible to order appropriate closures in the form of grip stoppers and screw caps.


- / With round or conical bottom
- / Available in polystyrene or polypropylene
- / Appropriate grip stoppers available separately

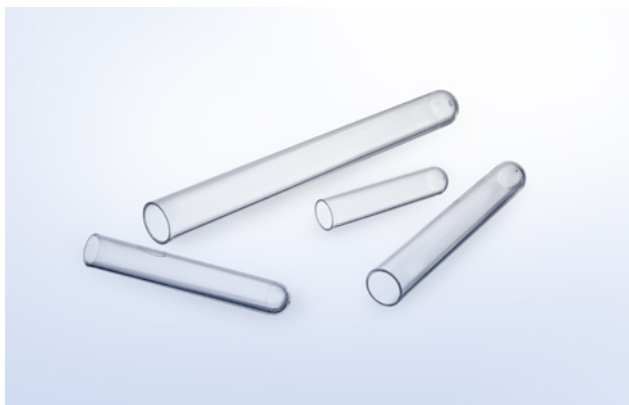**Tubes****Polystyrene**

- / High transparency
- / Available in different sizes and packaging units

Raw material: PS, Bottom shape: round

Item no.	Height	Ø	Working volume	Nominal volume	Qty. inner / outer
112101	55 mm	12 mm	≤3 ml	4 ml	240 / 3,600
115101	75 mm	12 mm	≤4 ml	5 ml	250 / 2,000
136101	100 mm	14 mm	≤8 ml	10 ml	1,400 / 1,400
160101	100 mm	16 mm	≤10.5 ml	12 ml	1,600 / 1,600
187101	100 mm	17 mm	≤12 ml	14 ml	1,500 / 1,500
169101	152 mm	16 mm	≤18 ml	20 ml	1,500 / 1,500

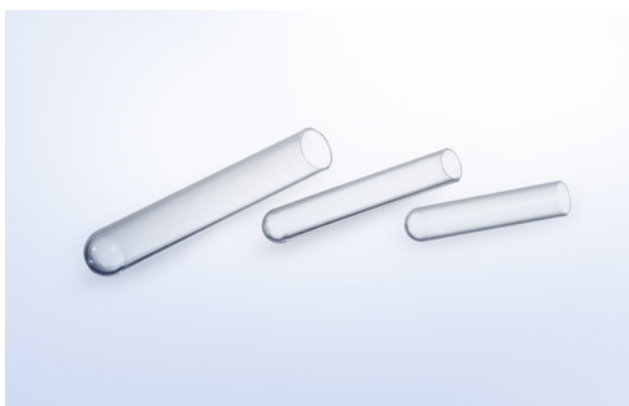**Tubes****Polypropylene**

- / Suited for sample storage
- / Thermal, chemical and mechanical stability
- / Available in different sizes and packaging units

Raw material: PP, Bottom shape: round

Item no.	Height	Ø	Working volume	Nominal volume	Qty. inner / outer
112201	55 mm	12 mm	≤3 ml	4 ml	240 / 3,600
115201	75 mm	12 mm	≤4 ml	5 ml	250 / 2,000
160201	100 mm	16 mm	≤10 ml	12 ml	- / 1,600
187201	100 mm	17 mm	≤12 ml	14 ml	1,500 / 1,500

Tubes / Multipurpose Beakers

Tubes without closures



Grip Stoppers

/ Made of high-grade polyethylene

Description: Grip stopper, Raw material: PE

Item no.	Product colour	Suitable for tubes Ø	Qty. inner / outer
303321	○ natural	12 mm	500 / 20,000
310321	○ natural	16 mm	100 / 8,000




Tubes Polypropylene for Storage Box

/ Tubes for storage box item no. 975502

Height: 44 mm, Ø: 8.5 mm, Raw material: PP, Bottom shape: round, Working volume: ≤1 ml, Nominal volume: 1.2 ml

Item no.	Feature	Qty. inner / outer
102201	Tube	1,000 / 5,000
102270	Tube chain	120 / 600

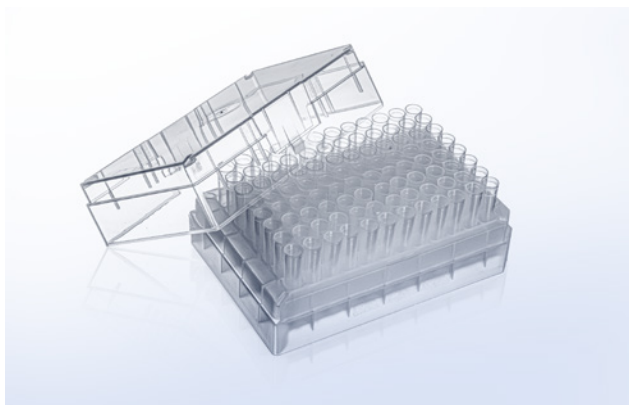**96 Well Storage Box**

- / Storage box made of polypropylen
- / With / without 96 polypropylene vessels
- / All components autoclavable

Item no.	Qty. inner / outer
975502	5 / 50
975570	5 / 50



/ Selected tubes are certified with ACT[®] Ecolabel. Learn more:



TUBES WITH CLOSURES

Greiner Bio-One tubes with closures are made of different materials: **Polystyrene (PS)** is ideally suited for optical measurements as a result of its high clarity. **Polypropylene (PP)** is recommended for the storage of chemical and biological samples due to its thermal, mechanical and chemical resistance. Greiner Bio-One tubes without closure are available with round or with conical bottom as well as with or without skirt. They are available in sterile or non-sterile versions.

The product range is completed by tubes with a two-position cap. This special cap enables ventilation or an airtight closure of the tube, depending on the position of the stopper.

To promote sustainability and transparency, many of our centrifugation tubes have already been certified by My Green Lab[®] with an excellent rating. The respective ACT[®] Ecolabel allows a clear and reliable assessment of the environmental impact of our products.



Centrifuge Tubes
Polypropylene - 15 ml

- / Optimal mechanical, thermal and chemical stability
- / Triple-packed options available
- / Selected product versions certified with ACT® Ecolabel (*)

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	--

Height: 120 mm, Ø: 17 mm, Support skirt: no, Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤14 ml, Nominal volume: 15 ml, Cap design: screw cap, Sterile: +

Item no.	Product colour	Cap colour	Colour writing field	Graduation colour	Type of packaging	Sterile	Qty. inner / outer
188261-N*	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	rack	+	50 / 500
188271-N*	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	bag	+	50 / 1,000
188271-TRI	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	triple packed	+	5 / 500

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



Centrifuge Tubes
Polypropylene - 50 ml

- / Optimal mechanical, thermal and chemical stability
- / Triple-packed options available
- / Selected product versions certified with ACT® Ecolabel (*)

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	--

Height: 115 mm, Ø: 30 mm, Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤50 ml, Nominal volume: 50 ml, Cap design: screw cap, Sterile: +

Item no.	Support skirt	Product colour	Cap colour	Colour writing field	Graduation colour	Type of packaging	Sterile	Qty. inner / outer
210650	yes	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	rack	+	25 / 300
210651	yes	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	bag	+	25 / 500
227261	no	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	bag	+	20 / 500
227261-N*	no	<input type="radio"/> natural	<input checked="" type="radio"/> blue	<input type="radio"/> white	<input checked="" type="radio"/> blue	bag	+	25 / 500

Tubes / Multipurpose Beakers

Tubes with closures

Item no.	Support skirt	Product colour	Cap colour	Colour writing field	Graduation colour	Type of packaging	Sterile	Qty. inner / outer
227270	no	○ natural	● blue	○ white	● blue	rack	+	25 / 300
227270-N*	no	○ natural	● blue	○ white	● blue	rack	+	25 / 250
227261-TRI	no	○ natural	● blue	○ white	● blue	triple packed	+	5 / 250

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

/ Please contact your local distribution partner for the availability of 227261-N / 227270-N or 227261 / 227270.



CELLreactor

15 ml and 50 ml polypropylene tube with filter screw cap

- / For cultivation of suspension cells and expansion of aerobic microorganisms
- / Facilitates a high number of parallel experiments
- / Maximal sterility and excellent gas exchange
- / Conical tube design and in-tube harvest

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Graduation: yes, Writing field: yes, Raw material: PP, Bottom shape: conical, Cap design: filter screw cap, Sterile: +

Item no.	Height	Ø	Cap colour	Working volume	Nominal volume	Sterile	Qty. inner / outer
188241	120 mm	17 mm	● blue	1 ml - 5 ml	15 ml	+	20 / 500
227245	115 mm	30 mm	● blue	1 ml - 35 ml	50 ml	+	20 / 500



Tubes with closures

Polystyrene

- / High transparency
- / In bulk or rack packaging

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------

Height: 120 mm, Ø: 17 mm, Support skirt: no, Graduation: yes, Writing field: yes, Raw material: PS, Bottom shape: conical, Working volume: ≤14 ml, Nominal volume: 15 ml, Cap design: screw cap, Sterile: +

Item no.	Cap colour	Sterile	Qty. inner / outer
188161	● blue	+	50 / 500
188171	● blue	+	100 / 1,000



Tubes

Two-position Cap

The two-position cap enables ventilation or an airtight closure of the tube, depending on the position of the stopper.

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------

Bottom shape: round, Cap design: Two-position Cap, Sterile: +

Item no.	Height	Ø	Graduation	Writing field	Raw material	Graduation colour	Working volume	Nominal volume	Sterile	Qty. inner / outer
115262	75 mm	12 mm	yes	no	PP		≤4 ml	5 ml	+	25 / 2,000
120180	75 mm	12.4 mm	-	no	PS	● blue	≤4 ml	4.5 ml	+	25 / 2,000
187262	95 mm	18 mm	yes	yes	PP		≤12 ml	14 ml	+	25 / 1,000
191180	95 mm	18 mm	-	no	PS		≤12.5 ml	14 ml	+	25 / 1,000



/ For further information, please refer to our website: www.gbo.com



LEUCOSEP SEPARATION TUBES

Different separation techniques can be used to enrich certain particles (DNA, RNA, proteins, cells etc.) specifically from complex biological mixtures such as cell and tissue homogenates, blood or urine, so that they can then be selectively investigated. The separation of cells by density gradient centrifugation has proven to be the most often used method.

Leucosep has been developed for optimal separation of lymphocytes and peripheral mononuclear cells (so-called PBMCs) from human whole blood and bone marrow by means of density gradient centrifugation. The key feature of Leucosep is the porous barrier incorporated into the centrifuge tube made of highly translucent polypropylene.

/ Leucosep for the separation of lymphocytes/ mononuclear cells from blood/bone marrow

**Leucosep****12 ml and 50 ml**

Efficient separation of lymphocytes and mononuclear cells from peripheral blood and bone marrow.

- / Easy to fill
- / No re-contamination with unwanted red blood cells
- / PBMC isolation from whole blood in 15 min

Type of sample: Blood

Item no.	Separation medium	Bottom shape	Sample volume	Sterile	Aseptic	Qty. inner / outer
163288	pre-filled with Leucosep separation medium	round	3 - 8 ml		+	50 / 500
163290		round	3 - 8 ml	+		50 / 500
227288	pre-filled with Leucosep separation medium	conical	15 - 30 ml		+	25 / 250
227290		conical	15 - 30 ml	+		25 / 300

/ For IFU(F071043) please refer to our website: www.gbo.com



/ For further information, please refer to our website: www.gbo.com



MULTIPURPOSE CONTAINERS / BEAKERS

Multipurpose containers or beakers can be used for a variety of different applications. They can be applied universally for academic and non-human medical purposes. Multipurpose containers are made of polystyrene or polypropylene and are available in different sizes. They feature either plastic or metal caps and can be supplied with or without plain or printed label.

Containers for plant culture are made of a very clear material which ensures maximum light transmission and thus rapid and successful growth of plant cul-

tures. These culture containers are not only suitable for the proliferation of plant cultures but can also be used as transport containers and are available with or without lids. Our **drosophila containers** are ideally suited for the cultivation of *Drosophila melanogaster*. The bottom part of the container is made of polystyrene. Drosophila containers and containers for plant culture can be supplied with a separately orderable ceaprene stopper. This stopper is gas-permeable and made of water-repellent material.

- / Multipurpose containers in polystyrene or polypropylene
- / Polystyrene containers for plant culture
- / Polystyrene drosophila containers



Multipurpose Containers

Polystyrene

- / Manufactured from crystal clear polystyrene
- / Available in different sizes
- / Can be used universally for academic and non-human medical purposes
- / With or without plain or printed label
- / Plastic or metal cap

Raw material: PS, Cap design: screw cap

Item no.	Feature	Height	Ø	Support skirt	Bottom shape	Cap colour	Working volume	Nominal volume	Cap material	Sterile	Qty. inner / outer
201150		91 mm	25 mm	yes	conical	○ white	≤23 ml	30 ml	plastics		400 / 400
201170		91 mm	25 mm	yes	conical	○ white	≤23 ml	30 ml	plastics	+	400 / 400
201152	with plain label	91 mm	25 mm	yes	conical	○ white	≤23 ml	30 ml	plastics		400 / 400
201172	with plain label	91 mm	25 mm	yes	conical	○ white	≤23 ml	30 ml	plastics	+	400 / 400
224170		77 mm	49 mm			● silver	≤80 ml	100 ml	metal		20 / 160
225170		115 mm	49 mm			● silver	≤150 ml	150 ml	metal		20 / 120
225180	with plain label	115 mm	49 mm			● silver	≤150 ml	150 ml	metal		20 / 120



Multipurpose Beakers

Polypropylene

- / Can be used universally for non human medical purposes
- / Sterile and non-sterile

Height: 70 mm, Ø: 55 mm, Graduation: yes, Working volume: ≤120 ml, Nominal volume: 120 ml, Cap design: screw cap

Item no.	Product colour	Cap colour	Sterile	Qty. inner / outer
724410	○ natural	● yellow	+	300 / 300
724411	○ natural	○ white		300 / 300
724412	○ natural	○ white	+	300 / 300

Tubes / Multipurpose Beakers

Multipurpose Containers / Beakers

Item no.	Product colour	Cap colour	Sterile	Qty. inner / outer
724414	○ natural	● red	+	300 / 300

/ Item No., lot number and expiry date are printed on the screw cap of the beaker to provide additional information and improve traceability



Drosophila Containers

- / Ideally suited for the cultivation of *Drosophila melanogaster*
- / Choice of different sizes
- / Bottom part of container made of polystyrene

Raw material: PS

Item no.	Height	Ø	Working volume	Nominal volume	Qty. inner / outer
205101	64 mm	27 mm		28 ml	1,500 / 1,500
217101	82 mm	36 mm		68 ml	605 / 605
960177	100 mm	53 mm	≤150 ml	175 ml	315 / 315

/ Can be supplied with a separately orderable ceaprene stopper. This stopper is gas-permeable and made of water-repellent material.

/ In combination with top part (Item no. 960178), 960177 can also be used as container for plant culture.



Ceaprene Stoppers for Drosophila Containers

- / Gas-permeable
- / Made of water-repellent material

Item no.	Suitable for tubes Ø	Qty. inner / outer
354070	27 mm	- / 1,500
330070	36 mm	- / 605
332070	53 mm	- / 315



For sample freezing and storage, Greiner Bio-One offers a comprehensive product portfolio including Cryo.s tubes, racks as well as electronic equipment for tube closure and sample tracking. With the brand Cryo.s Greiner Bio-One combines more than 30 years of experience in the manufacturing of cryogenic tubes with latest technology and innovation.

CRYOTECHNICS

/	Cryo.s Cryo Tubes	182
	Cryo.s - 1 ml	183
	Cryo.s - 2 ml	183
	Cryo.s - 4 ml	185
	Cryo.s - 5 ml	185
	Support Rack	186
	Cryo storage Boxes	186
/	Cryo.s with Barcode	188
	Cryo.s with Barcode – Preproduced unique codes.....	189
	Cryo.s with Barcode – Customised Codes	189
	48-way Datamatrix Cryo Rack	190
	81-way Datamatrix Cryo Rack	190
/	Cryo.s Biobanking Tubes	191
	Cryo.s Biobanking Tubes sterile 300 µl	192
	Cryo.s Biobanking Tubes non-sterile 300 µl...	193
	Cryo.s Biobanking Tubes sterile 600 µl	194
	Cryo.s Biobanking Tubes non-sterile 600 µl...	195
	Cryo.s Biobanking Tubes sterile 1000 µl.....	196
	Cryo.s Biobanking Tubes non-sterile 1000 µl ..	197
	Screw Caps Biobanking Tubes	198
	96-way Datamatrix Cryo Rack	198
	Cryo.s Decapper	199



- / For a freezing and thawing protocol as well as safety advisory for Cryo.s please refer to the technical appendix.

CRYO.S CRYO TUBES

Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

An outstanding feature of all Cryo.s is the utilisation of a USP class VI certified, medical grade polypropylene for tube manu-

facturing. This material does not release significant amounts of leachables, hence stored samples stay clean and free of contamination. Sterile product versions are sterilised applying an ISO 11137 validated irradiation procedure yielding an SAL level of 10^{-6} .

The portfolio of Cryo.s tubes includes cryogenic tubes with different volumes, different base forms as well as several cap colours. Cryo.s offer several features for the identification and labelling of individual samples, to include coloured screw caps, a white, scratch-resistant writing area and barcoding options.

- / CE-marked according to IVDR
- / USP class VI certified medical grade polypropylene
- / Sterilised applying an ISO 11137 validated irradiation procedure yielding an SAL level of 10^{-6}
- / Suited for storage in liquid nitrogen vapour phase (-196°C)



Cryo.s - 1 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene
- / ACT® Ecolabel certification



Height: 42 mm, Ø: 12.5 mm, Starfoot: yes, Raw material: PP, Bottom shape: conical, Working volume: ≤1.2 ml, Cap design: screw cap, Sterile: +

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Qty. inner / outer
123261	no	no	no	○ natural	internal	+	100 / 500
123263	no	yes	yes	○ natural	internal	+	100 / 500
123277	no	yes	yes	● green	internal	+	100 / 500
123278	no	yes	yes	● yellow	internal	+	100 / 500
123279	no	yes	yes	● blue	internal	+	100 / 500
123280	no	yes	yes	● red	internal	+	100 / 500

/ with 150 inserts per case

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



Cryo.s - 2 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene
- / Selected product versions certified with ACT® Ecolabel (*)



Raw material: PP, Bottom shape: round, Cap design: screw cap, Sterile: +

Item no.	Height	Ø	Starfoot	Barcode	Graduation	Writing field	Cap colour	Working volume	Thread type	Triple-packed	Sterile	Qty. inner / outer
121261*	48 mm	12.5 mm	no	no	-	no	○ natural	≤2 ml	internal		+	100 / 500

Item no.	Height	Ø	Starfoot	Barcode	Graduation	Writing field	Cap colour	Working volume	Thread type	Triple-packed	Sterile	Qty. inner / outer
121263*	48 mm	12.5 mm	no	no	yes	yes	○ natural	≤2 ml	internal		+	100 / 500
121277*	48 mm	12.5 mm	no	no	yes	yes	● green	≤2 ml	internal		+	100 / 500
121278*	48 mm	12.5 mm	no	no	yes	yes	● yellow	≤2 ml	internal		+	100 / 500
121279*	48 mm	12.5 mm	no	no	yes	yes	● blue	≤2 ml	internal		+	100 / 500
121280*	48 mm	12.5 mm	no	no	yes	yes	● red	≤2 ml	internal		+	100 / 500
122261*	48 mm	12.5 mm	yes	no	-	no	○ natural	≤2 ml	internal		+	100 / 500
122263*	48 mm	12.5 mm	yes	no	yes	yes	○ natural	≤2 ml	internal		+	100 / 500
122263-TRI	48 mm	12.5 mm	yes	no	yes	yes	○ natural	≤2 ml	internal	yes	+	10 / 200
122277*	48 mm	12.5 mm	yes	no	yes	yes	● green	≤2 ml	internal		+	100 / 500
122278*	48 mm	12.5 mm	yes	no	yes	yes	● yellow	≤2 ml	internal		+	100 / 500
122279*	48 mm	12.5 mm	yes	no	yes	yes	● blue	≤2 ml	internal		+	100 / 500
122280*	48 mm	12.5 mm	yes	no	yes	yes	● red	≤2 ml	internal		+	100 / 500
126261*	47 mm	12.4 mm	yes	no	-	no	○ natural	≤2.2 ml	external		+	100 / 500
126263*	47 mm	12.4 mm	yes	no	yes	yes	○ natural	≤2.2 ml	external		+	100 / 500
126263-TRI	47 mm	12.4 mm	yes	no	yes	yes	○ natural	≤2.2 ml	external	yes	+	10 / 200
126277*	47 mm	12.4 mm	yes	no	yes	yes	● green	≤2.2 ml	external		+	100 / 500
126278*	47 mm	12.4 mm	yes	no	yes	yes	● yellow	≤2.2 ml	external		+	100 / 500
126279*	47 mm	12.4 mm	yes	no	yes	yes	● blue	≤2.2 ml	external		+	100 / 500
126280*	47 mm	12.4 mm	yes	no	yes	yes	● red	≤2.2 ml	external		+	100 / 500

/ Item numbers marked with * are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc.
More details at <https://act.mygreenlab.org>

/ with 150 inserts per case



Cryo.s - 4 ml

- / High thermal resistance
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene
- / ACT® Ecolabel certification

CE 0197	IVD	STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	-----	---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------

Description: Cryo.s 4 ml, Height: 83 mm, Ø: 12.4 mm, Starfoot: yes, Raw material: PP, Bottom shape: round, Working volume: ≤4 ml, Cap design: screw cap, Sterile: +

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Qty. inner / outer
127261	no	no	no	○ natural	external	+	50 / 300
127263	no	yes	yes	○ natural	external	+	50 / 300
127277	no	yes	yes	● green	external	+	50 / 300
127278	no	yes	yes	● yellow	external	+	50 / 300
127279	no	yes	yes	● blue	external	+	50 / 300
127280	no	yes	yes	● red	external	+	50 / 300

/ with 100 inserts per case

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



Cryo.s - 5 ml

- / High thermal resistance
- / Cryo.s with internal thread have a screw cap with silicone gasket
- / Cap inserts item no. 304134 (50 pieces per bag)
- / USP class VI certified medical grade polypropylene
- / ACT® Ecolabel certification

CE 0197	IVD	STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	-----	---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------

Description: Cryo.s 5 ml, Height: 86 mm, Ø: 12.5 mm, Starfoot: no, Raw material: PP, Bottom shape: round, Working volume: ≤4.5 ml, Cap design: screw cap, Sterile: +

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Qty. inner / outer
124261	no	-	no	○ natural	internal	+	50 / 300

Item no.	Barcode	Graduation	Writing field	Cap colour	Thread type	Sterile	Qty. inner / outer
124263	no	yes	yes	○ natural	internal	+	50 / 300
124273	no	yes	yes	● red	internal	+	50 / 300
124275	no	yes	yes	● green	internal	+	50 / 300
124276	no	yes	yes	● yellow	internal	+	50 / 300
124274	no	yes	yes	● blue	internal	+	50 / 300

/ with 100 inserts per case

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>




Support Rack

- / Suitable for Cryo.s with starfoot base (Item No. 122XXX, 123XXX, 126XXX, 127XXX)
- / Improved handling since the tubes can be opened with one hand
- / Rubber base to prevent slipping
- / Offers space for up to 40 Cryo.s

Description: support rack for one-hand operation, Height: 22 mm, Length: 200 mm, Width: 100 mm, Raw material: PC

Item no.	Rack colour	Qty. inner / outer
802501	● blue	1 / 50

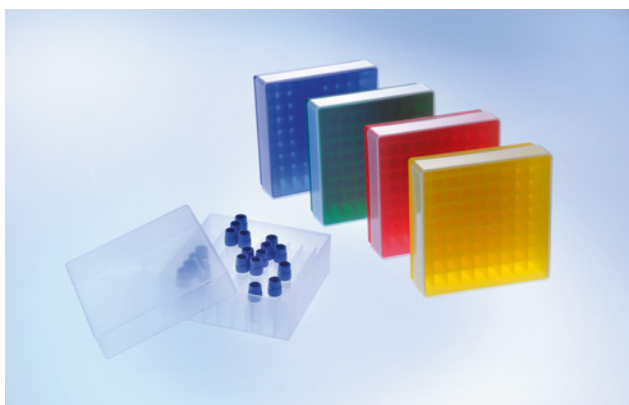

Cryo Storage Boxes

The box material polypropylene is very temperature and shock resistant, thus allowing storage temperatures as low as -90 °C.

- / Holds 81 Cryo.s sample tubes
- / Transparent lid for optimum visibility of box content

Description: Cryo storage box, Height: 51 mm, Length: 126.5 mm, Width: 126.5 mm, Raw material: PP

Item no.	Rack colour	Qty. inner / outer
802202	○ natural	5 / 20
802203	● red	5 / 20
802204	● blue	5 / 20
802206	● yellow	5 / 20

Item no.	Rack colour	Qty. inner / outer
802225	● green	5 / 20

/ The indicated height refers to the Cryo Storage Box filled with 2 ml Cryo.s and covered with a lid.



- / Selected Cryo.s are certified with ACT® Ecolabel. Learn more:



CRYO.S WITH BARCODE

Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the

gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

With Cryo.s with Datamatrix and the accessory racks in 48-way and 81-way formats, Greiner Bio-One expands its portfolio by optimum solutions for semi-automated and automated sample handling and storage.

- / Datamatrix code ECC200 on the tube bottom with reliable Reed-Solomon error correction
- / Linear barcode type 128 with plain text on the tube side
- / Application of barcodes and datamatrix codes by means of innovative laser technology
- / 100 % controlled code readability
- / Suited for storage in liquid nitrogen vapour phase (-196 °C)
- / Coded tubes are suitable for airfreight (IATA conformity)



Cryo.s with Barcode

– Preproduced unique codes

- / Pre-produced unique datamatrix code ECC 200 with 14 x 14 dots on tube bottom, additional linear barcode type 128 with human readable text on tube side
- / Highest code resistance against chemicals, mechanical stress and extremely low storage temperatures (down to -196 °C)
- / ACT® Ecolabel certification



Starfoot: yes, Barcode type: linear barcode and datamatrix code, Raw material: PP, Cap design: screw cap, Sterile: +

Item no.	Height	Ø	Barcode	Bottom shape	Cap colour	Working volume	Thread type	Sterile	Qty. inner / outer
123263-2DG	42 mm	12.5 mm	yes	conical	○ natural	≤1.2 ml	internal	+	100 / 500
122263-2DG	48 mm	12.5 mm	yes	round	○ natural	≤2 ml	internal	+	100 / 500
126263-2DG	47 mm	12.4 mm	yes	round	○ natural	≤2.2 ml	external	+	100 / 500
127263-2DG	83 mm	12.4 mm	yes	round	○ natural	≤4 ml	external	+	50 / 300

/ Other screw cap colours on request.

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



Cryo.s with Barcode

– Customised Codes

- / Customised datamatrix code ECC 200 with 14 x 14 dots on tube bottom, additional customised linear barcode type 128 with human readable text on tube side
- / Highest code resistance against chemicals, mechanical stress and extremely low storage temperatures (down to -196 °C)
- / ACT® Ecolabel certification



Starfoot: yes, Barcode type: linear barcode and datamatrix code, Raw material: PP, Cap design: screw cap, Sterile: +

Item no.	Height	Ø	Barcode	Bottom shape	Cap colour	Working volume	Nominal volume	Thread type	Sterile	Qty. inner / outer
123263-2D3	42 mm	12.5 mm	yes	conical	○ natural	≤1.2 ml	1 ml	internal	+	100 / 500
122263-2D3	48 mm	12.5 mm	yes	round	○ natural	≤2 ml	2 ml	internal	+	100 / 500
126263-2D1	47 mm	12.4 mm	yes	round	○ natural	≤2.2 ml	2 ml	external	+	100 / 500

Item no.	Height	Ø	Barcode	Bottom shape	Cap colour	Working volume	Nominal volume	Thread type	Sterile	Qty. inner / outer
127263-2D1	83 mm	12.4 mm	yes	round	○ natural	≤4 ml	4 ml	external	+	50 / 300

- / Custom barcode orders have a minimum order quantity of 10,000 tubes per order and request.
- / Other screw cap colours on request.
- / Barcodes to be specified in order form F071004.
- / All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>



48-way Datamatrix Cryo Rack

- / SLAS-standard format
- / Automation compatible
- / Suitable for storage in vapour phase above liquid nitrogen
- / Barcoded versions on request

Length: 85.5 mm, Width: 127.8 mm, Raw material: PP, Lid type: with lid

Item no.	Description	Height	Barcode	Rack colour	Qty. inner / outer
803202	low rack	52.5 mm	no	● black	- / 20
803270	high rack	88.5 mm	no	● black	- / 15



81-way Datamatrix Cryo Rack

- / Shock-proof and temperature-resistant polycarbonate
- / Stacking feature and rotation stoppers
- / 133 x 133 mm footprint for storage in classic freezer setups
- / Suitable for storage in vapour phase above liquid nitrogen
- / Barcoded versions on request

Description: low rack, Height: 52 mm, Length: 132.4 mm, Width: 132.4 mm, Raw material: PC, Lid type: with lid, Suitable for Cryos: 1/2 ml

Item no.	Rack colour	Qty. inner / outer
802576	● black	1 / 10



- / Forum No. 21: Sample Storage Tubes as Quality-Critical Components in Biobanking (F073072)
- / You can find more information about this product in our web-shop.



CRYO.S BIOBANKING TUBES

With Cryo.s biobanking tubes, Greiner Bio-One offers an ideal solution for the efficient storage of biological samples in large-scale biorepositories.

Cryo.s biobanking tubes are available with working volumes of 235 µl, 580 µl and 975 µl and offered in highly automation-friendly 96-way racks. The innovative design of tubes and racks allows for a very space-efficient storage with up to 30 % better utilisation of storage space in freezers or liquid nitrogen tanks.

Cryo.s with datamatrix and linear barcode and Cryo.s biobanking tubes are for storage of tissue, cells, fungi, bacteria, spores, cellular extracts or body fluids at ultra-low temperature for research and development purposes as well as in-vitro diagnostics. They must not be stored in the liquid phase of liquid nitrogen, but only in the gas phase above. Cryo.s tubes are not intended for any application in the context of reproductive medicine.

- / Application of barcodes and datamatrix codes by means of innovative laser technology
- / 100 % controlled code readability
- / Coded tubes are suitable for airfreight (IATA conformity)
- / Suited for storage at -20 °C, -80 °C and in vapour phase above liquid nitrogen



Cryo.s Biobanking Tubes sterile

300 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)
- / Sterile

CE 0197	IVD	STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	-----	---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Height: 18.7 mm, Total rack height: 21.6 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤235 µl, Cap design: screw cap, Sterile: +

Item no.	Barcode	Cap colour	Thread type	Type of packaging	Sterile	Qty. inner / outer
976561	yes	○ natural	internal	rack	+	480 / 960
976566	yes	● green	internal	rack	+	480 / 960
976565	yes	● yellow	internal	rack	+	480 / 960
976564	yes	● blue	internal	rack	+	480 / 960
976563	yes	● red	internal	rack	+	480 / 960
976568	yes	● pink	internal	rack	+	480 / 960
976569	yes	● brown	internal	rack	+	480 / 960
976567	yes	● black	internal	rack	+	480 / 960
131263	yes	○ natural	internal	bulk	+	480 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes non-sterile
300 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)



Height: 18.7 mm, Total rack height: 21.6 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤235 µl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Qty. inner / outer
976570	yes		without screw cap	internal	rack	480 / 960
976580	yes	○ natural	screw cap	internal	rack	480 / 960
976586	yes	● green	screw cap	internal	rack	480 / 960
976585	yes	● yellow	screw cap	internal	rack	480 / 960
976584	yes	● blue	screw cap	internal	rack	480 / 960
976583	yes	● red	screw cap	internal	rack	480 / 960
976588	yes	● pink	screw cap	internal	rack	480 / 960
976589	yes	● brown	screw cap	internal	rack	480 / 960
976587	yes	● black	screw cap	internal	rack	480 / 960
131202	yes	○ natural	screw cap	internal	bulk	480 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes sterile

600 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)
- / Sterile

CE 0197	IVD	STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	-----	---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Height: 33.3 mm, Total rack height: 36.2 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤580 µl, Cap design: screw cap, Sterile: +

Item no.	Barcode	Cap colour	Thread type	Type of packaging	Sterile	Qty. inner / outer
977561	yes	○ natural	internal	rack	+	192 / 960
977566	yes	● green	internal	rack	+	192 / 960
977565	yes	● yellow	internal	rack	+	192 / 960
977564	yes	● blue	internal	rack	+	192 / 960
977563	yes	● red	internal	rack	+	192 / 960
977568	yes	● pink	internal	rack	+	192 / 960
977569	yes	● brown	internal	rack	+	192 / 960
977567	yes	● black	internal	rack	+	192 / 960
132263	yes	○ natural	internal	bulk	+	192 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes non-sterile

600 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)



Height: 33.3 mm, Total rack height: 36.2 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤580 µl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Qty. inner / outer
977570	yes		without screw cap	internal	rack	192 / 960
977580	yes	○ natural	screw cap	internal	rack	192 / 960
977586	yes	● green	screw cap	internal	rack	192 / 960
977585	yes	● yellow	screw cap	internal	rack	192 / 960
977584	yes	● blue	screw cap	internal	rack	192 / 960
977583	yes	● red	screw cap	internal	rack	192 / 960
977588	yes	● pink	screw cap	internal	rack	192 / 960
977589	yes	● brown	screw cap	internal	rack	192 / 960
977587	yes	● black	screw cap	internal	rack	192 / 960
132202	yes	○ natural	screw cap	internal	bulk	192 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes sterile

1000 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)
- / USP class VI certified medical grade polypropylene
- / Sterile

CE 0197	IVD	STERILE	FREE OF detectable DNase	FREE OF detectable human DNA
FREE OF detectable RNase	non-cytotoxic	non-pyrogenic		

Height: 50.8 mm, Total rack height: 53.9 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤975 µl, Cap design: screw cap, Sterile: +

Item no.	Barcode	Cap colour	Thread type	Type of packaging	Sterile	Qty. inner / outer
978561	yes	○ natural	internal	rack	+	192 / 960
978566	yes	● green	internal	rack	+	192 / 960
978565	yes	● yellow	internal	rack	+	192 / 960
978564	yes	● blue	internal	rack	+	192 / 960
978563	yes	● red	internal	rack	+	192 / 960
978568	yes	● pink	internal	rack	+	192 / 960
978569	yes	● brown	internal	rack	+	192 / 960
978567	yes	● black	internal	rack	+	192 / 960
133263	yes	○ natural	internal	bulk	+	192 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Cryo.s Biobanking Tubes non-sterile
1000 µl

- / Pre-produced unique datamatrix code on tube, datamatrix code and linear barcode on rack
- / Height-reduced screw cap conserving up to 30 % freezer space
- / Customised code sequences on request (order form F071003)
- / USP class VI certified medical grade polypropylene



Height: 50.8 mm, Total rack height: 53.9 mm, Ø: 8.8 mm, Barcode type: datamatrix code, Raw material: PP, Working volume: ≤975 µl

Item no.	Barcode	Cap colour	Cap design	Thread type	Type of packaging	Qty. inner / outer
978570	yes		without screw cap	internal	rack	192 / 960
978580	yes	○ natural	screw cap	internal	rack	192 / 960
978586	yes	● green	screw cap	internal	rack	192 / 960
978585	yes	● yellow	screw cap	internal	rack	192 / 960
978584	yes	● blue	screw cap	internal	rack	192 / 960
978583	yes	● red	screw cap	internal	rack	192 / 960
978588	yes	● pink	screw cap	internal	rack	192 / 960
978589	yes	● brown	screw cap	internal	rack	192 / 960
978587	yes	● black	screw cap	internal	rack	192 / 960
133202	yes	○ natural	screw cap	internal	bulk	192 / 960

- / Follow the instructions of use provided within each box.
- / USP class VI certified medical grade polypropylene



Screw Caps
Biobanking Tubes

- / 96 screw caps in cap carrier
- / Non-sterile

CE	IVD	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
----	-----	--------------------------	------------------------------	--------------------------	---------------	---------------

Description: 96 screw caps in cap carrier, Raw material: PP, Cap design: screw cap

Item no.	Barcode	Cap colour	Qty. inner / outer
385270	no	○ natural	960 / 960
385276	no	● green	960 / 960
385275	no	● yellow	960 / 960
385274	no	● blue	960 / 960
385273	no	● red	960 / 960
385278	no	● pink	960 / 960
385279	no	● brown	960 / 960
385277	no	● black	960 / 960

96-way Datamatrix Cryo Rack



- / Pre-produced unique datamatrix code ECC 200 and linear barcode type 128
- / Made from polycarbonate

CE	IVD	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
----	-----	--------------------------	------------------------------	--------------------------	---------------	---------------

Barcode type: linear barcode and datamatrix code, Raw material: PC

Item no.	Height	Barcode	Rack colour	Qty. inner / outer
976501	19.1 mm	yes	● black	5 / 10
977501	33.7 mm	yes	● black	2 / 10

Item no.	Height	Barcode	Rack colour	Qty. inner / outer
978501	51.4 mm	yes	● black	2 / 10



Cryo.s Decapper

The Cryo.s Decapper enables the precise simultaneous de- and recapping of cryotubes.

- / Individual motors for precise application of defined torque values
- / 6 channel version for 1, 2 and 4 ml Cryo.s in 48 way cryo racks
- / 8 channel version for Cryo.s Biobanking Tubes
- / Docking station

Content kit: 1 Decapper, Power supply unit, Docking station with charging function, Quick start guide

Item no.	Description	Plug	Qty. inner / outer
852076-EU	6-channel decapper	Europe	- /1
852076-UK	6-channel decapper	UK	- /1
852076-NA	6-channel decapper	US	- /1
852076-CN	6-channel decapper	China	- /1
852076-AU	6-channel decapper	Australia / NZ	- /1
852076-JP	6-channel decapper	Japan	- /1
852078-EU	8-channel decapper	Europe	- /1
852078-NA	8-channel decapper	US	- /1
852078-UK	8-channel decapper	UK	- /1
852078-CN	8-channel decapper	China	- /1
852078-AU	8-channel decapper	Australia / NZ	- /1
852078-JP	8-channel decapper	Japan	- /1



For the handling and analysis of chemical and biological samples with small or medium volumes, Greiner Bio-One offers a wide range of different analyser cups and reaction vessels with corresponding closures and cuvettes.

REACTION TUBES / ANALYSER CUPS

/	Reaction Tubes Analyser Cups	202
	Reaction Tubes	203
	Reaction Tubes 5 ml	203
	Closures for Reaction Tubes	204
	Analyser Cups	204
/	Semi-micro / Macro Cuvette	206
	Semi-micro / Macro Cuvette	207



/ For the max. relative centrifugal force (RCF), please refer to the technical appendix.

REACTION TUBES / ANALYSER CUPS

Reaction tubes from Greiner Bio-One are available in different versions. They have a high chemical and temperature resistance and are available from 1.5 ml to 5 ml for small and medium volumes. The reaction tubes are made of high-quality polypropylene and available with screw cap or snap cap, depending on the version. The lids of the corresponding vessels are flat and frosted to ensure easy labelling.

For working with liquid volumes from 1 to 5 ml, Greiner Bio-One offers 5 ml reaction tubes with screw cap or snap cap. The tubes are also made of highly

transparent polypropylene and allow contamination-free access when pipetting medium sample volumes. They also have a frosted, writable surface. This ensures reliable sample identification in everyday laboratory work.

Analyser cups are available with conical or flat bottom design and with or without cap. All analyser cups are manufactured of highly transparent polystyrene and are available with volumes of 1.7 ml and 25 ml. We offer analyser cups that are suitable for the analytical systems of Hitachi or Coulter/Hycel.

- / Reaction tubes for volumes from 1.5 ml up to 5 ml
- / With screw cap or snap cap
- / With graduation

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
pyrogenic

Reaction Tubes

- / High chemical resistance and temperature tolerance
- / Flat, frosted lids and caps for easy labelling

Description: reaction tube, Graduation: yes

Item no.	Support skirt	Product colour	Volume range	Caps attached	Cap design	Sterile	Qty. inner / outer
616201		○ natural	≤1.5 ml	yes	snap cap		500 / 4,000
616261		○ natural	≤1.5 ml	yes	snap cap	+	500 / 4,000
623201		○ natural	≤2 ml	yes	snap cap		500 / 4,000
716201	no	○ natural	≤1.5 ml		screw cap		500 / 5,000
717201	yes	○ natural	≤1.5 ml		screw cap		500 / 5,000
717261	yes	○ natural	≤1.5 ml		screw cap	+	500 / 3,000
722201	yes	○ natural	≤2 ml		screw cap		500 / 5,000
722261	yes	○ natural	≤2 ml		screw cap	+	- / 3,000

/ 0.5 ml PCR reaction tubes can be found in the Molecular Biology section (Item No. 682201)

FREE OF
detectable
DNaseFREE OF
detectable
human DNAFREE OF
detectable
RNase
 non-
pyrogenic

Reaction Tubes

5 ml


5 ml reaction tubes with either a snap cap for fast opening and closing or a screw cap for secure sample storage.

- / Outer diameter and shape similar to 15 ml tube
- / Resealable packaging

Description: reaction tube, Support skirt: no, Graduation: yes, Raw material: PP

Item no.	Product colour	Cap colour	Caps attached	Cap design	Sterile	Qty. inner / outer
725201	○ natural	● blue		screw cap		100 / 500

Item no.	Product colour	Cap colour	Caps attached	Cap design	Sterile	Qty. inner / outer
725261	○ natural	● blue		screw cap	+	50 / 500
622201	○ natural		yes	snap cap		100 / 1,000
622261	○ natural		yes	snap cap	+	100 / 1,000

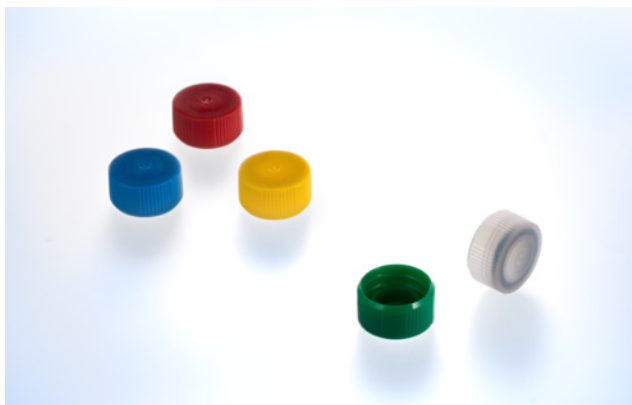

Closures for Reaction Tubes

- / Suitable for item no. 716XXX, 717XXX and 722XXX
- / Available in natural, red, blue, green and yellow
- / With gasket

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

non-
pyrogenic

Description: screw cap, Feature: with sealing ring, Suitable for tubes Ø: 12 mm, Cap design: screw cap

Item no.	Product colour	Qty. inner / outer
366380	○ natural	500 / 5,000
366383	● red	500 / 5,000
366384	● blue	500 / 5,000
366385	● green	500 / 5,000
366386	● yellow	500 / 5,000



Analyser Cups

- / For handling and analysis of chemical and biological samples
- / Available for different analytical systems

Description: analyser cup, Raw material: PS, Caps attached: no

Item no.	Feature	Support skirt	Bottom shape	Volume range	Lid type	Qty. inner / outer
668102	Coulter / Hycel	no	flat	≤25 ml	with lid	250 / 1,250

Item no.	Feature	Support skirt	Bottom shape	Volume range	Lid type	Qty. inner / outer
729101	Hitachi	yes	conical	≤1.7 ml	without lid	250 / 5,000



/ For further information, please refer to our website: www.gbo.com



SEMI-MICRO / MACRO CUVETTE

Cuvettes are used for UV/VIS spectroscopy. Greiner Bio-One offers semi-micro cuvettes with a total volume of 1.6 ml and macro cuvettes with a total volume of 4 ml. Both versions are made of crystal-clear polystyrene and are particularly suitable for enzymatic determinations, as the

very thin wall thickness enables fast and even temperature control. Our cuvettes are characterised by their low light scattering combined with a high transmission rate. Semi-micro and macro cuvettes can be used in a wavelength range from 340 nm to 900 nm.

- / Available as semi-micro or macro cuvettes
- / Very thin walls for fast and even temperature control
- / Low light scattering and high transmission rate
- / Suitable for a wavelength range from 340-900 nm



Semi-micro / Macro Cuvette

- / Ideal for enzymatic determinations
- / Manufactured from crystal clear polystyrene
- / Minimal light scatter but high transmission rate
- / Applicable wavelength from 340 to 900 nm

Height: 45 mm, Length: 12.5 mm, Width: 12.5 mm, Pathlength: 10 mm, Raw material: PS


Item no.	Description	Working volume	Total volume	Qty. inner / outer
613101	semi-micro cuvette	≥0.95 ml	1.6 ml	100 / 1,000
614101	macro cuvette	≥2.5 ml	4 ml	100 / 1,000



PCR (Polymerase Chain Reaction) is an important molecular biological method to determine processes in living organisms at the molecular level or genetic changes in the genetic material. It is used for the detection of diseases and parentage relationships as well as in forensics.

MOLECULAR BIOLOGY

/	Sapphire PCR Tubes.....	210
	PCR Tubes	211
	PCR 8-Tube Strips	212
	PCR 8-Cap Strips	213
/	Sapphire PCR Microplates.....	214
	96 Well PCR Microplates	215
	384 Well PCR Microplates	215

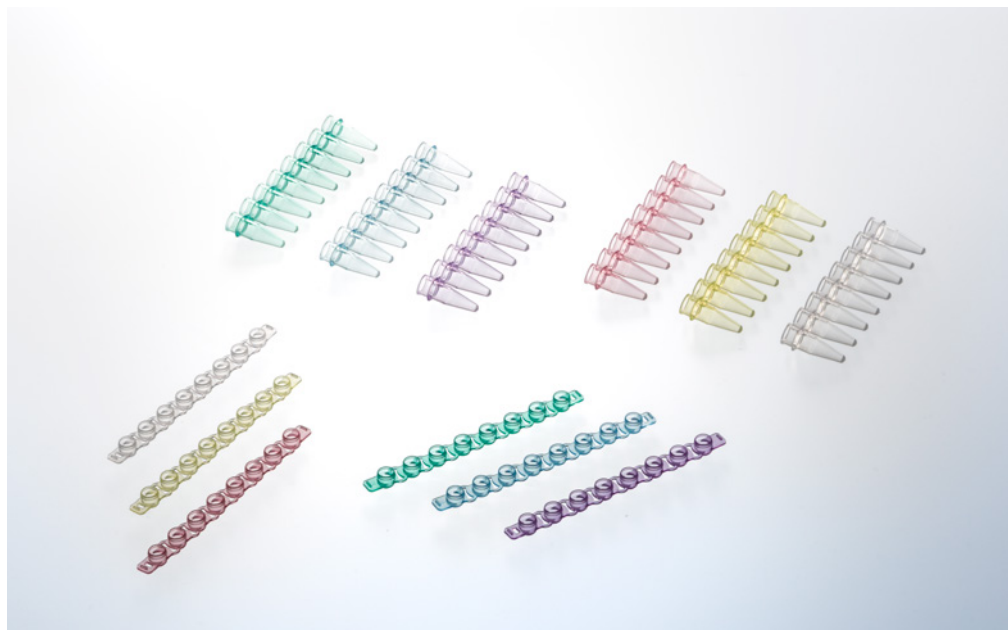

/ For standard reaction tubes please refer to chapter reaction tubes / analyser cups.

SAPPHIRE PCR TUBES

The Sapphire PCR tube portfolio comprises a full range of PCR reaction tubes (0.2 ml and 0.5 ml) as well as PCR 8-tube strips and the matching 8-cap strips.

The thin wall construction optimises the heat transfer from the block to the reaction solution. PCR tubes with flat caps are suitable for real-time / quantitative PCR and have frosted caps for easy labelling.

The classical Sapphire PCR 8-tube strips are also available with individually attached caps.

The angular attached domed or flat caps can be closed and opened individually minimising the risk of cross contamination. Moreover, individual closing and opening helps to avoid pipetting errors and facilitates sample processing.

PCR 8-tube strips are also available as **low profile tubes** with a well volume of 0.1 ml ideal for FAST, standard and real-time PCR supporting the reduction of cycle time without compromises in accuracy and efficiency.

- / Manufactured out of medical-grade polypropylene without interfering additives
- / Thin-walled for efficient heat transfer during amplification
- / Comprehensive thermal cycler compatibility



PCR Tubes

Tubes with flat caps are suitable for realtime/ quantitative PCR and have frosted caps for easy labelling.

/ Packaging labelled with product symbols



Description: PCR tube, Raw material: PP

Item no.	Product colour	Volume	Volume range	Caps attached	Cap design	Lid type	Qty. inner / outer
671201	○ natural	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671273	● red	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671274	● blue	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671275	● green	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671276	● yellow	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671277	● violet	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
671281	assorted	0.2 ml	≤0.2 ml	yes	individually attached	domed	1,000 / 10,000
683201	○ natural	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683273	● red	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683274	● blue	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683275	● green	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683276	● yellow	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683277	● violet	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
683271	assorted	0.2 ml	≤0.2 ml	yes	individually attached	flat	1,000 / 10,000
682201	○ natural	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682273	● red	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682274	● blue	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000

Item no.	Product colour	Volume	Volume range	Caps attached	Cap design	Lid type	Qty. inner / outer
682275	● green	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682276	● yellow	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682277	● violet	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
682281	assorted	0.5 ml	≤0.5 ml	yes	individually attached	flat	1,000 / 10,000
684201	○ natural	0.2 ml	≤0.2 ml	no		without lid	1,000 / 10,000

/ Assorted colour contains the following colours: red, blue, green, yellow, violet



PCR 8-Tube Strips

- / Thin wall for optimal heat transfer
- / Low evaporation rate in PCR
- / Also available with individually attached caps or as low-profile version
- / Packaging labelled with product symbols

Description: PCR 8-Tube Strips, Raw material: PP

Item no.	Feature	Product colour	Volume range	Caps attached	Cap design	Lid type	Qty. inner / outer
673210		○ natural	≤0.2 ml	no	separate cap strip		125 / 1,250
673273		● red	≤0.2 ml	no	separate cap strip		125 / 1,250
673274		● blue	≤0.2 ml	no	separate cap strip		125 / 1,250
673275		● green	≤0.2 ml	no	separate cap strip		125 / 1,250
673276		● yellow	≤0.2 ml	no	separate cap strip		125 / 1,250
673277		● violet	≤0.2 ml	no	separate cap strip		125 / 1,250
673271		assorted	≤0.2 ml	no	separate cap strip		125 / 1,250
671221	low profile	○ natural	≤0.1 ml	yes	individually attached	flat	120 / 1,200
608281		○ natural	≤0.2 ml	yes	individually attached	flat	120 / 1,200
673281	low profile	○ natural	≤0.1 ml	no	separate cap strip	flat	125 / 1,250
673283	breakable	○ natural	≤0.2 ml	yes	individually attached	domed	120 / 1,200

/ Assorted colour contains the following colours: red, blue, green, yellow, violet



PCR 8-Cap Strips


- / Real-time PCR cap strips (Item No. 373250) have a flat lid and are made of highly transparent polypropylene
- / These are ideal for real-time PCR applications and are compatible with most real-time PCR devices

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	-----

Description: 8-cap strip, Raw material: PP

Item no.	Feature	Product colour	Qty. inner / outer
373270		○ natural	125 / 1,250
373273		● red	125 / 1,250
373274		● blue	125 / 1,250
373275		● green	125 / 1,250
373276		● yellow	125 / 1,250
373277		● violet	125 / 1,250
373281		assorted	125 / 1,250
373250	for RT PCR	○ natural	125 / 1,250

/ Assorted colour contains the following colours: red, blue, green, yellow, violet

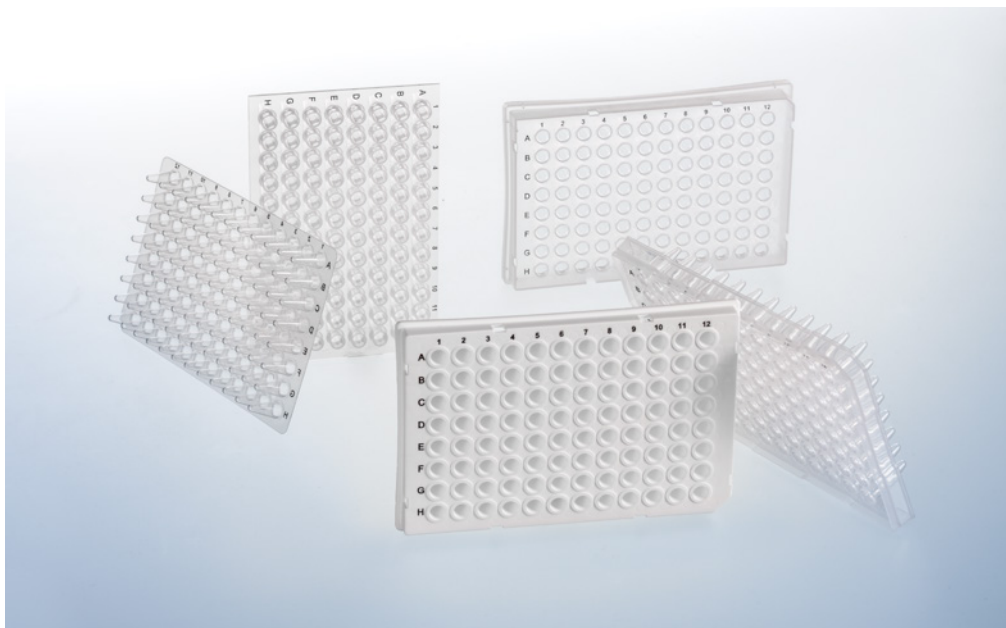

/ For details regarding plate geometry as well as compability with thermal cyclers, please refer to the technical appendix.


SAPPHIRE PCR MICROPLATES

The use of the 96 well format allows the scale up of basic PCR work, while the 384 well format is ideal for high-throughput screening projects.

All microplates are made of thin-walled polypropylene. This opti-

mises the heat transfer from the thermoblock to the reaction solution. Our heat-resistant sealers AMPLIseal, VIEWseal and SILVERseal are ideal for sealing the microplates during PCR, and the 96 well microplate may also be sealed with 8-cap strips.

- / Available in 96 well and 384 well format
- / Ultra-thin polypropylene for optimal heat transfer
- / Sealable with sealers SILVERseal, VIEWseal and AMPLIseal

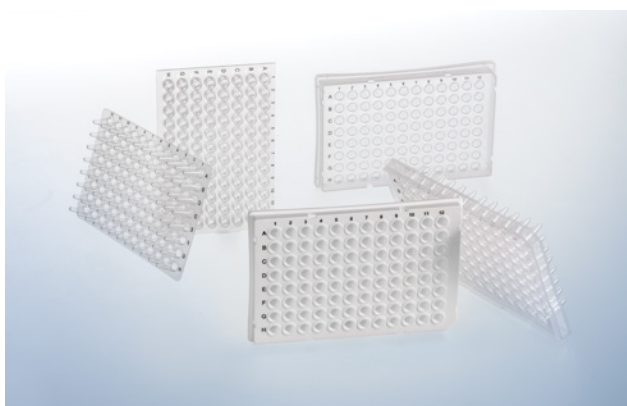


FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	-----

96 Well PCR Microplates

- / Without skirt, half-skirt and full-skirt
- / Alphanumeric printing for rapid sample identification
- / Sealable with sealers SILVERseal , VIEWseal and AMPLIseal or with compatible 8-cap strips
- / Low profile microplate for reduced sample volume and cycle time

Well format: 96, Raw material: PP

Item no.	Feature	Skirt	Product colour	Colour alphanumeric code	Well volume, max	Qty. inner / outer
652201		without skirt	○ natural	● black	0.2 ml	10 / 100
652250	flat, universal	without skirt	○ natural	● black	0.2 ml	10 / 100
652210	low profile	without skirt	○ natural	● black	0.1 ml	20 / 100
669285	suitable for LightCycler®	half-skirt	○ white	● black	0.2 ml	10 / 100
652290	suitable for ABI	half-skirt	○ natural	● black	0.2 ml	10 / 100
652260	ABI design	half-skirt	○ natural	● black	0.2 ml	10 / 100
652270		yes	○ natural	● black	0.2 ml	10 / 100

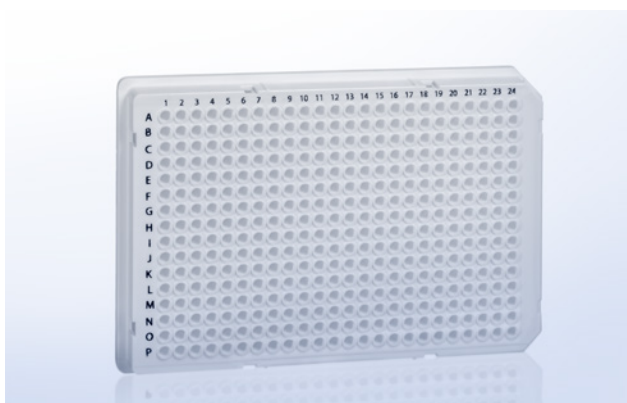

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	-----

384 Well PCR Microplates

- / Thin wall for optimal heat transfer
- / Alphanumeric well coding
- / Sealable with sealers SILVERseal , VIEWseal and AMPLIseal

Feature: suitable for LightCycler®, Well format: 384, Skirt: yes, Raw material: PP, Total volume (well): 25 µl

Item no.	Product colour	Colour alphanumeric code	Qty. inner / outer
785285	○ white	● black	50 / 100



To achieve consistent, reproducible results when handling liquids in the laboratory, reliable dispensers and pipette tips are essential. With high quality materials and pipetting solutions, maximum precision can be achieved while reducing the consumption of reagents and time. Comfortable and fatigue-free working is also crucial for everyday laboratory life.

LIQUID HANDLING

/	Sapphire Pipettes	218
	Single-Channel Pipettes	219
	Multi-Channel Pipettes	219
	Pipette Carrousel	220
/	Sapphire Pipette Tips	221
	Pipette Tips 10 µl	222
	Pipette Tips 10 µl extended.....	223
	Filter Tips 20 µl	224
	Filter Tips 100 µl	224
	Pipette Tips 200 µl.....	225
	Pipette Tips 300 µl.....	226
	Pipette Tips 1000 µl	227
	Pipette Tips 1250 µl.....	227
	Racks (unfilled)	228
	Macro Pipette Tips 1-5 ml.....	229
	Gel-Load Pipette Tips 1-200 µl.....	229
/	CELLSTAR® Serological Pipettes	230
	Serological Pipettes	231
	Serological Pipettes Triple-packed	232
	Aspirating Pipette	232
	MaxiPette	233
	Pasteur / Serum Pipettes	233



/ Watch our video:



SAPPHIRE PIPETTES

Sapphire pipettes are a variable volume range of air-displacement pipettes designed and built to deliver optimal performance, ergonomics, and robustness in day-to-day use.

The lightweight design coupled with a state-of-the-art mechanism reduces pipetting forces and makes our pipettes both comfortable and easy-to-use. The contoured body shape allows for a relaxed grip whatever size your hand, and the 3-position tip ejector button makes

pipetting comfortable for both left-handed and right-handed users. The colour-coded push button allows for easier identification of the relevant tips.

Sapphire pipettes provide an excellent performance within ISO 8655 which is considered the basis for determining the accuracy and precision of air-displacement pipettes. In combination with Sapphire pipette tips they show one of the lowest error tolerances on the market. All Sapphire pipettes are fully autoclavable.

- / Outstanding accuracy and precision
- / Ergonomic design and minimal pipetting forces for a new level of comfort
- / Lightweight, robust and fully autoclavable
- / Digital volume setting
- / Colour-coded for tip identification



Single-Channel Pipettes

- / Light and comfortable design
- / Low pipetting forces for ease of use
- / Digital volume setting
- / Colour-coded for tip identification
- / 5 ml & 10 ml pipettes without ejector lever
- / Fully autoclavable

Description: single-channel pipette

Item no.	Colour code	Volume range	Qty. inner / outer
89000002	● orange	0.2 - 2 µl	1 / 1
89000010	● red	1 - 10 µl	1 / 1
89000020	● light yellow	2 - 20 µl	1 / 1
89000100	● light orange	10 - 100 µl	1 / 1
89000200	● yellow	20 - 200 µl	1 / 1
89001000	● blue	100 - 1,000 µl	1 / 1
89000500	● violet	500 - 5,000 µl	1 / 1
89010000	● light blue	1,000 - 10,000 µl	1 / 1

/ Spare parts and other accessories available on request



Multi-Channel Pipettes

- / Light and comfortable design
- / Low pipetting forces for ease of use
- / Digital volume setting
- / Colour-coded for tip identification
- / Optimised for use with Greiner Bio-One tips
- / Fully autoclavable

Item no.	Description	Colour code	Volume range	Qty. inner / outer
89000810	8-channel pipette	● red	0.5 - 10 µl	1 / 1
89000820	8-channel pipette	● light yellow	2 - 20 µl	1 / 1
89008200	8-channel pipette	● yellow	20 - 200 µl	1 / 1
89008300	8-channel pipette	● green	20 - 300 µl	1 / 1
89001210	12-channel pipette	● red	0.5 - 10 µl	1 / 1
89001220	12-channel pipette	● light yellow	2 - 20 µl	1 / 1

Item no.	Description	Colour code	Volume range	Qty. inner / outer
89012200	12-channel pipette	● yellow	20 - 200 µl	1 / 1
89012300	12-channel pipette	● green	20 - 300 µl	1 / 1

/ Spare parts and other accessories available on request



Pipette Carrousel

- / Holds up to 7 pipettes
- / Robust and space-saving
- / Holds single and multi-channel pipettes

Item no.	Qty. inner / outer
89000099	1 / 1



/ Watch our video:



SAPPHIRE PIPETTE TIPS

- / Fully automated production and packaging
- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Coloured box inserts for easy volume identification
- / Low retention surface properties for high recovery rate and maximum precision

We understand that flexibility is critical in terms of modern liquid handling solutions. This is why the Sapphire tips are manufactured in eight different sizes in a volume range from 10 μ l to 1250 μ l, including an extended length 10 μ l tip for the recovery of small sample volumes.

All our tips are made of medical grade polypropylene and are

available as standard, low-retention, filter or low-retention filter tips.

Depending on the format, Sapphire pipette tips are packed in racks, bulk or refill units. Sapphire tips and pipettes when used together provide a fully optimised and harmonised liquid handling solution.



Pipette Tips

10 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	---------------	-----

Raw material: PP, Volume range: 0.2 - 10 µl, Nominal volume: 10 µl, Matching rack: 770310

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
771350	standard	○ natural	● red	bag		1,000 / 10,000
771351	standard	○ natural	● red	rack		960 / 5,760
771352	standard	○ natural	● red	rack	+	960 / 5,760
771354	standard	○ natural	● red	refill		960 / 4,800
771355	standard	○ natural	● red	refill	+	960 / 4,800
771361	low retention	○ natural	● red	rack		960 / 5,760
771362	low retention	○ natural	● red	rack	+	960 / 5,760
771364	low retention	○ natural	● red	refill		960 / 4,800
771365	low retention	○ natural	● red	refill	+	960 / 4,800
771353	filter	○ natural	● red	rack	+	960 / 5,760
771363	low retention filter	○ natural	● red	rack	+	960 / 5,760



Pipette Tips
10 µl extended

- / Extended 10 µl tip for better recovery of small sample volumes
- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	---------------	-----

Raw material: PP, Volume range: 0.5 - 10 µl, Nominal volume: 10 µl, Matching rack: 770310

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
772350	standard	○ natural	● red	bag		1,000 / 5,000
772351	standard	○ natural	● red	rack		960 / 5,760
772352	standard	○ natural	● red	rack	+	960 / 5,760
772354	standard	○ natural	● red	refill		960 / 4,800
772355	standard	○ natural	● red	refill	+	960 / 4,800
772361	low retention	○ natural	● red	rack		960 / 5,760
772362	low retention	○ natural	● red	rack	+	960 / 5,760
772364	low retention	○ natural	● red	refill		960 / 4,800
772365	low retention	○ natural	● red	refill	+	960 / 4,800
772353	filter	○ natural	● red	rack	+	960 / 5,760
772363	low retention filter	○ natural	● red	rack	+	960 / 5,760



Filter Tips

20 µl

- / Prevent contamination with liquids and aerosols during pipetting
- / Filters manufactured without additives
- / Easy-to-read graduation
- / Colour-coded for tip identification



Raw material: PP, Volume range: 2 - 20 µl, Nominal volume: 20 µl, Type of packaging: rack, Sterile: +

Item no.	Feature	Product colour	Box insert colour	Sterile	Qty. inner / outer
773353	filter	<input type="radio"/> natural	<input checked="" type="radio"/> orange	+	960 / 5,760
773363	low retention filter	<input type="radio"/> natural	<input checked="" type="radio"/> orange	+	960 / 5,760



Filter Tips

100 µl

- / Prevent contamination with liquids and aerosols during pipetting
- / Filters manufactured without additives
- / Easy-to-read graduation
- / Colour-coded for tip identification



Raw material: PP, Volume range: 5 - 100 µl, Nominal volume: 100 µl, Type of packaging: rack, Sterile: +

Item no.	Feature	Product colour	Box insert colour	Sterile	Qty. inner / outer
774353	filter	<input type="radio"/> natural	<input checked="" type="radio"/> dark yellow	+	960 / 5,760
774363	low retention filter	<input type="radio"/> natural	<input checked="" type="radio"/> dark yellow	+	960 / 5,760



Pipette Tips

200 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification

FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	PCR
--------------------------	------------------------------	--------------------------	---------------	---------------	-----

Raw material: PP, Volume range: 5 - 200 µl, Nominal volume: 200 µl, Matching rack: 770320

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
775350	standard	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	bag		1,000 / 15,000
775351	standard	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack		960 / 5,760
775352	standard	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack	+	960 / 5,760
775354	standard	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	refill		960 / 4,800
775355	standard	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	refill	+	960 / 4,800
775361	low retention	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack		960 / 5,760
775362	low retention	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack	+	960 / 5,760
775364	low retention	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	refill		960 / 4,800
775365	low retention	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	refill	+	960 / 4,800
775353	filter	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack	+	960 / 5,760
775363	low retention filter	<input type="radio"/> natural	<input checked="" type="radio"/> yellow	rack	+	960 / 5,760



Pipette Tips

300 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification

FREE OF
detectable
DNase

FREE OF
detectable
human DNA

FREE OF
detectable
RNase

 non-
cytotoxic

 non-
pyrogenic

 PCR

Raw material: PP, Volume range: 10 - 300 µl, Nominal volume: 300 µl, Matching rack: 770330

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
776350	standard	○ natural	● green	bag		1,000 / 10,000
776351	standard	○ natural	● green	rack		960 / 5,760
776352	standard	○ natural	● green	rack	+	960 / 5,760
776354	standard	○ natural	● green	refill		960 / 4,800
776355	standard	○ natural	● green	refill	+	960 / 4,800
776361	low retention	○ natural	● green	rack		960 / 5,760
776362	low retention	○ natural	● green	rack	+	960 / 5,760
776364	low retention	○ natural	● green	refill		960 / 4,800
776365	low retention	○ natural	● green	refill	+	960 / 4,800
776353	filter	○ natural	● green	rack	+	960 / 5,760
776363	low retention filter	○ natural	● green	rack	+	960 / 5,760



Pipette Tips

1000 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification



Raw material: PP, Volume range: 50 - 1,000 µl, Nominal volume: 1,000 µl, Matching rack: 770340

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
777350	standard	○ natural	● blue	bag		1,000 / 5,000
777351	standard	○ natural	● blue	rack		960 / 3,840
777352	standard	○ natural	● blue	rack	+	960 / 3,840
777354	standard	○ natural	● blue	refill		960 / 4,800
777355	standard	○ natural	● blue	refill	+	960 / 4,800
777361	low retention	○ natural	● blue	rack		960 / 3,840
777362	low retention	○ natural	● blue	rack	+	960 / 3,840
777364	low retention	○ natural	● blue	refill		960 / 4,800
777365	low retention	○ natural	● blue	refill	+	960 / 4,800



Pipette Tips

1250 µl

- / Universal fit for all common pipettes
- / Tips (without filter) and racks are autoclavable
- / Easy-to-read graduation
- / Colour-coded for tip identification



Raw material: PP, Volume range: 50 - 1,250 µl, Nominal volume: 1,250 µl, Matching rack: 770340

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
778350	standard	○ natural	● blue	bag		1,000 / 5,000

Item no.	Feature	Product colour	Box insert colour	Type of packaging	Sterile	Qty. inner / outer
778351	standard	○ natural	● blue	rack		960 / 3,840
778352	standard	○ natural	● blue	rack	+	960 / 3,840
778354	standard	○ natural	● blue	refill		960 / 4,800
778355	standard	○ natural	● blue	refill	+	960 / 4,800
778361	low retention	○ natural	● blue	rack		960 / 3,840
778362	low retention	○ natural	● blue	rack	+	960 / 3,840
778364	low retention	○ natural	● blue	refill		960 / 4,800
778365	low retention	○ natural	● blue	refill	+	960 / 4,800
778353	filter	○ natural	● blue	rack	+	960 / 3,840
778363	low retention filter	○ natural	● blue	rack	+	960 / 3,840



FREE OF
detectable
DNase


FREE OF
detectable
human DNA

FREE OF
detectable
RNase

Racks (unfilled)

- / Highly transparent and stable
- / Can be autoclaved and reused up to 50 times
- / Stackable and automation-friendly
- / Colour-coded rack inserts for quick volume identification

Item no.	Description	Feature	Box insert colour	Rack colour	Qty. inner / outer
770310	suitable for 10 µl / 10 µl ext.		● red	○ natural	10 / 60
770320	suitable for 200 µl		● yellow	○ natural	10 / 60
770330	suitable for 300 µl		● green	○ natural	10 / 60
770340	suitable for 1000 µl / 1250 µl		● blue	○ natural	10 / 40
770370	suitable for 10 - 300 µl	without rack insert		○ natural	10 / 60
770380	suitable for 1000 - 1250 µl	without rack insert		○ natural	10 / 40

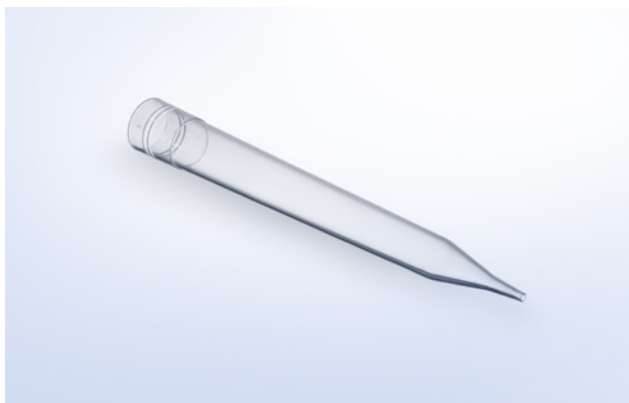


Macro Pipette Tips

1-5 ml

- / Volume range from 1 - 5 ml
- / Suitable for Gilson® P5000

Feature: Macro, Graduation: -, Volume range: 1 - 5 ml

Item no.	Product colour	Qty. inner / outer
745290	○ natural	250 / 2,500

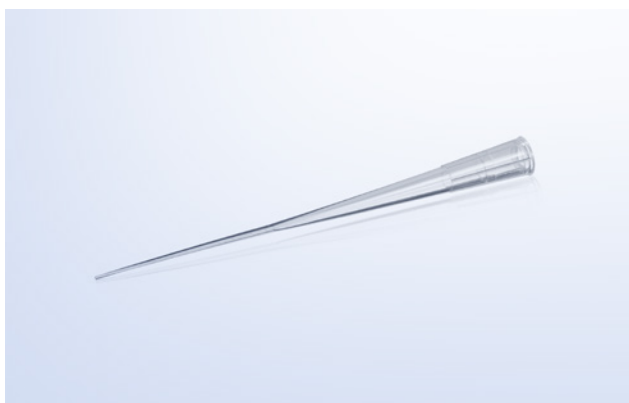

Gel-Load Pipette Tips

1-200 µl

- / Made from high-quality polypropylene
- / Universal fit for all common pipettes
- / 5 µl graduation for easy control of the pipetted volume
- / Tip with 0.57 mm outer diameter suitable for loading gel pockets
- / 83 mm total length to reach even the bottom of narrow vessels

FREE OF
detectable
DNase

FREE OF
detectable
human DNA


FREE OF
detectable
RNase

non-
pyrogenic

PCR

Volume: 200 µl, Volume range: 1 - 200 µl, Type of packaging: bag

Item no.	Product colour	Qty. inner / outer
775390	○ natural	1,000 / 10,000

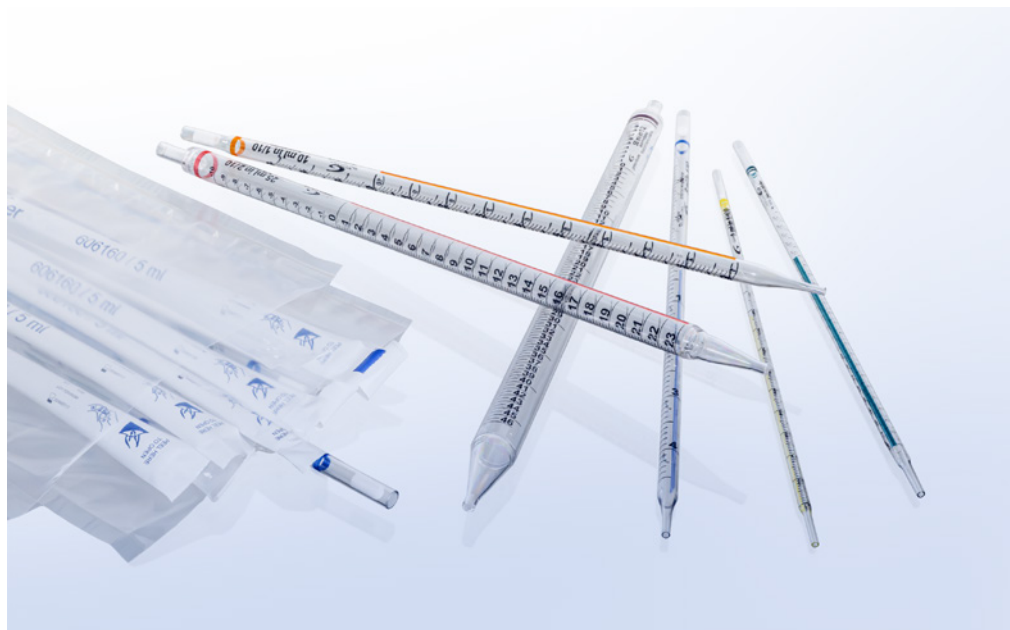

/ Selected pipettes are certified with ACT® Ecolabel. Learn more:



CELLSTAR® SEROLOGICAL PIPETTES


Greiner Bio-One offers a wide range of different serological pipettes and transfer pipettes.

The volume capacity of the serological pipettes is increased by a negative graduation. The pipettes are provided with colour codes according to international standards. The use of high-grade polystyrene guarantees maximum clarity. Expiry date and LOT number are printed on each single packaging unit. All pipettes are supplied with a filter for protection against the suction of liquid into the pipetting device.

CELLSTAR® serological pipettes are available in three different packaging options: in plastic bulk packs as well as individually wrapped in plastic or paper/plastic, both with peel-off function and additional break-through function.

To promote sustainability and transparency, many of our serological pipettes have already been **certified by My Green Lab®** with an excellent rating. The respective ACT® Ecolabel allows a clear and reliable assessment of the environmental impact of our products.

- / Tip design guarantees drop-free pipetting
- / Maximum accuracy
- / High optical clarity
- / Clear, easy to read graduation
- / A coloured vertical Schellbach stripe on 1, 2, 5, 10 and 25 ml pipettes makes it considerably easier to read off the volume
- / Colour code according to international standards

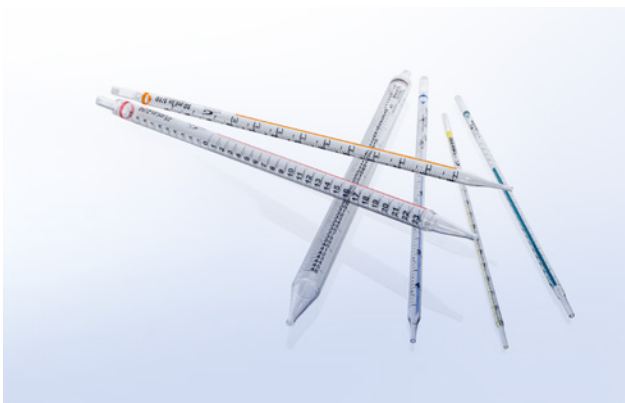

Serological Pipettes


- / Increased volume range with negative graduations
- / Lot number and expiry date on each bag
- / Filter for protection against suction of liquids into the pipetting device
- / Various packaging options
- / ACT® Ecolabel certification

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic	my green lab. ACT Ecolabel
---------	--------------------------	------------------------------	--------------------------	---------------	---------------	----------------------------

Raw material: PS

Item no.	Description	Graduation	Nominal volume	Packaging	Type of packaging	Sterile	Oty. inner / outer
604107	1 ml pipette	1/100	1 ml		bulk	+	25 / 1,000
604160	1 ml pipette	1/100	1 ml	plastic / plastic	single-packed	+	100 / 1,000
604181	1 ml pipette	1/100	1 ml	paper / plastic	single-packed	+	100 / 1,000
710107	2 ml pipette	1/100	2 ml		bulk	+	25 / 1,000
710160	2 ml pipette	1/100	2 ml	plastic / plastic	single-packed	+	100 / 1,000
710180	2 ml pipette	1/100	2 ml	paper / plastic	single-packed	+	100 / 1,000
606107	5 ml pipette	1/10	5 ml		bulk	+	25 / 500
606160	5 ml pipette	1/10	5 ml	plastic / plastic	single-packed	+	50 / 200
606180	5 ml pipette	1/10	5 ml	paper / plastic	single-packed	+	50 / 200
607107N	10 ml pipette	1/10	10 ml		bulk	+	25 / 450
607180N	10 ml pipette	1/10	10 ml	paper / plastic	single-packed	+	50 / 200
607160N	10 ml pipette	1/10	10 ml	plastic / plastic	single-packed	+	50 / 200
760107	25 ml pipette	2/10	25 ml		bulk	+	25 / 200
760160	25 ml pipette	2/10	25 ml	plastic / plastic	single-packed	+	50 / 200
760180	25 ml pipette	2/10	25 ml	paper / plastic	single-packed	+	50 / 200
768160	50 ml pipette	1/2	50 ml	plastic / plastic	single-packed	+	20 / 100
768180	50 ml pipette	1/2	50 ml	paper / plastic	single-packed	+	20 / 100

/ All products in this overview are certified with ACT® Ecolabel. The ACT® registered trademark is owned and controlled by My Green Lab Inc. More details at <https://act.mygreenlab.org>

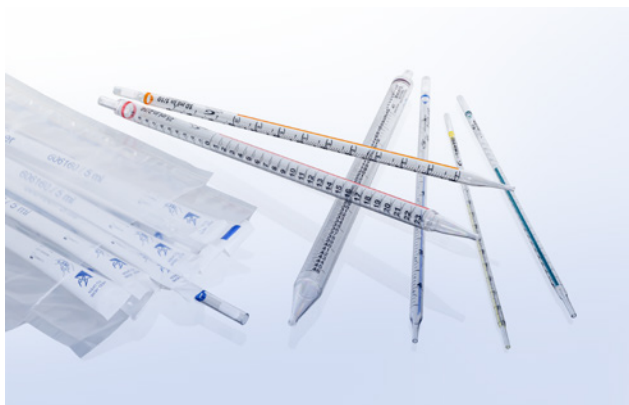

Serological Pipettes

Triple-packed

- / Sterile
- / Increased volume range with negative graduations
- / Lot number and expiry date on each bag
- / Filter for protection against suction of liquids into the pipetting device
- / Triple-packed

STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Raw material: PS, Packaging: plastic / plastic, Triple-packed: yes, Sterile: +

Item no.	Description	Graduation	Nominal volume	Sterile	Qty. inner / outer
604160-TRI	1 ml pipette	1/100	1 ml	+	10 / 100
710160-TRI	2 ml pipette	1/100	2 ml	+	10 / 100
606160-TRI	5 ml pipette	1/10	5 ml	+	10 / 100
607160N-TRI	10 ml pipette	1/10	10 ml	+	10 / 100
760160-TRI	25 ml pipette	2/10	25 ml	+	10 / 100
768160-TRI	50 ml pipette	1/2	50 ml	+	10 / 100

Aspirating Pipette

- / 2 ml aspirating pipette without plug and graduation
- / Single-packed



STERILE	FREE OF detectable DNase	FREE OF detectable human DNA	FREE OF detectable RNase	non-cytotoxic	non-pyrogenic
---------	--------------------------	------------------------------	--------------------------	---------------	---------------

Description: 2 ml aspirating pipette, Feature: no plug, Graduation: -, Raw material: PS, Nominal volume: 2 ml, Packaging: paper / plastic, Sterile: +

Item no.	Sterile	Qty. inner / outer
710183	+	100 / 1,000



MaxiPette

- / Ergonomic design for easy pipetting
- / Variable speed control
- / Cordless and rechargeable
- / Up to eight hours of continuous operation
- / Compatible with Greiner Bio-One serological pipettes

Description: MaxiPette, Content kit: pipette controller charger benchtop stand extra filter replaceable lithium battery

Item no.	Qty. inner / outer
847070	1 / 1

/ Device with US or UK plug available on request



Pasteur / Serum Pipettes

- / Ideally suited for the rapid transfer of liquids
- / Sterile / non-sterile

Length: 153 mm

Item no.	Description	Graduation	Working volume	Sterile	Qty. inner / outer
700370	pasteur pipette	-	≤0.1 ml		500 / 1,500
612301	serum pipette	yes	≤2.5 ml		500 / 1,500
612362	serum pipette	yes	≤2.5 ml	+	25 / 1,000



The Greiner Bio-One portfolio includes small laboratory equipment to match the consumables such as PCR tubes, plates and tubes, in order to provide our customers with comprehensive service, compatibility and highest quality: Mini-centrifuges, Vortex Mixers, Block Heaters, Microplate Centrifuges and ergonomic pipettes are used for modern handling in everyday laboratory work. All from one source - for highest quality requirements.

LAB EQUIPMENT

/ Lab Equipment	236
Mini Centrifuge	237
Microplate Centrifuge	237
Mini Vortex Mixer	238
Vortex Mixer	238
Mini Block Heater	239
Inserts for Mini Block Heater	239



- / For Sapphire Maxi-Pette pipette controller please refer to chapter Liquid Handling
- / For further information, please refer to our website: www.gbo.com

LAB EQUIPMENT

Greiner Bio-One, a leading provider of labware, completes its product portfolio by compatible lab equipment.

Mini centrifuges are ideal for quick spin-downs of microtubes and PCR tubes. Braking system is activated when the lid is opened, smoothly decelerating to a complete stop in just one second.

The microplate centrifuge allows centrifugation of two pla-

tes, even without a lid or seal.

Our mini vortex mixer can vortex even the largest samples instantly with its 4 mm orbit and speed of 2800 rpm. Despite its powerful motor, it has a size smaller than 10 x 10 cm, allowing it to fit on even the most crowded bench.

The Mini Bock Heater is the perfect tool for laboratory incubations, with its compact construction, broad temperature range and multiple block options.

- / Centrifuges
- / Vortex Mixer
- / Mini Block Heater



Mini Centrifuge

- / Ideal for quick spin-downs of microtubes and PCR tubes
- / Rotor instantly reaches 6,000 rpm
- / Storage compartment for PCR tube rotor
- / Decelerates in just one second
- / Compatible with all Greiner Bio-One PCR and reaction tubes

Feature: capacity 8 x 1.5 / 2.0 ml tubes, 4 x PCR strips (8 x 0.2 ml), 32 x 0.2 ml PCR tubes, Height: 11.4 cm, Length: 11.4 cm, Width: 15 cm, Speed: 6,000 RPM, Operating temperature: 4 - 45 °C, Max. radius: 4.9 cm

Item no.	Plug	Qty. inner / outer
843070	Europe	- / 1



Microplate Centrifuge

- / Rapid spin-down of microplates with 2,550 rpm
- / Unique rotor design holds two plates without spillage
- / 50 % smaller than traditional centrifuges
- / Compatible with most Greiner Bio-One microplates

Description: Microplate Centrifuge, Feature: capacity x 96 well microplates, 2 x 96 well PCR microplates, Height: 19.7 cm, Length: 23 cm, Width: 26 cm, Speed: 2,550 RPM, Operating temperature: 4 - 45 °C, Max. radius: 8 cm

Item no.	Plug	Qty. inner / outer
846070	Europe	- / 1

/ Device with US or UK plug available on request



Mini Vortex Mixer

- / Powerful, quiet vortexing for tubes up to 50 ml
- / Instant pressure-activated operation
- / Compact, fits in the palm of your hand
- / Compatible with all Greiner Bio-One tubes

Description: Mini Vortex Mixer, Feature: head: Standard rubber lid for mixing single tubes, Height: 6.6 cm, Length: 9.4 cm, Width: 9.9 cm, Mode: TOUCH Mix, Orbit: 4 mm, Operating temperature: 4 - 45 °C

Item no.	Plug	Qty. inner / outer
845070	Europe	- /1

/ Device with US or UK plug available on request



Vortex Mixer

- / Powerful motor for instant vortexing of up to 50 ml tubes
- / Dynamic balance system to minimise noise and vibration
- / Touch or continuous operation
- / Variable speed control from 200 to 3,200 rpm
- / Compatible with all Greiner Bio-One tubes

Description: Vortex Mixer, Feature: head: standard rubber cap single tube, base with four suction cups, Height: 17 cm, Length: 13 cm, Width: 16 cm, Mode: ON (continuous), OFF and TOUCH mix, Orbit: 3 mm, Operating temperature: 4 - 45 °C

Item no.	Plug	Qty. inner / outer
844070	Europe	- /1

/ Device with US or UK plug available on request

Mini Block Heater




- / Simple touchpad operation
- / Easy-to-read digital display
- / Compact, fits in the palm of your hand
- / Exchangeable blocks for tubes from 0.2 to 50 ml

Description: Mini block heater (without insert), block options, high-grade aluminium, Feature: ambient temperature from 5 - 100 °C, Height: 10 cm, Length: 11.2 cm, Width: 15 cm, Temperature increments: +/- 0.5 °C, Temperature uniformity: 0.2 °C, Temperature incements: 0.1 °C, Operating temperature: 4 - 45 °C

Item no.	Plug	Qty. inner / outer
848070	Europe	- /1

/ Device with US or UK plug available on request

Inserts for Mini Block Heater

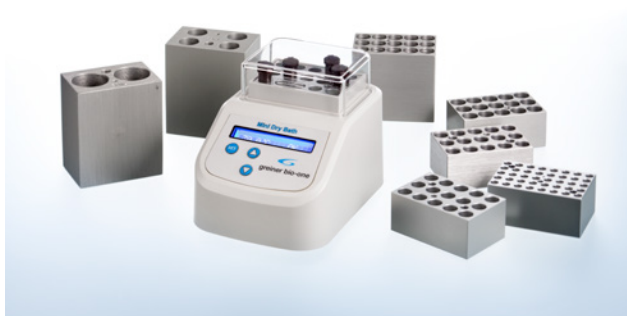

- / Exchangeable blocks for tubes from 0.2 to 50 ml

Item no.	Feature	Qty. inner / outer
848916	suitable for 15 pcs. 1.5 ml reaction tube, 0.5 ml PCR tube	- /1
848923	suitable for 15 pcs. 1.5 / 2 ml reaction tube, 0.5 ml PCR tube	- /1
848902	suitable for 40 pcs. 0.2 ml PCR tube, PCR 8-tube strip	- /1
848913	suitable for 15 pcs. 4 / 5 ml Cryo.s, 4.5 / 5 / 7 ml tube	- /1
848921	suitable for 15 pcs. 1 / 2 ml Cryo.s, 4 ml tube	- /1
848915	suitable for 4 pcs. 15 ml conical tube, 12 / 14 / 20 ml tube	- / 1
848950	suitable for 2 pcs. 50 ml conical tube	- / 1



DIN EN ISO 9001 Certification



EN ISO 13485 Certification



DIN EN ISO 50001 Certification

QUALITY STANDARDS AT GREINER BIO ONE

Greiner Bio-One is certified according to the international standards DIN EN ISO 9001 and EN ISO 13485 for Medical Devices. Since 2013,

Greiner Bio-One Frickenhausen (Germany) is also certified according to DIN EN ISO 50001 (systematic energy management).

TECHNICAL APPENDIX

<ul style="list-style-type: none"> / General Information for the Lab 242 <ul style="list-style-type: none"> Chemical Resistance of Various Materials 242 Chemical Resistance of Cycloolefins (COC / COP)..... 244 Chemical Resistance of Polyethylene Terephthalate (PET) Capillary Pore Membranes (ThinCert® Cell Culture Inserts)..... 245 Chemical Resistance of Sealers 246 Temperature Stability of Sealers 246 Physical Properties of Various Materials..... 247 Manual Calculation 248 Overview Metric Prefixes 249 / Laboratory Information for Microplates..... 250 <ul style="list-style-type: none"> Well Profiles of 96 Well Microplates 250 Well Profiles of 96 Well ELISA Strip Plates 252 / Laboratory Information for Liquid Handling..... 254 <ul style="list-style-type: none"> Compatibility for Sapphire Pipette Tips and Pipettes..... 254 / Laboratory Information for PCR 258 <ul style="list-style-type: none"> Overview PCR Microplates 258 Compatibility for PCR Microplates 260 Compatibility for Mini Block Heater Inserts..... 262 	<ul style="list-style-type: none"> / Laboratory Information for Centrifugation 262 <ul style="list-style-type: none"> Centrifugation - Principle and Calculation of the RCF (Relative Centrifugal Force) 262 Alignment Chart 264 Maximal Centrifuge Capacity of Tubes, Reaction Tubes and Microplates 264 Reaction Tubes..... 264 Polystyrene Tubes 265 Multiwell Plates..... 266 Microplates..... 266 PCR Plates..... 267 / Laboratory Information for Cryo.s Sample Storage 268 <ul style="list-style-type: none"> Freezing Protocol 268 Thawing Protocol..... 269 / Laboratory Information for Immunology..... 270 <ul style="list-style-type: none"> Volume-Dependent Wetting of Immunological Products..... 270 / Glossary 272
---	---

GENERAL INFORMATION FOR THE LAB

CHEMICAL RESISTANCE OF VARIOUS MATERIALS

	PS 20°C	PS 50°C	PP 20°C	PP 50°C	HDPE 20°C	HDPE 50°C	LDPE 20°C	LDPE 50°C	PC 20°C	PC 50°C
Acetic acid 10 %	1	1	1	1	1	1	1	1	1	2
Acetic acid 50 %	2	2	1	1	1	1	1	1	1	2
Acetic acid 90 %	4	4	1	2	1	1	1	2	4	4
Acetone	4	4	1	3	1	1	3	3	4	4
Acetonitrile	4	4	3	4	1	1	1	1	4	4
Ammonia 25 %	2	2	1	1	1	1	1	1	4	4
Ammonium acetate	1	1	1	1	1	1	1	1	1	1
Amyl alcohol	1	1	1	1	1	1	1	2	1	
Ascorbic acid			1	1	1		1		2	2
Benzene	4	4	4	4	4	4	4	4	4	4
Benzyl alcohol	4	4	4	4	3	4	4	4	4	4
Boric acid 10 %	1	1	1	1	1	1	1	1	1	1
Carbon tetrachloride	4	4	4	4	3	4	4	4	4	4
Carbonic acid	1	1	1	1	1	3	1	1	1	
Chloroform 100 %	4	4	3	4	3		3		4	4
Citric acid 10 %	1	1	1	1	1	1	1	1	1	2
Cyclohexanol	3	3	1	3	1	1	1	1	3	
Detergents			1	1						
Dichloroacetic acid			1	1	1	1			4	4
Diethyl ether	4	4	4	4	3	4	4	4	4	4
Dimethyl acetamide	4	4	1	1	1	1	3	4		
Dimethylsulfox. (DMSO)	1	2	1	1	1	1	1	1	4	4
Emulsifier			1	1						
Ethanol 50 %	1	1	1		1	1	1	2	1	1
Ethanol 96 %	1	1	1	1	1		1		1	3
Ether	4	4	4	4	3	4	4	4	4	4
Formaldehyde 10 %	3	4	1	1	1	1	1	1	1	2
Formaldehyde 40 %	4	4	1	2	1	2	2	3	1	2
Formamide	1	1	1	1	1	1	1	1	3	3
Formic acid 50 %	3	3	1	2	1	1	1	2	3	3
Glucose	1	1	1	1	1	1	1	1	1	1
Glycerine	1	1	1	1	1	1	1	1	3	3
Heptane	4	4	3	3	2	3	3	4	1	2
Hexanol			1		1		1		2	2
Hydrochloric acid 20 %	1	1	1	1	1	1	1	1	2	3
Hydrochloric acid conc.	3	3	1	1	1	1	1	1	4	4
Hydrogen peroxide 3 %	1	1	1	1	1	1	1	1	3	3
Hydroquinone	4	4	1				1	3	3	3
Isoamyl alcohol	1	1							3	0
Isobutanol	2	2	1	1	1	1	1	1	1	2
Isopropanol	2	2	1	1	1	1	1	1	1	2
Isopropyl acetate	4	4	2	3	1	2	2	3	4	4
Isopropyl benzene	4	4	3	4	2	3	3	4	4	4
Isopropyl ether	4	4	4	4	4	4	4	4	4	4
Lactic acid 3 %	2	2	1	2	1	1	1	2	1	2
Lactic acid 85 %	2	2	1	2	1	1	1	1	1	2
Liquid paraffin	1	1	1	3	1	1	1	3	1	

CHEMICAL RESISTANCE OF VARIOUS MATERIALS

	PS 20°C	PS 50°C	PP 20°C	PP 50°C	HDPE 20°C	HDPE 50°C	LDPE 20°C	LDPE 50°C	PC 20°C	PC 50°C
Methanol	3	4	1	1	1	1	1	1	4	4
Methyl acetate	4	4	2	3	3	3	3	4	4	4
Methyl phenyl ether 100 %	4	4	3				3		4	4
Methyl propyl ketone	4	4	2	3	1	2	2	3	4	4
Methylamine 32 %			1		1		1		4	4
Methylene chloride	4	4	3	4	4	4	4	4	4	4
Naphthalene			1		1	3			3	3
Nitric acid 70%	4	4	4	4	4	4	4	4	4	4
Nitrobenzene	4	4	4	4	3	4	4	4	4	4
Oxalic acid	1	1	1	1	1	1	1	1	3	4
Ozone	3	3	1	2	1	1	1	2	1	2
Palmitic acid	1	1	3	4	3		2		2	2
Phenol 10 %	4	4	1	1	1	1	1	1	4	4
Phenol 100 %	4	4	1	1	2	3	3	3	4	4
Phosphoric acid 1-5 %	2	2	1	1	1	1	1	1	1	1
Phosphoric acid 85 %	1	1	1	2	1	1	1	1	1	2
Phthalic acid	1	1	1	1	1	1	1	1	3	3
Potassium carbonate	1	1	1	1	1	1	1	1	3	3
Potassium chromate	1	1	1	1	1	1	1		2	2
Potass. permanganate	2	3	1	1	1	1	1	1	1	
Propanol	3	3	1	1	1	1	1	1	1	
Sodium acetate	2	2	1	1	1	1	1	1	1	2
Sodium chloride	1	1	1	1	1	1	1	1	1	1
Sodium hydroxide 30 %	1	1	1	1	1	1	1	1	4	4
Sodium hydroxide 45 %	1	1	1	1	1	1	1	1	4	4
Sodium hydroxide 60 %	1	1	1	1					4	4
Sodium hypochloride	1	1	2	3	2	3	2	3	2	3
Sodium permanganate	2	3	1	1	1	1	1	1		
Sodium thiosulfate	1	1	1	1	1	1	1	1	2	2
Stearic acid	1	2	1	1	1	1	1	1	1	2
Sulphuric acid 1-6 %	1	2	1	1	1	1	1	1	1	1
Sulphuric acid 60 %	2	4	1	3	1	3	1	3	3	3
Sulphuric acid conc.	4	4	4	4	4	4	4	4	4	4
Tannin acid	1	1	1	1					3	3
Terpentine oil					3	4	3	4	4	4
Tetrahydrofuran	4	4	3	4	3	4	4	4	4	4
Toluene	4	4	3	4	3	4	3	4	4	4
Trichloroacetic acid	4	4	3	4	3	3	3	4	4	4
Urea	1	2	1	1	1	1	1	1	1	1
Uric acid			1		1		1		1	
Urine	3	3	1	1	1	1	1	1	1	
Xylene	4	4	4	4	2	3	2	4	4	4

1 = resistant 2 = limited resistant 3 = moderate resistant 4 = no resistance

These tables are a general guide only. As many factors can affect the chemical resistance of a given product, its suitability for a specific application should be tested.

CHEMICAL RESISTANCE OF CYCLOOLEFINS (COC / COP)

Cycloolefin		Cycloolefin		Cycloolefin	
Acetic acid 99 %	1	Detergents	1	Hydrogen peroxide water 30 %	1
Acetone	1	Dibutyl ether	4	Isopropanol	1
Acetonitrile	1	Dichloroethane	4	Methanol	1
Acrylonitrile	1	Dichloromethane	4	Methylene chloride	4
Ammonia 33 %	1	Diethyl ether	4	Nitric acid (HNO ₃)	1
Benzaldehyde	3	Dimethyl sulfoxide	1	Octane	4
Benzene	4	DMSO	1	Pentane	4
Benzine	4	Ethanol 50 %	1	Sodium hydrox. (NaOH) 50 %	1
Butanon	1	Ethanol 96 %	1	Sulphuric acid (H ₂ SO ₄) 40 %	1
Carbon tetrachloride	4	Fatty acid	4	Sulphuric acid (H ₂ SO ₄) 95 %	1
Chloroform	4	Heptane (n-Heptane)	4	Toluene	4
Cyclohexane	4	Hexane	4	Xylene	4
Cyclohexanone	4	Hydrochloric acid (HCl) 36 %	1		

1 = resistant 2 = limited resistant 3 = moderate resistant 4 = no resistance

This table is a general guide only. As many factors can affect the chemical resistance of a given product, its suitability for a specific application should be tested.

CHEMICAL RESISTANCE OF POLYETHYLENE TEREPHTHALATE (PET) CAPILLARY PORE MEMBRANES (THINCERT® CELL CULTURE INSERTS)

PET		PET		PET	
Acetaldehyde	1	Ethanol	1	Monochlorbenzene	1
Acetic acid (10 %)	1	Ethyl acetate	1	Nitric acid (30 %)	1
Acetic acid (100 %)	3	Ethyl ether	1	Nitrobenzene	1
Acetone	1	Ethylendichloride	1	Nitropropane	1
Ammonium hydroxide (5 %)	1	Ethylene glycol	1	n-Propanol	1
Amyl acetate	1	Fluoric acid (35 %)	1	Pentane	1
Amyl alcohol	1	Formaldehyde	1	Perchlorethylene	1
Aniline	1	Formic acid (50 %)	1	Petroleum ether	1
Benzene	3	Freon	1	Phosphoric acid (85 %)	3
Benzyl alcohol	1	Glutaraldehyde	1	Potassium hydroxide	4
Benzyl benzoate	1	Glycerol	1	Propyl acetate	1
Boric acid (5 %)	1	H ₂ O ₂ (30 %)	1	Pyridine	1
Butanol	1	Halogenated phenoles	4	Silicon oil	1
Butyl acetate	1	Hexane	1	Sodium hydroxide	4
Butyl cellulose	1	Hydrochloric acid (20 %)	1	Sulphuric acid (25 %)	1
Carbon tetrachloride	1	i-Propanol	1	Terpentine oil	1
Chloroform	1	Isopropyl myristate	1	Tetrahydrofurane	1
Concentrated strong acids	4	Methanol	1	Tetraline	1
Cyclohexane	1	Methyl acetate	1	Toluene	3
Cyclohexanone	3	Methyl cellulose	1	Trichlorbenzene	1
Dekaline	1	Methylenchloride	3	Trichlorethylene	1
Dimethylacetamide	1	Methylethylketone	1	Triethanolamin	1
Dimethylformamide	1	Methylglycol acetate	1	Trikresyl phosphate	1
Dimethylsulfoxide	1	Methylisobutylketone	1	Xylene	3
Dioxane	1	Mineral oils	1		

For the solvents effecting slight changes the user should test the compatibility under the specific application conditions. All tests have been performed at RT. Please be aware that ThinCert® cell culture inserts are made of PET membranes sealed on polystyrene housings. Therefore, solvents shown compatible with PET membranes in the above table might be incompatible with the polystyrene housing. Please check solvent compatibility with polystyrene.

Resistance scale from 1 to 4

1 = resistant	i.e. the plastics may be treated with the chemical compound at mentioned temperature over several years without any significant alterations in its physical, optical and chemical properties
2 = limited resistant	i.e. the plastics may be treated with the chemical compound at mentioned temperature over several weeks without any significant alterations in its physical, optical and chemical properties
3 = moderate resistant	i.e. the plastics may be treated with the chemical compound at mentioned temperature for short time only (several minutes to one hour) without any alterations in physical, optical and chemical properties (mixing and measuring is possible)
4 = no resistance	i.e. treating the plastics with the substance named may cause alterations in physical, optical and chemical properties within seconds

CHEMICAL RESISTANCE OF SEALERS

	EASYseal (Art. No. 676001)	VIEWseal (Art. No. 676070)	AMPLiseal (Art. No. 676040)	SILVERseal (Art. No. 676090)
Acetone	4	4	4	3
Acetonitrile	3	3	4	1
Acetic acid 1 %	3	1	4	3
Glacial acetic acid	1	3	4	3
Chloroform	4	4	4	4
DMSO	3	3	3	1
Ethanol	3	1	1	1
Hydrochloric acid 32 %	3	1	3	4
Isopropanol	3	1	1	1
Methanol	3	1	4	1
Phenol	3	3	4	3
Sulphuric acid 0.5 M	1	1	1	1

1 = Stable
 3 = Moderately stable
 4 = Unstable

no visible change in the sealer after one week's incubation
 after one week, optical and physical changes in the sealer (clouding tears on removal)
 adhesive and foil are dissolved, wells not leak-tight

This table can only be used as an orientation aid for the suitability of the respective sealers, since their behaviour against chemicals depends on the respective application. Tests under practical conditions are absolutely essential in many cases.

TEMPERATURE STABILITY OF SEALERS

	Temperature Stability
EASYseal	-40 °C to + 120 °C
VIEWseal	-70 °C to + 100 °C
AMPLiseal	-80 °C to + 110 °C
SILVERseal	-70 °C to + 100 °C
BREATHseal	n.a. / Evaporation rate 4200 g H ₂ O/m ² in 24 h

This table can basically be used as an orientation aid for the temperature stability of the respective sealers, since the behaviour of the product depends on the respective application. Tests under practical conditions are absolutely essential in many cases.

PHYSICAL PROPERTIES OF VARIOUS MATERIALS

Material	Sterilisation by				Auto-clavability	Thermal Stability [°C]	Transparency	Gas Permeability ²			WVTR ³
	gamma irradiation	chemicals (formalin, ethanol)	dry heat	gas ¹				O ₂	N ₂	CO ₂	
Polystyrene	•	•		•		-20 to +60	clear	4.7	853	17.8	108 - 155
Polypropylene	•	•		•	•	-196 to +121	translucent	3.7	744	12.4	3.9
HDPE	•	•		•		-50 to +100	translucent	2.9	651	9	4.6 - 6.2
LDPE	•	•		•		-50 to +80	translucent	7.8	2.8	41.9	15.5 - 23.3
UV-Star®	•			•		-20 to +40	clear				
PETG	•	•		•		-40 to +60	clear	388	155	1.2	62
PET	•	some		•		-40 to +60	clear	46.5	10.9	236	15 - 20
Cycloolefin	•			•		-80 to +100	clear				

Exemptions are mentioned in the respective product data sheets.

¹ Ethylene oxide, formaldehyde

² [cc x mm/m² x 24 h x Bar]

³ at 37°C, 90 % humidity [g x mm/m² x 24 h x Bar]

Material	Refractive Index
Polystyrene	1.59
UV-Star®	1.53
Cycloolefin	1.53
Glass	1.53

These tables are a general guide only. As many factors can affect the resistance of a given product, its suitability for a specific application should be tested.

MANUAL CALCULATION

COEFFICIENT OF VARIATION (CV)


The coefficient of variation compares the variability of several random samples with different means, taking into account the different dimensions of means:

$$CV \% = \frac{S}{|\bar{X}|} \cdot 100 \%$$

/ where S is the standard deviation and $|\bar{X}|$ is the absolute value of the arithmetic mean.

VOLUME OF DIVERSE BODIES

VOLUME OF A CYLINDER:

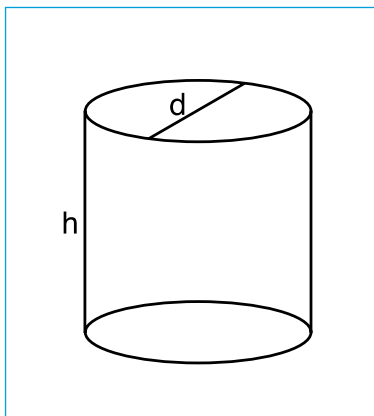


$$V = \frac{\pi \cdot d^2 \cdot h}{4}$$

$$h = \frac{4 \cdot V}{\pi}$$

/ This formula can be used for calculating the filling level in relation to the filling volume in a 96 well microplate with cylindrical wells.

VOLUME OF A CUBOID:

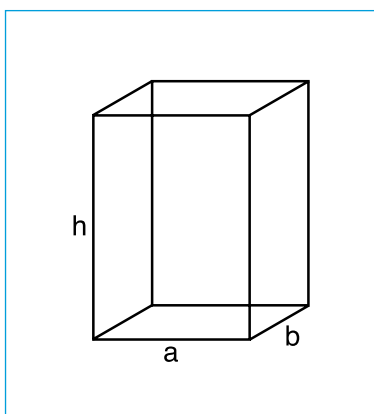

$$V = a \cdot b \cdot h$$

$$h = \frac{V}{a \cdot b}$$


/ This formula can be used for calculating the filling level in relation to the filling volume in 384 and 1536 well microplates with rectangular wells.

OVERVIEW METRIC PREFIXES

G	= giga	= 10^9
M	= mega	= 10^6
k	= kilo	= 10^3
c	= centi	= 10^{-2}
m	= milli	= 10^{-3}
μ	= micro	= 10^{-6}
n	= nano	= 10^{-9}
p	= pico	= 10^{-12}
f	= femto	= 10^{-15}
a	= atto	= 10^{-18}
z	= zepto	= 10^{-21}

LABORATORY INFORMATION FOR MICROPLATES

WELL PROFILES OF 96 WELL MICROPLATES

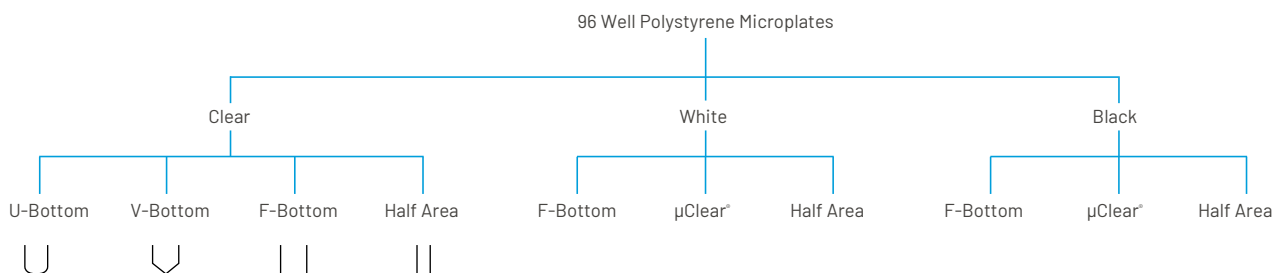

WELL PROFILE

The well profile is a critical aspect in a 96 well microplate. Different well shapes are available for each application (Fig. 1 – Fig. 4):

U-BOTTOM

The “U” describes the round bottom shape (Fig. 1). U-bottom microplates are ideally suited for agglutination tests.

- / No sharp corners to facilitate easy and residue-free pipetting
- / Suitable for +/- analyses

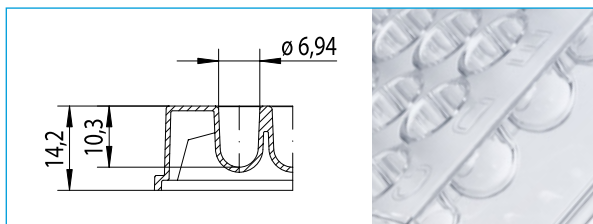


Figure 1:
Well profile: 96 well U-bottom, polystyrene
Total volume: 323 µl
Working volume: 40 – 280 µl

V-BOTTOM

The “V” stands for the conically tapered well bottom (Fig. 2). These microplates are ideally suited for applications in which the entire sample volume must be pipetted off.

- / For precise pipetting
- / Ideally suited for the storage of samples
- / Suitable for +/- analyses

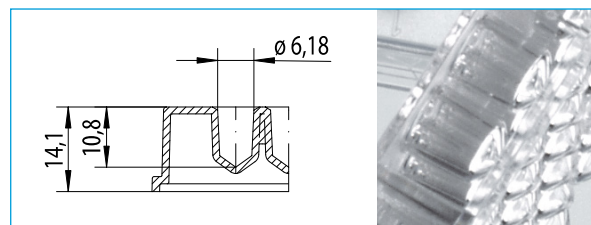


Figure 2:
Well profile: 96 well V-bottom, polystyrene
Total volume: 234 µl
Working volume: 40 – 200 µl

F-BOTTOM/STANDARD (ST)

The "F" refers to the flat well bottom (Fig. 3). This well type is ideal for precise optical measurements. The measuring light source is not deflected by the well profile.

- / Excellent optical properties
- / For precise optical measurements
- / For microscopic applications (bottom reading)

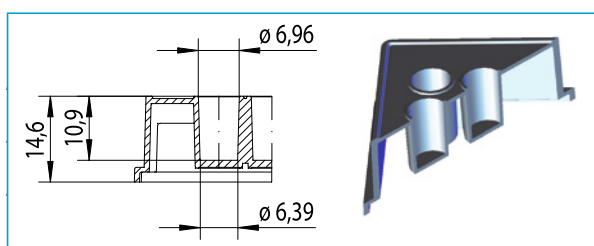


Figure 3:
Well profile: 96 well F-bottom / ST, polystyrene
Total volume: 382 μ l
Working volume: 25–340 μ l
Growth area: 32 mm²

HALF AREA

Half Area microplates (Fig. 5) offer an interesting alternative to conventional 96 well microplates. They can be pipetted manually without any problem but at the same time allow a reduction of the sample volume by up to 50 %.

- / Reduction of sample volume of up to 50 %
- / Standardised pathlength (1 cm=170 μ l, 0.5 cm=80 μ l)

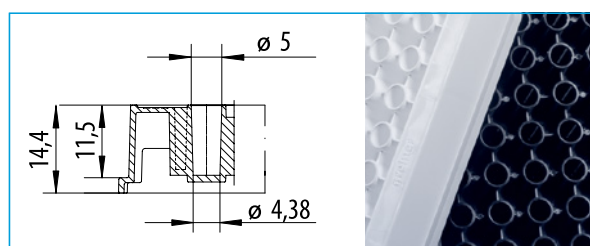


Figure 5:
Well profile: 96 well Half Area
Total volume: 199 μ l
Working volume: 15–175 μ l
Growth area: 15 mm²

F-BOTTOM/CHIMNEY WELL

The standard flat bottom microplate (Fig. 3) has the same well profile as the chimney well microplate (Fig. 4). The difference from the standard plate is the chimney-like arrangement of the wells. Each well stands on its own. Therefore the risk of sample carryover and cross contamination is minimised.


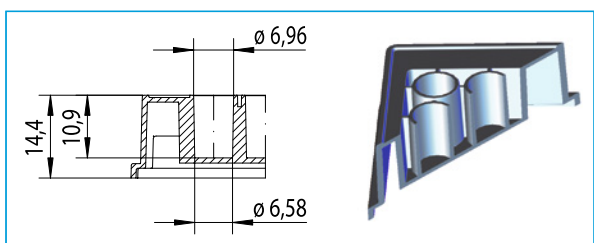


Figure 4:
Well profile: 96 well F-bottom / chimney well, polystyrene
Total volume: 392 μ l
Working volume: 25–340 μ l
Growth area: 34 mm²

WELL PROFILES OF 96 WELL ELISA STRIP PLATES

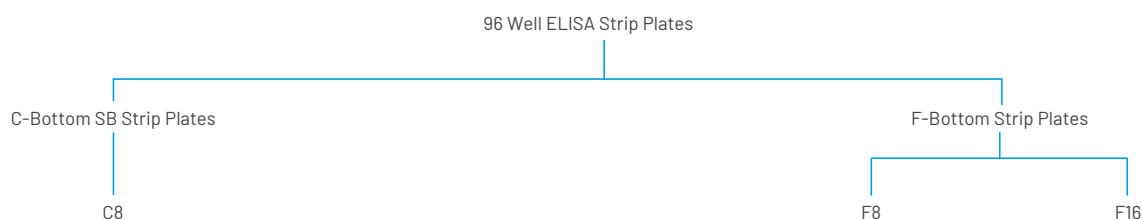

WELL PROFILE

Microplates in strip format offer the advantage of greater flexibility in diagnostics. Individual strips can be removed from the support frame so that the number of tests to be performed can be adjusted to the number of samples, and is not predetermined by the microplate format used. In addition, the individual strips of a microplate can be subjected to a wide variety of different test conditions.

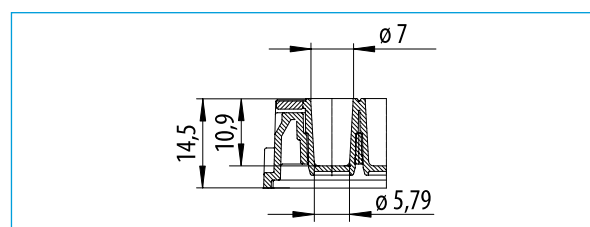


Figure 1:
 Well profile: C-bottom
 Total volume: 346 μ l
 Working volume: 20–300 μ l

C-BOTTOM SB (SINGLE-BREAK) STRIP PLATES

The “C” describes a flat-bottom profile with rounded corners (Fig. 1). The rounded corners mean that the individual wells can be pipetted without leaving a residue, and the flat bottom still enables precise optical measurements. C-bottom SB strips are supplied as twelve 8 well strips, in a support frame with 96 spaces (12 x 8 matrix). The individual wells can be broken off separately (“single-break” option) and the number of tests performed can thus be precisely adjusted to the number of samples. C-bottom SB strip plates are available in clear polystyrene and additionally with a red, green, blue or yellow colour coding on the well rim.

F-BOTTOM STRIP PLATES

96 well strip plates with a flat bottom (F-bottom, Fig. 2) are available as 8 well strips and as 16 well strips. The “F” stands for the flat bottom of the wells. This well type is ideal for the most precise optical measurements as the measuring light beam is not deflected by the well profile. F-bottom strip plates are available in clear polystyrene (MICROLON) and additionally in black (FLUOTRAC) and white (LUMITRAC).

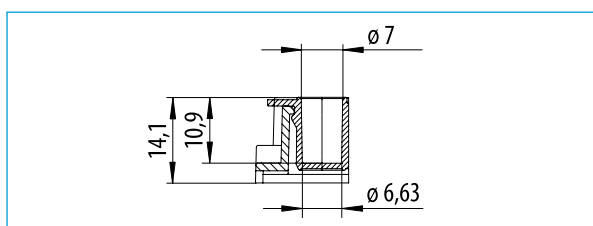


Figure 2:

Well profile: F-bottom

Total volume: 388 μ l

Working volume: 20-350 μ l

LABORATORY INFORMATION FOR LIQUID HANDLING

COMPATIBILITY FOR SAPPHIRE PIPETTE TIPS AND PIPETTES

	PIPETTE TIPS (STANDARD / LOW RETENTION)							
Volume [μ l]	10	10 extended	20	100	200	300	1000	1250
Volume range [μ l]	0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version	Tip/ Filter tip	Tip/ Filter tip	Filter tip	Filter tip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip
Article no.	7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX

PIPETTE MODELS

GREINER BIO-ONE

		Single-channel pipettes							
Sapphire	0.2-2 μ l	●	●						
	1-10 μ l	●	●						
	2-20 μ l			●		T ● F -			
	10-100 μ l			■	●	T ● F -			
	20-200 μ l			■	■	●	●		
	100-1000 μ l							●	●
	500-5000 μ l								
1000-10000 μ l									
		Multi-channel pipettes							
Sapphire	8/12 CH 0.5-10 μ l	●	●						
	8/12 CH 2-20 μ l			●		T ● F -			
	8/12 CH 20-200 μ l			■	■	●	●		
	8/12 CH 20-300 μ l			■	■	■	●		

BRAND

		Single-channel pipettes							
Transferpette® S	0.5-10 μ l	●	●						
	2-20 μ l	■	■	●		T ● F -			
	10-100 μ l			■		T ● F -			
	20-200 μ l			■	■	●	●		
	100-1000 μ l							●	T ● F -

EPPENDORF

		Single-channel pipettes							
Reference®	0.5-10 μ l	●	●						
	2-20 μ l			●	●	T ● F -			
	10-100 μ l			■	●	T ● F -			
	50-200 μ l				■	●	●		
	100-1000 μ l							●	●
Research®	0.5-10 μ l	●	●						
	2-20 μ l			●	●	T ● F -			
	10-100 μ l			■	●	T ● F -			
	20-200 μ l			■	■	●	●		
	100-1000 μ l							●	●
Research® plus	0.5-10 μ l	●	●						
	10-100 μ l			■	●	T ● F -			
	20-200 μ l				■	●	●		
	100-1000 μ l							●	●

PIPETTE TIPS (STANDARD / LOW RETENTION)

Volume [μl]	10	10 extended	20	100	200	300	1000	1250
Volume range [μl]	0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version	Tip/ Filter tip	Tip/ Filter tip	Filter tip	Filter tip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip
Article no.	7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX

PIPETTE MODELS

Research [®] pro	0.5-10 μl	●	●					
	5-100 μl				●	T ● F -		
	20-300 μl						●	
Fixed-volume pipettes								
Reference [®]	50 μl				●	T ● F -		
	100 μl				●	T ● F -		
	500 μl						●	●
	1000 μl						●	●
Research [®]	500 μl						●	●
Multi-channel pipettes								
Eppendorf	8CH 10-200 μl			■	■	●	●	
Research [®]	12CH 0.5-10 μl	●	●					
Research [®] plus	8CH 0.5-10 μl	●	●					
	8CH 10-100 μl			■	●	●	●	
	8CH 20-300 μl			■	■	■	●	
Electronic pipettes								
Xplorer [®]	50-1000 μl						●	●
	8CH 15-300 μl			■	■	■	●	

GILSON

Single-channel pipettes								
Pipetman [®]	P2 (0.5-2 μl)	●	●					
	P10 (1-10 μl)	●	●					
	P20 (2-20 μl)			●		T ● F -		
	P100 (20-100 μl)			■	●	●		
	P200 (50-200 μl)			■	■	●	●	
	P1000 (200-1000 μl)						●	●
Pipetman [®] L	P2L (0.2-2 μl)	●	●					
	P10L (1-10 μl)	●	●					
	P20L (2-20 μl)			●		T ● F -		
	P200L (20-200 μl)				■	●	●	
	P1000L (100-1000 μl)						●	●
Fixed-volume pipettes								
Pipetman [®]	F5 (5 μl)			●		T ● F -		
	F10 (10 μl)			●		T ● F -		
	F20 (20 μl)			●		T ● F -		
	F25 (25 μl)				●	T ● F -		

T = Pipette tip
F = Filter tip

● compatible - not compatible
■ Pipetting volume limited to the maximum volume of the pipette tips

PIPETTE TIPS (STANDARD / LOW RETENTION)

Volume [µl]	10	10 extended	20	100	200	300	1000	1250
Volume range [µl]	0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version	Tip/ Filter tip	Tip/ Filter tip	Filter tip	Filter tip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip
Article no.	7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX

PIPETTE MODELS

GILSON (continued)

Fixed-volume pipettes

Pipetman®	F100 (100 µl)						T ● F -	
	F200 (200 µl)				●	●		
	F250 (250 µl)						●	
	F300 (300 µl)						●	●
	F500 (500 µl)						●	
	F1000 (1000 µl)						●	●

Multi-channel pipettes

Pipetman® L	P8X 10L (0.5-10 µl)	●	●					
	P8X 200L (20-200 µl)			■	■	●	●	
	P8X 300L (20-300 µl)			■	■	■	●	

IKA

Single-channel pipettes

IKA Pette vario	100-1000 µl						●	
-----------------	-------------	--	--	--	--	--	---	--

METTLER TOLEDO

Single-channel pipettes

Rainin Pipette-Lite™	SL200 (20-200 µl)			■	■	●	●	
Rainin Pipette-Lite™	SL1000 (100-1000 µl)						●	●

NICHIRYO

Single-channel pipettes

Nichipet	0.5-10 µl	●	●					
	2-20 µl						T ● F -	
	10-100 µl			■	●	●	●	
	20-200 µl			■	■	●	●	
	100-1000 µl						●	●

Multi-channel pipettes

Nichipet	8K 30-300 µl				■	■	●	
----------	--------------	--	--	--	---	---	---	--

SARTORIUS

Single-channel pipettes

Biohit Proline®	0.5-10 µl	●	●					
Biohit Proline® plus	0.5-10 µl	●	●					
	2-20 µl			●			T ● F -	
	10-100 µl			■	●	●	●	
	20-200 µl			■	■	●	●	
	100-1000 µl						●	●

Multi-channel pipettes

Biohit Proline®	8CH 0.5-10 µl	●	●					
	8CH 25-250 µl					■	●	
	8CH 50-300 µl				■	■	●	

SOCOREX

Single-channel pipettes

Acura® 825	0.5-10 µl	●	●					
------------	-----------	---	---	--	--	--	--	--

PIPETTE TIPS (STANDARD / LOW RETENTION)

Volume [µl]	10	10 extended	20	100	200	300	1000	1250
Volume range [µl]	0.2-10	0.5-10	2-20	2-100	5-200	10-300	50-1000	50-1250
Tip version	Tip/ Filter tip	Tip/ Filter tip	Filter tip	Filter tip	Tip/ Filter tip	Tip/ Filter tip	Tip	Tip/ Filter tip
Article no.	7713XX	7723XX	7733XX	7743XX	7753XX	7763XX	7773XX	7783XX

PIPETTE MODELS

	2-20 µl			●		T ● F -		
	5-50 µl				●	T ● F -		
	10-100 µl			■	●	T ● F -		
	20-200 µl			■	■	●	●	
	100-1000 µl						●	●
Calibra® 822	1-10 µl	●						
	2-20 µl			●		T ● F -		
	10-100 µl			■		T ● F -		
	20-200 µl			■	■	●	●	
	100-1000 µl						●	●
Multi-channel pipettes								
Acura® 855	8CH 5-50µl			■	●	T ● F -		
Calibra® 852	8CH 1-10 µl	●	●					
	8CH 20-200 µl			■	■	T ● F -		
	12CH 10-100 µl			■		T ● F -		

STARLAB

Single-channel pipettes								
ErgoOne®	0.1-2.5 µl		●					
	0.5-10 µl	●	●					
	10-100 µl			■	●	T ● F -		
	100-1000 µl						●	●
Multi-channel pipettes								
ErgoOne®	8 CH 30-300 µl				■	■	●	

THERMO FISHER SCIENTIFIC

Single-channel pipettes								
Finnpipette™ F1	1-10 µl	●	●					
	2-20 µl			●		T ● F -	●	
	20-200 µl			■	■	●	●	
	30-300 µl						●	
	100-1000 µl						●	●
Finnpipette™ 4500	200-1000 µl						●	●
Multi-channel pipettes								
Finnpipette™ F1 MCP8	8 CH 1-10 µl	●	●					
	8 CH 10-100 µl			■	●	●	●	
	8CH 30-300 µl				■	■	●	
Finnpipette™ 4510	8CH 50-300 µl				■	■	●	
Finnpipette™ 4500	8CH 200-1000 µl						●	●

VWR

Single-channel pipettes								
Erg. High Perform.	0.5-10 µl	●	●					
Erg. High Perform.	2-20 µl			●		T ● F -		
Erg. High Perform.	10-100 µl			■	●	●		
Erg. High Perform.	20-200 µl			■	■	●	●	
Erg. High Perform.	100-1000 µl						●	●

LABORATORY INFORMATION FOR PCR

OVERVIEW PCR MICROPLATES

The use of the 96 well format allows the scale up of basic PCR work, while the 384 well format is ideal for high-throughput screening projects. All microplates are made of thin-walled polypropylene. This optimises the heat transfer from the thermoblock to

the reaction solution. Our heat-resistant sealers AMPLIseal, VIEWseal and SILVERseal are ideal for sealing the microplates during PCR, and the 96 well microplate may also be sealed with 8-cap strips.

96 WELL POLYPROPYLENE MICROPLATES FOR PCR

1/ NON-SKIRTED MICROPLATES

Non-skirted microplates may be used in all commonly available thermocyclers with a 96 well block.

1a) Non-skirted microplate with raised well rims (Art. No. 652201)

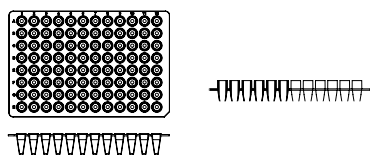


Figure 1a: View of a non-skirted microplate with raised well rims

1b) Non-skirted microplate with flat surface (Art. No. 652250)

Black alphanumeric coding enables a quick identification of samples

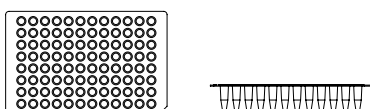


Figure 1b: View of a non-skirted microplate with flat surface

1c) Non-skirted microplate, low profile (Art. No. 652210)

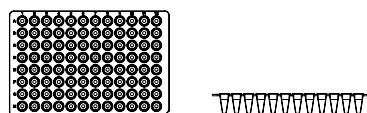


Figure 1c: View of a non-skirted low profile microplate

2/ HALF-SKIRTED MICROPLATES

2a) Half-skirted microplate with one notch suitable for real-time PCR systems such as LightCycler® 480 (Art. No. 669285)

- / Maximal pigmented white polypropylene and therefore most suitable for sensitive real-time PCR reactions
- / Black alphanumeric coding enables a quick identification of samples
- / Notches in the rim facilitate automation due to better gripping in robotic systems

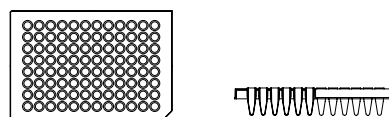


Figure 2a: View of a half-skirted microplate with one notch suitable for real-time PCR systems such as LightCycler® 480 from Roche

The registered trademarks of the mentioned manufacturers belong to the above mentioned companies.

2b) Half-skirted microplate with one notch suitable for ABI (Art. No. 652290)

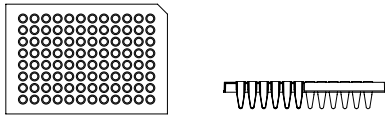


Figure 2b: View of a half-skirted microplate with one notch suitable for ABI

2c) Half-skirted microplate, recessed rim, ABI design with one notch (Art. No. 652260)

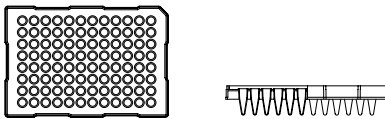


Figure 2c: View of a half-skirted microplate, recessed rim, ABI design with one notch

3/ FULL-SKIRTED MICROPLATE WITH ONE NOTCH (ART. NO. 652270)

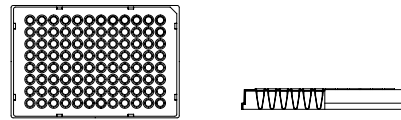


Figure 3: View of full-skirted microplate with one notch

384 WELL POLYPROPYLENE MICROPLATES FOR PCR

The 384 well PCR microplates from Greiner Bio-One are manufactured in an advanced injection moulding process following stringent quality criteria. Minimal distortion and sagging curvature, homogeneous heat transfer and sealing of the individual wells are essential quality criteria here. The footprint of all 384 well PCR microplates is compatible with automated systems.

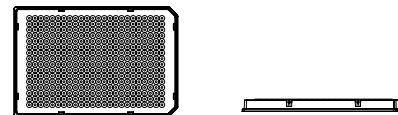


Figure 4: Full-skirted 384 well microplate with two notches for real-time PCR systems such as LightCycler® 480 from Roche

1/ FULL-SKIRTED 384 WELL MICROPLATE WITH TWO NOTCHES FOR REAL-TIME PCR SYSTEMS (ART. NO. 785285)

- / White pigmentation boosts real-time PCR signal
- / Black alphanumeric coding enables a quick identification of samples

The registered trademarks of the mentioned manufacturers belong to the above mentioned companies.

COMPATIBILITY FOR PCR MICROPLATES

▲ Optimal Fit	Model	652201 unskirted, chimney top	652250 unskirted, universal	652210 unskirted, low-profile	652260 semi-skirted, ABI-Design	652270 skirted	652290 semi-skirted, suitable for ABI	669265 semi-skirted, RT-PCR	785285 skirted, RT-PCR
		96 well							384 well
Applied Biosystems	Thermal Cyclers	2700	•	•		•		•	
		2720				▲			
		6100				▲			
		9600	•	•		•		•	
		9700	•	•		▲		•	
		Veriti 0.2 ml				▲		▲	
	Real Time Thermal Cyclers	5700		•		•		•	
		PRISM 7000		•		▲		▲	
		7300		•		▲		▲	
		7500		•		▲		▲	
		7700		•		•		▲	
		7900HT		•		▲			
		ViiA 7				▲			
		Quant Studio 12K Flex				•			
	Sequencers	PRISM 310	•	•		•		▲	
		PRISM 3100	•	•		▲		•	
		3130 (XL)	•	•		▲		▲	
		3700	•	•		▲		▲	
PRISM 3730 (XL)		•	•		▲		▲		
3500 (XL)					•				
Bio-Rad / MJ Research	Thermal Cyclers	Gene Cyclers	•	•		•			
		PTC-100	•	•	•		▲	•	
		PTC-200	•	•	•		▲	•	
		PTC-225 Tetrad	•	•	•		▲	•	
		Dyad/Dyad Disciple	•	•	•		▲	•	
		iCycler	•	•			•	▲	
		MyCycler	•	•					
		Mini Gradient		•	•				
		Personal		•				•	
		T100		▲			▲		
	Real Time Thermal Cyclers	DNA Engine family		•	•		•	•	
		C1000/S1000		•	▲		•	•	
		Opticon/Opticon2		•	•		▲		
		Chromo-4			•		•		
		iCycler	•	•			•	▲	
		MyiQ	•	•			•	▲	
		iQ5	•	•			•	▲	
		CFX Connect			•		•	•	
Sequencers	CFX96					▲			
	CFX384								
	BaseStation					•			
Eppendorf	Thermal Cyclers	Mastercycler	•	•	•		•	•	
		Mastercycler ep Gradient	•	•		•	▲	•	
		Mastercycler M384							
		Mastercycler Nexus		•	•		•	•	
		Mastercycler Nexus Gradient		•	•		▲	•	
		Mastercycler Nexus Eco		•	•		•	•	
	Mastercycler Pro		•			•	•		
Real Time T.C.	Mastercycler ep Realplex	•	•			▲	•		

The registered trademarks of the mentioned manufacturers belong to the above mentioned companies. This compatibility chart is a general guide only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

▲ Optimal Fit		Model	652201 unskirted, chimney top	652260 unskirted, universal	652210 unskirted, low-profile	652260 semi-skirted, ABI-Design	652270 skirted	652290 semi-skirted, suitable for ABI	669285 semi-skirted, RT-PCR	785285 skirted, RT-PCR
Amers- ham	Sequencers	MegaBACE 500								
		MegaBACE 1000								
		MegaBACE 4000								
Beck- mann	Sequencers	CEQ	•							
Biometra	Thermal Cyclers	Uno	•	•	•		•	•		
		Uno II	•	•	•		•	•		
		T1 Thermal Cycler	•	•	•		•	•		
		Tgradient	•	•	•		•	•		
		Trobot	•	•	•		•	•		
		T3000		•	•		•			
	T Professional		•	•		•				
Real Time T.C.	Toptical		•	•		•				
Cor- bett	Thermal Cyclers	PalmCycler 96					•	•		
		PalmCycler 384								
Ericom	Thermal Cyclers	Power Block I	•	•	•	•				
		Deltacycler I	•	•	•			•		
		Deltacycler II	•	•	•	•				
		Single Block	•	•	•				•	
		Twin Block	•	•	•					•
Esco	Thermal Cyclers	Swift	•	•					•	
		Gene	•	•			•	•		
		Genius	•	•			•	•		
G-Storm	Thermal Cyclers	GS1	•	•	•	•			•	
		GS2	•	•	•	•			•	
		GS4	•	•	•	•			•	
		GSX	•	•	•	•			•	
		GSXs	•	•	•	•			•	
MWG	Thermal Cyclers	Primus 96	•	•	•	•	•	•		
		Primus 384								
Strate- gene	Thermal Cyclers	Robocycler 96		•					•	
		Robocycler Gradient	•	•	•	•	•	•		
Real Time T.C.	Mx4000 and Mx3005P	•	•	•	•					
TaKa- Ra	Thermal Cyclers	TP240					•			
		TP3000	•	•	•	•	•	•		
Techne	Thermal Cyclers	Touchgene	•	•	•	•			•	
		Cyclogene	•	•	•	•			•	
		Genius	•	•	•	•			•	
		Genius Quad	•	•	•	•			•	
		Genius (TC412)	•	•	•	•			•	
		Flexigene	•	•	•	•			•	
	Touchgene X		•	•	•	•	•	•		
Real Time T.C.	Touchgene Gradient (TC512)	•	•	•	•	•	•			
Thermo Hybaid	Thermal Cyclers	PCR Sprint	•	•	•	•	•	•		
		MBS Satellite (Multiblock)	•	•	•	•	•	•		
		Px2 and PxE	•	•	•	•	•	•		
		PCR Express and Omni-E	•	•	•	•	•	•		
		Touchdown	•	•	•	•	•	•		
Omnigene	•	•	•	•	•	•				
Trans- gen- omic	Sequencers	WAVE System					•			
Ro- che	Real Time T.C.	LightCycler 96							•	
		LightCycler 480							•	•

The registered trademarks of the mentioned manufacturers belong to the above mentioned companies. This compatibility chart is a general guide only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

COMPATIBILITY FOR MINI BLOCK HEATER INSERTS

Art. no.	848916	848923	848902	848913	848921	848915	848950
No. of Tubes per Insert	15x	15x	40x	15x	15x	4x	2x
1ml Cryo.s					•		
2ml Cryo.s					•		
4ml Cryo.s				•			
5ml Cryo.s				•			
1.5ml Reaction Tube	•	•					
2ml Reaction Tube		•					
0.5ml PCR Tube	•	•					
0.2ml PCR Tube			•				
1x8 PCR Tube Strip			•				
15ml Conical Tube						•	
50ml Conical Tube							•
4ml, 12x55mm Tube					•		
5ml, 12x75mm Tube				•			
4.5ml, 12x75mm Tube				•			
7ml, 13x100mm Tube				•			
12ml, 16x100mm Tube						•	
12ml, 17x100mm Tube						•	
20ml, 16x152mm Tube						•	
14ml, 17x95mm Tube						•	
14ml, 18x95mm Tube						•	

This compatibility chart is a general guide only and subject to error and modifications. We cannot accept any liability or responsibility for the above information.

LABORATORY INFORMATION FOR CENTRIFUGATION

CENTRIFUGATION - PRINCIPLE AND CALCULATION OF THE RCF (RELATIVE CENTRIFUGAL FORCE)

Sedimentation of particles in a gravitational field

If a mixture of sand and water is shaken thoroughly and then left to stand, the sedimentation of the solid particles takes place according to their size. As a result of gravitational acceleration ($g = 9.81\text{m/s}^2$), all of the particles are located in a gravitational field under the influence of which the coarse

grains of sand collect at the bottom first and the smaller grains of sand are deposited later. After around 10–20 minutes, the following layering is produced (from bottom to top): coarse grains of sand – fine grains of sand – water. However, other particles (proteins, nucleic acids, viruses, pro- or eucaryotic cells) do not necessarily precipitate or only sediment out after they have been

exposed to higher forces than the force of gravity resulting from the gravitational acceleration. If these forces exceed the counter-forces resulting from convection (heat circulation) and Brownian molecular motion, both of which cause constant mixing of solutions and suspensions, sedimentation takes place.

The sedimentation rate can be calculated on the basis of Stoke's law as follows:

$$V = \frac{d^2(p_p - p_L)g}{18\mu}$$

However, a particle will only sediment out if $p_p > p_L$. If $p_p < p_L$, V becomes negative, consequently the particle floats rather than sedimenting out.

- / V = sedimentation rate
- p_p = density of the particle
- p_L = density of the liquid
- g = 9.81m/s^2
- μ = viscosity of the liquid

Influence of the Centrifugation and Calculation of the RCF respectively RPM

A centrifuge can be used to create a transient gravitational field under the influence of which the sedimentation of cells, cellular components and macro-molecules takes place. In a centrifuge, a suspension located in a centrifuge tube

rotates around a rotational axis. Each particle of the suspension is subject to centrifugal force, which moves it radially away from the rotational axis.

The centrifugal force F_c is calculated as follows:

$$F_c = m_p w^2 r$$

- / where m_p = mass of the particle,
- w = angular velocity (s^{-1}) and
- r = distance of the particle from the rotational axis

The force acting on a particle in a centrifugal field is stated relative to gravitational acceleration, usually as so-called relative centrifugal force (**RCF**) or g-force (**x g**). It is calculated as follows:


$$RCF = 11.18r \left(\frac{R}{1000} \right)^2$$

- / where R = rotor revolutions per min and r = distance of the particle from the rotational axis (cm)

For easier conversion of the **RCF** or g-force into revolutions per min, the equation can be transformed as follows:

$$R = 299 \sqrt{\frac{RCF}{r}}$$

ALIGNMENT CHART

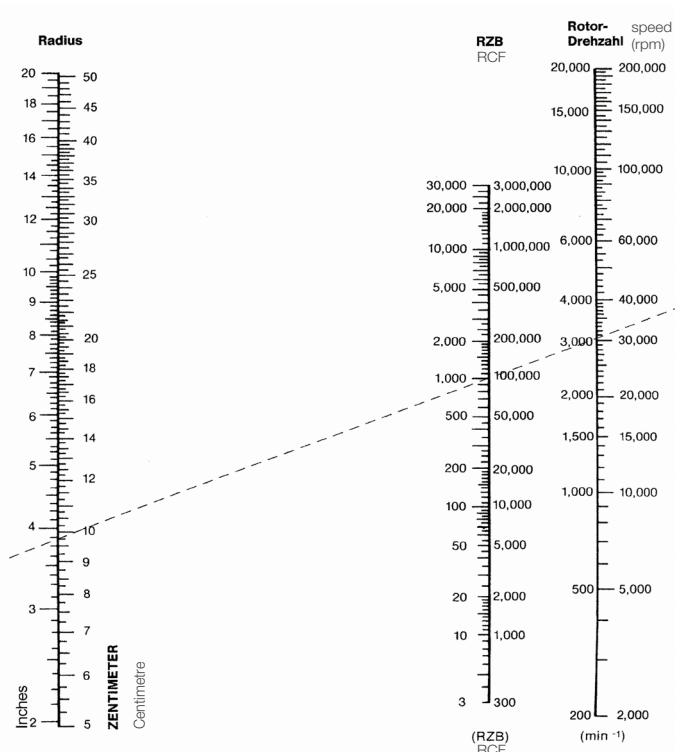

/ By use of a ruler, the third value relating to two known scale values can be read from the alignment chart.

MAXIMAL CENTRIFUGE CAPACITY OF TUBES, REACTION TUBES AND MICROPLATES

The maximum centrifuge capacity for Greiner Bio-One tubes, reaction tubes and microplates is listed in the form of the RCF in the tables below.

For centrifugation, all products were filled with water up to their maximum filling volume. Determination of the maximum RCF in a swinging-bucket rotor was conducted in a Thermo Scientific Centrifuge (Heraeus Multifuge 3 SR).

Determination of the maximum RCF in a fixed angle rotor was conducted in a Sorvall Centrifuge (Evolution RC). Therefore special rotor inserts for different vessel shapes and sizes were used for a stable fit.

REACTION TUBES

Art. no.	Volume [ml]	max. RCF [g] fixed-angle rotor
6162XX	1.5	18000
6232XX	2.0	16000
7162XX	1.5	20000
7172XX	1.5	20000
7222XX	2.0	22000

The stated maximum RCF values are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes for a specific application using high centrifugation forces has to be tested.

POLYSTYRENE TUBES

Art. no.	Dimensions \emptyset [mm] x height [mm]	max. RCF [g] swinging-bucket rotor	max. RCF [g] fixed-angle rotor
1121XX	12 x 55	5800	6200
1151XX	12 x 75	5800	5800
1201XX	12.4 x 75	4800	5000
1361XX	14 x 100	4000	5200
1601XX	16 x 100	3500	5000
1871XX	17 x 100	5200	6600
1881XX	17 x 120	2500	4500
1911XX	18 x 95	4000	5500
2011XX	24 x 90	1000	3500

POLYPROPYLENE TUBES

Art. no.	Dimensions \emptyset [mm] x height [mm]	max. RCF [g] swinging-bucket rotor	max. RCF [g] fixed-angle rotor
1122XX	12 x 55	5800	30000
1152XX	12 x 75	5800	34000
1212XX	12.5 x 48	5800	34000
1222XX	12.5 x 48	5800	34000
1232XX	12.5 x 42	5800	26000
1242XX	12.5 x 86	5800	34000
1262XX	12.4 x 47	5800	26000
1272XX	12.4 x 83	5800	34000
1602XX	16 x 100	5800	33000
187201	17 x 100	4800	34000
1882XX	17 x 120	4000	15000
1912XX	18 x 95	4800	34000
2102XX	30 x 115	2800	9500
2106XX	30 x 115	2800	3500
2272XX	30 x 115	3200	9500
227261	30 x 115	3200	17000
227270	30 x 115	3200	17000

The stated maximum RCF values are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes for a specific application using high centrifugation forces has to be tested.

MULTIWELL PLATES

Art. no.	Multiwell Plate	max. RCF [g] swinging-bucket rotor
657160	6 well, PS, clear	4800
665102	12 well, PS, clear	4800
662160	24 well, PS, clear	4800
677180	48 well, PS, clear	4800

MICROPLATES

Art. no.	Microplate	max. RCF [g] swinging-bucket rotor
650101	96 well, PS, U-bottom, clear	1000
651101	96 well, PS, V-bottom, clear	1000
655101	96 well, PS, F-bottom, clear	4800
650201	96 well, PP, U-bottom, natural	4800
651201	96 well, PP, V-bottom, natural	4800
655201	96 well, PP, F-bottom, natural	4800
655209	96 well, PP, U-bottom, black	4800
655074	96 well, PS, F-bottom, white	4800
655076	96 well, PS, F-bottom, black	4800
655094	96 well, PS, μ Clear [®] , white	4800
655096	96 well, PS, μ Clear [®] , black	4800
655801	96 well, PS, UV-Star [®]	4800
780201	96 well, PP, MASTERBLOCK [®] 1 ml	4800
780270	96 well, PP, MASTERBLOCK [®] 2 ml	4800
786201	96 well, PP, MASTERBLOCK [®] 0.5 ml	4800
781101	384 well, PS, clear	4800
781077	384 well, PS, black	4800
781096	384 well, PS, μ Clear [®] , black	3000
781201	384 well, PP, F-bottom, natural	4800
781280	384 well, PP, V-bottom, natural	4800
781270	384 well, PP, V-bottom, Deep Well, natural	4800
781801	384 well, PS, UV-Star [®]	4800

The stated maximum RCF values in these tables are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes and plates for a specific application using high centrifugation forces has to be tested.

MICROPLATES

Art. no.	Microplate	max. RCF [g] swinging-bucket rotor
784101	384 well, PS, Small Volume, clear	800
784075	384 well, PS, Small Volume, white	800
784076	384 well, PS, Small Volume, black	800
784201	384 well, PP, Small Volume, natural	4800
782101	1536 well, PS, HiBase, clear	1800
782074	1536 well, PS, HiBase, white	1500
782096	1536 well, PS, μ Clear [®] , HiBase, black	1500
782270	1536 well, PP, V-bottom, Deep Well, natural	4800
783076	1536 well, PS, LoBase, black	4800
783096	1536 well, PS, μ Clear [®] , LoBase, black	4800

PCR PLATES

Art. no.	PCR Plate	max. RCF [g] swinging-bucket rotor
652270	96 well, PP, natural, full-skirt	4800
652290	96 well, PP, natural, half-skirt, suitable for ABI	4800

For centrifugation the plates were filled with water as follows: 96 well = 300 μ l | 384 well = 50 μ l | 1536 well = 5 μ l

The stated maximum RCF values in these tables are guidelines only, depending on a variety of factors such as rotor, temperature, density, pH and type of liquid. The suitability of tubes and plates for a specific application using high centrifugation forces has to be tested.

LABORATORY INFORMATION FOR CRYO.S SAMPLE STORAGE

FREEZING PROTOCOL

1

Wash the cells with warm PBS solution, aspirate the solution and cover the cells with a solution containing trypsin and EDTA (a thin liquid film is enough; the concentration should be evaluated for each cell line).

2

Incubate the cells for max. 3–5 min at 37°C.

3

Once the cells detach from the bottom, stop incubation by adding cell culture medium supplemented with serum and slightly suspend cells using a pipette.

4

Spin down the suspension (500xg, 5 min) and re-suspend the pellet with medium containing serum.

5

Determine the cell number (using a Neubauer chamber).

6

Spin down the cells for 5 min at 500xg and discard the supernatant. Resuspend the pellet with an adequate volume of cell culture medium containing serum.

7

Mix the cell suspension 1:1 with freezing medium (60 % medium, 20 % FCS, 20 % DMSO) and transfer it in Cryo.s. For freezing in Cryo.s, the concentration of cells should be $1-5 \times 10^6$ cells/ml.

8

Cryo.s containing cells should be frozen at a cooling rate of -1K/min . This can be achieved by placing them into an isopropanol-filled chamber at -70°C . If other types of samples are contained, Cryo.s may be frozen directly at -20°C , -70°C or in the gas phase of liquid nitrogen. In order to assure even freezing of the sample, 4 and 5 ml Cryo.s should be frozen at -20°C overnight before transferring them to -70°C or to the gas phase of liquid nitrogen.

9

Then transfer the Cryo.s into the nitrogen tank. To avoid contamination (e.g. mycoplasma) and due to safety precautions Cryo.s must only be stored in the gas phase above and not in the liquid nitrogen.

THAWING PROTOCOL

1

Immediately after removing them out of the nitrogen tank the frozen cells are thawed in about 1-2 min brandishing the Cryo.s in a water bath at 37°C. The thawing process should be performed as fast as possible.

2

Transfer the thawed cell suspension into a 15ml tube and mix it immediately with copious amounts of cell culture medium containing serum.

3

After spinning down the cells (500xg, 5min) discard the supernatant and resuspend the pellet in an appropriate cell culture medium supplemented with serum and transfer it into one or more cell culture flasks.

4

Follow the recommended cell concentration for seeding.

5

During the next 12 hours cells should rest.

6

A change of medium is recommended after 24 resp. 48 hours.

SAFETY ADVISORY FOR WORKING WITH CRYO.S

Cryo.s tubes are intended for sample storage exclusively in the gas phase over liquid nitrogen or in freezers! If Cryo.s are stored in the liquid phase, nitrogen can seep into the tubes. Then upon thawing the vaporising nitrogen can generate high pressure, ultimately resulting in an explosion, as well as the release of any infectious material.

Always take appropriate personal safety measures when working with Cryo.s, including wearing safety clothing, using goggles

and working at a safety laboratory bench. When undertaking cryogenic preservation, Cryo.s must be evenly exposed to freezing temperatures. Uneven temperature exposures can cause formation of ice plugs (i. e. at tube top) that inhibit the expansion of freezing liquid (i. e. at tube bottom), resulting in dangerous high pressure and subsequent harm or damage of tubes.

Never exceed maximum working volumes as specified.

LABORATORY INFORMATION FOR IMMUNOLOGY

VOLUME-DEPENDENT WETTING OF IMMUNOLOGICAL PRODUCTS

Liquid volume [μl]	Covered area [mm ²]	Liquid height [mm]	Area / volume ratio [cm ² / cm ³]
96 Well ELISA Microplate, U-Bottom			
25	34	1.7	13.6
50	52	2.6	10.4
75	68	3.4	9.1
100	84.6	4.2	8.5
125	99	4.9	7.9
150	115.5	5.7	7.7
175	130	6.4	7.4
200	145	7.1	7.3
225	160	7.8	7.1
250	174.7	8.5	7.0
275	190	9.2	6.9
300	205	9.9	6.8
96 Well ELISA Microplate, V-Bottom			
25	35	2.3	14.0
50	54.6	3.4	10.9
75	72.4	4.4	9.7
100	88.6	5.3	8.9
125	105	6.2	8.4
150	123.8	7.2	8.3
175	140.8	8.1	8.0
200	156	8.9	7.8
96 Well ELISA Microplate, F-Bottom / Standard			
25	47	0.8	18.8
50	62	1.55	12.4
75	77.5	2.3	10.3
100	92	3.0	9.2
125	108	3.8	8.6
150	123	4.5	8.2
175	137.6	5.2	7.9
200	152.3	5.9	7.6
225	168	6.65	7.5
250	183	7.35	7.3
275	197	8.0	7.2
300	212	8.7	7.1
96 Well ELISA Microplate, F-Bottom / Chimney Well			
25	47	0.7	18.8
50	64	1.5	12.8
75	78.5	2.2	10.5
100	93	2.9	9.3
125	108	3.6	8.6
150	122.6	4.3	8.2
175	137.5	5.0	7.9
200	152	5.7	7.6
225	167	6.4	7.4
250	182	7.1	7.3
275	197	7.8	7.2
300	212	8.4	7.1

Liquid volume [μl]	Covered area [mm ²]	Liquid height [mm]	Area / volume ratio [cm ² / cm ³]
96 Well ELISA Microplate, F-Bottom, Half Area			
25	38	1.65	15.2
50	60	3.2	12.0
75	81.5	4.7	10.9
100	103.6	6.2	10.4
125	124.5	7.6	10.0
150	144	8.9	9.6
175	165.8	10.3	9.5
200	181.7	11.5	9.1
C8 Strip Plate			
25	39	1.0	15.6
50	56	1.9	11.2
75	73	2.8	9.7
100	88.6	3.6	8.9
125	104.3	4.4	8.3
150	120	5.2	8.0
175	136.5	6.0	7.8
200	150.8	6.7	7.5
225	165.4	7.4	7.4
250	181	8.15	7.2
275	196	8.85	7.1
300	211	9.55	7.0
F8 Strip Plate			
25	50.4	0.8	20.2
50	64	1.45	12.8
75	79.7	2.2	10.6
100	93.5	2.85	9.4
125	108.3	3.55	8.7
150	123	4.25	8.2
175	138	4.95	7.9
200	153	5.65	7.7
225	167	6.3	7.4
250	182	7.0	7.3
275	196	7.65	7.1
300	211.5	8.35	7.1
F16 Strip Plate			
25	49	0.8	19.6
50	63	1.5	12.6
75	79.8	2.3	10.6
100	94.3	3.0	9.4
125	108	3.7	8.6
150	123.5	4.4	8.2
175	138	5.1	7.9
200	153	5.8	7.7
225	168	6.5	7.5
250	183	7.2	7.3
275	198	7.9	7.2
300	213	8.6	7.1

Liquid volume [μl]	Covered area [mm ²]	Liquid height [mm]	Area / volume ratio [cm ² / cm ³]
384 Well Microplate, F-Bottom			
25	39.07	2.50	15.6
50	66.60	4.8	13.3
75	94.03	7.00	12.5
100	119.63	9.05	12.0
125	145.6	11.05	11.6
132	152.6	11.50	11.6

Abbreviations

ANSI	American National Standards Institute
COC	Cycloolefin co-polymer
COP	Cycloolefin polymer
CV	Coefficient of Variation
DMSO	Dimethyl Sulphoxide
DNA	Deoxyribonucleic Acid
DNase	Deoxyribonuclease
ECM	Extracellular Matrix
ELISA	Enzyme Linked Immuno Sorbent Assay
EVA	Ethyl Vinyl Acetate
FDA	Food and Drug Administration
FIA	Fluorescence Immuno Assay
F-Rack	Filter Tip Rack
HDPE	High Density Polyethylene
HLA	Human Leucocyte Antigen
HTS	High-Throughput Screening
IgG	Immunoglobulin G
ID-Card	Identity Card
LAL	Limulus Amoebocyte Lysate
LIA	Luminescence Immuno Assay
med.	Medium
NMWCO	Nominal Molecular Weight Cut-Offs
PC	Polycarbonate
PCR	Polymerase Chain Reaction
PDL	Poly-D-Lysine
PET	Polyethylene Terephthalate
PETG	Polyethylene Terephthalate Copolymer
pH	pH Value
PLA	Poly lactate
PLL	Poly-L-Lysine
PP	Polypropylene
PS	Polystyrene
PTFE	Polytetrafluoroethylene
RNA	Ribonucleic Acid
RNase	Ribonuclease
rRNA	Ribosomal RNA
RT	Room Temperature
SBS	Society for Biomolecular Sciences
SPA	Scintillation Proximity Assays
ST-Rack	Standard Rack
TC	Tissue Culture
USP	United States Pharmacopoeia
UV Spectrum	Ultraviolet Spectrum
VIS Spectrum	Visible Spectrum

Units

°C	Degree Centigrade
Da	Dalton, the unit of molecular mass
g	Gram or Gravitational Acceleration (about 9.81 m/s ²)
Gy	Gray, Radiation Unit
h	Hour
l	Liter
M	Molarity, moles of solute per litre of solution
m	Meter
min	Minute
Mol	Absolute Amount of Substance
s	Second

GLOSSARY

Advanced TC is a polymer modification increasing the cellular primary and long-term adhesion of Greiner Bio-One cell culture vessels. Based on the innovative technique the surface of the cell culture vessels is modified to positively influence cellular features and functions. Enhanced cellular adhesion and higher proliferation rates improve cell expansion and cultivation of sensitive cells or cells under restricted growth conditions.

Biobanking Tubes are 300 µl, 600 µl and 1000 µl Cryo.s tubes for the efficient storage of biological samples in large-scale biorepositories. The design of tubes and racks allows for a very space-efficient storage with up to 30 % better utilisation of storage space in freezers or liquid nitrogen tanks. In addition, they are optimised for sample storage at extremely low temperatures over long periods of time.

Bioburden is used to describe the colonisation of viable microorganisms on a material or product and is the basis for determining the necessary radiation dose for sterilisation.

C-bottom stands for the well profile of a flat well bottom with rounded corners.

CELLCOAT® is the Greiner Bio-One brand name for all protein-coated cell culture vessels for adherent cell culture.

CELLMASTER is a quality term that refers to all roller bottles.

CELLreactor is a 15ml/50ml polypropylene tube with filter screw cap for the cultivation of suspension and spheroid cells, expansion of aerobic bacteria, yeast or other microorganisms as well as storage of components and liquids requiring gas exchange.

Cell-repellent surface reliably prevents cell attachment in suspension cultures of semi-adherent and adherent cell lines where standard hydrophobic surfaces generally used for suspension culture are insufficient.

CELLSTAR® is a Greiner Bio-One brand name and includes culture vessels with physically modified surfaces for adherent or suspension cell cultures.

CELLview is a quality term for cell culture products with glass bottom for high-resolution microscopic applications.

Datamatrix Code is a 2D barcode which can also be used for tracking biological and medical reagents and samples. Its small footprint provides nearly infinite scalability and large data capacity. Datamatrix codes can be scanned independent of their orientation and are very accurate due to the Reed-Solomon error correction method.

Deep Well microplates have conical bottom wells and are ideally suited for the storage of non-human samples.

EASYstrainer are cell strainers for the fast and safe filtration of cell suspensions such as those from tissue dissociation or for flow cytometry.

F-bottom stands for a flat bottom well profile.

F-bottom/ chimney well stands for the well profile of a flat well bottom in a chimney-like arrangement. In other words, each well stands on its own. The risk of contamination from sample material being carried over is minimised.

FLUOTRAC is a quality term for immunological products, referring to black microplates (fluorescence measurement).

FourWell Plate is a subdivided plate for microscopic applications facilitating the cultivation of cells and the storage of microscopic slides in an HTS-compatible plate complying with ANSI standards.

Hanging Drop is a technique for protein crystallisation based on vapour diffusion, where droplets literally hang from the top of an upper substrate.

HiBase is a special plate profile of 384 well > Small Volume and 1536 well microplates. In contrast to the > LoBase profile, the HiBase profile is particularly well suited for top-reading systems, since the measuring optic has a minimal separation from the upper edge of the well in this plate profile.

High binding microplates (= MICROLON 600, FLUOTRAC 600 and LUMITRAC 600) are immunological microplates with a high-binding polystyrene surface. Hydrophilic groups are introduced to the polystyrene surface by physical treatment. The high binding surface contains more hydrophilic groups than the less hydrophilic > medium binding surface.

LoBase is a special plate profile in 1536 well microplates. In contrast to the > HiBase profile, the LoBase profile is particularly well suited for bottom-reading systems, since the measuring optic has a minimal separation from the well bottom in this plate profile.

LUMITRAC is a quality term for immunological products, referring to white microplates (luminescence measurement).

MASTERBLOCK® is a brand name that stands for polypropylene microplates that are suitable for the storage of non-human sample material. They are also ideally suited for cultivating bacteria or yeast.

Med. binding (medium binding) microplates (= MICROLON 200, FLUOTRAC 200 and LUMITRAC 200) are immunological microplates with a less hydrophobic surface than > high binding microplates.

µClear® (Micro-Clear) microplates, in contrast to standard microplates with a solid bottom, have a very thin foil bottom. µClear® microplates are ideal for cell-based test systems, microscopic analyses, as well as for bottom-reading systems.

MICROLON is a quality term for immunological products, referring to clear microplates (transmission measurement).

Non-binding microplates are characterised by low protein, DNA, RNA and peptide binding properties.

OneWell Plate is a non-divided HTS plate for tissue culture applications complying with the ANSI standards. The plate is also available in a non-TC-treated version for bacteriology.

Sapphire is a quality term for pipette and filter tips. The product family comprises standard pipette tips, standard filter tips as well as a low-retention version of both. All tips are transparent, graduated and allow precise pipetting with maximal recovery. They can be used with all common pipettes.

SCREENSTAR is a quality term for microplates manufactured out of high-quality cycloolefin with an ultra-clear film bottom for high-content and high-throughput screening.

Sitting Drop is a technique for protein crystallisation based on vapour diffusion, where droplets sit on the bottom of a substrate.

Small Volume is a well profile that was developed in 384 well format for reducing the sample volume. In contrast to the 384 well standard microplate, the sample volume can be considerably reduced, while the detection limit remains the same or is even improved.

TC surface treatment stands for a special physical procedure with which the surfaces of CELLSTAR® products for adherent cell culture are treated. This treatment leads to the incorporation of polar groups, such as carboxy and hydroxy groups, into the plastic surface making it hydrophilic. This enables the adhesion of cells to the plastic surface.




U-bottom stands for well profile of wells with round bottom.

UV-Star® microplates are made of polyolefin and have a film bottom. In contrast to standard microplates with a solid bottom, they are characterised by an extended transparency range to as low as 200 nm.

Vapour diffusion is the most commonly used method for protein crystallisation. In this method a crystallisation droplet, formed by combining a protein solution with a reagent solution, is incubated together with a larger volume of the same reagent solution within a closed system. The reagent solution can contain a wide range of chemicals, e.g. buffers, salts or precipitating agents. Due to mixing the reagent and protein solutions, the concentration of reagents within the crystallisation droplet becomes lower than the concentration of the reagent solution itself. This causes water to evaporate out of the droplet until equilibrium is reached. During this process, the concentration of protein and chemicals in the crystallisation droplet is continuously rising, and, if optimal conditions have been chosen, protein crystals will begin to form. Vapour diffusion experiments are most often set up as hanging or sitting drop.

V-bottom stands for the well profile of wells with a conically tapered well bottom.

GLOSSARY OF SYMBOLS

	Manufacturer		Keep dry		Use-by date
	In vitro diagnostic medical device		Caution		This way up
	Batch code		Do not re-use		Fragile, handle with care
	Article number		Consult instructions for use		Temperature limit
	Sterilised by irradiation		Do not use if package is damaged		

NUMERICAL INDEX

102201.....	168	163290	175	541080	43	639160	45	655096.....	117
102270	168	169101.....	167	542000	71	639160-TRI.....	45	655097.....	117
112101.....	167	187101.....	167	542040	71	639161.....	159	655098.....	62
112201	167	187201.....	167	542070.....	71	639960.....	45	655101.....	117
115001	152	187262.....	173	542100	71	643201.....	162	655160	61
115061.....	152	188161.....	173	542120	71	643203.....	162	655161.....	117
115101.....	167	188171.....	173	542140	71	643401.....	162	655162.....	61
115201.....	167	188241	69	542170	71	644201.....	162	655167.....	58
115262	173	188241	172	543078.....	107	644203.....	162	655169.....	58
120180.....	173	188261-N.....	171	543079.....	107	644401.....	162	655180	61
121261.....	183	188271-N.....	171	543978.....	107	646201.....	162	655182.....	61
121263	184	188271-TRI	171	543979	107	646203.....	162	655185.....	64
121277	184	191180.....	173	604107	231	646401.....	162	655201	119
121278	184	201150.....	177	604160	231	649201.....	162	655209.....	119
121279	184	201152	177	604160-TRI	232	649203.....	162	655630.....	57
121280.....	184	201170	177	604181.....	231	649401.....	162	655640.....	57
122261.....	184	201172	177	606107	231	650001.....	147	655641.....	57
122263	184	205101.....	178	606160	231	650061.....	147	655642.....	58
122263-2D3....	189	210650.....	171	606160-TRI	232	650101.....	117	655643.....	58
122263-2DG ...	189	210651.....	171	606180	231	650160	61	655644.....	58
122263-TRI.....	184	217101.....	178	607107N.....	231	650161.....	117	655680.....	57
122277	184	224170.....	177	607160N.....	231	650180	61	655801	137
122278	184	225170	177	607160N-TRI..	232	650185	64	655809.....	137
122279	184	225180	177	607180N.....	231	650201	119	655830.....	93
122280	184	227245.....	69	608281	212	650209.....	119	655840.....	93
123261.....	183	227245.....	172	609120	142	650261	119	655841.....	93
123263.....	183	227261	171	609171.....	142	650901	132	655846.....	95
123263-2D3....	189	227261-N	171	609820	142	650970.....	90	655850.....	93
123263-2DG ...	189	227261-TRI.....	172	609871	142	650979.....	90	655866.....	111
123277.....	183	227270.....	172	612301.....	233	651001.....	147	655891	109
123278.....	183	227270-N.....	172	612362	233	651061.....	147	655892	113
123279.....	183	227288.....	175	613101.....	207	651101.....	117	655900.....	132
123280	183	227290.....	175	614101.....	207	651160	61	655901	132
124261.....	185	303321	168	616201.....	203	651161.....	117	655903.....	132
124263.....	186	310321.....	168	616261	203	651180	61	655904.....	132
124273.....	186	330070.....	178	622201	204	651201.....	119	655906.....	132
124274.....	186	332070.....	178	622261	204	651209	119	655930.....	63
124275.....	186	354070.....	178	623201	203	651524	95	655936.....	63
124276.....	186	366380.....	204	627102	159	651596.....	96	655940.....	63
126261.....	184	366383.....	204	627160	45	651901.....	132	655944.....	63
126263	184	366384.....	204	627860.....	47	651970.....	90	655946.....	63
126263-2D1	189	366385.....	204	627860.....	105	652201	215	655948.....	63
126263-2DG ...	189	366386.....	204	627870.....	47	652210	215	655950.....	63
126263-TRI.....	184	373250.....	213	627870.....	105	652250.....	215	655956.....	63
126277.....	184	373270.....	213	627965.....	47	652260.....	215	655970.....	90
126278.....	184	373273	213	627965.....	105	652270.....	215	655976.....	90
126279.....	184	373274	213	627975.....	47	652290.....	215	655976-SIN....	90
126280	184	373275	213	627975.....	105	653180	154	655980.....	64
127261.....	185	373276	213	627979.....	46	653190	154	655983.....	64
127263.....	185	373277	213	627979.....	89	655001	147	655986.....	64
127263-2D1	190	373281.....	213	628102	159	655061	147	655990.....	135
127263-2DG ...	189	381061.....	140	628160	45	655073.....	61	655994.....	135
127277.....	185	381070	140	628160-TRI.....	45	655074	117	655996.....	135
127278.....	185	381080	140	628161	159	655075.....	117	655997.....	135
127279.....	185	381081	140	628910.....	46	655076.....	117	656101.....	67
127280	185	385270.....	198	628920.....	46	655077.....	117	656101.....	139
131202.....	193	385273.....	198	628930.....	46	655079.....	61	656161.....	67
131263	192	385274	198	628940.....	46	655080.....	147	656161.....	139
132202	195	385275	198	628950.....	46	655081	147	656170	67
132263	194	385276	198	628979.....	46	655083.....	61	656170	139
133202	197	385277	198	628979.....	89	655086.....	61	656171	67
133263.....	196	385278	198	632180	159	655087.....	62	656171.....	139
136101.....	167	385279.....	198	633180	159	655088.....	62	656190	67
160101	167	420161.....	161	633181.....	159	655090.....	62	656190	139
160201.....	167	421161.....	161	635102	160	655094.....	117	656191.....	67
163288	175	541070	43	639102	159	655095.....	117	656191.....	139

657110.....	55	662630.....	55	673273.....	212	680648.....	77	724410.....	177
657110.....	100	662630.....	100	673274.....	212	680658.....	77	724411.....	177
657160.....	49	662631.....	55	673275.....	212	680660.....	77	724412.....	177
657185.....	49	662631.....	100	673276.....	212	680665.....	77	724414.....	178
657610.....	53	662638.....	55	673277.....	212	680668.....	77	725201.....	203
657610.....	98	662638.....	100	673281.....	212	681670.....	77	725261.....	204
657630.....	53	662640.....	55	673283.....	212	681672.....	77	729101.....	205
657630.....	98	662640.....	100	675001.....	147	681675.....	77	731101.....	161
657631.....	53	662641.....	55	675061.....	147	682201.....	211	731161.....	161
657631.....	98	662641.....	100	675074.....	118	682273.....	211	731165.....	161
657638.....	53	662825.....	94	675075.....	118	682274.....	211	731170.....	161
657638.....	98	662830.....	94	675076.....	118	682275.....	212	731171.....	161
657640.....	53	662840.....	94	675077.....	118	682276.....	212	731175.....	161
657640.....	98	662892.....	113	675083.....	62	682277.....	212	731180.....	161
657641.....	53	662930.....	50	675086.....	62	682281.....	212	731181.....	161
657641.....	98	662940.....	50	675090.....	63	682612.....	77	731185.....	162
657824.....	95	662950.....	50	675096.....	118	682615.....	77	745290.....	229
657825.....	94	662960.....	50	675101.....	118	682624.....	77	756070.....	149
657830.....	94	662970.....	51	675161.....	118	682625.....	77	756071.....	149
657840.....	94	662970.....	90	675180.....	62	682660.....	77	760107.....	231
657841.....	96	664102.....	159	675801.....	137	682670.....	77	760160.....	231
657843.....	96	664160.....	45	676001.....	139	682672.....	77	760160-TRI.....	232
657846.....	96	664160-TRI.....	45	676040.....	139	682678.....	77	760180.....	231
657850.....	95	664161.....	159	676050.....	139	683201.....	211	762070.....	149
657852.....	96	664910.....	46	676051.....	140	683271.....	211	762071.....	149
657896.....	95	664920.....	46	676070.....	139	683273.....	211	762074.....	149
657930.....	50	664940.....	46	676090.....	139	683274.....	211	762075.....	149
657940.....	50	664950.....	46	677102.....	49	683275.....	211	762076.....	149
657950.....	50	664960.....	45	677180.....	49	683276.....	211	762077.....	149
657960.....	49	664970.....	46	677970.....	51	683277.....	211	768160.....	231
657970.....	51	664970.....	89	677970.....	90	684201.....	212	768160-TRI.....	232
657970.....	90	665102.....	49	678101.....	80	688102.....	160	768180.....	231
658170.....	39	665110.....	55	678101-CF1.....	82	690170.....	39	770310.....	228
658175.....	39	665110.....	101	678101-EXF.....	81	690175.....	39	770320.....	228
658175-TRI.....	39	665180.....	49	678104.....	80	690175-TRI.....	39	770330.....	228
658190.....	41	665610.....	54	678104-CF1.....	82	690190.....	41	770340.....	228
658195.....	41	665610.....	99	678104-CF2.....	82	690195.....	41	770370.....	228
658195-TRI.....	41	665630.....	54	678104-EXF.....	81	690195-TRI.....	41	770380.....	228
658910.....	41	665630.....	99	678108.....	80	690910.....	41	771350.....	222
658920.....	41	665631.....	54	678108-CF1.....	82	690920.....	41	771351.....	222
658940.....	41	665631.....	99	678108-CF2.....	82	690940.....	40	771352.....	222
658950.....	40	665638.....	54	678108-EXF.....	81	690950.....	40	771353.....	222
658970.....	40	665638.....	99	678112.....	80	690985.....	42	771354.....	222
658975.....	40	665640.....	54	678112-CF1.....	82	690985.....	89	771355.....	222
658985.....	42	665640.....	99	678112-CF2.....	82	691101.....	67	771361.....	222
658985.....	89	665641.....	54	678112-EXF.....	81	691101.....	139	771362.....	222
660160.....	39	665641.....	99	678116.....	80	691161.....	67	771363.....	222
660175.....	39	665970.....	51	678116-CF1.....	82	691161.....	139	771364.....	222
660175-TRI.....	39	665970.....	90	678116-CF2.....	82	700370.....	233	771365.....	222
660190.....	41	665980.....	49	678116-EXF.....	81	705063.....	150	772350.....	223
660960.....	40	668102.....	204	678124.....	80	705065.....	150	772351.....	223
660975.....	40	669285.....	215	678124-CF1.....	82	705070.....	149	772352.....	223
660985.....	42	670102.....	160	678124-CF2.....	82	705071.....	150	772353.....	223
660985.....	89	670180.....	47	678124-EXF.....	81	705073.....	150	772354.....	223
661160.....	39	670190.....	47	678140.....	80	705074.....	150	772355.....	223
661175.....	39	670640.....	59	678140-CF1.....	82	705075.....	150	772361.....	223
661190.....	41	670641.....	59	678140-CF2.....	82	705076.....	150	772362.....	223
661195.....	41	671201.....	211	678140-EXF.....	81	710107.....	231	772363.....	223
661195-TRI.....	41	671221.....	212	678904.....	80	710160.....	231	772364.....	223
661910.....	41	671273.....	211	678908.....	80	710160-TRI.....	232	772365.....	223
661920.....	41	671274.....	211	678912.....	80	710180.....	231	773353.....	224
661940.....	41	671275.....	211	678916.....	80	710183.....	232	773363.....	224
661950.....	40	671276.....	211	678924.....	80	716201.....	203	774353.....	224
662102.....	49	671277.....	211	678940.....	80	717201.....	203	774363.....	224
662160.....	49	671281.....	211	680190.....	78	717261.....	203	775350.....	225
662610.....	55	673210.....	212	680195.....	78	722201.....	203	775351.....	225
662610.....	100	673271.....	212	680645.....	77	722261.....	203	775352.....	225

775353.....	225	781201-906	122	787979	91	977501	198
775354.....	225	781201-906	130	788860-906 ...	130	977561.....	194
775355.....	225	781209	123	788876	137	977563.....	194
775361.....	225	781270.....	127	789866.....	111	977564.....	194
775362.....	225	781271	127	789979.....	91	977565.....	194
775363.....	225	781280	123	792870-906 ...	130	977566.....	194
775364.....	225	781801	137	793855.....	130	977567.....	194
775365.....	225	781830	94	802202	186	977568.....	194
775390.....	229	781840	93	802203	186	977569.....	194
776350.....	226	781841	93	802204	186	977570.....	195
776351.....	226	781846.....	95	802206	186	977580.....	195
776352.....	226	781850	93	802225.....	187	977583.....	195
776353.....	226	781856.....	113	802501.....	186	977584.....	195
776354.....	226	781866.....	111	802576.....	190	977585.....	195
776355.....	226	781892.....	113	803202	190	977586.....	195
776361.....	226	781900.....	133	803270.....	190	977587.....	195
776362.....	226	781901	132	843070.....	237	977588.....	195
776363.....	226	781903	133	844070	238	977589.....	195
776364.....	226	781904	132	845070.....	238	978501	199
776365.....	226	781906	133	846070.....	237	978561.....	196
777350.....	227	781936.....	65	847070	233	978563.....	196
777351.....	227	781940.....	65	848070.....	239	978564.....	196
777352.....	227	781944	65	848902	239	978565.....	196
777354.....	227	781946.....	65	848913	239	978566.....	196
777355.....	227	781948.....	65	848915	239	978567.....	196
777361.....	227	781950	65	848916	239	978568.....	196
777362.....	227	781956.....	65	848921	239	978569.....	196
777364.....	227	781970.....	90	848923.....	239	978570.....	197
777365.....	227	781974.....	90	848950.....	239	978580.....	197
778350.....	227	781976.....	90	852076-AU....	199	978583.....	197
778351.....	228	781976-SIN	90	852076-CN	199	978584.....	197
778352.....	228	781983.....	65	852076-EU....	199	978585.....	197
778353.....	228	781986.....	65	852076-JP....	199	978586.....	197
778354.....	228	781990.....	135	852076-NA	199	978587.....	197
778355.....	228	781995.....	135	852076-UK....	199	978588.....	197
778361.....	228	781997.....	135	852078-AU....	199	978589.....	197
778362.....	228	782061	125	852078-CN	199	89000002	219
778363.....	228	782073.....	66	852078-EU....	199	89000010.....	219
778364.....	228	782074.....	125	852078-JP....	199	89000020.....	219
778365.....	228	782075.....	125	852078-NA	199	89000099.....	220
779160.....	42	782076.....	125	852078-UK....	199	89000100.....	219
779190.....	42	782078	66	878071	83	89000200.....	219
780201.....	127	782080.....	66	878072.....	83	89000500.....	219
780215.....	127	782086.....	66	878073.....	83	89000810.....	219
780261.....	127	782092.....	66	878075.....	84	89000820.....	219
780270.....	127	782093.....	66	950710	73	89001000.....	219
780271.....	127	782095.....	125	950720.....	73	89001210.....	219
780285.....	127	782096.....	125	951710	73	89001220.....	219
781061.....	121	782101.....	125	951720	73	89008200.....	219
781073.....	65	782180	66	952710	73	89008300.....	219
781074.....	121	782261.....	128	952720.....	73	89010000.....	219
781075.....	121	782270.....	128	960177.....	178	89012200.....	220
781076.....	121	782855.....	130	975502.....	169	89012300.....	220
781077.....	121	782865.....	130	975570.....	169	96077307.....	47
781079.....	65	782900.....	133	976501	198		
781080.....	65	782904.....	133	976561.....	192		
781086.....	65	782946.....	66	976563.....	192		
781090.....	65	782974	91	976564.....	192		
781091.....	65	783892.....	113	976565.....	192		
781092.....	65	784075.....	122	976566.....	192		
781093.....	65	784075-25	122	976567.....	192		
781095.....	121	784076.....	122	976568.....	192		
781096.....	121	784076-25	122	976569.....	192		
781097.....	121	784080.....	66	976570.....	193		
781098.....	65	784101.....	122	976580.....	193		
781101.....	121	784201.....	122	976583.....	193		
781162.....	121	784900.....	133	976584.....	193		
781165.....	65	784904.....	133	976585.....	193		
781182.....	65	784946.....	66	976586.....	193		
781185.....	121	785285.....	215	976587.....	193		
781186.....	121	786201.....	127	976588.....	193		
781201.....	122	786261.....	127	976589.....	193		

GLOSSARY

1536 Well Microplates	125,128,130,132,133	Gel-Load Pipette Tips.....	229
384 Well Microplates.....	57,62,64,121,122,123,127,130,132,133,135,137	Glass Bottom Microplates.....	113
384 Well PCR Microplates	215	Grip stopper.....	168
3D Cell Culture	89,93,94,95,96		
48-way Datamatrix Cryo Rack	190	HTS-Microplates	57,62,64,117,118,119,121,122,123,125,127
81-way Datamatrix Cryo Rack	190		
96 Well ELISA Microplates	147	Immuno Tubes	152
96 Well ELISA Strip Plates	149,150	Inserts for Mini Block Heater	239
96 Well Microplates.....	117,118,119,127,132,135,137		
96 Well PCR Microplates	215	Lab Equipment	237,238,239
96 Well Storage Box.....	169	Leucosep.....	174,175
96-way Datamatrix Cryo Rack	198,199	Leucosep Separation Tubes.....	174,175
		Lids.....	139
Advanced TC.....	49,50	Lids / Sealers / CapMats	139,140
Analyser Cups.....	204,205		
Aspirating Pipette	232	MASTERBLOCK®	127,128
AutoFlask - Cell Culture Flask.....	42	Macro Pipette Tip	229
		MagPen	95,96
Bacteriology	159,160,161	Magnetic 3D Cell Culture.....	93,94,95,96
Biobanking Tubes	198,199	Magnetic Levitation.....	94
		Mass Cell Culture	77
CELLCOAT®	40,41,50	MaxiPette	231,233
CELLdisc.....	80,81,82	Media Bottles.....	73
CELLlevator.....	83	Microbiology	159,160,161
CELLreactor	69,172,173	Microplate Centrifuge	237
CELLring	84	Microplates.....	57,62,64,117,118,119,121,122,123,125,127
CELLstage.....	83	Microplates for Compound Storage.....	130
CELLview Dish	47,105	Mini Block Heater	239
CELLview Plate.....	109	Mini Centrifuge	237
CELLview Slide.....	105,107	Mini Vortex Mixer	238
CapMats.....	140	Molecular Biology	211,212,213,215
Ceaprene stopper	178	Multi-Channel Pipettes.....	219,220
Cell Culture	36,37	Multipurpose Containers / Beakers.....	177,178
Cell Culture Dishes	45	Multiwell Plates.....	49,50,51,90,91
Cell Culture Dishes - Advanced TC.....	45,46,50		
Cell Culture Dishes - CELLCOAT®	46	Non-binding Microplates	132,133
Cell Culture Dishes - Cell-Repellent Surface	46,89		
Cell Culture Flasks	39,42,89	PCR 8-Tube Strips.....	212
Cell Culture Flasks - Advanced TC	40	PCR Microplates	215
Cell Culture Flasks - CELLCOAT®	40,41	PCR Tubes.....	211,212,213
Cell Culture Inserts.....	53,54,55,56,98,99,100,101	Pasteur / Serum Pipettes	233
Cell Culture Microplates	65,66,90,91	Petri Dishes	159,160
Cell Culture Multiwell Plates	49,50,51	Pipette Carousel	220
Cell Culture Roller Bottles	76,77,78	Pipette Tips	222,223,224,225,226,227,228,229
Cell Culture Tubes.....	69	Pipettes.....	219,220
Cell Scraper	43	Polypropylene Tubes	167,168,172,173
Cell Strainer.....	71,73	Polystyrene Tubes	167,173
Cell-Repellent Surface	42,51,89,90,91,96	Polystyrene drosophila containers	178
Cryo Tubes.....	183,185,186	Products for Microscopy	105,107,109,111,113
Cryo storage box.....	186,187	Protein Crystallisation Plates.....	142
Cryo.s	183,185,186,187		
Cryo.s Biobanking Tubes sterile	198,199	Rack for pipette tips	228
Cryo.s with Barcode	189,190	Reaction Tubes.....	203,204,211,212
CrystalQuick	142	Receiver Plates for ThinCert®.....	58
CrystalQuick Plus.....	142	Roller Bottles.....	76,77,78
Datamatrix coded Cryo.s tubes	189,190	SCREENSTAR Microplates.....	91,96,111
Datamatrix rack for cryo tubes.....	190	Sapphire PCR 8-Cap Strips.....	213
Disposable Inoculation Loops / Needles	161,162	Sapphire Pipette Tips.....	219,222,223,224,225,226,227,228
Disposal Bags	162	Sapphire Pipettes.....	219,220
Drosophila Containers / Containers for Plant Culture	178	Screening	95
		Screw Caps.....	198,204
EASYstrainer Cell Strainers.....	71,73	Sealers	139,140
ELISA Microplates	147	Semi-micro / Macro Cuvette.....	207
ELISA Strip Plates	149,150	SensoPlate	113
		Serological Pipettes.....	231,232
Feeder Plates for ThinCert®	59	Single-Channel Pipettes	219
FourWell Plate	47	Single-break Strip Plates	149,150

Support Rack.....	186
Suspension Culture Flasks.....	41
Swab Tube	161
Terasaki Plates.....	154
ThinCert®.....	56,57,58,59,98,99,100
ThinCert® Cell Culture Inserts	53,54,55
ThinCert® Plate.....	55,56,100,101
Transfer Pipettes.....	233
Tube.....	69
Tubes / Multipurpose Beakers	167,172,173,177,178
Two-position Cap	173
UV-Star® Microplates	137
Vortex Mixer.....	238
inoculation loop.....	161,162
inoculation needle.....	161,162

making a difference

making a difference

www.gbo.com

GREINER BIO-ONE GMBH
FRICKENHAUSEN, GERMANY

PHONE +49 7022 948-0
FAX +49 7022 948-514
E-MAIL info@de.gbo.com



GREINER BIO-ONE IS A GLOBAL PLAYER.
FIND THE CONTACT DETAILS OF YOUR
LOCAL PARTNER ON OUR WEBSITE.



Devices of Greiner Bio-One are to be used by properly qualified persons only in accordance with the relevant Instructions for Use (IFU), where applicable.
For more information contact your local Greiner Bio-One sales representative or visit our website (www.gbo.com).

All information is provided without guarantee despite careful processing. Any liability, warranty or guarantee of Greiner Bio-One GmbH is excluded. All rights, errors and changes are reserved. If not stated otherwise, Greiner Bio-One GmbH has all copyrights and/or other (user-)rights in this documents, in particular to signs such as the mentioned (word-picture-)brands and logos. Any use, duplication or any other use of the rights of Greiner Bio-One GmbH is expressly prohibited.

Media owner: Greiner Bio-One GmbH / Represented by Managing Director Bernd Klingel.
The company is registered in the Commercial Register at the first instance court in Stuttgart, HRB 224604 / VAT Number: DE812585719.

F071075 EN Rev. 12/2025


greiner
BIO-ONE